Know what happens

View Your Dissolution Vessels in Real Time — at Half the Cost



With the Merel dissoGUARD Dissolution Surveillance System — specially designed for the Agilent 708-DS Dissolution Apparatus — you can monitor and review videos from tests already performed. And now, for a limited time, you can save 50 % on this real-time surveillance solution.





Limited time offer

Receive a 50 % discount on a dissoGUARD Surveillance System with PRO software with the purchase of an Agilent 708-DS Dissolution Apparatus!

Promotional period ends May 31, 2018. **Use Code 9959**

Don't wait! Talk to your Agilent representative now. www.agilent.com/chem/contactus

Promotion Conditions

Promotional discount cannot be combined with the purchase agreement discount. This promotion cannot be combined with any other current promotions or special offers. Customers must reference the promotion code upon order placement. The order quantities for dissoGUARD systems must be equivalent to the order quantities of the 708-DS dissolution apparatus. Agilent reserves the right to modify or discontinue this offering, during the promotional period.



About dissoGUARD

Using a camera placed beneath each vessel, dissoGUARD requires no additional bench space. In addition, an external camera is employed for an alternate view, determined at your discretion.

The dissoGUARD software gives you a real-time view as well as preset recording options. The videos or snapshots can be marked with key events such as dosage introduction, sampling, or unusual behavior, and then easily exported for sharing or documentation purposes. Each camera does more than simply record events, but can also provide information regarding key physical parameters such as RPM, vessel-to-shaft centering, and wobble. The dissoGUARD PRO software alerts you of any abnormal status of these parameters during a test, preventing lost time and aiding in any failure investigation.

An innovative lighting system using white/red LEDs that are manually or automatically adjusted to red lighting allow you to observe light-sensitive products as well.

Overview software features

Real time preview of 6/7 vessels	\checkmark
Acquire video for 6/7 vessels	\checkmark
Preview and export of video for 6/7 vessels	\checkmark
Dynamic LED (white/red) illumination for 6/7 vessels	\checkmark
Motion detector	\checkmark
Light Shield around the bath	Optional
RPM measurements	\checkmark
Wobble detection	\checkmark
Detection of vessel/shaft centering	\checkmark
Vibration/level sensor	Optional

dissoGUARD PRO

Have you seen this?

Merel dissoGUARD was featured in a recent issue of Dissolution Technologies:

www.dissolutiontech.com/issues/ 201611/DT201611_A03.pdf

Don't wait

Learn more about dissoGUARD at www.dissoguard.com Contact your Agilent representative for your individual offer at www.agilent.com/chem/contactus

Information, descriptions and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc. 2017 Published in USA, August 1, 2017 5991-8293EN

Agilent Trusted Answers