

LC-UV-MS-based Synthetic Peptide Identification and Impurity Profiling Using the ACQUITY QDa Detector with ProMass Software

Ximo Zhang, Brooke Koshel, Jing Fang, William Alley, Robert Birdsall, and Yinq Qing Yu
Waters Corporation, Milford, MA, USA

APPLICATION BENEFITS

- Automated workflow with MassLynx® and ProMass for synthetic peptide target mass confirmation and impurity profiling
- A cost-effective solution for quality control of synthetic peptides using the ACQUITY® QDa® Detector
- Quantitative determination and confirmation of peptide impurities by online UV and MS detection

WATERS SOLUTIONS

[ACQUITY UPLC® H-Class Bio System](#)

[ACQUITY QDa Detector](#)

[ACQUITY UPLC Tunable Ultraviolet \(TUV\) Detector](#)

[Waters® CSH column](#)

[MassLynx 4.1 Software](#)

KEYWORDS

Synthetic peptide, impurity profiling, batch processing, mass detection, ACQUITY QDa, ProMass

INTRODUCTION

The market for peptide-based drugs is growing due to the broad range of activity and low toxicity of peptides.¹ The use of solid-phase synthesis to produce the majority of peptide drugs often introduces process impurities such as incomplete deprotection of peptides and side reaction products.^{2,3} In addition, product-related impurities, such as oxidation and deamidation, add to the complexity of the impurity profile of the drug product.^{2,3} To ensure product efficacy and safety, the International Council for Harmonisation (ICH) guidelines require impurities to be identified based on toxicity and dosage levels of a new product.⁴ Optical thresholding based on LC-UV detection is commonly used to determine the impurity levels. However, optical detection can only be applied to known or pre-characterized samples, which typically require a laborious process of fractionation and enrichment followed by analysis using MS-based techniques for identification. The continuous growth in peptide-based therapeutics requires efficient workflows that can bring products to market in a timely manner while preserving product quality. Techniques that incorporate orthogonal detection methods, such as mass spectrometry (MS), offer the ability to address the challenges of impurity profiling in synthetic peptide production.

The ACQUITY QDa Detector offers an efficient and cost-effective solution to incorporating MS detection to LC-UV-based workflows.⁵ As demonstrated by previous work, the software ProMass (Novatia, LLC) works with MassLynx to provide an automated processing method for MS data generated by the ACQUITY QDa Detector and therefore enables high throughput screening for synthetic peptides.⁶

The purpose of this application note is to demonstrate a cost-effective LC-UV-MS-based workflow using ProMass with MassLynx for mass confirmation and impurity monitoring of synthetic peptides using the ACQUITY QDa Detector as an in-line orthogonal detector. To demonstrate this workflow, a clinically relevant synthetic peptide, eledoisin, is used for analysis.

EXPERIMENTAL

Chemical and reagents

The synthetic peptide eledoisin (pGlu-Pro-Ser-Lys-Asp-Ala-Phe-Ile-Gly-Leu-Met-NH₂) was purchased from New England Peptide. HPLC grade water, acetonitrile, and formic acid were purchased from Fisher Scientific and used as received. A stock solution of the peptide was prepared at a concentration of 2 mg/mL in water and diluted to the desired concentration. Assay optimization was performed by varying the mass load of the peptide from 0.05 µg to 5.00 µg, while the injection volume was kept constant at 5 µL. A 2 µg sample load was found to be the optimal mass load and was used for all analyses unless otherwise noted.

LC conditions

LC system:	ACQUITY UPLC H-Class Bio
Detectors:	ACQUITY UPLC TUV Detector, 5 mm flow cell, $\lambda = 215$ nm ACQUITY QDa Detector (Performance model)
LC column:	ACQUITY UPLC CSH C ₁₈ , 1.7 µm, 130 Å, 2.1 mm × 100 mm (P/N 186006937)
Column temp.:	60 °C
Sample vial:	12 mm × 32 mm glass vial, Total recovery (P/N 600000750cv)
Injection volume:	5 µL
Mobile phase A:	H ₂ O, 0.1% formic acid
Mobile phase B:	Acetonitrile, 0.1% formic acid

Gradient table:

Time (min)	Flow rate (min)	%A	%B
Initial	0.200	82.0	18.0
2.00	0.200	82.0	18.0
2.01	0.200	82.0	18.0
22.01	0.200	72.0	28.0
22.02	0.200	20.0	80.0
24.02	0.200	20.0	80.0
24.03	0.200	82.0	18.0
30.00	0.200	82.0	18.0

ACQUITY QDa Detector settings

Mass range:	350–1,250 <i>m/z</i>
Mode:	ESI Positive
Collection mode:	Continuum
Sample rate:	2 points/sec
Cone voltage:	10 V
Probe temp.:	500 °C
Capillary voltage:	1.5 kV

Data management

MassLynx v 4.1
ProMass for MassLynx

RESULTS AND DISCUSSION

METHOD OPTIMIZATION

Synthetic peptide impurity profiling via LC-UV-based techniques requires optimization of peak resolution and detector response. For this study, a 20 minute gradient was found to resolve 11 impurity peaks from the main peak as shown in Figure 1. ICH guidelines recommend impurity identification to be based on total daily intake. For this study, a cut-off of 0.2% relative abundance was used along with $S/N \geq 10$ to determine the broadest applicability of the method. Eight of the 11 peaks shown in Figure 1 were identified as peaks meeting this criteria based on automatic peak integration performed by MassLynx and are annotated in blue. Peaks above the optical cut-off of S/N ratio were all observed in the MS spectrum (Figure 1B).

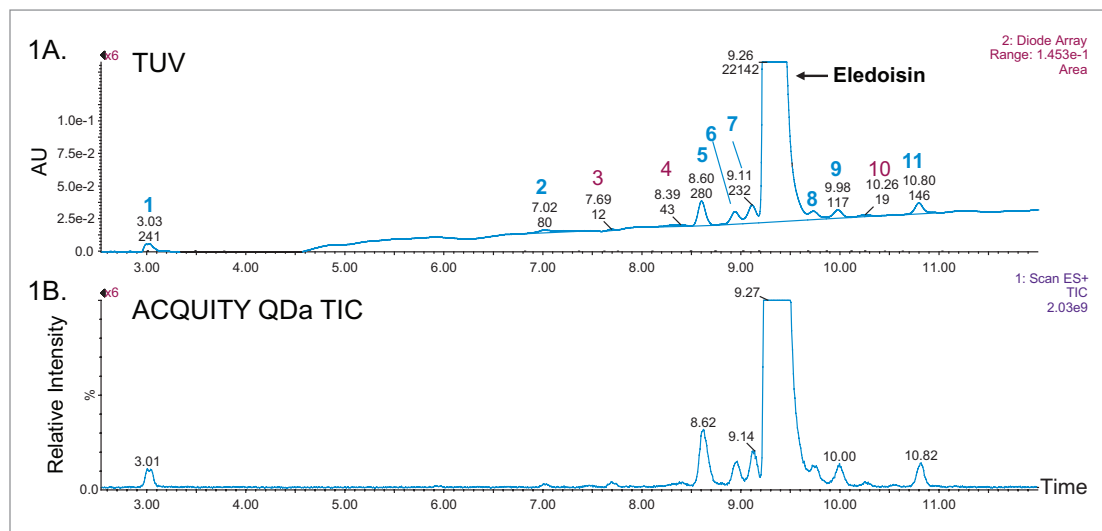


Figure 1. Optimized separation of synthetic peptide eleudoisin with orthogonal mass detection. 1A) UV detection at 215 nm. Using the MassLynx integration tool, 11 impurity peaks were identified and annotated with their retention time and area. Peaks above 0.2% area are labeled in blue, otherwise in red. 1B) MS detection by ACQUITY QDa Detector.

Percent purity determined by mass load was used in lieu of a spiking study to determine the working range of the assay. Ten samples were prepared with increasing concentration (from 0.01 mg/mL to 1.0 mg/mL) and evaluated with the optimized gradient. Purity assessment based on total peaks detected in the optical trace was determined to be stable above mass loads of 0.5 μg at 92.74% \pm 0.45%. To evaluate assay precision, three replicate runs were acquired to measure the %RSD of purity. Results showed the %RSD was below 5% at all mass loads. Using the combined data it was determined that a 2 μg mass load offered optimal chromatographic performance while maintaining an $S/N \geq 10$ for eleudoisin and its impurities as shown in Figure 2.

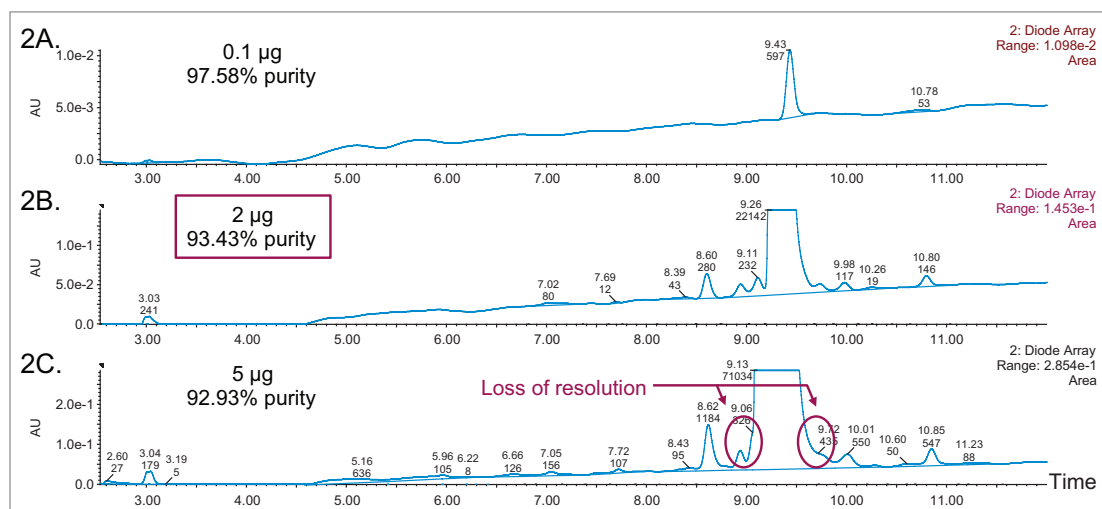


Figure 2. Mass load optimization. Comparison of mass load at 2A) 0.1 μg , 2B) 2 μg , and 2C) 5 μg for synthetic peptide separations are performed using a 20 minute gradient from 18% B to 28% B. Purity was calculated and reported as area percent of target peptide relative to total peak area detected. The optimal mass load was determined to be 2 μg , as higher mass loads resulted in a loss of resolution.

PARAMETERS OF SPECTRUM DECONVOLUTION

With the optimized method determined, the benefit of mass detection can now be fully utilized. The ACQUITY QDa Detector enables the software to determine mass difference between impurities and the target peptide for impurity identification. Automated data acquisition and processing of synthetic peptides can be performed by MassLynx when used in conjunction with ProMass. Figure 3A is an example of the modified Sample List in MassLynx, which contains the instrument methods for data acquisition, as well as information needed for target mass calculation, and parameter files for data deconvolution of synthetic peptides using ProMass. For peptide impurity profiling, the information of the target peptide needs to be entered in the Sequence column as well as the desired Target Info as shown in the zoom-in image of Figure 3A. In this example, 'Ladder = internal' was used to detect all possible fragments with either termini. The target synthetic peptide, eledoisin, has a primary amine group (NH₂) at the C-terminus rather than an OH, thus requiring the termini to be specified as such for correct mass determination. Parameters for MS spectra analysis are defined by the files entered into the ProMassBridge Parameter File column and the Znova Parameter file column. The ProMassBridge Parameter File enables communication between ProMass and MassLynx, and contains user definable parameters such as the retention time window to process, smooth functions and baseline subtraction. As part of the ProMass software, the Znova™ Parameter file provides parameters for spectrum deconvolution, and thresholding for streamlining reporting. The Znova processing file contains a customizable database for target identification and impurity profiling as shown in Figure 3B. Modified and derivatized amino acids can be entered along with naturally occurring amino acids as part of the "building-blocks" for calculating target masses. In the case of eledoisin, Pyroglutamic Acid (/Pyr/=111.1) was entered to account for the amino acid derivative at the terminus of the synthetic peptide. Similarly, previously characterized impurities can be manually entered as their calculated mass differences from the target peak. As shown in the blue rectangle of Figure 3B, deletion of leucine from the oxidized target peptide (loss of Leu from OxiMet = -97.2), possible peptide fragment (Impurity A = -313.92), and C-terminal substitution (OH terminal group instead of NH₂ = 1) were previously identified by high resolution MS prior to this study. Matching mass differences will be shown in deconvolution reports along with their identities.

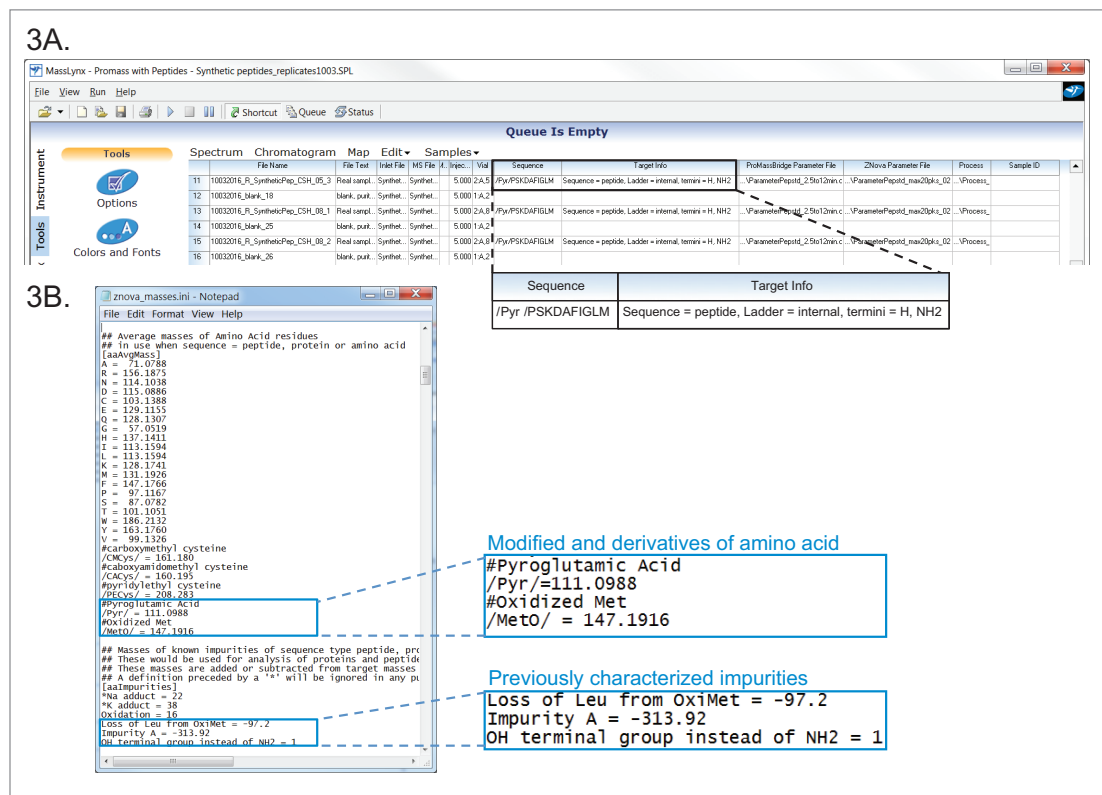


Figure 3. Parameters for data deconvolution. 3A) An example of the modified Sample List in MassLynx. Zoom-in image shows the format for peptide target mass calculation. 3B) The customizable database in Znova Processing File. Naturally occurring amino acids and selective impurities were already included in the database, while entries in the blue rectangle were manually added for target identification and impurity profiling.

RESULTS OF IMPURITY PROFILING

Using the method described above, an analysis of eledoisin was performed for target mass confirmation and impurity screening. Purity of the sample was reported to be 88.87%, which is estimated based on percent abundance of the target mass (including adducts) in the MS spectrum and normalized by chromatographic peak area percent. Percent purity can be found in the Target Mass Summary table, along with other information of the target peptide such as mass error and percent abundance (Figure 4). The sequence Ladder Summary table shows the sequence of identified peptide fragments. The Chromatogram Summary table lists all detected peaks above threshold (0.2%) across the chromatogram. The Spectral Quality, which is based on abundance and the number of charge states detected, was reported "ok" for all impurities with the exception of peaks 4 and 5 which were below the cutoff value (Score ≥ 4). In this example, the Low Score warning at 8.40 min (peak 4) results from a high degree of spectral noise, and the one at 8.62 min (peak 5) is from the misidentification of the base peak charge due to only one charge state being present in the spectrum. The Spectrum Quality threshold, which can be set by the user, enhances the confidence of impurity identification results by indicating the quality of both input and output spectra for deconvolution.

ProMass results:

Target Mass Summary								
RT (min)	Target Mass (Da)	Observed Mass (Da)	Mass Error	Intensity	% Abundance (in Spectrum)	% Purity (Estimate)	Identity	Result Code
9.28	1188.4	1188.2	-0.2 Da (-0.017 %)	2.42E+009	91.67	88.87	Target Mass	OK

Sequence Ladder Summary							
RT (min)	Calculated Mass (Da)	Observed Mass (Da)	Mass Error	Intensity	Sequence		
9.28	1188.4	1188.2	-0.2 Da (-0.017 %)	2.42E+009	/Pyr/1-M11		
8.40	1058.2	1057.2	-1.0 Da (-0.095 %)	9.47E+005	/Pyr/1-L10		
8.95	566.6	566.4	-0.2 Da (-0.035 %)	3.67E+005	S3-F7		

Chromatogram Summary						
Peak	RT (min)	Base Peak Mass (Da)	Intensity	Spectral Quality	LC/MS Peak Area	LC/MS Area Percent
1	3.02	1204.9	1.43E+007	ok	2.58E+006	0.68
2	7.02	874.8	3.78E+006	ok	4.62E+005	0.12
3	7.70	1091.2	4.08E+006	ok	6.98E+005	0.18
4	8.40	1003.6	2.83E+006	low score	7.03E+005	0.19
5	8.62	637.4	2.42E+007	low score	9.02E+006	2.38
6	8.95	1131.1	2.04E+007	ok	2.77E+006	0.73
7	9.12	1116.9	2.79E+007	ok	2.92E+006	0.77
Elledoisin	9.28	1188.2	2.42E+009	ok	3.52E+008	92.80
8	9.76	1259.5	1.70E+007	ok	1.21E+006	0.32
9	9.99	1101.0	1.27E+007	ok	3.02E+006	0.80
10	10.26	1188.5	2.71E+006	ok	5.25E+005	0.14
11	10.82	1189.1	2.17E+007	ok	3.38E+006	0.89

Hyperlinks

Figure 4. Summary tables of peptide impurity profiling. Target mass and percent purity are shown in the Target Mass Summary table. Peaks above 0.2% with UV detection are labeled in blue numbers in the Chromatogram Summary table. All color coded entries in the tables are hyperlinks that lead to further details of deconvolution.

As part of the ProMass software, additional details can be obtained for each analysis by clicking the hyperlinks in Chromatogram Summary table. An example for the peaks at 9.99 min is shown in Figure 5. The top panel shows a list of peaks with their scores and presumed identities. The bottom panel shows the deconvoluted mass spectrum for all detected peaks at 9.99 min. In this list, it is suggested that the mass 1101.0 Da could be a deletion of Serine (S), from the target peptide. Data processing results obtained from ProMass are summarized in Table 1, including retention time, percent area and its %RSD, mass, mass difference between impurity and the target peptide, and presumed identity of detected impurities. The identification of impurities were highly reproducible for all monitored impurities in three replicate runs demonstrating the reliability of LC-MS based workflows in impurity profiling of synthetic peptides.

Hyperlinked embedded data:

Deconvolution Peak Report

Mass Peak List Sorted by Intensity:							
Mass(Da)	+- Std. Dev.	Intensity	Score	Delta Mass	%Relative	%Total	Presumed Identity
1101.0	0.2	1.27E+007	4.62	0.0	100.00	55.48	1188.4 (Minus S)
1154.3	0.5	3.65E+006	4.94	53.3	28.61	15.87	
1241.1	0.4	2.42E+006	3.98	140.1	19.02	10.55	
1242.6	0.7	1.75E+006	3.01	141.6	13.75	7.63	
1123.7	0.9	1.41E+006	3.87	22.7	11.03	6.12	
1189.6	2.0	3.44E+005	3.64	88.6	2.70	1.50	
1170.7	0.8	3.35E+005	3.26	69.7	2.63	1.46	
759.9	1.5	1.77E+005	3.10	-341.1	1.39	0.77	
769.6	0.7	1.42E+005	3.21	-331.4	1.11	0.62	

Mass Peak List Sorted by Mass:							
Mass(Da)	+- Std. Dev.	Intensity	Score	Delta Mass	%Relative	%Total	Presumed Identity
759.9	1.5	1.77E+005	3.10	-341.1	1.39	0.77	
769.6	0.7	1.42E+005	3.21	-331.4	1.11	0.62	
1101.0	0.2	1.27E+007	4.62	0.0	100.00	55.48	1188.4 (Minus S)
1123.7	0.9	1.41E+006	3.87	22.7	11.03	6.12	
1154.3	0.5	3.65E+006	4.94	53.3	28.61	15.87	
1170.7	0.8	3.35E+005	3.26	69.7	2.63	1.46	
1189.6	2.0	3.44E+005	3.64	88.6	2.70	1.50	
1241.1	0.4	2.42E+006	3.98	140.1	19.02	10.55	
1242.6	0.7	1.75E+006	3.01	141.6	13.75	7.63	

Deconvoluted Mass Spectrum

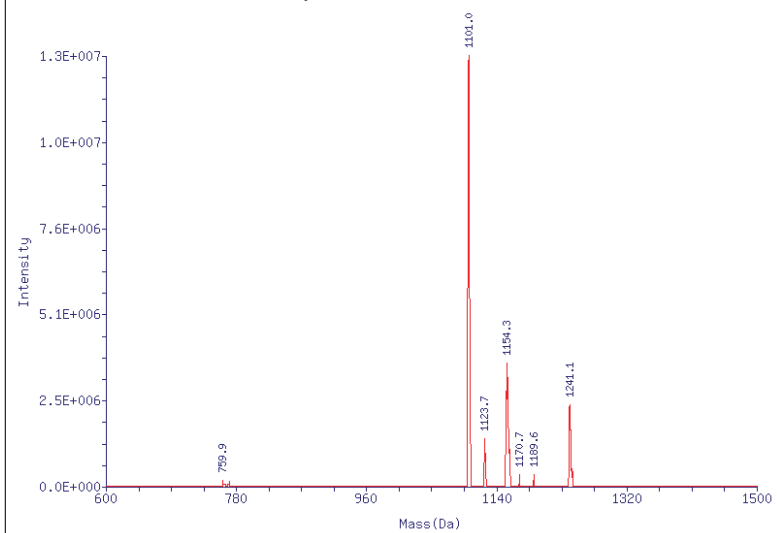


Figure 5. Hyperlinked embedded details of impurity profiling results, such as deconvoluted mass spectra and peak reports, can be accessed through the hyperlinks in the summary tables in Figure 4.

RT (min)	MS %area	%RSD	Mass (Da)	Δ Mass (Da)	Score	Identity
3.02	0.68	3.53	1204.9	16.7	4.86	Oxidation
7.02	0.12	5.09	874.8	-313.4	5.01	Impurity A
7.70	0.18	9.35	1091.2	-97	4.09	Minus P or L
8.40	0.19	12.01	1003.6	-184.6	3.77	pE-KDAFIGLM-NH ₂
8.62	2.38	2.73	637.4	-550.8	3.04	Not identified
			1188.9	0.7	5.11	pE-PSKDAFIGLM-OH
			1245.0	56.8	5.34	Plus G
8.95	0.73	1.57	1131.1	-57.1	4.95	Minus G
9.12	0.77	2.95	1116.9	-71.3	5.12	Minus A
9.28	92.80	0.08	1188.2	0	6.66	pE-PSKDAFIGLM-NH ₂
9.76	0.32	8.17	1259.5	71.3	4.59	Plus A
9.99	0.80	4.56	1099.5	-88.7	4.62	Minus S
10.26	0.14	13.32	1188.4	0.3	4.34	pE-PSKDAFIGLM-NH ₂
10.82	0.89	2.90	1189.1	0.9	4.82	pE-PSKDAFIGLM-OH

Table 1. Summary of impurities with their percent peak area, %RSD of percent peak area based on three replicate runs, mass, mass difference between impurity and the target peptide, score, and presumed identity.

CONCLUSION

An automated workflow of synthetic peptide mass confirmation and impurities profiling was developed using the ACQUITY QDa Detector with MassLynx and ProMass. Impurities above 0.2% optical threshold were successfully detected and identified for eledoisin at 2 µg mass load. Retention time, mass, and presumed identity of both the target peptide and impurities were summarized in ProMass deconvolution summary tables and reports. Confidence of deconvolution was represented as a Score for each peak. The addition of the ACQUITY QDa Detector to an optical workflow significantly reduces the time of impurity analysis. With ProMass for MassLynx, the automated workflow allows high throughput screening and therefore improves the productivity of synthetic peptide impurity profiling in development and quality control environments.

References

1. Vlieghe, M., Lisowski, V., Martinez, J., Khrestchartisky, M. Synthetic therapeutic peptides: science and market. *Drug Discovery Today*. 2010, 15:40–56.
2. Sanz-Nebot, V., Toro, I., Barbosa, J. Fractionation and characterization of a crude peptide mixture for the synthesis of eledoisin by liquid chromatography-electrospray ionization mass spectrometry. *Journal of Chromatography A*. 1999, 846:25–38.
3. Zeng, K., Geerlof-Vidavisky, I., Gucinski, A., Jiang, X., Boyne, M. T. Liquid chromatography-high resolution mass spectrometry for peptide drug quality control. *The AAPS Journal*. 2015, 17:643–651.
4. Guidance for industry Q3B (R2) impurities in new drug products. 2006, Revision 2.
5. Birdsall, R., E., McCarthy, S., M. Increasing specificity and sensitivity in routine peptide analyses using mass detection with the ACQUITY QDa Detector. Waters Application Note. 2015 ([p/n 720005377EN](#)).
6. Birdsall, R., E., Yu, Y., Q. High-throughput screening of oligonucleotides for identity and purity assessment using the ACQUITY QDa Detector and ProMass for MassLynx. Waters Application Note. 2016 ([p/n 720005681EN](#)).

Waters

THE SCIENCE OF WHAT'S POSSIBLE.®

Waters, The Science of What's Possible, MassLynx, ACQUITY, QDa, and ACQUITY UPLC are registered trademarks of Waters Corporation. All other trademarks are the property of their respective owners.

©2017 Waters Corporation. Produced in the U.S.A. February 2017 720005921EN AG-PDF

Waters Corporation
34 Maple Street
Milford, MA 01757 U.S.A.
T: 1 508 478 2000
F: 1 508 872 1990
www.waters.com