

916 Ti-Touch



Manual

8.916.8005EN / 2022-04-30



Metrohm AG
CH-9100 Herisau
Switzerland
+41 71 353 85 85
info@metrohm.com
www.metrohm.com

916 Ti-Touch

Program version 5.916.0045

Manual

Technical Communication
Metrohm AG
CH-9100 Herisau

This documentation is protected by copyright. All rights reserved.

This documentation has been prepared with great care. However, errors can never be entirely ruled out. Please send comments regarding possible errors to the address above.

Disclaimer

Deficiencies arising from circumstances that are not the responsibility of Metrohm, such as improper storage or improper use, etc., are expressly excluded from the warranty. Unauthorized modifications to the product (e.g. conversions or attachments) exclude any liability on the part of the manufacturer for resulting damage and its consequences. Instructions and notes in the Metrohm product documentation must be strictly followed. Otherwise, Metrohm's liability is excluded.

Melody for the BEEP command: excerpt from "En Altfrentsche", with kind permission of the Laseyer Quartett, Appenzell.

Table of contents

1	Introduction	1
1.1	Instrument description	1
1.1.1	Titration and measuring modes	1
1.1.2	Connectors	3
1.1.3	Intended use	3
1.2	About the documentation	4
1.2.1	Symbols and conventions	4
2	Safety instructions	6
2.1	General notes on safety	6
2.2	Electrical safety	6
2.3	Tubing and capillary connections	7
2.4	Flammable solvents and chemicals	7
2.5	Recycling and disposal	8
3	Overview of the instrument	9
3.1	Front of the instrument	9
3.2	Rear of the instrument	10
4	Installation	11
4.1	Setting up the instrument	11
4.1.1	Packaging	11
4.1.2	Checks	11
4.1.3	Location	11
4.2	Connecting the power supply unit	11
4.3	Mounting the electrode holder	13
4.4	Connecting the propeller stirrer (depending on the product version)	14
4.5	Connecting MSB devices	15
4.5.1	Connecting a dosing device	16
4.5.2	Connecting an additional stirrer or titration stand	17
4.5.3	Connecting a Remote Box	17
4.6	Connecting USB devices	18
4.6.1	General	18
4.6.2	Connecting a USB hub	18
4.6.3	Connecting a printer	18
4.6.4	Connecting a balance	19
4.6.5	Connecting a PC keyboard	20
4.6.6	Connecting a barcode reader	21



4.6.7	Connecting a Sample Processor	21
4.7	Connecting sensors	22
4.7.1	General	22
4.7.2	Connecting a pH, metal or ion-selective electrode	23
4.7.3	Connecting a reference electrode	23
4.7.4	Connecting a polarizable electrode	24
4.7.5	Connecting the temperature sensor or electrode with integrated temperature sensor	25
4.7.6	Connecting the iConnect	26
4.7.7	Differential potentiometry	29
4.8	Connecting the Ti-Touch to a network	30
5	Titration	31
5.1	Dynamic equivalence point titration (DET)	31
5.2	Monotonic equivalence point titration (MET)	31
5.3	Endpoint titration (SET)	32
5.4	Titration with constant maintenance of the measured value (STAT)	32
5.5	Manual titration (MAT)	33
6	Operation	35
6.1	Switching the instrument on and off	35
6.2	Fundamentals of operation	36
6.2.1	Touch-sensitive screen	36
6.2.2	Display elements and controls	37
6.2.3	Status display	38
6.2.4	Entering text and numbers	39
7	System settings	41
7.1	General system settings	41
7.1.1	Selecting the dialog language	41
7.1.2	Setting the date, time and local time	42
7.2	System-specific dialog options	43
7.3	User administration	46
7.3.1	Editing the user configuration	48
7.3.2	Creating an identification profile	51
7.3.3	Defining login options	52
7.3.4	Password options	54
7.3.5	Modification options	56
7.3.6	Reasons	56
7.3.7	Audit Trail	57
7.4	Measured value display	59
7.5	Acoustic signals	59

8	Titrants	60
8.1	Adding a new titrant	61
8.2	Editing titrant data	62
8.3	Monitoring the working life	64
8.4	Dosing unit	65
8.4.1	Parameters for preparing (PREP) and emptying (EMPTY)	66
8.4.2	Tubing parameters	68
8.4.3	Shift direction of the valve disk	71
8.5	Exchange unit	71
8.5.1	Parameters for the preparation (PREP)	73
8.5.2	Tubing parameters	74
8.6	GLP test for exchange unit and dosing unit	75
8.7	Titer determination options and data	77
8.7.1	Titer validity	77
8.7.2	Properties of the previous titer determinations	78
9	Reagents	80
9.1	Editing reagent data	80
9.2	Reagent monitoring	81
10	Sensors	84
10.1	Adding a new sensor	85
10.2	Editing the sensor data	86
10.3	Monitoring the working life	87
10.4	Calibration data (for pH electrodes only)	88
10.4.1	Properties of the previous calibrations	90
10.5	Limit values for the calibration data	92
10.6	Monitoring the calibration interval (for pH electrodes only)	93
11	Device manager	95
11.1	Adding a new device	96
11.2	Configuring the instrument	96
11.3	Ti-Touch	97
11.3.1	E-mail	98
11.3.2	PC/LIMS report	99
11.3.3	Shared memory	100
11.3.4	TCP/IP settings	102
11.4	Metrohm control instruments	103
11.4.1	Properties – Measuring input	104
11.4.2	Properties – MSB connector	105



11.4.3	Properties – Peripheral devices	106
11.5	Sample Processor	107
11.5.1	Properties – Sample Processor	108
11.5.2	Properties – Tower	108
11.5.3	Properties – Swing Head	110
11.6	Sample racks	115
11.6.1	Editing rack data	117
11.6.2	Rack adjustment	123
11.7	Printer	124
11.7.1	PDF settings	125
11.7.2	Network printer	126
11.7.3	More options	128
11.8	Balance	128
11.9	USB/RS-232 adapter	130
11.10	PC keyboard	132
11.11	Barcode reader	133
12	File manager	136
12.1	Managing files	136
12.1.1	Copying a file	138
12.1.2	Renaming a file	138
12.1.3	File properties	139
12.2	External storage medium	140
12.3	Creating backups / Restoring data	142
12.3.1	Restoring data	142
13	GLP manager	144
13.1	Automatic system test	145
13.2	Test tools	145
13.3	GLP tests for measurement and titration	146
13.3.1	Parameter description	147
13.4	System validation	150
13.4.1	Parameter description	151
13.5	System monitoring	154
13.5.1	Service interval	154
13.5.2	Backup interval	154
14	Common variables	156
14.1	Editing common variables	157
14.2	Properties of common variables	158
14.3	Monitoring validity	159

14.4	Assigning a result automatically to a common variable	160
15	Templates	162
15.1	Sample data	162
15.1.1	Sample identification list	163
15.1.2	Sample assignment table	164
15.2	Custom result templates	166
15.2.1	Editing result templates	167
15.3	Input lines	170
15.3.1	Editing the input signal	171
15.4	Output lines	172
15.4.1	Editing the output signal	174
15.5	Custom calibration buffers	175
15.5.1	Defining calibration buffers	175
15.6	Report header	177
15.7	Custom electrode type	178
15.7.1	Limit values for the electrode rating	179
16	Methods	182
16.1	Creating a new method	182
16.2	Saving a method	183
16.3	Loading a method	184
16.4	Editing a method	185
16.4.1	Inserting a command	186
16.5	Method options	187
16.5.1	Start options	188
16.5.2	Stop options	190
16.5.3	Sample data	191
16.5.4	Method properties	195
16.5.5	Note	196
16.5.6	Saving a determination automatically	197
17	Electronic signatures	199
17.1	Signing methods/determinations electronically	199
17.2	Deleting electronic signatures	200
18	Control	201
19	Favorites	205
19.1	Creating favorites	206
19.1.1	Editing favorites	206

20 Sample data	209
20.1 Entering sample data in the main dialog	209
20.2 Requesting sample data at the start of the determination	210
21 Sample table	212
21.1 General	212
21.2 Editing sample data	215
21.3 Properties	217
22 Determination run	220
22.1 Carrying out a single determination	220
22.2 Performing a sample series	221
22.3 Canceling determinations manually	222
23 Live modifications	223
23.1 Editing the sample data of the running determination	223
23.2 Editing the sample table while a determination is running	224
23.3 Live display	225
23.4 Live parameters	227
24 Results and more determination data	229
24.1 More determination data	230
24.1.1 Details	231
24.2 Messages	233
24.3 Local common variables	234
24.4 Determination properties	234
24.5 Loading a determination	237
24.5.1 Determination list	238
24.6 Saving a determination	242
24.7 Curves	242
24.8 Recalculation and reevaluation	244
25 Statistics	246
25.1 Displaying details for a result	248
25.2 Deleting statistical data	249
25.3 Adding a determination to a determination series	249

26 Result table	251
26.1 Properties	252
26.2 Saving the result table	256
26.3 Loading the result table	256
27 Printing	257
27.1 General report options	259
27.2 Settings of the individual reports	260
27.3 List of all printable reports	261
28 Manual control	265
28.1 Opening and closing the manual control	266
28.2 Measuring	267
28.2.1 Parameter description	268
28.3 Dosing	269
28.3.1 Continuous dosing	272
28.3.2 Dosing fixed volumes	273
28.3.3 Preparing	276
28.3.4 Emptying	277
28.3.5 Filling	277
28.3.6 Replacing reagent	277
28.4 Stirring	278
28.5 Remote	280
28.6 Sample Processor	281
28.6.1 Moving the lift	283
28.6.2 Moving to a rack position	285
28.6.3 External positions	286
29 Parameters	291
29.1 Dynamic equivalence point titrations (DET)	291
29.1.1 Start conditions	291
29.1.2 Titration parameters	294
29.1.3 Stop conditions	298
29.1.4 Potentiometric evaluation	299
29.1.5 Control instrument	304
29.1.6 Sensor	304
29.1.7 Dosing device	306
29.1.8 Stirrer	306
29.2 Monotonic equivalence point titrations (MET)	307
29.2.1 Start conditions	307
29.2.2 Titration parameters	310
29.2.3 Stop conditions	314
29.2.4 Potentiometric evaluation	315



29.2.5	Control instrument	320
29.2.6	Sensor	320
29.2.7	Dosing device	322
29.2.8	Stirrer	323
29.3	Endpoint titrations (SET)	324
29.3.1	Start conditions	324
29.3.2	Control parameters	325
29.3.3	Titration parameters	329
29.3.4	Stop conditions	330
29.3.5	Conditioning	331
29.3.6	Cell	333
29.3.7	Control instrument	333
29.3.8	Sensor	333
29.3.9	Dosing device	335
29.3.10	Stirrer	336
29.4	STAT titrations (STAT)	337
29.4.1	Start conditions	337
29.4.2	Control parameters	338
29.4.3	Titration parameters	341
29.4.4	Stop conditions	342
29.4.5	Monitoring	344
29.4.6	Control instrument	350
29.4.7	Sensor	350
29.4.8	Dosing device	351
29.4.9	Stirrer	354
29.5	Manual titrations (MAT)	355
29.5.1	Dosing parameters	356
29.5.2	Control instrument	357
29.5.3	Sensor	357
29.5.4	Dosing device	358
29.5.5	Stirrer	361
29.6	Measurements (MEAS)	362
29.6.1	Measuring parameters	362
29.6.2	Control instrument	364
29.6.3	Sensor	365
29.6.4	Stirrer	366
29.7	Calibration of sensors (CAL)	367
29.7.1	Calibration parameters (CAL pH)	367
29.7.2	Control instrument	371
29.7.3	Sensor	371
29.7.4	Stirrer	372
29.8	Electrode test for pH electrodes (ELT)	373
29.8.1	Electrode test parameters	374
29.8.2	Control instrument	375
29.8.3	Sensor	375
29.8.4	Stirrer	376
29.8.5	Additional information	377

29.9 Evaluations (EVAL)	379
29.9.1 Fixed endpoint evaluation (EVAL FIX-EP)	380
29.9.2 pK value and half neutralization potential evaluation (EVAL pK/HNP)	381
29.9.3 Minimum and maximum evaluation (EVAL MIN/MAX)	382
29.9.4 Break point evaluation (EVAL BREAK)	384
29.9.5 Rate evaluation (EVAL RATE)	386
29.10 Calculations	387
29.10.1 Calculations (CALC)	387
29.10.2 Calculations (CALC LIVE)	395
29.10.3 Formula editor	397
29.11 Reports (REPORT)	403
29.11.1 General report options	404
29.11.2 Settings of the individual reports	405
29.11.3 List of reports	405
29.12 Dosing and Liquid Handling	408
29.12.1 Preparing an exchange or dosing unit (PREP)	408
29.12.2 Emptying a dosing unit (EMPTY)	409
29.12.3 Dosing a specified volume (ADD)	410
29.12.4 Liquid Handling (LQH)	415
29.13 Communication	418
29.13.1 Scanning remote lines (SCAN)	418
29.13.2 Setting remote lines (CTRL)	420
29.13.3 Scanning the RS-232 interface (SCAN RS)	421
29.13.4 Defining RS-232 commands (CONTROL RS)	422
29.14 Automation	423
29.14.1 Rotating sample rack (MOVE)	423
29.14.2 Moving the lift (LIFT)	425
29.14.3 Controlling pumps (PUMP)	426
29.14.4 Resetting the rack (RACK)	427
29.14.5 Defining the sample variable (SAMPLE)	427
29.14.6 Creating a subsequence (SUBSEQ)	428
29.15 Miscellaneous commands	431
29.15.1 Controlling a stirrer (STIR)	431
29.15.2 Pausing the method run (WAIT)	432
29.15.3 Scan data (REQUEST)	433
29.15.4 Defining an acoustic signal (BEEP)	434
29.15.5 Signing a determination (SIGN)	434
29.15.6 Canceling the method run (END)	434
30 Operation and maintenance	435
30.1 System initialization	435
30.2 Maintenance	436
31 Troubleshooting	437
31.1 Editing methods	437

31.2	Sample table	437
31.3	Results/Statistics	438
31.4	Printing	438
31.5	Manual control	439
31.6	Miscellaneous	439
31.7	SET titration	439
31.8	STAT titration	440
32	Appendix	442
32.1	Manual Dosing Controller, optional accessory for manual titration	442
32.2	Dosing unit	443
32.2.1	Maximum dosing and filling rate	443
32.2.2	Default parameters for preparing (PREP) and emptying (EMPTY)	443
32.3	Exchange unit	444
32.3.1	Maximum dosing and filling rate	444
32.3.2	Default parameters for preparing (PREP)	444
32.4	Stirring rate	444
32.5	Balance	445
32.6	Remote interface	446
32.6.1	Pin assignment of the remote interface	446
32.7	Result variables as parameter setting	449
32.8	Electrode calibration with Sample Processors	450
32.9	Stored buffer series for pH calibration	452
32.9.1	Metrohm	452
32.9.2	NIST (according to DIN standard 19266, 2015)	453
32.9.3	DIN (according to DIN standard 19267, 2012)	454
32.9.4	Fisher	455
32.9.5	Honeywell Fluka	457
32.9.6	Mettler Toledo	458
32.9.7	Beckmann	459
32.9.8	Radiometer Analytical	460
32.9.9	Baker	462
32.9.10	Hamilton DURACAL	463
32.9.11	Precisa	464
32.9.12	Merck CertiPUR 20 / Titrisol	465
32.9.13	Merck CertiPUR 25	466
32.10	Using AuditTrailViewer	468
32.10.1	Installing AuditTrailViewer	468
32.10.2	Opening the Audit Trail	469
32.10.3	Contents of the Audit Trail	470

32.10.4	Filtering the Audit Trail	471
32.10.5	Exporting the Audit Trail	472
32.11	Diagnosis	472
32.11.1	LCD test	473
32.11.2	Formatting an external storage medium	473
32.11.3	Removing an external storage medium	474
32.11.4	Adjusting the touch screen	474
32.11.5	Testing the touch screen	475
32.11.6	Software update (loading program versions and language files)	476
32.11.7	Service	479
32.12	Arithmetic algorithms in the Ti-Touch	479
33	Technical specifications	482
33.1	Touch screen	482
33.2	Measuring inputs	483
33.2.1	Potentiometry	483
33.2.2	Polarizer	483
33.2.3	Temperature	484
33.3	Built-in magnetic stirrer (depending on the product version)	485
33.4	Interfaces	485
33.5	Power supply	486
33.6	Declarations of Conformity	486
33.7	Ambient temperature	486
33.8	Reference conditions	486
33.9	Dimensions	487
33.10	Storage capacity	487
34	Accessories	488
	Index	489

Table of figures

Figure 1	Front 916 Ti-Touch	9
Figure 2	Rear 916 Ti-Touch	10
Figure 3	Connect power supply unit	12
Figure 4	Connecting the propeller stirrer	14
Figure 5	MSB connections	15
Figure 6	MSB connector	16
Figure 7	Connecting a pH, metal or ion-selective electrode	23
Figure 8	Connecting a reference electrode	24
Figure 9	Connecting a polarizable electrode (stirrer connector; not available, depending on the product version)	24
Figure 10	Connecting the temperature sensor or electrode with integrated temperature sensor	25
Figure 11	Plugging in the mini USB adapter cable	26
Figure 12	Plugging in the 854 iConnect	27
Figure 13	Removing the protective cap	27
Figure 14	Aligning the guide pin	28
Figure 15	Attaching the electrode	28
Figure 16	Unplugging the 854 iConnect	29
Figure 17	Connecting the Ti-Touch to a network	30
Figure 18	Reagent dosing for DET	31
Figure 19	Reagent dosing for MET	31
Figure 20	Reagent dosing for SET	32
Figure 21	Reagent dosing for STAT	33
Figure 22	Dosing unit – port assignment	70
Figure 23	Exchange unit – tubing connections	75
Figure 24	Directory structure on the external storage medium	141
Figure 25	Live display "Preparing the dosing unit"	276
Figure 26	Live display "Preparing the exchange unit"	277
Figure 27	Equivalence point recognition and equivalence point numbering within windows	303
Figure 28	Tubbs method for determining the equivalence point	303
Figure 29	Equivalence point recognition and equivalence point numbering within windows	319
Figure 30	Action "Exit method" or "Cancel command"	347
Figure 31	Action "Pause"	347
Figure 32	Action "Wait"	348
Figure 33	Evaluation of pK value / half neutralization potential	382
Figure 34	Evaluation of minimum and maximum	383
Figure 35	Evaluation of a break point	384
Figure 36	Manual Dosing Controller 6.2107.120 – Overview	442
Figure 37	Rotational speed depending on the stirring rate	445
Figure 38	Connectors of the Remote Box	446
Figure 39	Pin assignment of remote socket and plug	446

1 Introduction

1.1 Instrument description

The 916 Ti-Touch is a compact titration system for volumetric titration. This titrator combines in a single instrument the touch-sensitive color monitor for convenient and efficient operation, the titration unit and as well as the magnetic stirrer or an integrated stirrer interface for a rod stirrer, depending on the product version. The upper side of the housing offers space for the titrant and the titration vessel. The titrator is standard-equipped for operation with an external dosing drive of the *800 Dosino* type with a dosing unit. You can, however, also use a *805 Dosimat* with an exchange unit. Thanks to its compact construction, you can use the 916 Ti-Touch in a small space as a stand-alone titrator.

You manage titrants, sensors, methods, etc. conveniently in the internal memory of the 916 Ti-Touch. You can also save your files externally through the USB connector, e.g. on a USB flash drive. On this storage medium you can not only store your methods and determinations, but also create a backup together with all of the data and settings of your system.

The integrated Ethernet connector is available to you should you wish to connect your 916 Ti-Touch to a network. The network connection offers you the following advantages:

- Saving data to a computer within the network
- Printing reports on a network printer
- Sending displayed messages as e-mails

1.1.1 Titration and measuring modes

The 916 Ti-Touch supports the following titration and measuring modes.

- **DET**
Dynamic equivalence point titration. The reagent is added in variable volume steps.
Measuring modes:
 - **pH** (pH measurement)
 - **U** (potentiometric voltage measurement)
 - **Ipol** (voltametric measurement with selectable polarization current)
 - **Upol** (amperometric measurement with selectable polarization voltage)



- **MET**
Monotonic equivalence point titration. The reagent is added in constant volume steps.
Measuring modes:
 - **pH** (pH measurement)
 - **U** (potentiometric voltage measurement)
 - **Ipol** (voltametric measurement with selectable polarization current)
 - **Upol** (amperometric measurement with selectable polarization voltage)
- **SET**
Endpoint titration at one or two specified endpoints.
Measuring modes:
 - **pH** (pH measurement)
 - **U** (potentiometric voltage measurement)
 - **Ipol** (voltametric measurement with selectable polarization current)
 - **Upol** (amperometric measurement with selectable polarization voltage)
- **STAT**
Titration with constant maintenance of the measured value.
Measuring modes:
 - **pH** (pH measurement)
 - **U** (potentiometric voltage measurement)
- **MAT**
Manual titration with manual dosing and manual ending.
Measuring modes (optional):
 - **pH** (pH measurement)
 - **U** (potentiometric voltage measurement)
- **MEAS**
Measuring modes:
 - **pH** (pH measurement)
 - **U** (potentiometric voltage measurement)
 - **Ipol** (voltametric measurement with selectable polarization current)
 - **Upol** (amperometric measurement with selectable polarization voltage)
 - **T** (temperature measurement)
- **CAL**
pH electrode calibration.
Measuring mode:
 - **pH** (calibration of pH electrodes)
 - **ELT** (Electrode test for pH electrodes)

1.1.2 Connectors

The 916 Ti-Touch is equipped with the following connectors:

- **Power connection**
For connecting to the power grid using the power supply unit provided.
- **Two MSB connectors (Metrohm Serial Bus)**
For connecting dosing devices, one additional stirrer or a Remote Box.
- **USB connector**
For connecting peripheral devices (printer, PC keyboard, etc.), a USB flash drive, a Sample Processor or a USB hub.
- **Sensor connectors**
One connector each for:
 - Potentiometric electrodes (pH, ISE, metal)
 - Reference electrodes
 - Polarizable electrodes
 - Intelligent electrodes (iTrodes)
 - Temperature sensor (Pt1000 or NTC)
- **Stirrer connector (depending on the product version)**
For connecting a propeller stirrer (only available for the version without magnetic stirrer).
- **Ethernet connector**
For connecting the Ti-Touch to a network.

1.1.3 Intended use

The 916 Ti-Touch is designed for usage as a titrator in analytical laboratories. Its main application field is volumetric titration.

This instrument is suitable for processing chemicals and flammable samples. Therefore, the use of the instrument requires the user to have basic knowledge and experience in handling toxic and caustic substances. Knowledge regarding the application of fire prevention measures prescribed for laboratories is also mandatory.

1.2 About the documentation



CAUTION

Please read through this documentation carefully before putting the instrument into operation. The documentation contains information and warnings which the user must follow in order to ensure safe operation of the instrument.

1.2.1 Symbols and conventions

The following symbols and formatting may appear in this documentation:

(5-12)

Cross-reference to figure legend

The first number refers to the figure number, the second to the instrument part in the figure.

1

Instruction step

Perform the steps one after the other.

Method

Dialog text, parameter in the software

File ► New

Menu or menu item

[Continue]

Button or **key**



WARNING

This symbol draws attention to a possible life-threatening hazard or risk of injury.



WARNING

This symbol draws attention to a possible hazard due to electrical current.



WARNING

This symbol draws attention to a possible hazard due to heat or hot instrument parts.



WARNING

This symbol draws attention to a possible biological hazard.



WARNING

Warning of optical radiation



CAUTION

This symbol draws attention to possible damage to instruments or instrument parts.



NOTICE

This symbol highlights additional information and tips.



2 Safety instructions

2.1 General notes on safety



WARNING

Operate this instrument only according to the information contained in this documentation.

This instrument left the factory in a flawless state in terms of technical safety. To maintain this state and ensure non-hazardous operation of the instrument, the following instructions must be observed carefully.

2.2 Electrical safety

The electrical safety when working with the instrument is ensured as part of the international standard IEC 61010.



WARNING

Only personnel qualified by Metrohm are authorized to carry out service work on electronic components.



WARNING

Never open the housing of the instrument. The instrument could be damaged by this. There is also a risk of serious injury if live components are touched.

There are no parts inside the housing which can be serviced or replaced by the user.

Supply voltage



WARNING

An incorrect supply voltage can damage the instrument.

Only operate this instrument with a supply voltage specified for it (see rear panel of the instrument).

Protection against electrostatic charges



WARNING

Electronic components are sensitive to electrostatic charges and can be destroyed by discharges.

Do not fail to pull the power cord out of the power socket before you set up or disconnect electrical plug connections at the rear of the instrument.

2.3 Tubing and capillary connections



CAUTION

Leaks in tubing and capillary connections are a safety risk. Tighten all connections well by hand. Avoid applying excessive force to tubing connections. Damaged tubing ends lead to leakage. Appropriate tools can be used to loosen connections.

Check the connections regularly for leakage. If the instrument is used mainly in unattended operation, then weekly inspections are mandatory.

2.4 Flammable solvents and chemicals




WARNING

All relevant safety measures are to be observed when working with flammable solvents and chemicals.

- Set up the instrument in a well-ventilated location (e.g. fume cupboard).
- Keep all sources of flame far from the workplace.
- Clean up spilled liquids and solids immediately.
- Follow the safety instructions of the chemical manufacturer.



2.5 Recycling and disposal

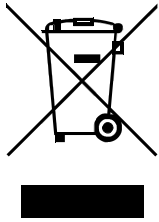

This product is covered by European Directive 2012/19/EU, WEEE – Waste Electrical and Electronic Equipment.

The correct disposal of your old instrument will help to prevent negative effects on the environment and public health.

More details about the disposal of your old instrument can be obtained from your local authorities, from waste disposal companies or from your local dealer.

3 Overview of the instrument

3.1 Front of the instrument

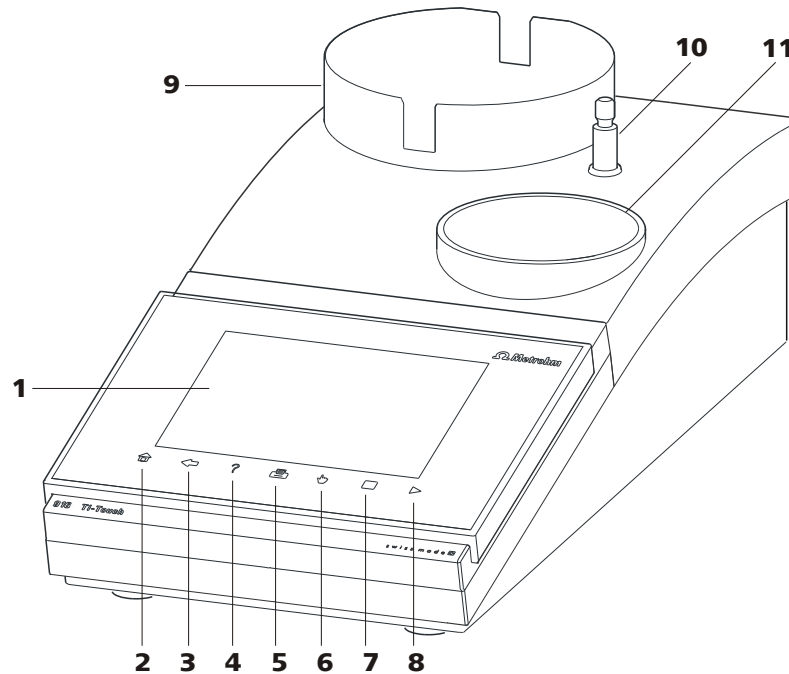


Figure 1 Front 916 Ti-Touch

<p>1 Display Touch screen.</p>	<p>2 Fixed key [Home] Opens the main dialog.</p>
<p>3 Fixed key [Back] Saves the entry and opens the next-higher dialog page.</p>	<p>4 Fixed key [Help] Opens the online help for the dialog displayed.</p>
<p>5 Fixed key [Print] Opens the print dialog.</p>	<p>6 Fixed key [Manual] Opens the manual control.</p>
<p>7 Fixed key [STOP] Cancels the running determination.</p>	<p>8 Fixed key [START] Starts a determination.</p>
<p>9 Bottle holder With holding clips, for reagent bottle.</p>	<p>10 Support rod (lower part) For mounting the support rod (upper part).</p>
<p>11 Titration stand For placing the titration vessel (with built-in magnetic stirrer, depending on the product version).</p>	



3.2 Rear of the instrument

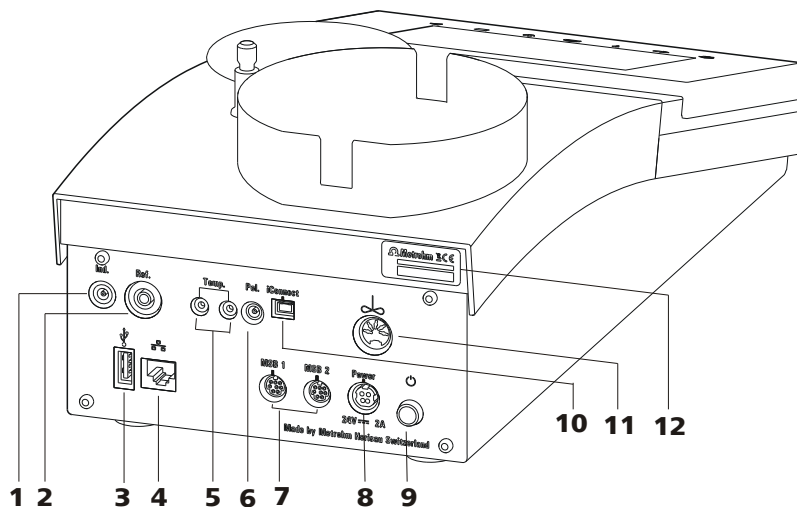


Figure 2 Rear 916 Ti-Touch

1 Electrode connector (Ind.)

For connecting pH, metal or ion-selective electrodes with integrated or separate reference electrode. Socket F.

2 Electrode connector (Ref.)

For connecting reference electrodes, e.g. Ag/AgCl reference electrode. Socket B, 4 mm.

3 USB connector (type A)

For connecting a printer, USB flash drive, USB hub, Sample Processor etc.

4 Ethernet connector (RJ-45)

For connecting to a network.

5 Temperature sensor connector (Temp.)

For connecting temperature sensors (Pt1000 or NTC). Two B sockets, 2 mm.

6 Electrode connector (Pol.)

For connecting polarizable electrodes, e.g. Pt wire electrodes. Socket F.

7 MSB connector (MSB 1 and MSB 2)

Metrohm Serial Bus. For connecting external dosing devices, one additional stirrer or a Remote Box. Mini DIN, 8-pin.

8 Power socket (Power)

For connecting the external power supply unit.

9 Power switch

Switches the instrument on/off.

10 Electrode connector (iConnect)

For connecting electrodes with integrated data chip (iTrodes).

11 Stirrer connector (present, depending on the product version)

For connecting the propeller stirrer (802 Stirrer) to instruments without built-in magnetic stirrer.

12 Type plate

Contains the serial number.

4 Installation

4.1 Setting up the instrument

4.1.1 Packaging

The instrument is supplied in protective packaging together with the separately packed accessories. Keep this packaging, as only this ensures safe transportation of the instrument.

4.1.2 Checks

Immediately after receipt, check whether the shipment has arrived complete and without damage by comparing it with the delivery note.

4.1.3 Location

The instrument has been developed for operation indoors and may not be used in explosive environments.

Place the instrument in a location of the laboratory which is suitable for operation and free of vibrations and which provides protection against corrosive atmosphere and contamination by chemicals.

The instrument should be protected against excessive temperature fluctuations and direct sunlight.

4.2 Connecting the power supply unit

The 916 Ti-Touch has an external power supply unit for a 24 V power supply (DC). This is connected to the power socket of the Ti-Touch.



WARNING

An incorrect supply voltage can damage the instrument.

Operate the instrument only with the supply voltage specified for it. Use the supplied power supply unit exclusively, see chapter *Technical specifications*.

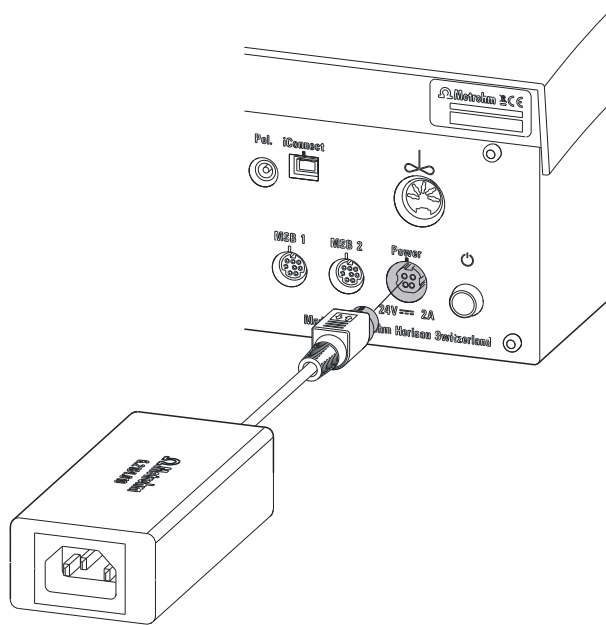


Figure 3 Connect power supply unit

Proceed as follows:

- 1 Connect the plug of the external power supply unit with the power socket of the Ti-Touch (see figure 3, page 12).



NOTICE

The plug of the power supply unit is protected against accidental disconnection of the cable by means of a pull-out protection feature. If you wish to pull out the plug, you first need to pull back the outer plug sleeve marked with arrows.

- 2 Connect the power cord with the external power supply unit of the Ti-Touch and with the power grid.



CAUTION

Switch off the Ti-Touch correctly by pressing the power switch before you disconnect the electricity supply. If this is not done, then there is a danger of data loss.

4.3 Mounting the electrode holder

An electrode or, depending on the product version, a propeller stirrer can be mounted onto the 916 Ti-Touch with the aid of the electrode holder.



NOTICE

To prevent damage to the electrode or rod stirrer, the permissible immersion depth of the products applied can be fixed with the aid of the clamping ring on the electrode holder.

Proceed as follows:

- 1** Push the clamping ring (6.2013.010) down as far as possible onto the mounted support rod (6.2016.050). Make sure while doing so that the indent is facing upwards.
- 2** Attach the electrode holder (6.2021.020) on the support rod.
- 3** Keep the locking lever pressed down and push the electrode holder down.
- 4** Release the locking lever at the desired position. The electrode holder is fixed.
- 5** Readjust the clamping ring to the electrode holder position and screw tight to fix the position of the electrode holder. Make sure that the nose of the electrode holder is placed in the opening of the clamping ring.

4.4 Connecting the propeller stirrer (depending on the product version)

The 916 Ti-Touch is equipped with a built-in magnetic stirrer or a built-in stirrer connector, depending on the product version. You can connect the 802 Stirrer propeller stirrer to the stirrer connector.



NOTICE

Before the rod stirrer is connected, the electrode holder with the clamping ring must be mounted on the support rod, see (see chapter 4.3, page 13).

Proceed as follows:

- 1 Plug in the connection cable of the 802 Stirrer propeller stirrer to the stirrer connector of the Ti-Touch (2-11).

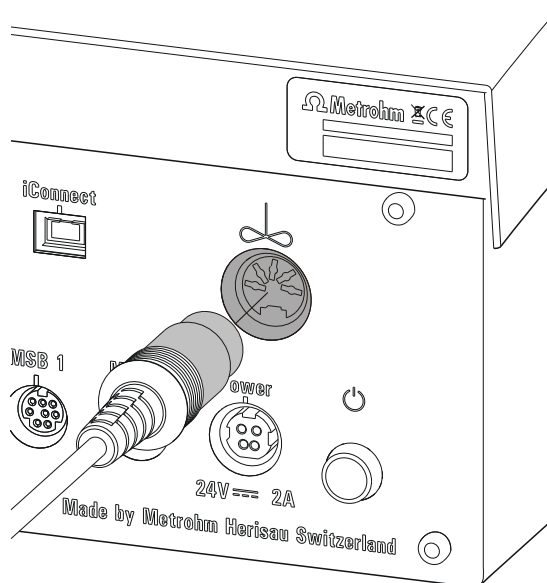


Figure 4 Connecting the propeller stirrer

- 2 Insert the propeller stirrer 802 Stirrer without the stirring propeller from above into the center opening of the electrode holder.
- 3 Plug the stirring propeller (6.1909.010) from below to the propeller stirrer.

- 4 Check whether the height of the electrode holder on the support rod is correctly selected and fixed to prevent damages on the propeller.

4.5 Connecting MSB devices

In order to connect MSB devices, e.g. dosing device or Remote Box, the Ti-Touch has two connectors at what is referred to as the *Metrohm Serial Bus* (MSB). Various peripheral devices can be connected in sequence (daisy chain) at a single MSB connector (8-pin Mini DIN socket) and be controlled simultaneously by the Ti-Touch. In addition to the connection cable, stirrers and the Remote Box are each equipped with their own MSB socket for this purpose.



NOTICE

The additional stirrer must not be connected to the MSB 1!

The following figure provides an overview of the instruments that can be connected to an MSB socket, along with a number of different cabling variations.

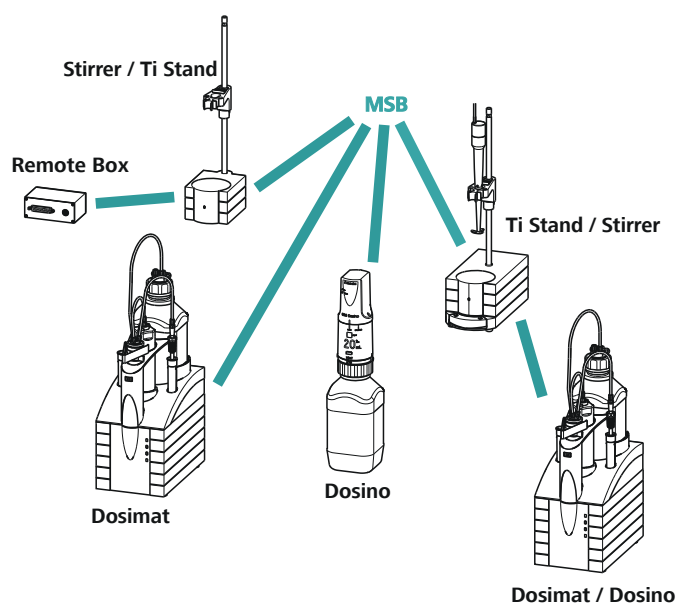


Figure 5 MSB connections



NOTICE

When connecting MSB devices together, the following must be observed:

- No additional stirrer may be connected to the MSB 1!
- Only one device of the same type may be used at a single MSB connector at one time.
- When making the connection, take care to ensure that the flat part of the MSB plug marked with arrows is pointing in the direction of the marking on the MSB connector (see figure 6, page 16).

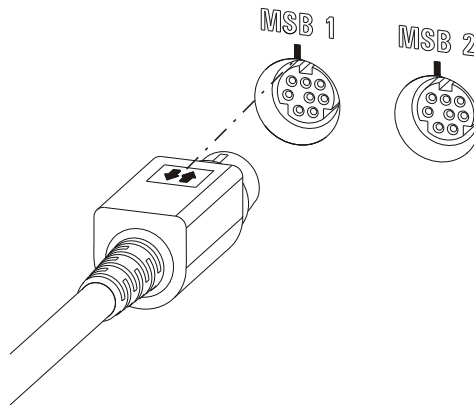


Figure 6 MSB connector



CAUTION

Switch off the Ti-Touch before you plug in MSB devices. When it is switched on, the Ti-Touch automatically recognizes which device is connected to which MSB connector. The connected MSB devices are entered automatically in the device manager.

MSB connections can be extended with the 6.2151.010 cable. The maximum connection length permitted is 6 m.

4.5.1 Connecting a dosing device

You can connect two dosing devices to the Ti-Touch.

The types of dosing devices that are supported are:

- 800 Dosino
- 805 Dosimat

Proceed as follows:

- 1 Switch off the Ti-Touch.
- 2 Connect the dosing device connection cable to an MSB connector (2-7) on the rear side of the Ti-Touch.
- 3 Switch on the Ti-Touch.

4.5.2 Connecting an additional stirrer or titration stand

Depending on the product version, you can use a further stirrer in addition to the built-in magnetic stirrer or the built-in stirrer connector for the propeller stirrer 802 Stirrer, namely the magnetic stirrer 801 Stirrer, the 803 Ti Stand or the 804 Ti Stand with the propeller stirrer 802 Stirrer .

Proceed as follows:

- 1 Switch off the Ti-Touch.
- 2 Connect the connection cable of the magnetic stirrer or of the titration stand to MSB 2 (2-7) on the rear of the Ti-Touch.
- 3 Switch on the Ti-Touch.

4.5.3 Connecting a Remote Box

Instruments that are controlled via remote lines and/or that send control signals via remote lines can be connected via the 6.2148.010 Remote Box. In addition to Metrohm, other instrument manufacturers also use similar connectors that make it possible to connect different instruments together. These interfaces are also frequently given the designations "TTL Logic", "I/O Control" or "Relay Control" and they generally have a signal level of 5 volts.

Control signals are understood to be electrical line statuses or brief (> 200 ms) electrical pulses which display the operational state of an instrument or which trigger or report an event. Sequences on a variety of instruments can thus be coordinated in a single complex automation system. However, no exchange of data is possible.

Proceed as follows:

- 1 Switch off the Ti-Touch.

2 Connect the Remote Box connection cable to an MSB connector (2-7) on the rear side of the Ti-Touch.

3 Switch on the Ti-Touch.

You can connect an 869 Compact Sample Changer. The Remote Box also has an MSB socket at which a further MSB device, e.g. a dosing device, can be connected.

You will find precise information concerning the pin assignment of the interface on the Remote Box in the *appendix*.

4.6 Connecting USB devices

4.6.1 General

The 916 Ti-Touch has a USB connector (type A socket) for peripheral devices with USB interface and for Sample Processors. If you wish to connect more than one device to the USB, you can use a commercially available USB hub.



NOTICE

We recommend that the Ti-Touch be switched off while you set up or disconnect connections between the devices.

4.6.2 Connecting a USB hub

Use a USB hub with its own power supply.

Connect the USB hub as follows:

1 With the help of the 6.2151.030 cable (length 0.6 m) or the 6.2151.020 cable (length 1.8 m), connect the USB connector of the Ti-Touch (Type A) with the USB connector of the hub (Type B, see manual for the USB hub).

The USB hub is recognized automatically.

4.6.3 Connecting a printer

Printers connected to the 916 Ti-Touch must meet the following requirements:

- Printer languages: HP-PCL, HP-PCL-GUI, Canon BJT Commands or Epson ESC P/2
- Paper size: A4 or Letter.

Connect the printer as follows:

- 1 With the aid of the 6.2151.020 cable, connect the USB connector of the Ti-Touch (type A) with the USB connector of the printer (type B, see manual for the printer).
- 2 Configure the printer in the device manager of the Ti-Touch (*see chapter 11.7, page 124*).

4.6.4 Connecting a balance

If you want to connect a balance to the Ti-Touch you need a USB/RS-232 adapter (6.2148.050).

The following table offers an overview of the balances that you can use together with the Ti-Touch and of which cable you will need for connection to the RS-232 interface:

Balance	Cables
AND ER, FR, FX with RS-232 interface (OP-03)	6.2125.020 + 6.2125.010
Mettler AB, AG, PR (LC-RS9)	In the scope of delivery for the balance
Mettler AM, PM, PE with interface option 016 or Mettler AJ, PJ with interface option 018	6.2146.020 + 6.2125.010 Also from Mettler: ME 47473 adapter and either ME 42500 hand switch or ME 46278 foot switch
Mettler AT	6.2146.020 + 6.2125.010 Also from Mettler: ME 42500 hand switch or ME 46278 foot switch
Mettler AX, MX, UMX, PG, AB-S, PB-S, XP, XS	6.2134.120
Mettler AE with interface option 011 or 012	6.2125.020 + 6.2125.010 Also from Mettler: ME 42500 hand switch or ME 46278 foot switch
Ohaus Voyager, Explorer, Analytical Plus	Cable AS017-09 from Ohaus
Precisa balances with RS-232-C interface	6.2125.080 + 6.2125.010



Balance	Cables
Sartorius MP8, MC, LA, Genius, Cubis	6.2134.060
Shimadzu BX, BW	6.2125.080 + 6.2125.010

Connect the balance as follows:

- 1 Connect the USB plug of the USB/RS-232 adapter with the USB connector of the Ti-Touch (Type A).
The USB/RS-232 adapter will be recognized automatically and entered in the device manager of the Ti-Touch.
- 2 Connect the RS-232 interface of the USB/RS-232 adapter with the RS-232 interface of the balance (see table for cable).
- 3 Switch on the balance.
- 4 If necessary, switch on the RS-232 interface of the balance.
- 5 Configure the RS-232 interface of the USB/RS-232 adapter in the device manager of the Ti-Touch (*see chapter 11.9, page 130*).
- 6 Enter and configure the balance in the device manager of the Ti-Touch (*see chapter 11.8, page 128*).
- 7 Make sure that the parameters of the USB/RS-232 adapter configured in the device manager match those of the balance.

4.6.5 Connecting a PC keyboard

The PC keyboard is used as an aid for text and numerical input.

Connect the PC keyboard as follows:

- 1 Connect the USB plug of the keyboard with the USB connector of the Ti-Touch (Type A).
- 2 Enter and configure the keyboard in the device manager of the Ti-Touch (*see chapter 11.10, page 132*).

4.6.6 Connecting a barcode reader

The barcode reader is used as an aid for text and numerical input. You can connect a barcode reader with USB interface.

Connect the barcode reader as follows:

- 1 Connect the USB plug of the barcode reader with the USB connector of the Ti-Touch (Type A).
- 2 Enter and configure the barcode reader in the device manager of the (see chapter 11.11, page 133).

Settings on the barcode reader:

Program the barcode reader as follows (also see manual for the barcode reader):

- 1 Switch the barcode reader to programming mode.
- 2 Specify the desired layout for the keyboard (USA, Germany, France, Spain, German-speaking Switzerland).
This setting must match the setting in the device manager.
- 3 Make sure that the barcode reader is set in such a way that Ctrl characters (ASCII 00 to 31) can be sent.
- 4 Program the barcode reader in such a way that the ASCII character 02 (STX or Ctrl B) is sent as the first character. This first character is normally referred to as the "Preamble" or "Prefix Code".
- 5 Program the barcode reader in such a way that the ASCII character 04 (EOT or Ctrl D) is sent as the last character. This last character is normally referred to as the "Postamble", "Record Suffix" or "Postfix Code".
- 6 Exit the programming mode.

4.6.7 Connecting a Sample Processor

If you wish to integrate your Ti-Touch in an automation system, then you can connect the following Sample Processors to the USB connector:

- 814 USB Sample Processor
- 815 Robotic USB Sample Processor XL
- 810 Sample Processor (from firmware version 5.916.0040)

To connect a Sample Processor, you need the 6.2151.000 controller cable.

Connect the Sample Processor as follows:

- 1 Connect the Sample Processor to the power grid.
- 2 Connect the Sample Processor to the Ti-Touch with the controller cable.

The Sample Processor will be recognized automatically and entered in the device manager of the Ti-Touch.



NOTICE

The plug on the controller cable is protected against accidental disconnection by means of a pull-out protection feature. If you wish to pull out the plug, you first need to pull back the outer plug sleeve marked with arrows.

- 3 Configure the Sample Processor in the device manager (*see chapter 11.5, page 107*).

4.7 Connecting sensors

4.7.1 General

The measuring interface includes one high-ohm measuring input (**Ind.**) for a pH, metal or ion-selective electrode, one input (**Ref.**) for a separate reference electrode, one measuring input (**Temp.**) for a temperature sensor (Pt1000 or NTC), one measuring input (**Pol.**) for a polarizable electrode and one connector (**iConnect**) for an iConnect. This contains a measuring interface for electrodes with integrated data chip, referred to as iTrodes.



NOTICE

Before the sensor is connected, the electrode holder must be mounted on the support rod with the clamping ring, see (*see chapter 4.3, page 13*).

4.7.2 Connecting a pH, metal or ion-selective electrode

Connect the pH, metal or ion-selective electrode as follows:

- 1 Plug the electrode plug into the **Ind.** socket of the Ti-Touch.

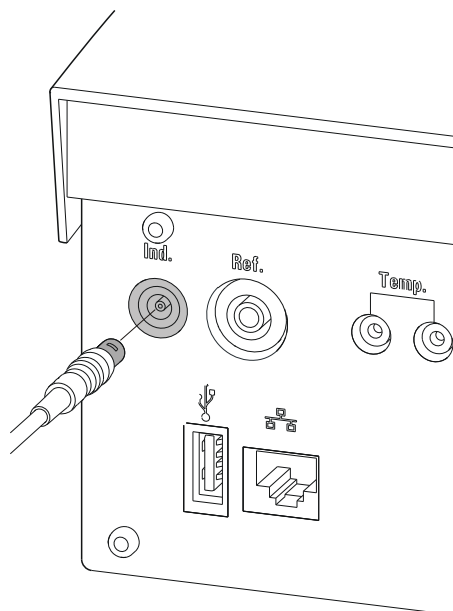


Figure 7 Connecting a pH, metal or ion-selective electrode



NOTICE

The electrode cable is protected against accidental disconnection of the cable by means of a pull-out protection. If you wish to pull out the plug again, you first need to pull back the outer plug sleeve.

4.7.3 Connecting a reference electrode

Connect the reference electrode as follows:

- 1 Plug the electrode plug into the **Ref.** socket of the Ti-Touch.

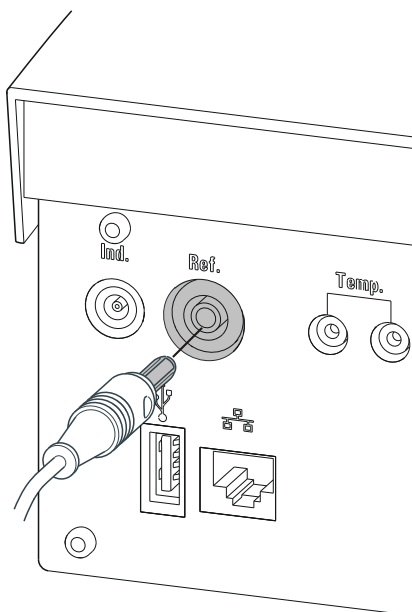


Figure 8 Connecting a reference electrode

4.7.4 Connecting a polarizable electrode

Connect the polarizable electrode as follows:

- 1 Plug the electrode plug into the **Pol.** socket of the Ti-Touch.

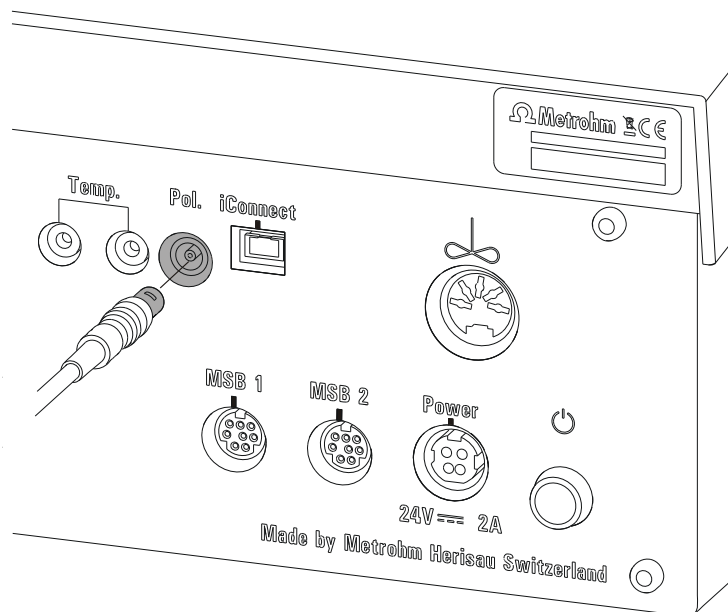


Figure 9 Connecting a polarizable electrode (stirrer connector; not available, depending on the product version)

**NOTICE**

The electrode cable is protected against accidental disconnection of the cable by means of a pull-out protection. If you wish to pull out the plug again, you first need to pull back the outer plug sleeve.

4.7.5 Connecting the temperature sensor or electrode with integrated temperature sensor

A temperature sensor of the Pt1000 or NTC type can be connected to the **Temp.** connector.

Connect the temperature sensor or the electrode with integrated temperature sensor as follows:

- 1 Plug the temperature sensor plugs into the **Temp.** sockets of the Ti-Touch.

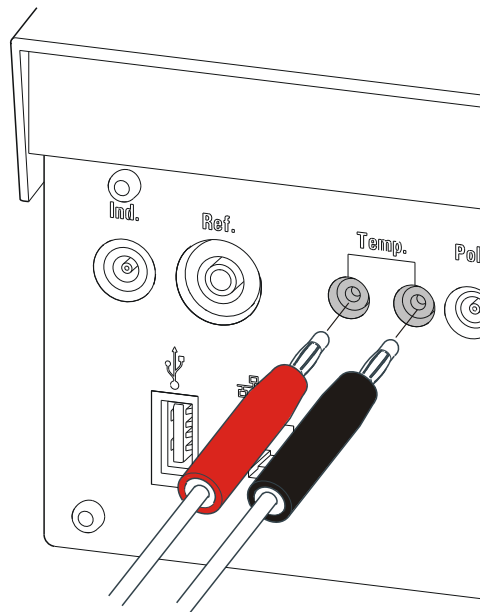


Figure 10 Connecting the temperature sensor or electrode with integrated temperature sensor



NOTICE

The red plug must always be plugged into the red socket for the purpose of shielding against disruptions.

If you use an electrode with an integrated NTC sensor, then you must plug the red plug into the red socket.

4.7.6 Connecting the iConnect

The external measuring interface 854 iConnect can be connected to the **iConnect** connector.

Connecting the 6.2168.000 mini USB adapter cable to an instrument

If the mini USB adapter cable is not connected to the instrument yet, proceed as follows:

- 1 Plug the mini USB adapter cable (2) into the **iConnect** connector of the instrument (1). Observe the correct orientation (markings).

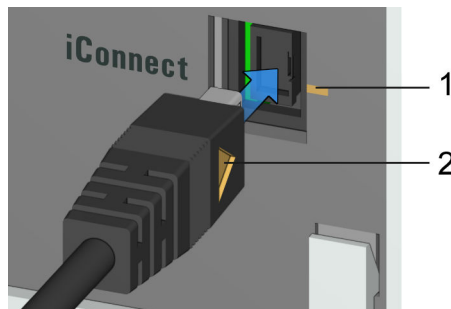


Figure 11 Plugging in the mini USB adapter cable

Depending on the instrument's model version, there may be several **iConnect** connectors.

- 2 Leave the adapter cable plugged in to protect the connector inside the instrument (1) against mechanical impact.

Connecting the 854 iConnect to the adapter cable

Make sure that the mini USB adapter cable is connected to the instrument.

The 854 iConnect can also be connected while the instrument is switched on.

- 1 Plug the plug of the 854 iConnect (3) into the socket of the mini USB adapter cable (2). Observe the correct orientation (markings).

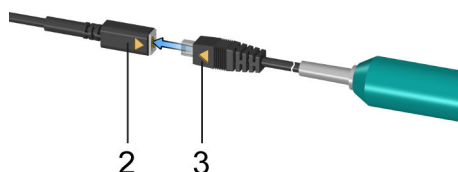


Figure 12 Plugging in the 854 iConnect

As soon as the instrument is switched on, the 854 iConnect is detected automatically and entered as measuring input into the device properties.

Connecting the electrode

The 854 iConnect is used as measuring input for iTrodes (electrodes with integrated memory chip).

Connect the electrode as follows:

- 1 Remove the protective cap of the 854 iConnect.

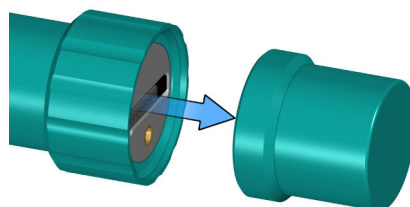


Figure 13 Removing the protective cap

- 2 Align the guide pin (5) of the electrode with the recess in the 854 iConnect (4).

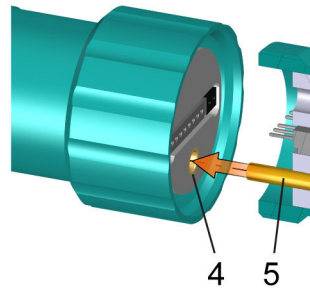


Figure 14 Aligning the guide pin

- 3 Attach the electrode to the 854 iConnect.

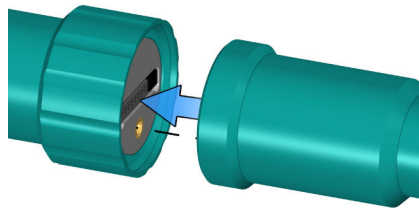


Figure 15 Attaching the electrode

The guide pin guarantees correct connection in such a way that the contact pins cannot be damaged.

- 4 Tighten the screw cap by hand.

If there is an electrode in the sensor list of the firmware or software, the electrode is detected automatically when connecting it.



NOTICE

Mount the protective cap again as soon as the 854 iConnect is not in use anymore and no electrode is connected.

Removing the 854 iConnect

The 854 iConnect can also be removed while the instrument is switched on.

- 1 Unplug the 854 iConnect (3) from the socket of the mini USB adapter cable (2).

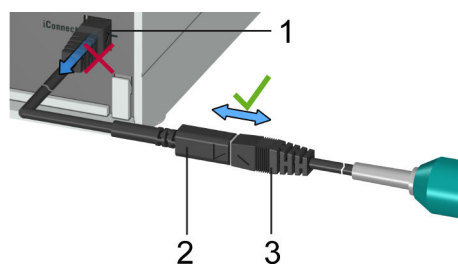


Figure 16 Unplugging the 854 iConnect

- 2 Leave the mini USB adapter cable (2) plugged into the socket of the instrument (1).



NOTICE

Position the mini USB adapter cable in such a way that it cannot be removed by accident.



NOTICE

For further information on the 854 iConnect, see *Manual iConnect 854*, article number 8.854.8002.

4.7.7 Differential potentiometry

Potentiometric measurements with high-ohm measuring chains can be disrupted by electrostatic and electromagnetic fields in media with low conductivity. Use our 6.0229.100 Solvotrode or other special electrodes for pH measurements in organic solvents. If no reliable measurements are possible with these, then a 6.5104.030 (230 V) or 6.5104.040 (115 V) differential amplifier can be used. The differential amplifier is connected to the high-ohm measuring input (**Ind.**).



4.8 Connecting the Ti-Touch to a network

The 916 Ti-Touch has a network connection (Ethernet). This can be used to integrate your Ti-Touch in your network. You can, for example, store data on a PC within the network or print reports on a network printer. In *Chapter 11, page 95*, you will find information as to which settings are necessary for the network connection.

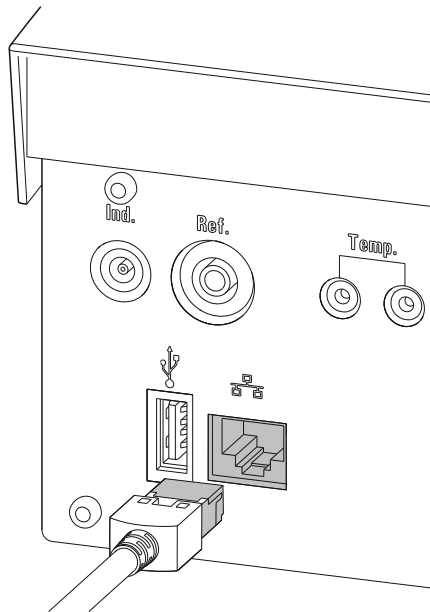


Figure 17 Connecting the Ti-Touch to a network

5 Titrations

5.1 Dynamic equivalence point titration (DET)

Dynamic equivalence point titration is a titration mode for all standard titrations with an s-shaped curve progression. The reagent is added in variable volume steps. The volume steps vary as a function of the slope of the curve. An attempt is made to reach constant measured value changes with each dosing. The optimal volume for dosing is determined from the measured value changes of the previous dosings. Measured value acceptance is measured value drift-controlled (equilibrium titration) or after a waiting time. Equivalence points are evaluated automatically.

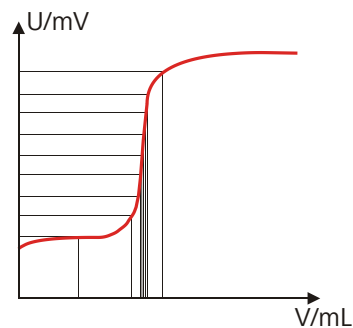


Figure 18 Reagent dosing for DET

5.2 Monotonic equivalence point titration (MET)

Monotonic equivalence point titration is a robust titration mode for titrations with any curve shape and for slow titrations or slow-response electrodes. The reagent is added in constant volume steps. Measured value acceptance is measured value drift-controlled (equilibrium titration) or after a waiting time. Equivalence points are evaluated automatically.

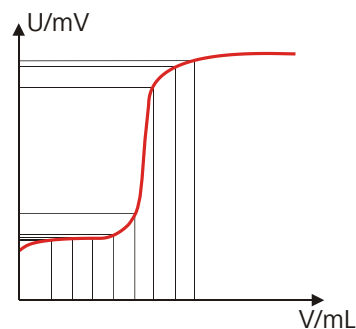


Figure 19 Reagent dosing for MET

5.3 Endpoint titration (SET)

Endpoint titration is a titration mode for rapid routine determinations to a preset endpoint (e.g. titrations in accordance with special standards) and titrations for which reagent overflow must be avoided. The titration termination at the endpoint takes place either volume drift-controlled or after a waiting time. The volume dosed until the endpoint can be used for further calculations (e.g. the content of the sample).

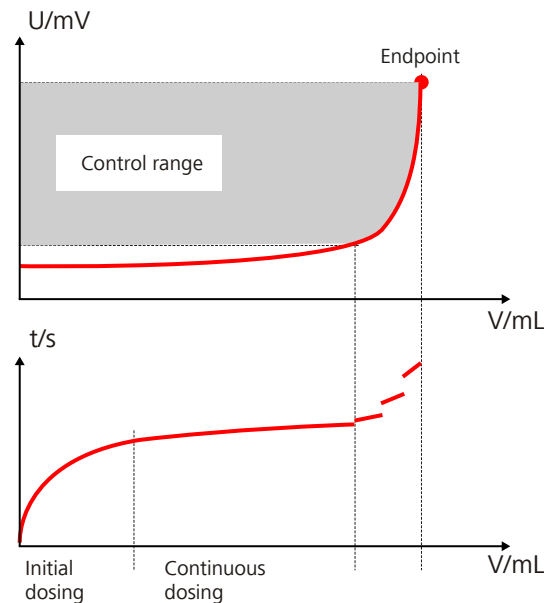


Figure 20 Reagent dosing for SET

5.4 Titration with constant maintenance of the measured value (STAT)

In STAT titrations, reagent is added until a preset measured value (control point) has been reached, which is then maintained at a constant level. Because the substance that is released as a result of the reaction is immediately titrated off with the titrant, the STAT controller has to be able to maintain a preset control point at a constant level until the preset stop condition has been reached, i.e. without oscillating around the control point.

STAT titrations are used, for example, in enzyme analysis where the enzyme activity can be determined from the resulting dosing rate (in mL/min).

Reagent dosing takes place in two phases during the titration:

- **Dosing outside the control range**
The dosing rate increases continuously during this phase. It starts with the Min. rate and increases to the Max. rate. The dosing continues to be carried out at the Max. rate until the Control range is reached. In this range, STAT mode and SET mode control equally.
- **Dosing inside the control range**
The STAT controller compensates minimal measured value deviations from the control point quickly by dosing in the control range faster than for the SET controller.

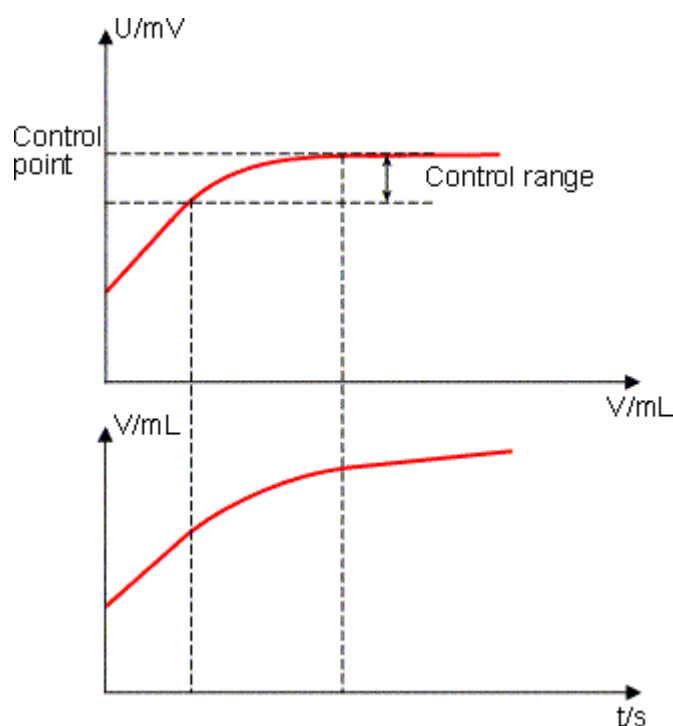


Figure 21 Reagent dosing for STAT

5.5 Manual titration (MAT)

Manual titration is a titration mode that is suitable in particular for titrations that are carried out conventionally manually with a Schellbach buret or a Dosimat.

The reagent is added by manual dosing. The endpoint is detected by a color change or with an electrode connected over a measuring signal. The command always has to be canceled manually with the **Cancel command** button or with the **[FILL]** key on the Manual Dosing Controller.

An advantage of the MAT titration mode is that the manual reagent addition as well as the measuring signal can be recorded optionally. These data points (time, volume, measured value) can be evaluated after the titration, e.g. to assess the titration quality, especially the accuracy of reaching the transition point.



In addition, the method can be equipped with further commands, e.g. calculating a result, creating a report or scanning sample data.

The system is FDA compliant and enables a manual titration according to the FDA guidelines with this mode.

6 Operation

6.1 Switching the instrument on and off

Switching on the instrument



CAUTION

Peripheral devices (e.g., printers) must be connected and switched on before you switch on the 916 Ti-Touch.



NOTICE

English is set as the default dialog language when the instrument is switched on for the first time.


Proceed as follows:

- 1 ▪ Press the power switch on the left-hand side of the back panel of the 916 Ti-Touch.
The 916 Ti-Touch is initialized. A system test is performed. This process takes some time.



NOTICE

If a buret unit is connected, then a request appears to carry out the **Prepare** function:

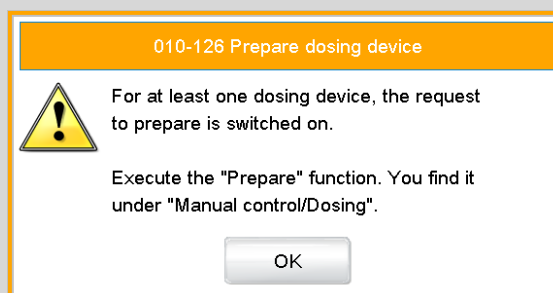

All tubings and the cylinder are rinsed with the **Prepare** function. The preparing of the buret unit is described in *chapter 28.3.3, page 276*.



- Confirm the message with **[OK]**.

The main dialog is displayed:

Switching off the instrument



CAUTION

The 916 Ti-Touch must be switched off by pressing the power switch on the rear of the instrument before the electricity supply is interrupted. If this is not done, then there is a danger of data loss.


Proceed as follows:

- 1 Press the power switch on the left-hand side of the back panel of the 916 Ti-Touch.

The current data is saved and the system is shut down. This process takes just a short time. At the same time, all other instruments connected to the 916 Ti-Touch via a USB cable are also being switched off.

6.2 Fundamentals of operation

6.2.1 Touch-sensitive screen

The entire 916 Ti-Touch user interface is touch-sensitive. Simply touch a few of the buttons on the interface to learn how a touch-sensitive screen reacts. You can always return to the main dialog by touching [].

In order to enable an element on the 916 Ti-Touch user interface, just touch the screen with your fingertip, the eraser of a pencil or a stylus (special tool for operating instruments with touch-sensitive screens).




CAUTION

Never touch the touch screen with a pointed or sharp object such as a ballpoint pen.





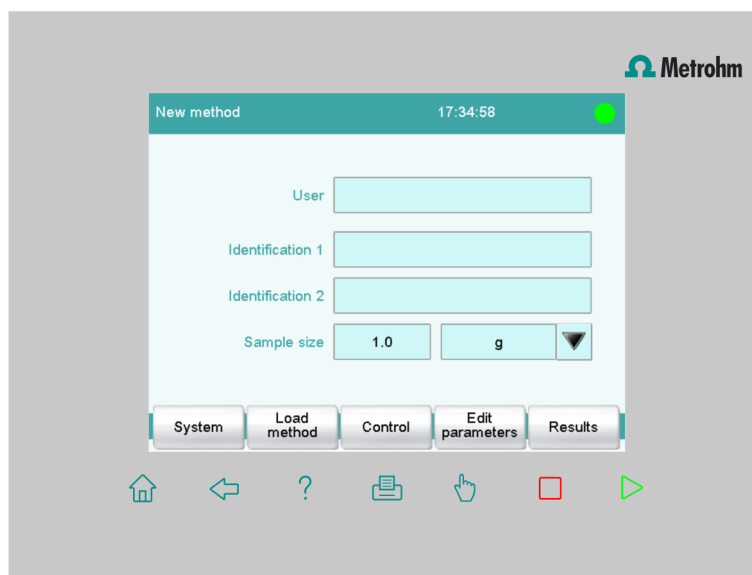


In the default setting, the software is configured in such a way that an acoustic signal will be generated every time an active control is touched. This setting can be deactivated in the system settings (*see chapter 7.5, page 59*).

6.2.2 Display elements and controls



The following display elements and controls are available:



Table 1 Fixed keys which are always available


	[Home] always opens the main dialog.
	[Back] saves the entry and opens the next-higher dialog page.
	[Help] opens the online help for the dialog displayed.
	[Print] opens the printing dialog.
	[Manual] opens the manual control.
	[Stop] cancels the running determination.
	[Start] starts a determination.

The file name of the loaded method, the time and the system status are displayed in the main dialog in the **Title bar**.


In the other dialogs, the title bar shows the headings of the next upper level and of the displayed dialog. This is an aid for orientation during navigation through the user dialog.

Table 2 Screen elements

	Buttons open a new dialog when they are tapped.
	

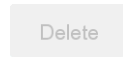

Inactive buttons with gray lettering indicate that the respective function is not available at the moment.

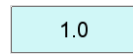

Input fields open an input dialog when tapped.



Tapping on the **selection symbol** opens a selection list.



A **check box** can also be activated or deactivated by tapping on it.

6.2.3 Status display

The current status of the system is displayed in the upper right-hand corner of the title bar.

Table 3 Status displays



The instrument is in normal status.



The working medium is being conditioned.



Conditioning has been paused.



The working medium is conditioned.



A method has been started.



A method has been paused.



An action has been started in manual control.

6.2.4 Entering text and numbers

In the editing dialog for text or numerical input, enter the individual characters by tapping in the input field. The following functions are available:

Text editor

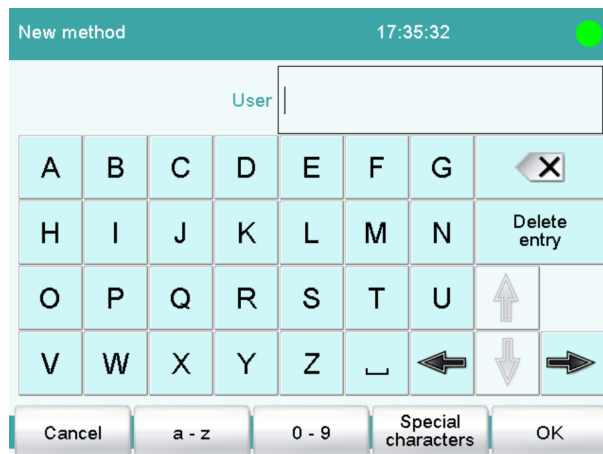


Table 4 Editing functions

Editing function	Description
[OK]	The modification is applied and the editing dialog is exited.
[Cancel]	The editing dialog is exited without applying the modification.
[Delete entry]	The content of the input field is deleted completely.
[X]	The character in front of the cursor is deleted.
[←]	The cursor within the input field is shifted to the left by one character.
[⇒]	The cursor within the input field is shifted to the right by one character.
[a...z]	The lower-case letters are displayed. The label changes to [A...Z]. The upper-case letters are displayed again by tapping.
[0-9]	Numbers and mathematical characters are displayed.
[Special characters]	Special characters are displayed. You can use the button [More] to navigate through all available characters.

Number editor

Edit command / Stop conditions

Stop time s

Input:
1 ... 999999

Default value:
off

7	8	9	off
4	5	6	R1 ▼
1	2	3	
0	+/-	.	

Cancel Delete entry OK

Table 5 Editing functions

Editing function	Description
[OK]	The modification is applied and the editing dialog is exited.
[Cancel]	The editing dialog is exited without applying the modification.
[Delete entry]	The content of the input field is deleted completely.
[off]	If not only numbers but also special values (e.g., off) can be entered, then the corresponding buttons will be shown to the right of the numerical keypad.
[R1]	For many parameters, a result previously defined in the method can also be entered in place of a number (see chapter 32.7, page 449). You can select the result variable by touching [R1] .




NOTICE

A commercially available USB keyboard can be connected to make text and numerical input easier. The key assignment is described in chapter 11.10, page 132.

7 System settings

Main dialog: **System ► System settings**

This chapter describes the various system settings and configurations.

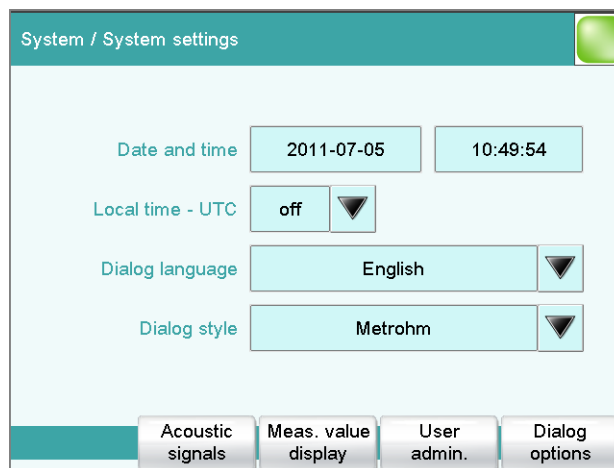

- Selecting the dialog language (see chapter 7.1.1, page 41).
- Setting the date and time (see chapter 7.1.2, page 42).
- Defining system-specific dialog options (see chapter 7.2, page 43).
- User administration (see chapter 7.3, page 46).
- Defining settings for the measured value display (see chapter 7.4, page 59).
- Configuring settings for acoustic signals (see chapter 7.5, page 59).

7.1 General system settings

7.1.1 Selecting the dialog language

The user interface is available in several languages. In addition to the two default dialog languages *English* and *German*, additional languages can be selected.

Proceed as follows to select the dialog language:

1 Opening the system settings

- In the main dialog, tap on **[System]**.
- Tap on **[System settings]**.

The **System / System settings** dialog is displayed.



2 Selecting the dialog language

- Tap on the **Dialog language** list box and select the desired language.

3 Saving the settings

Tap on the fixed keys [] or [].

The main dialog is displayed in the respective dialog language.

7.1.2 Setting the date, time and local time

The Ti-Touch displays the date and time in accordance with ISO standard 8601.

Proceed as follows to set the date and time:

1 Opening the system settings

- In the main dialog, tap on **[System]**.
- Tap on **[System settings]**.

The **System / System settings** dialog is displayed.

2 Entering the date

- Tap on the input field for the date.
The editor opens.
- Enter the current date in the format **YYYY-MM-DD** and confirm with **[OK]**.
The arrow keys [] and [] are used to move the cursor to the left or to the right by one character.

The input is saved and the editor is closed.

3 Entering the time

- Tap on the input field for the time.
The editor opens.
- Enter the current time in the format **hh:mm:ss** (24-hour format) and confirm with **[OK]**.
The arrow keys [] and [] are used to move the cursor to the left or to the right by one character.

The input is saved and the editor is closed.

4 Enter the local time

- Tap on the list box **Local time - UTC** and select the difference from the UTC (Coordinated Universal Time).
The selection **off** means that the time is saved with no difference from the UTC.


5 Saving the settings

Tap on the fixed keys [] or [].

The time settings are saved.

7.2 System-specific dialog options

Main dialog: **System ► System settings ► Dialog options**

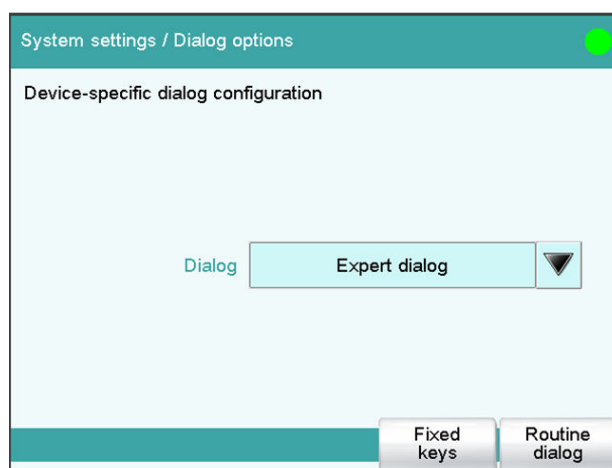

If you work without a login function, in this dialog you can define whether the system should generally be operated in expert mode or in routine mode. If you work with the login function activated, you must define this setting separately for each user (*see chapter 7.3.1, page 48*).

Dialog

Dialog mode in which the user may operate the system.

Selection	Expert dialog Routine dialog
Expert dialog	All functions that are supported by the system are available.
Routine dialog	The user dialog can be limited for routine operations. Only selected functions are available (<i>see Configuring the routine dialog, page 45</i>).

[Fixed keys]

Block unneeded fixed keys (see "Blocking unneeded fixed keys", page 44).

[Routine dialog]

Configure functions for the routine dialog (see "Configuring the routine dialog", page 45).

Selecting the dialog mode

Proceed as follows to change the dialog mode:

1 Select the dialog mode

Open the selection list **Dialog** and select either **Expert dialog** or **Routine dialog**.

2 Saving the settings

Tap on the fixed keys [] or [].

The setting will apply to all dialogs.



NOTICE

If you have selected **Routine dialog** and if the routine dialog was configured in such a way that the dialog **System settings / Dialog options** is blocked, then you can switch back over to the expert dialog as follows:

- Operation without login function:
In the main dialog, enter **User = Metrohm**.
- Operation with login function:
A user who works with expert dialog must log in.

Blocking unneeded fixed keys

This following configurations apply for **both** dialog modes: routine dialog and expert dialog.

Blocking fixed keys

Proceed as follows to block unneeded fixed keys:

1 Display fixed keys which can be blocked

Tap on the **[Fixed keys]** button.

All fixed keys which can be blocked are displayed.

2 Deactivate fixed keys

Deactivate those fixed keys which are not permitted to be used.

3 Saving the settings

Tap on the fixed keys [] or [].

Deactivated fixed keys cannot be used.

Configuring the routine dialog

A suitable **Standard configuration** has already been saved for routine operations:

- Methods can only be loaded, but not modified or created.
- Determinations cannot be recalculated.

You can readjust this standard configuration by disabling additional functions or re-enabling disabled functions.



NOTICE


The configuration of the routine dialog applies for all routine users. You also have the option of defining user-specific routine settings. To do this, you must work with the login function activated and create an identification profile for each user (see chapter 7.3.2, page 51).

Proceed as follows to modify the configuration for the routine dialog:

1 Opening the dialog

Tap on the [**Routine dialog**] button.

The list of all buttons in the main dialog, in the manual control, etc. is displayed:

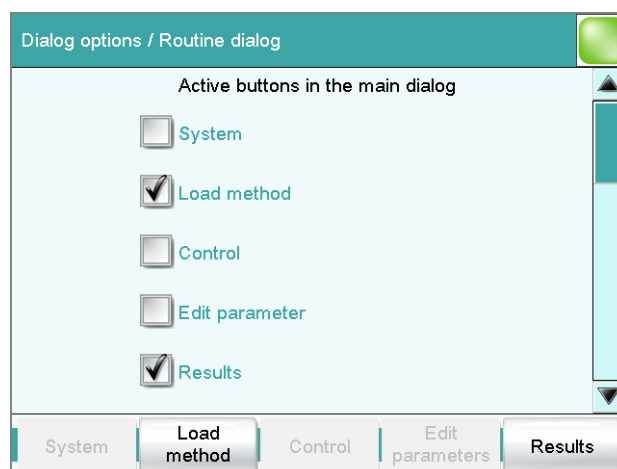

2 Deactivate buttons

Deactivate those buttons which are not permitted to be used.

All deactivated buttons will appear grayed-out, i.e. they are inactive.

3 Deactivate other functions

Many additional buttons and parameters can be disabled in the dialogs of **[System]**, **[Load method]**, **[Control]**, **[Edit parameters]** and **[Results]**. The corresponding option must be activated in order to enable these buttons.

4 Saving the settings

Tap on the fixed keys [] or [].

All deactivated functions will appear grayed-out, i.e. they are inactive.

7.3 User administration

Main dialog: **System ► System settings ► User admin.**

All of the functions of the user administration are described in this chapter. No user administration is mandatory for the simple operation of a titration system. If, however, you wish to be in compliance with *FDA Guidance 21 CFR Part 11*, then you must use the functions of the user administration. Additional information regarding the requirements of the FDA Guidance can be found in the document *Compliance Guide Touch Control*.



NOTICE

If you work with the login function activated, then the user administration is accessible only for users with administrator rights. This means that you must ensure that at least two users have administrator rights so at least one of them will be available. Keep the access rights for a user with administrator rights in a safe place so that they are accessible in an emergency.

System settings / User administration		
User	Dialog	Status
Administrator	Expert dialog	active
ak	Routine dialog	active
ga	Routine dialog	active
jb	Expert dialog	active
wl	Expert dialog	active

The following data is displayed in the user list for each user:

- Name
- Dialog mode in which the user may operate the system.
- Status

You can use the user list two different ways:

- Operation with login function:
If you work with the login function activated, i.e. if each user must log into the system before starting work, then only those users entered in the list can log in. The user who is currently logged in is shown in the main dialog.
- Operation without login function:
If you work without the login function, the users whose names are entered in the list can be selected in the main dialog or a user name can be entered. This makes it possible to document who has operated the titration system.

[Login options]

Define the settings for the login, Audit Trail, etc. (see chapter 7.3.3, page 52).

[Create ID profile]

Create an identification profile for the selected user on a storage medium (see chapter 7.3.2, page 51).

[New]

Adding a new user to the list (see chapter 7.3.1, page 48).

[Delete]

Delete the selected user from the list.



NOTICE

Once users have worked with the login function and password protection, they can no longer be deleted, even if the password protection is disabled again. The status of these users must be set to **inactive** (requirement of FDA Guidance 21 CFR Part 11).

The last user with administrator rights cannot be deleted.

[Edit]

Editing the data of the selected user (see chapter 7.3.1, page 48).

7.3.1 Editing the user configuration

User list: **User** ▶ **New / Edit**

User

The designation of the user is used for unambiguous identification, e.g. the company internal shorthand symbol or the personal number. The user name is printed out in all reports containing determination data and stored in the determination file. Each file always contains the name of the user who created it and the name of the last user to edit it.

Entry	24 characters maximum
Default value	empty

Full name

Complete name of the user.

Entry	24 characters maximum
Default value	empty

Dialog

Dialog mode in which the user may operate the system.

Selection	Expert dialog Routine dialog
-----------	--

Expert dialog

All functions that are supported by the system are available.

Routine dialog

The user dialog can be limited for routine operations. Only selected functions are available (*see Configuring the routine dialog, page 45*).

Status

Status of the user. Users can be deactivated. This function is useful, for instance, if the user is no longer authorized to operate the system or no longer works for the company.

Selection	active inactive
Default value	active

active

The user is authorized to operate the system.

inactive

The user is not authorized to operate the system and can no longer log in.

Admin. rights

on | **off** (Default value: **off**)

If this option is activated, then the user has administrator rights.



NOTICE

The last user with administrator rights cannot be deleted anymore.

[Favorites]

Saving methods and sample tables as user-specific favorites (*see chapter 19, page 205*).

[Signature method]

Assign the permission to use and sign methods (*see "Edit user / Signature method" dialog, page 49*).

[Signature determ.]

Assign the permission to use and sign determinations (*see "Edit user / Signature determination" dialog, page 50*).

"Edit user / Signature method" dialog

User: **Edit** ► **Signature method**



Authorizations for using and signing methods can be defined in this dialog. These settings are only effective if you are working with the login function enabled and password protection.

Use only released methods

on | off (Default value: **off**)

If this option is activated, then the user may only start methods that have been released.

Review methods (signature level 1)

on | off (Default value: **off**)

If this option is activated, then the user may only sign methods which have the status **saved**. The method is assigned the status **reviewed**.

Release methods (signature level 2)

on | off (Default value: **off**)

If this option is activated, then the user may only sign methods which have the status **reviewed**. The method is assigned the status **released**.



NOTICE

If the two options **Review methods** and **Release methods** are selected for a user, then this user can sign different methods at either level 1 or level 2, but not the same method at both level 1 and level 2.

Delete signatures

on | off (Default value: **off**)

If this option is activated, then the user may delete the signatures of a method which has been released. The method is assigned the status **saved**. The signatures of a method can only be deleted if it has the status **released**.

"Edit user / Signature determination" dialog

User: **Edit ▶ Signature determ.**

Authorizations for signing determinations can be defined in this dialog. These settings are only effective if you are working with the login function enabled and password protection.

Review determinations (signature level 1)**on | off** (Default value: **off**)

If this option is activated, then the user may only sign determinations at the first level. The determination is assigned the status **reviewed**.

Release determinations (signature level 2)**on | off** (Default value: **off**)

If this option is activated, then the user may only sign determinations at the second level. The determination is assigned the status **released**.

Delete signatures**on | off** (Default value: **off**)

If this option is activated, then the user may delete the signatures of a determination which has been released. The signatures of a determination can only be deleted if it has the status **released**.

7.3.2 Creating an identification profileUser list: **User ► Create ID profile**

If you plan to carry out the login with an identification profile (*see chapter 7.3.3, page 52*), then you must first create an identification profile on a storage medium for each user. A check can then be made at the time of login as to whether or not the user does exist and whether or not he or she is working in the expert dialog or in the routine dialog. After a successful login the routine dialog settings stored on the card are loaded.

**NOTICE**

In addition to the user name, the current routine dialog settings are also stored in this identification profile. This means you can define user-specific routine dialog settings for each user. However, you must configure them (*see "Configuring the routine dialog", page 45*), **before** you create the identification profile.

Before you create the identification profile, check whether the desired routine dialog settings are active.

1 Plug in a storage medium

Plug in the USB storage medium.

2 Select a user

In the user list, select the user for whom the profile is to be created.

3 Create the identification profile

Tap on **[Create ID profile]**.

The user configuration and the current routine dialog settings are saved.

7.3.3 Defining login options

Main dialog: **System ► System settings ► User admin. ► Login options**

There are a number of different ways to log onto the system:

- Without login
- Login via user name
- Login via user name and password
- Login via identification profile
- Login via identification profile and password



NOTICE


If you work with the login function activated, then the user administration is accessible only for users with administrator rights. This means that you must ensure that at least two users have administrator rights so at least one of them will be available. Keep the access rights for a user with administrator rights in a safe place so that they are accessible in an emergency.



NOTICE

If you exit this dialog with [] or [], and if you have selected one of the login variants **Login via user name** or **Login via identification profile**, then the login dialog will open automatically and you must also log in to the system.

This means that you must make sure that you have first defined all the users and created the identification profiles before you activate the login function.

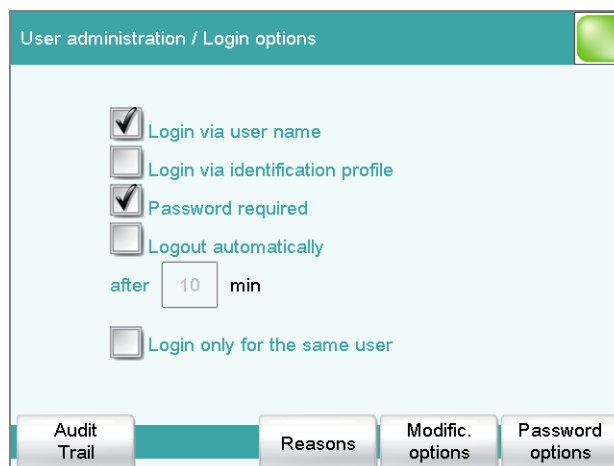

Login via user name

on | off (Default value: **off**)

If this option is activated, then the user must log in with his or her unambiguous identification.

Login via identification profile

on | off (Default value: **off**)

If this option is activated, then the login will take place via USB storage medium with the identification profile stored on it.

Password required

on | off (Default value: **off**)

If this option is activated, then the user must enter a password in addition to his or her user name or identification profile.

Logout automatically

on | off (Default value: **off**)

If this option is activated, then the user will be logged out automatically after the specified time.

Input range	1 to 60 min
-------------	--------------------

Login only for the same user

on | off (Default value: **off**)

If this option is activated, then only the same user may log in again after he or she has logged out. Users with administrator rights can, however, log in at any time.

[Audit Trail]

Define the settings for recording an Audit Trail (*see chapter 7.3.7, page 57*).



[Reasons]

Predefining a list of reasons which can be selected for the modification/ signature of a method or determination (see chapter 7.3.6, page 56).

[Modific. options]

Defining the modifications for which a reason is required (see chapter 7.3.5, page 56).


[Password options]

Define the settings for the password, see following chapter.

7.3.4 Password options

Main dialog: **System ► System settings ► User admin. ► Login options ► Password options**

You can make various settings for password entry in the password options.

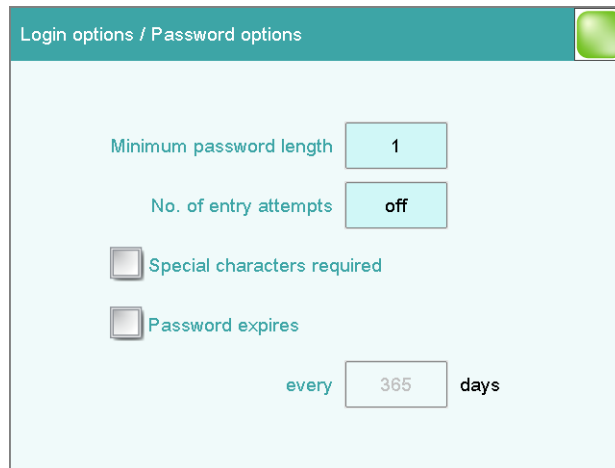

Minimum password length

Minimum number of characters of the passwords.

Input range	1 to 10
Default value	1

No. of entry attempts

If the user has logged in incorrectly this many times, then it will automatically be deactivated. It can only be reactivated by a user with administrator rights.

Input range	2 to 5
Selection	off
Default value	off

Special characters required

on | off (Default value: **off**)

If this option is activated, then the password must contain one of the following special characters: ° § + | @ * # ç % & ¬ () = ' ^ ` ~] [] { - _ : . ; , > < £ !

Password expires

on | off (Default value: **off**)

If this option is activated, then the user must define a new password after the time specified. A password that has already been used cannot be used again.

Input range	1 to 999 days
Default value	365 days

Forgotten password



CAUTION

If a user has forgotten his password, a new user name must be defined. The same user name can only be used again after a re-installation and re-creating the user list.

Proceed as follows:

1 Create a backup

Create a backup (see chapter 12.3, page 142).

2 Executing a system initialization

Switch off the Ti-Touch and execute a system initialization (see chapter 30.1, page 435).

3 Restore backed-up data

Use the function **Restore** to reload the data from the backup into your system (see chapter 12.3.1, page 142).

Deactivate the options **User list** and **System settings / User admin..**

4 Switch on the Ti-Touch

Switch the Ti-Touch back on again after a few seconds.

5 Restore the user list

Create the user list again and redefine the login options.

7.3.5 Modification options

Main dialog: **System ► System settings ► User admin. ► Login options ► Modific. options**

In the dialog **Login options / Modification options**, you can define for which actions a reason must be entered. These reasons are documented in the Audit Trail (*see chapter 7.3.7, page 57*) together with the modification. The reason for the last modification is shown in the properties of the method or determination.



NOTICE

The reasons are only requested when working with activated login function and password.

Saving modified method

on | off (Default value: **off**)

If this option is activated, then a reason must be given when saving a method modification.

Recalculating determination

on | off (Default value: **off**)

If this option is activated, then a reason must be given when determinations are being recalculated.

7.3.6 Reasons

Main dialog: **System ► System settings ► User admin. ► Login options ► Reasons**

In the dialog **Login options / Reasons**, you can create a selection list containing reasons from which a selection can be made when signing and modifying methods and determinations. Some reasons have already been provided.

**[New]**

Add a new reason to the list.

[Delete]

Delete the selected reason from the list.

[Edit]

Change the designation of the selected reason.

7.3.7 Audit Trail

Main dialog: **System ► System settings ► User admin. ► Login options ► Audit Trail**

An Audit Trail is an automatically generated log of all user activities. An Audit Trail contains precise logs of user actions (date, time, user, action, etc.). Recording an Audit Trail is important for compliance with *FDA Guidance 21 CFR Part 11* when using PC programs. A step-by-step description of how you must proceed to remain in compliance with the FDA Guidance is included in the document *Compliance Guide Touch Control*. The Audit Trail is saved to the internal memory.

You can also use the Audit Trail function specifically to record the data which are of interest to you.

You can view, filter and export the Audit Trail on a PC with the supplied software program *AuditTrailViewer*. Details concerning utilization of the *AuditTrailViewer* can be found in the Appendix (*chapter 32.10, page 468*).

If necessary, you can delete all of the entries in the Audit Trail with **[Delete Audit Trail]**. Do not fail to create a backup before doing so, however.

7.4 Measured value display

Main dialog: **System ► System settings ► Meas. value display**

The number of decimal places can be defined for pH values and voltages.

This setting refers exclusively to the display of the measured values in the live display and in the manual control. The values will, however, always be stored with their full accuracy.

7.5 Acoustic signals

Main dialog: **System ► System settings ► Acoustic signals**

You can define acoustic signals in order to direct attention to particular events. You can define signals for the following events:

- **Wrong manipulation**
An acoustic signal will sound each time an invalid action is attempted (e.g. pressing [?] again when Help is open).
- **Display a message**
A short beep will sound each time a message appears on the display. This informs the user that the message must be confirmed.
- **Button contact**
Each time a button is touched on the touch screen, this will be confirmed by an acoustic signal.
- **External data input**
An acoustic signal will provide confirmation each time data is received from external devices (e.g. balance, barcode reader).

8 Titrants

Main dialog: **System ► Titrants**

This chapter describes how you can create a list of titrants used in the system. Titrants can be used in intelligent buret units or in nonintelligent buret units. Intelligent buret units have a built-in data chip on which the data for the titrant is stored. This data is automatically read out during attachment and entered in the titrant list.

System / Titrants			
Titrant	Cyl.	Type	Dos.device
c(AgNO ₃) = 0.1 mol/L	10 mL	IDU	
c(HCl) = 0.1 mol/L	10 mL	IDU	D1/Ti-Touch
c(NaOH) = 0.1 mol/L	5 mL	IDU	

New Delete Edit

The titrant list can contain a maximum of 30 titrants. The following data is specified for each titrant:

- Designation
- Cylinder volume
- Type
 - **IDU**: dosing unit with integrated data chip
 - **IEU**: exchange unit with integrated data chip
 - **EU**: exchange unit without data chip
- MSB connector of the dosing device/control instrument (only when exchange unit/dosing unit is attached)

Titrants in exchange units/dosing units with integrated data chips are depicted in green lettering.

The following titrant data is stored in the titrant list:

- Name
 - Each titrant in the system is identified by its unambiguous name.
- Concentration
- Current titer
- Working life

- Data on exchange/dosing unit:
 - Parameters for the function **PREP**
 - Length and diameter of the tubings
 - Port assignment of the dosing unit
 - Cylinder volume
 - Serial number
 - etc.
- etc.



NOTICE

If data is read out from the data chip, then a check is made whether the titrant list already contains a titrant of the same type with the identical serial number. If this is the case, then the older data set will **always** be overwritten by the new data set, no matter whether the data set in the titrant list or the data set on the data chip is the most recent one.

[New]

Adding a new titrant to the list (see chapter 8.1, page 61).

[Delete]

Delete the selected titrant from the list.

[Edit]

Editing the data of the selected titrant (see chapter 8.2, page 62).

8.1 Adding a new titrant

Before you can use a titrant, you must add it to the titrant list. To do this, use the button **[New]**.

- Exchange unit/dosing unit with data chip:
The exchange unit or dosing unit must be attached. All dosing devices on which non-configured exchange or dosing units have been detected are included in a selection list. Tapping on the button **[Edit]** opens the properties dialog, see following chapter.
- Exchange unit without data chip:
The properties dialog is opened after the dosing unit type has been selected, see following chapter.

8.2 Editing titrant data

Titrant list: **Titrants** ▶ **New / Edit**

All of the data for the selected titrant is displayed in the dialog **Titrants / Edit**.

Titrant

The designation of the titrant is used for unambiguous identification.

Entry	24 characters maximum
Selection	Selection of frequently used titrants

Concentration

Concentration of the titrant.

Input range	-999999999 to 999999999
Default value	1.000

Unit of the concentration.

Entry	10 characters maximum
Selection	μmol/mL mmol/L mol/L g/L mg/L mg/mL μg/L ppm % mEq/L
Default value	mol/L

Comment

Entry	24 characters maximum
-------	------------------------------

Titer

Titer of the titrant.

Input range	-999999999 to 999999999
Default value	1.000

Unit of the titer.

Entry	10 characters maximum
Default value	empty
Selection	$\mu\text{mol/mL}$ mmol/L mol/L g/L mg/L mg/mL $\mu\text{g/L}$ ppm $\%$ mEq/L



NOTICE

If you modify the titer or the concentration of the titrant in a loaded determination at a later date and would then like to recalculate the determination with the corrected value, then you must modify the value in the determination data under **View data / Titrant data** (see "Calibration and titrant data", page 233).

Date titer det.

Date and time of the last titer determination. For new titrants, the time the preparation was made is specified until after the first time a titer determination has been carried out.

[Working life]

Defining the working life of the titrant (see chapter 8.3, page 64).

[Dosing unit]

This button is only displayed for **Type = IDU**.

Defining the properties of the dosing unit used (see chapter 8.4, page 65).

[Exchange unit]

This button is only displayed for **Type = IEU** or **EU**.

Defining the properties of the exchange unit used (see chapter 8.5, page 71).

[Titer options]

Displaying the properties for titer determination (see chapter 8.7, page 77).

8.3 Monitoring the working life

Titrant: **Edit ▶ Working life**

In the dialog **Edit titrant / Working life**, you can define the time interval after which the titrant must be replaced. This is particularly important if your titrant has a limited working life. If you do not wish to monitor the working life, then you can enter only the date of manufacture for documentation purposes.

Preparation date

Date on which the reagent was manufactured or the bottle was opened. For new titrants, the time the preparation was made will be specified.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the working life will be monitored.

Working life

If you define a time interval for the working life, then the **Expiry date** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Expiry date

If you define an expiry date, then the **Working life** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message

For all three options it is documented in the determination data (see dialog **More determination data / Messages**), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

8.4 Dosing unit

Titrant: **Edit ▶ Dosing unit**

Edit titrant / Dosing unit
 Titrant: c(I2) = 0.05 mol/L
 Name
 Order number 6.3032.250
 Serial number 10170552
 Cylinder volume 50 mL
 Serial no. cyl.
 Valve disk GLP test PREP param. Tubing param.

You can edit data for the dosing unit of the titrant in this dialog.

Name

Designation of the exchange or dosing unit.

Entry	24 characters maximum
Default value	empty

Order number

Order number of the exchange or dosing unit. It will be read out automatically on units with integrated data chips.

Entry	24 characters maximum
Default value	empty

Serial number

Serial number of the exchange or dosing unit. It will be read out automatically on units with integrated data chips.

Entry	8 digits maximum
-------	-------------------------



Cylinder volume

Cylinder volume of the dosing unit. It will be read out automatically on dosing units with integrated data chips.

Selection	2 5 10 20 50
Default value	20

Serial no. cyl.

Serial number of the dosing cylinder. It will be read out automatically on exchange or dosing units with integrated data chips. The number can be changed manually at any time, e.g. when a cylinder is replaced.

Entry	8 digits maximum
-------	-------------------------

[Valve disk]

Specifying the shift direction of the valve disk (*see chapter 8.4.3, page 71*).

[GLP test]

Defining the time interval for the GLP test (*see chapter 8.6, page 75*).

[PREP param.]

Entering the parameters for the preparation (*see chapter 8.4.1, page 66*).

[Tubing param.]

Entering the parameters for the connected tubing (*see chapter 8.4.2, page 68*).

8.4.1 Parameters for preparing (PREP) and emptying (EMPTY)

Titrant: **Edit ► Dosing unit ► PREP param.**

In the dialog **Dosing unit / PREP parameters**, you can adjust the parameters for the execution of the **Prepare** (command PREP) and **Empty** (command EMPTY) functions. The **Prepare** function is used to rinse the cylinder and tubing of the dosing unit and fill it air bubble-free. You should carry out this function before the first determination or once a day. The **EMPTY** function empties the cylinder and the tubings of the dosing unit.

Dosing port PREP/EMPTY

Dosing port through which the cylinder contents are ejected.

Selection	Dosing port 1 Dosing port 2 Fill port Special port
Default value	Dosing port 1

Dosing rate Dos. port 1

Rate used for the aspiration and ejection of the reagent via dosing port 1.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32.2, page 443*).

Dosing rate Dos. port 2

Rate used for the aspiration and ejection of the reagent via dosing port 2.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32.2, page 443*).

Dosing rate Fill port

Rate used for the aspiration and ejection of the reagent via fill port.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32.2, page 443*).

Dosing rate Spec.port

Rate used for the aspiration and ejection of the reagent via the special port.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32.2, page 443*).

8.4.2 Tubing parameters

Titrant: **Edit ► Dosing unit ► Tubing param.**

You can enter the length and diameter of the connected tubings in the dialog **Dosing unit / Tubing parameters**. The values which have already been entered correspond to the dimensions of the supplied standard tubings. In addition, the port assignment can be modified.



NOTICE

These parameters are important for the correct execution of the **Pre-paring** (PREP command) and **Emptying** (EMPTY command) functions, because the volumes of the tubing connections are taken into account.

Dosing port 1

Port

Port to be used as dosing port 1 for the **PREP** and **EMPTY** (see figure 22, page 70) functions.

Selection	Port 1 Port 2 Port 3 Port 4
Default value	Port 1

Length

Length of the tubing.

Input range	0.0 to 999.9 cm
Default value	40.0 cm
	The setting 0.0 means that this tubing will neither be rinsed nor emptied.

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

Dosing port 2

Port

Port to be used as dosing port 2 for the **PREP** and **EMPTY** functions (see figure 22, page 70).

Selection	Port 1 Port 2 Port 3 Port 4
Default value	Port 3

Length

Length of the tubing.

Input range	0.0 to 999.9 cm
Default value	0.0 cm
	The setting 0.0 means that this tubing will neither be rinsed nor emptied.

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

Fill port**Port**

Port to be used as fill port for the **PREP** and **EMPTY** functions (*see figure 22, page 70*).

Selection	Port 1 Port 2 Port 3 Port 4
Default value	Port 2

Length

Length of the tubing.

Input range	0.0 to 999.9 cm
Default value	25.0 cm
	The setting 0.0 means that this tubing will neither be rinsed nor emptied.

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

Special port**Port**

Port to be used as special port for the **PREP** and **EMPTY** functions (*see figure 22, page 70*).

Selection	Port 1 Port 2 Port 3 Port 4
Default value	Port 4

Length

Length of the tubing.



Input range	0.0 to 999.9 cm
Default value	0.0 cm
	The setting 0.0 means that this tubing will neither be rinsed nor emptied.

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

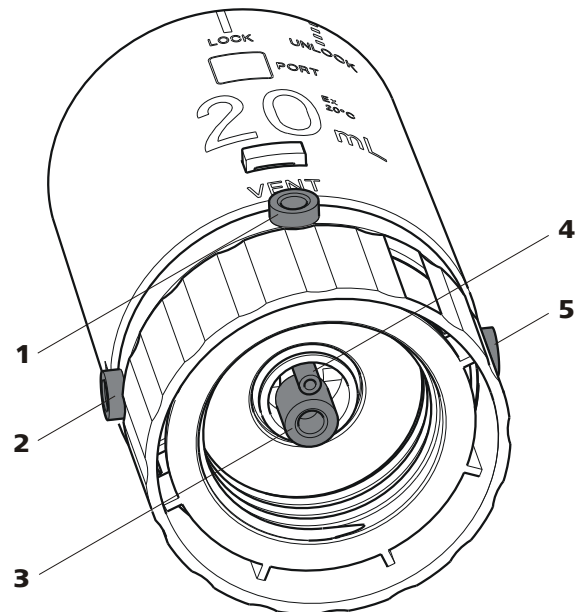


Figure 22 Dosing unit – port assignment

1 VENT

This port is set up for the deaeration of the reagent bottle. An adsorber tube (filled with desiccant) is usually mounted here.

3 Port 2

The default definition of this port is the filling port. A riser tube is usually mounted to it.

5 Port 3

The default definition of this port is dosing port 2.

2 Port 1

The default definition of this port is dosing port 1.

4 Port 4

Air is suctioned through this port during the **Emptying** function.

8.4.3 Shift direction of the valve disk

Titrant: **Edit ▶ Dosing unit ▶ Valve disk**

In this dialog, you can specify the shift direction of the valve disk.

Direction

Shift direction of the valve disc.

Selection	descending ascending automatic not over
Default value	automatic

descending

The ports are moved to in descending order.

ascending

The ports are moved to in ascending order.

automatic

The ports are moved to by the shortest path.

not over

Define a protected port.

Not over port

This parameter can only be edited with **Direction = not over**.

Define a protected port if the valve disc is not to be rotated over this port (useful with pipetting functions). The protected port can, however, be moved to directly.

Selection	1 2 3 4
Default value	4

8.5 Exchange unit

Titrant: **Edit ▶ Exchange unit**

Edit titrant / Exchange unit 🟢

Titrant: c(HCl) = 0.1 mol/L

Name

Order number 6.3026.150

Serial number 10950001

Cylinder volume 5 mL

Serial no. cyl.

GLP test
PREP param.
Tubing param.



You can edit data for the exchange unit of the titrant in this dialog.

Name

Designation of the exchange or dosing unit.

Entry	24 characters maximum
Default value	empty

Order number

Order number of the exchange or dosing unit. It will be read out automatically on units with integrated data chips.

Entry	24 characters maximum
Default value	empty

Serial number

Serial number of the exchange or dosing unit. It will be read out automatically on units with integrated data chips.

Entry	8 digits maximum
-------	-------------------------

Cylinder volume

Cylinder volume of the exchange unit. It will be read out automatically on exchange units with integrated data chips.

Selection	1 5 10 20 50
Default value	20

Serial no. cyl.

Serial number of the dosing cylinder. It will be read out automatically on exchange or dosing units with integrated data chips. The number can be changed manually at any time, e.g. when a cylinder is replaced.

Entry	8 digits maximum
-------	-------------------------

[GLP test]

Defining the time interval for the GLP test (*see chapter 8.6, page 75*).

[PREP param.]

Entering the parameters for the preparation (*see chapter 8.5.1, page 73*).

[Tubing param.]

Entering the parameters for the connected tubing (*see chapter 8.5.2, page 74*).

8.5.1 Parameters for the preparation (PREP)

Titrant: **Edit ► Exchange unit ► PREP param.**

In the dialog **Exchange unit / PREP parameters**, you can adjust the parameters for the execution of the **Prepare** function (command PREP). This function is used to rinse the cylinder and tubings of the exchange unit and fill it air bubble-free. You should carry out this function before the first determination or once a day.

Volume

Volume of titrant dosed during a rinsing cycle.

Input range	0.00000 to 99999.9 mL
Selection	Cylinder volume
Default value	Cylinder volume

Cylinder volume

The entire cylinder volume is being dosed.

Cycles

Number of rinsing cycles. We recommend carrying out at least two rinsing cycles in order to remove all air bubbles.

Selection	1 2 3 4 5 6 7 8 9
Default value	2

Dosing rate

The rate at which dosing takes place.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 32, page 442*).



8.5.2 Tubing parameters

Titrant: **Edit ▶ Exchange unit ▶ Tubing param.**

You can enter the length and diameter of the connected tubings in the dialog **Exchange unit / Tubing parameters**. The values which have already been entered correspond to the dimensions of the supplied standard tubings.

Dosing tip

Tubing to the dosing tip (23-2).

Length

Length of the tubing.

Input range	0.0 to 999.9 cm
Default value	40.0 cm

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

Dosing cylinder

Tubing to the dosing cylinder (23-3).

Length

Length of the tubing.

Input range	0.0 to 999.9 cm
Default value	13.0 cm

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

Reagent bottle

Tubing to the reagent bottle (23-1).

Length

Length of the tubing.

Input range	0.0 to 999.9 cm
Default value	25.0 cm

Diameter

Diameter of the tubing.

Input range	0.0 to 9.9 mm
Default value	2.0 mm

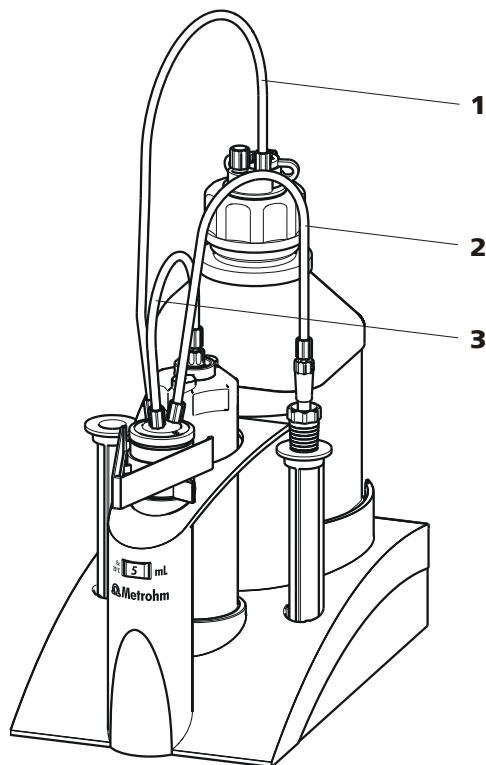


Figure 23 Exchange unit – tubing connections

1 Tubing to the reagent bottle

2 Tubing to the dosing tip

3 Tubing to the dosing cylinder

8.6 GLP test for exchange unit and dosing unit

Titrant: **Edit ► Dosing unit / Exchange unit ► GLP test**

In the dialog **Exchange unit / GLP test** or **Dosing unit / GLP test**, respectively, you can define the time interval after which a GLP test must be carried out again for the exchange unit or dosing unit.

GLP test date

Date on which the last GLP test was carried out.

Format: YYYY:MM:DD



Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which a GLP test has to be carried out again will be monitored.

GLP test interval

If you define a time interval for the GLP test, then the date in **Next GLP test** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next GLP test

If you define a date for the next GLP test, then the **GLP test interval** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

8.7 Titer determination options and data

Titration: **Edit ► Titer options**

Detailed information concerning the titer determination is displayed in the **Edit titrant / Titer options** dialog:

- **Titer method**
Method by which the titer was determined. If the titer was entered manually, then **manual** will be displayed.
- **User**
User who carried out the titer determination.
- **Statistical data**
The following information is also displayed for automatically assigned titers if the mean value of the results has been saved as the titer (*see "Save as titer", page 168*):
 - **n (titer det.)**
Number of titer determinations.
 - **s abs**
Absolute standard deviation
 - **s rel**
Relative standard deviation

[Validity]

Defining the time interval for the titer validity (*see chapter 8.7.1, page 77*).

[History]

Displaying information about the last ten titer determinations (*see chapter 8.7.2, page 78*).

8.7.1 Titer validity

Titration: **Edit ► Titer options ► Validity**

In the dialog **Titer options / Validity**, you can define the time interval after which the titer must be determined again.

Date titer det.

Date and time of the last titer determination. For new titrants, the time the preparation was made is specified until after the first time a titer determination has been carried out.

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the titer validity will be monitored.



Titer validity

If you define a time interval for the validity of the titer, then the date in **Next titer determ.** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next titer determ.

If you define a date for the next titer determination, then the time interval for the **Titer validity** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

8.7.2 Properties of the previous titer determinations

Dialog "Titer options / History"

Titant: **Edit ► Titer options ► History**

The date, time and titer of the last ten titer determinations are displayed in tabular form in the dialog **Titer options / History**. Titters that were determined automatically will be displayed in green; manually entered titer values will be displayed in black with the designation **(m)**. You can delete these entries, e.g. if you have opened a new bottle.

**NOTICE**

If multiple determinations have been carried out in order to determine the titer, then only one entry will be made in the history.

[Delete History]


Delete the entire history.

[Graph]

Open the diagram of the titer values, see following chapter.

Dialog "History / Graph"

Titrant: **Edit ▶ Titer options ▶ History ▶ Graph**

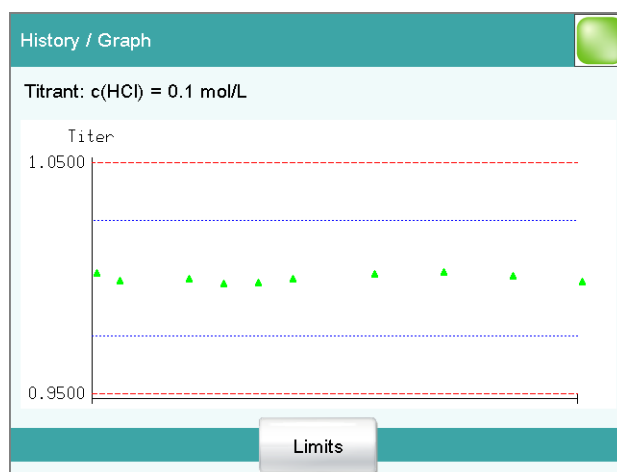

In this diagram, titer values are plotted against the date of the titer determination. You can define warning limits (blue dashed lines) and intervention limits (red dashed lines). These limits will not, however, be monitored.


[Limits]

Define warning and intervention limits.

9 Reagents

Main dialog: **System ► Reagents**

This chapter describes how you can create a list of the reagents used in the system.

System / Reagents 	
Reagent	Type
Reagent A	Volumetric
Reagent B	Volumetric
Reagent C	Volumetric
<div style="display: flex; justify-content: space-around; margin-top: 10px;"> New Delete Edit </div>	

The list of reagents gives the designation and type of each reagent.

[New]

Add a new reagent to the list, see following chapter.

[Delete]

Delete the selected reagent from the list.

[Edit]

Edit the data of the selected reagent, see following chapter.

9.1 Editing reagent data

Main dialog: **System ► Reagents ► New / Edit**

Reagent

The designation of the reagent is used for unambiguous identification.

Entry **24 characters maximum**

Comment

Entry **24 characters maximum**

[Reagent monitoring]

Set the parameters for the reagent monitoring, see following chapter.

9.2 Reagent monitoring

The conditions for the monitoring of the reagent are defined in the dialog **Edit reagent / Reagent monitoring**.

If one of the following values is reached, then the reagent must be replaced. The values are checked in the following cases:

- at the start of the determination.



- at the end of the determination.

Number of determ.

The number of determinations to be carried out with a certain amount of reagent depends on the type of sample and its amount.

Input range	1 to 999
Selection	off
Default value	off

Working life

Working life of the reagent.

Input range	1 to 999 days
Selection	off
Default value	off

Volume

Volume of titrant dosed.

Input range	1.0 to 999.9 mL
Selection	off
Default value	off

[Status]

Display the status overview of the current values of the reagent monitoring.

[Reagent replacement]

Edit the parameters for the reagent exchange.

"Reagent monitoring / Status" dialog

The current reagent monitoring values are displayed in this dialog.

[Reset]

Reset the values to zero.

Dialog "Reagent monitoring / Reagent replacement"

The parameters for the reagent replacement are defined in this dialog.

Reagent replacement

The reagent can either be exchanged manually or automatically.

Selection	manual auto
Default value	manual

manual

If a monitored parameter has reached the limit set, a message is being displayed. Then the reagent has to be exchanged manually.

auto

If a monitored parameter has reached the limit set, the method defined below is started automatically.

Memory

This parameter can only be edited with **Reagent replacement = auto**.

Memory location the method is loaded from. All memory locations are selectable, even if they are currently not accessible.

Selection	Internal memory External memory 1 External memory 2 Shared memory
Default value	Internal memory

Shared memory

Shared directory in the network.

Method

This parameter can only be edited with **Reagent replacement = auto**.

Method used for emptying the titration cell.

**NOTICE**

Make sure that the memory is accessible.

Entry	32 characters maximum
Selection	Selection of stored methods

10 Sensors

Main dialog: **System ► Sensors**

This chapter describes how you can create a list of the sensors used in the system.

System / Sensors		
Sensor	Sensor type	Meas. input
Ecotrode Plus	pH	
iUnitrode with Pt1000	pH IS	I1/Ti-Touch
pH electrode	pH	
Ag Titrode	Metal	
Metal electrode	Metal	
Temperature sensor	Temp.	

Three standard sensors are defined in the sensor list: **pH electrode**, **Metal electrode** and **Temperature sensor**. These sensors cannot be deleted or renamed. A maximum of 25 additional sensors can be added to these sensors.

For each sensor, the following data is displayed in the sensor list:

- Designation
- Type
 - **pH**: pH electrode
 - **Metal**: Metal electrode
 - **Temp.**: Temperature sensor
 - **Other**: Other sensor
- Measuring input/control instrument (only for intelligent sensors if they are connected)

Intelligent sensors are also indicated by **IS** and are depicted in green lettering.

The following sensor data is stored in the list of sensors:

- Name
 - Each sensor in the system is identified by its unambiguous name.
- Calibration data (for pH sensors only)
- Calibration interval (for pH sensors only)
- Working life
- etc.



NOTICE

If data is read out from the data chip of an intelligent sensor, then a check is made whether the sensor list already contains a sensor of the same serial number. If this is the case, then the older data set will **always** be overwritten by the new data set, no matter whether the data set in the sensor list or the data set on the data chip is the most recent one.

[New]

Adding a new sensor to the list (*see chapter 10.1, page 85*).

[Delete]

Delete the selected sensor from the list.

[Edit]

Editing the data of the selected sensor (*see chapter 10.2, page 86*).

10.1 Adding a new sensor


Before you can use a sensor, you must add it to the sensor list. To do this, use the button **[New]**.

- Conventional sensors:
The properties dialog is opened after the sensor type has been selected, see following chapter. The following sensor types can be selected:
 - pH electrode
 - Metal electrode (Pt electrode, Ag Titrode, Ag/AgCl reference electrode, etc.)
 - Other sensor, e.g. Spectrosense
 - Temperature sensor
- Intelligent sensors (also known as iTrodes):
If the 854 iConnect with iTrode is connected to the Ti-Touch, then the sensor is automatically entered in the sensor list and can be configured, see following chapter.



10.2 Editing the sensor data

Sensor list: **Sensor ▶ New / Edit**

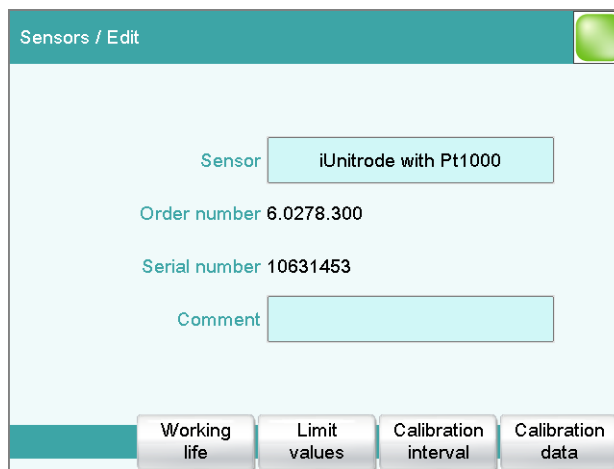

All of the data for the selected sensor is displayed in the dialog **Sensors / Edit**.

Sensor

The designation of the sensor is used for unambiguous identification.

Entry	24 characters maximum
-------	------------------------------

Order number

Order number of the sensor. With intelligent sensors it is read out automatically.

Entry	24 characters maximum
Default value	empty

Serial number

Serial number of the sensor. With intelligent sensors it is read out automatically.

Entry	8 digits maximum
-------	-------------------------

Comment

Entry	24 characters maximum
-------	------------------------------

[Working life]

Defining the working life of the sensor (*see chapter 10.3, page 87*).

[Limit values]

This button is displayed only for pH electrodes.

Defining the limit values for monitoring the slope and the electrode zero point (see chapter 10.5, page 92).

[Calibration interval]

This button is displayed only for pH electrodes.

Define the time interval for the next calibration (see chapter 10.6, page 93).

[Calibration data]

This button is displayed only for pH electrodes.

Display the properties for titer determination (see chapter 10.4, page 88).

10.3 Monitoring the working life

Sensor: **Edit ► Working life**

In the dialog **Edit sensor / Working life**, you can define the time interval after which the sensor must be replaced. If you do not wish to monitor the working life, then you can enter only the date of manufacture for documentation purposes.

Start-up

Date on which the sensor was used for the first time.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the working life will be monitored.

Working life

If you define a time interval for the working life, then the **Expiry date** will be tracked automatically.

Input range **1 to 999 days**

Default value **999 days**

Expiry date

If you define an expiry date, then the **Working life** will be tracked automatically.

Format: YYYY:MM:DD



Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message


In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

10.4 Calibration data (for pH electrodes only)

Sensor: **Edit ► Calibration data**

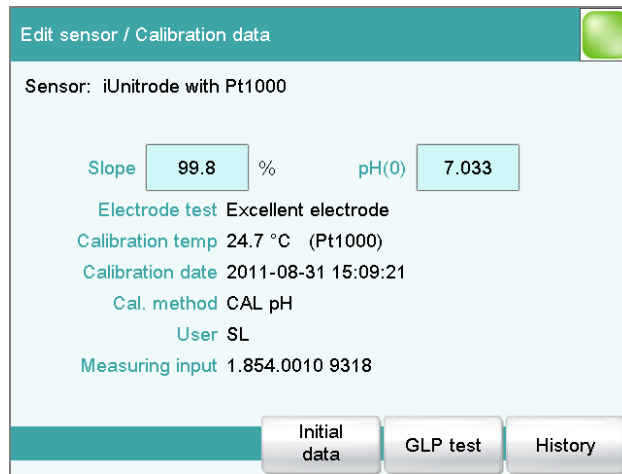

Detailed information concerning calibration is displayed in this dialog **Edit sensor / Calibration data**.

Slope

Slope of the electrode.

pH electrodes:

Input range	-999.9 to 999.9 %
Default value	100.0 %

pH(0)

This parameter is only visible with pH electrodes.

pH value of the electrode at 0 mV. Apart from the slope, pH(0) is the second characteristic of the calibration curve.

Input range	-20.000 to 20.000
Default value	7.000

The following data cannot be edited:

- **Electrode test** (only with intelligent sensors)
Result of the electrode test.
- **Calibration temp**
Temperature at which the calibration was carried out.
If the temperature was measured manually during the calibration, then **(manual)** will also be displayed. If the temperature was measured with a connected temperature sensor, then the sensor type (**(Pt1000)** or **(NTC)**) will be displayed.
- **Calibration date**
Date and time of the last calibration. For new sensors, the time the preparation was made is specified until after the first time a calibration has been carried out.
- **Cal. method**
Method with which the sensor was calibrated. If the calibration data was entered manually, then **manual** will be displayed.
- **User**
User who carried out the calibration.
- **Measuring input** (only with intelligent sensors)
The type and the serial number of the measuring input with which the calibration was carried out.

[Initial data]

This button is only displayed for intelligent sensors.

Display the initial calibration data determined at the time of the Metrohm quality control.

[GLP test]

Define the time interval for the GLP test (see "Dialog "Calibration data / GLP test"", page 90).

[History]

Display information about the last ten calibrations (see "Dialog "Calibration data / History"", page 90).

10.4.1 Properties of the previous calibrations

Dialog "Calibration data / History"

Sensor: **Edit ▶ Calibration data ▶ History**

The date, time and calibration data of the last ten calibrations are displayed in tabular form in the dialog **Calibration data / History**. Calibrations that were carried out automatically will be displayed in green; manually entered calibration data will be displayed in black with the designation **(m)**.

[Delete History]

Delete the entire history.

[Graph slope]


Open the diagram of the electrode slopes, see following chapter.

[Graph pH(0)]

Open the diagram of the electrode zero points, see following chapter.

Dialog "History / Graph"

Sensor: **Edit ▶ Calibration data ▶ History ▶ Graph Slope / pH(0)**.

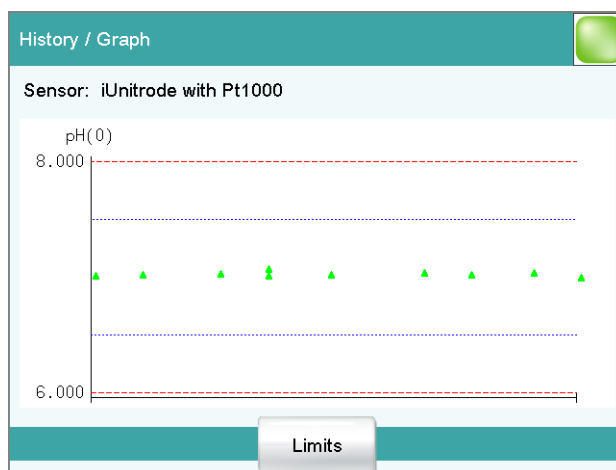

In this diagram, either slope or pH(0) is plotted against the date of the calibration. You can define warning limits (blue dashed lines) and intervention limits (red dashed lines). These limits will not, however, be monitored.

[Limits]

Define warning and intervention limits.

Dialog "Calibration data / GLP test"

Sensor: **Edit ▶ Calibration data ▶ GLP test**

In the dialog **Calibration data / GLP test**, you can define the time interval after which a GLP test must be carried out again for the sensor.

GLP test date

Date on which the last GLP test was carried out. After you have carried out an electrode test (ELT command), the date of the electrode test is automatically entered into this field. However, you also can enter the date manually.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which a GLP test has to be carried out again will be monitored.

GLP test interval

If you define a time interval for the GLP test, then the date in **Next GLP test** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next GLP test

If you define a date for the next GLP test, then the **GLP test interval** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.


Cancel determination

The determination is stopped.



10.5 Limit values for the calibration data

Sensor: **Edit ▶ Limit values**

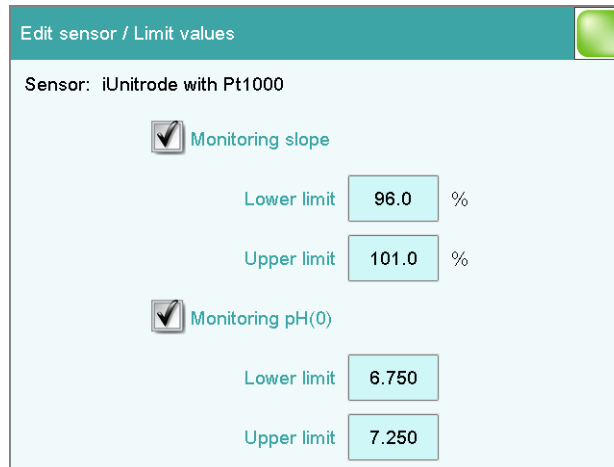

You can define the following limit values in the dialog **Edit sensor / Limit values**:

- Slope (pH electrodes)
- Electrode zero point (pH electrodes)

These values are monitored during the calibration. If these limits are infringed, then a message will be displayed and you can decide whether to accept the calibration data or not.

Monitoring slope

on | off (Default value: **off**)

If this parameter is activated, then the slope will be monitored.

Lower limit

pH electrodes:

Input range	-999.9 to 999.9 %
Default value	96.0 %

Upper limit

pH electrodes:

Input range	-999.9 to 999.9 %
Default value	101.0 %

Monitoring pH(0)**on | off** (Default value: **off**)

This parameter is only available with pH electrodes.

If this parameter is activated, then the electrode zero point pH(0) will be monitored.

Lower limit

Input range	-20.000 to 20.000
Default value	6.750

Upper limit

Input range	-20.000 to 20.000
Default value	7.250

10.6 Monitoring the calibration interval (for pH electrodes only)Sensor: **Edit ► Calibration interval**

In the dialog **Edit sensor / Calibration interval**, you can define the time interval after which the sensor must be recalibrated.

Calibration date

Date of the last calibration.

Monitoring**on | off** (Default value: **off**)

If this parameter is activated, then the validity of the calibration will be monitored.

Calibration interval

If you define a time interval for the validity of the calibration, then the date in **Next calibration** will be tracked automatically.

Input range	1 to 999 days
Default value	7 days

Next calibration

If you define a date for the next calibration, then the **Calibration interval** will be tracked automatically.

Format: YYYY:MM:DD



Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

11 Device manager

Main dialog: **System ► Device manager**

This chapter describes how you can configure the Ti-Touch, a Sample Processor and the connected peripheral devices. A detailed description of the hardware installation of the Sample Processor can be found in the corresponding installation manual.

System / Device manager	
Device name	Device type
916 Ti-Touch	Ti-Touch operation
Ti-Touch	Ti-Touch titrator
Sample Processor	815 Sample Processor
Printer	Printer
Balance	Balance
USB/RS-232 adapter	USB/RS-232 adapter

New Delete Edit

The instrument list gives the name and type of every instrument. The **Ti-Touch titrator** is a component of the 916 Ti-Touch and therefore permanently entered as control instrument in the device manager. A connected Sample Processor with the peripheral devices (dosing devices, stirrers, etc.) connected to the MSB connector is entered in the list automatically. A USB/RS-232 adapter will also be automatically recognized and entered in the instrument list with default settings. A printer is entered in the instrument list in the default settings. You must enter a PC keyboard, a barcode reader or a balance in the instrument list yourself.

[New]

Add a new device to the list (*see chapter 11.1, page 96*).

[Delete]

Delete the selected device from the list.



NOTICE

Devices which are recognized automatically cannot be deleted from the list while they are still connected.

[Edit]

Configure the selected device (*see chapter 11.2, page 96*).



11.1 Adding a new device

Of the following device types, you can enter one device each in the device list, even if it is not yet connected:

- Sample Processor
- Balance
- Barcode reader
- USB/RS-232 adapter
- PC keyboard

Proceed as follows:

1 Displaying the device selection

Tap on **[New]**.

2 Selecting the instrument

Tap on the button for the desired device.

The new device is entered in the list.

11.2 Configuring the instrument

Instrument list: **Instrument ► Edit**

The data stored for an instrument depends on the type of instrument. You can define an instrument name and a comment for each instrument. The instrument name of the Ti-Touch is printed out in the standard report header.

The description of the individual instruments can be found in the following specific chapters:

- Ti-Touch (*see chapter 11.3, page 97*)
- Metrohm control instruments (*see chapter 11.4, page 103*)
- Sample Processor (*see chapter 11.5, page 107*)
- Printer (*see chapter 11.7, page 124*)
- Balance (*see chapter 11.8, page 128*)
- PC keyboard (*see chapter 11.10, page 132*)
- USB/RS-232 adapter (*see chapter 11.9, page 130*)
- Barcode reader (*see chapter 11.11, page 133*)

11.3 Ti-Touch

Instrument list: **916 Ti-Touch** ► **Edit**

Device manager / Edit

Device type: Ti-Touch operation

Device name: 916 Ti-Touch

Comment:

Switch off display: off min

Program version: 5.916.0010

Serial number: 395612

Control Remote Box: Not available

E-mail PC/LIMS report Shared memory TCP/IP settings More information

Device name

This designation is used for identification purposes when selecting control devices (command, manual control).

Entry **24 characters maximum**

Comment

Entry **24 characters maximum**

Switch off display

If this time interval has expired without the Ti-Touch having been operated, then the display will be switched off. Touching the display will switch it back on again at any time.

Input range	1 to 999 min
Selection	off
Default value	off

Program version

Program version of the software.

Serial number

Shows the serial number of the device.

Control Remote Box

Shows to which control device and MSB connector the Remote Box is connected.



The "Control Remote Box" is the interface via which the system can be started and stopped externally. If multiple Remote Boxes are connected, then the one that is recognized first when the program starts will be used as the "Control Remote Box."

Selection	Name of the control device / Number of the MSB connector Not available
-----------	--

11.3.1 E-mail

916 Ti-Touch: **Edit ► E-mail**

The system allows you to send displayed messages as e-mails. The Ti-Touch must be connected to a network for this to function. The following types of messages can be sent:

- : general warning messages
- : error messages

Configuring e-mail dispatch

Proceed as follows so that messages can be sent as e-mails:

1 Activating the option

- In the instrument properties of the 916 Ti-Touch, tap on the **[E-mail]** button.
- In the **Edit device / E-mail** dialog, activate the **Send the following messages as e-mail:** option.

2 Configuring e-mail addresses



- Tap on the **[E-mail settings]** button.
The **E-mail / Settings** dialog is displayed.
- Enter the addresses of the mail server, the sender and the desired recipient.

Parameter description

Send the following messages as e-mail:

on | off (Default value: **off**)

If this parameter is activated, then messages with the following symbols will be sent as e-mails:

- : General warning messages
- : Error messages

Send only during running determination

on | off (Default value: **on**)

If this parameter is activated, then messages will be sent as e-mails only if a determination is running. Deactivate this parameter if messages are also to be sent in normal status.

Mail server

Address of the mail server for outgoing mail, e.g. mail.metrohm.ch. You can find the address of the mail server either in your e-mail program settings or obtain it from your IT department.

Entry	50 characters maximum
Default value	empty

Sender

E-mail address of the sender. This address must be formatted as an e-mail address, but need not necessarily correspond to an existing e-mail account, e.g. ti-touch@metrohm.com.

Entry	50 characters maximum
Default value	empty

Recipient


The messages will be sent to this e-mail address.

Entry	50 characters maximum
Default value	empty

11.3.2 PC/LIMS report

916 Ti-Touch: **Edit ► PC/LIMS report**

You can generate a machine-readable report with all of the important data concerning a determination, which is referred to as a PC/LIMS report. This report can be saved as a TXT file (according to ISO/IEC 8859-1) or as a UTF-8 file or sent to a terminal program or a LIMS via an RS-232 interface:

- manually with the [] fixed key (see chapter 27, page 257).
- automatically at the end of a determination (see chapter 16.5.6, page 197).

The file name of the report file is constructed as follows: *PC_LIMS_Report-ID1-YYYYMMDD-hhmmss.txt*. A detailed description of the contents of the PC/LIMS report can be found in the *PC/LIMS Report Guide*.

Memory

Memory location where the PC/LIMS report is stored as a file. The report will be saved in the directory *pc_lims_report*. This directory will be created



the first time a PC/LIMS report is generated. All three memory locations are listed as possible selections, even if they cannot be accessed at the moment.

Selection	off External memory 1 External memory 2 Shared memory
Default value	off

off

The report will not be saved as a file.

Shared memory

The report will be saved in a shared directory on the network. The shared directory is selected in the **Edit device / Shared memory** dialog (see chapter 11.3.3, page 100).

RS-232

The RS-232 interface via which the PC/LIMS report is sent. The interface parameters are adjusted in the **Edit device / Port parameters** dialog (see chapter 11.9, page 130).

Selection	off COM 1 COM 2
Default value	off

off

The report will not be sent via an RS-232 interface.

COM 2

This interface is inactive.

Coding

Format in which the PC/LIMS report is coded and stored.

Selection	ISO 8859-1 UTF-8
Default value	ISO 8859-1

ISO 8859-1

This format is recommended for all languages that use the extended ASCII code (e.g. German, English, Spanish, etc.).

UTF-8

This format is required for all languages that do not use the extended ASCII code (e.g. Russian, Chinese, Korean, etc.).

11.3.3 Shared memory

916 Ti-Touch: **Edit ► Shared memory**

If you have your Ti-Touch connected to your network, then you can define in this dialog a shared memory location on a computer within your network for the purpose of saving data (methods, determinations, etc.).

**CAUTION**

If the computer on which you share a memory location does not have the same subnet as the Ti-Touch, then a WINS server must be present. The computer must be entered in this server. In order for the communication to work, the SMB protocol version 1 (SMBv1) must be activated. Since this function is not standard anymore in Windows 10, it has to be activated manually. From firmware version 5.916.0044, the SMB protocol version 2 and 3 is also supported.

Computer

Host name of the computer on which a memory location is to be shared. If you are working with Windows, then you will find the host name of the computer as follows: In the Windows **command prompt** window, enter the command **ipconfig -all**. The host name is listed together with other parameters of the computer.

Do not under any circumstances enter an IP address in this input field.

Entry	max. 32 characters
Default value	empty

Share name

Share name of the shared memory location (file directory) on the above-specified computer. Please note that the Share name of a file directory often does not match the name of the file directory. The share name can be found on the Release tab in the Properties dialog of the shared file directory.

Entry	max. 32 characters
Default value	empty

Domain

Network domain in which the above-specified computer is located. If you are working with Windows, then you will find the domain name as follows: In the Windows **command window**, enter the command **ipconfig -all**. The domain is listed together with other parameters of the computer. Leave the field empty if the computer is not located within a domain.

Entry	max. 32 characters
Default value	empty

User

User name of the user authorized to access the shared memory location.



Entry	max. 32 characters
Default value	empty

Password

Password of the user configured on the computer.

Entry	max. 32 characters
Default value	empty

[Connect]

Establish the network connection. If the connection has been set up correctly, then all of the input fields will become inactive and the label switches to **[Disconnect]**. The network connection can be disconnected with this.

11.3.4 TCP/IP settings

916 Ti-Touch: **Edit ► TCP/IP settings**

If you have connected your Ti-Touch to your network, then you will have to define the network-relevant settings in this dialog. The Ti-Touch requires an IP address so that it is identifiable on the network. The Ti-Touch can acquire the IP address either dynamically from a DHCP server or you can enter the address directly.

Get IP address automatically (DHCP)

on | off (Default value: **on**)

If this parameter is activated, then the Ti-Touch will obtain its IP address directly from a DHCP server. In this case, the remaining parameters can no longer be edited.

IP address

IP address for the Ti-Touch. IP addresses are 32-bit numbers and are written as sequences of four decimals, each separated by a period, e.g. "10.157.212.8".

Entry	x.x.x.x
Default value	192.168.10.11
	"x" is a decimal between 0 and 255.

Subnet mask

The net mask or subnet mask, together with the IP address, indicates to which network the device to be connected belongs. Subnet masks are 32-bit numbers and are written as sequences of four decimals, each separated by a period.

Entry	x.x.x.x
Default value	255.255.255.0

Default gateway

IP address for the standard gateway. A gateway sets up connections to several networks. It is located in the same subnet as the device to be configured.

Entry	x.x.x.x
Default value	192.168.10.10

MAC address

Unambiguous address of the network adaptor for identification within a network. This parameter cannot be edited.

NetBIOS name

Unambiguous designation of the device within the network. This parameter cannot be edited.


11.4 Metrohm control instruments

The following Metrohm devices can be used as control instruments:

- Ti-Touch Titrator (permanently entered as control instrument in the device manager)
- Sample Processors

The following specifications apply for all of the control instruments mentioned above.

Additional device configurations for the Sample Processor are described in a separate chapter (*see chapter 11.5, page 107*).

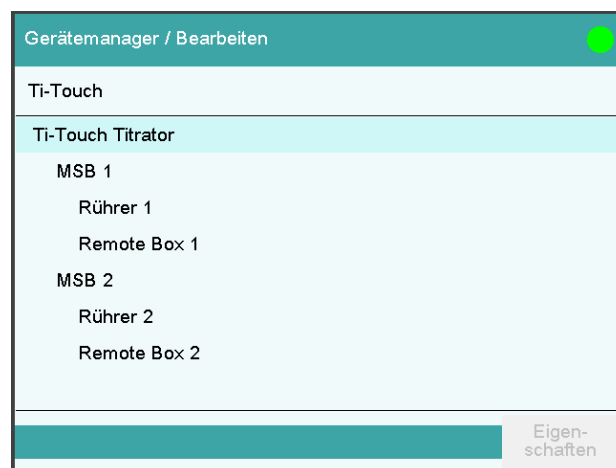

The dialog shows the control instrument with its measuring inputs, the MSB connectors and connected peripheral devices (dosing device, stirrer, Remote Box).

[Properties]

Open the properties dialog of the highlighted entry.

The following properties can be displayed and edited to a certain extent:

- Properties of the measuring inputs (*see chapter 11.4.1, page 104*)
- Properties of the MSB connectors (*see chapter 11.4.2, page 105*)
- Properties of the peripheral devices on the MSB connector (*see chapter 11.4.3, page 106*)

11.4.1 Properties – Measuring input

Instrument list: **Control instrument** ► **Edit** ► **Measuring input**
1 ► **Properties**

Device manager / Edit

Ti-Touch / Measuring input 1

ADC type 3.848.1210

Serial number 46180

Thick film type 3.680.0743

Temperature sensor

Type Pt1000 ▼ R (25 °C) 30000

B value 4100 K

In this dialog, you define the type of temperature sensor you have connected to the selected measuring input.

ADC type

Type of analog-digital converter.

Serial number

Serial number of the measuring interface.

Thick film type

Type of the thick film.

Temperature sensor

Type

The instrument supports the use of two different temperature measurement techniques:

- NTC (Negative Temperature Coefficient)
- Pt1000 (Platinum resistance)

Select here the type that has been connected to the instrument. If an NTC sensor is used, then it is also necessary to enter two sensor characteristics. These characteristics are listed in the specifications of the sensor.

Selection	Pt1000 NTC
Default value	Pt1000

R (25 °C)

This parameter can only be edited with **Type = NTC**.

Nominal resistance of the NTC sensor at 25 °C.

Input range	1000 to 99999 ohm
Default value	30000 ohm The default value applies to Metrohm sensors with an NTC sensor.

B value

This parameter can only be edited with **Type = NTC**.

Material constant of the NTC sensor. B values of NTC sensors are frequently based on different reference temperatures (usually 25 °C and 50...100 °C). When entering the B value the influence of the second reference temperature is negligible in comparison with the measuring accuracy of an NTC sensor.

Input range	1000 to 9999 K
Default value	4100 K The default value applies to Metrohm sensors with an NTC sensor. If no B value is given for your sensor then you can retain the default value.

11.4.2 Properties – MSB connector

Instrument list: **Control instrument** ► **Edit** ► **MSB connector**
1/2 ► **Properties**

Device manager / Edit

Ti-Touch / MSB 1

When would you like to be asked to prepare the dosing device on MSB 1?

Switch on

Attach an exchange or dosing unit

Time interval h



In this dialog, you can define when the request to carry out the **Prepare** function for connected dosing devices is to be displayed. This setting applies for all dosing devices of the selected MSB connector. The manual control contains a description of how to prepare the exchange unit and/or the dosing unit (*see chapter 28.3.3, page 276*).

Switch on

on | off (Default value: **on**)

If this parameter is activated, then you will be requested to prepare the dosing device when the Ti-Touch is switched on.

Attach an exchange or dosing unit

on | off (Default value: **on**)

If this parameter is activated, then you will be requested to prepare the dosing device when the exchange/dosing unit is attached.

Time interval

on | off (Default value: **off**)

Activate this parameter if you wish to receive a regular request to prepare the dosing unit.

Input range	0.1 to 999.9 h
Default value	12.0 h


11.4.3 Properties – Peripheral devices

Instrument list: **Control instrument ► Edit ► Peripheral device ► Properties**

The following data is displayed for the connected dosing devices and stirrers:

- Type
The display includes the device number, e.g. "800" (Dosino).
- Serial number

11.5 Sample Processor

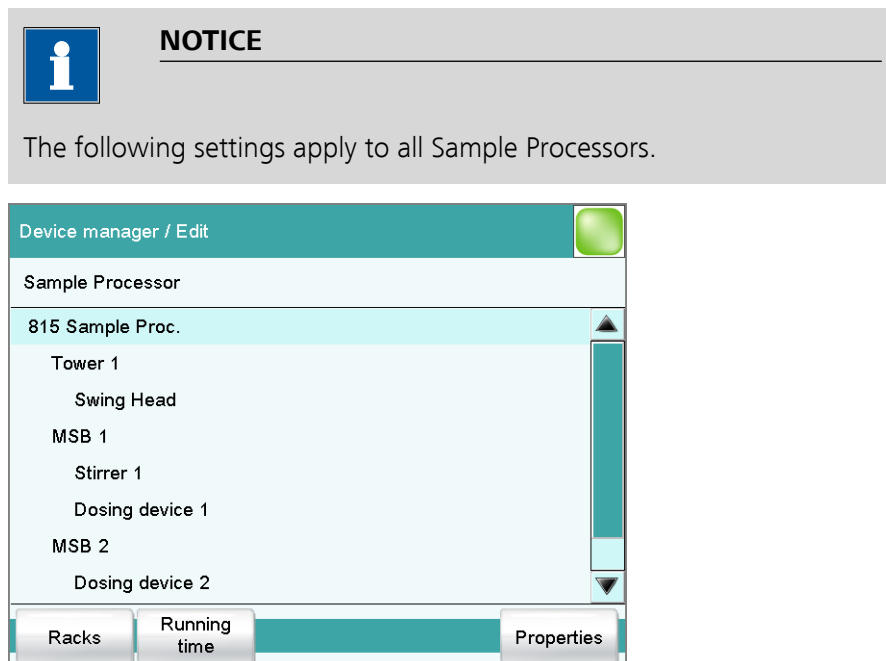

The dialog shows the Sample Processor with its tower (or towers), the MSB connectors and connected peripheral devices (dosing device, stirrer, Remote Box). If the control instrument is not connected, then the properties dialog of the control instrument will be displayed directly (*see chapter 11.5.1, page 108*).

The settings of the MSB connectors and the connected peripheral devices are described in *Chapter "Properties – MSB connector", page 105*.

[Racks]

Configure sample racks (*see chapter 11.6, page 115*).

[Running time]

Configure the running time meter.

The running time meter adds together the time while the Sample Processor is "busy" i.e. when an action is being carried out. A message with the prompt to service the device is displayed after expiry of the time limit set here. We recommend to carry out a service after 1,000 operating hours. A reset of the running time meter can only be carried out by a Metrohm service engineer.

[Properties]

Open the properties dialog of the highlighted entry.

11.5.1 Properties – Sample Processor

Instrument list: **Sample Processor ▶ Edit ▶ Properties**

Device name

This designation is used for identification purposes when selecting control devices (command, manual control).

Entry **24 characters maximum**

Comment

Entry **24 characters maximum**

Program version

Program version of the instrument software.

Serial number

Shows the serial number of the device.

Rack name

Name of the currently attached sample rack.

[Adjustment data]

Display the internal adjustment data (EEPROM data) of the Sample Processor.



CAUTION

Do not modify and save these settings under any circumstances. This data is required by the service engineer in case of positioning problems.

11.5.2 Properties – Tower

Instrument list: **Sample Processor ▶ Edit ▶ Tower 1/2 ▶ Properties**

Device manager / Edit		
Sample Processor / Tower 1		
Maximum stroke path	235	mm
Min. beaker radius	off	mm
Rotation increment	5.00	°
Axial distance	196.00	mm



NOTICE

In the case of a Sample Processor with 2 towers, Tower 1 is the tower on the right and Tower 2 is the tower on the left, when viewed from the front.

Maximum stroke path

Lowest permissible lift position for automatic and manual operation.

This is a safety setting. A correct specification of this value can prevent the glass of an electrode from breaking, because a titration head can not be moved below the position specified.

Input range	1 to 235 mm
Default value	235 mm

Min. beaker radius

Smallest permissible radius of the sample vessels for automatic and manual operation.

This is a safety setting. It is intended to prevent a completely loaded, wide titration head from being moved into a narrow sample vessel.

Input range	1.0 to 100.0 mm (Increment: 0.1) If the lift is to be moved to the work position, a check is made to see whether the beaker radius defined for the attached rack (see "Beaker radius samples", page 117) is less than this value.
Selection	off
Default value	off

off

No check takes place.



Rotation increment

Amount by which the sample rack can be rotated forward or backward relative to its current position. This parameter is required in the command **MOVE** for the settings **Destination = rotate +** and **rotate -**.

Input range	0 to 353.99 °
Default value	5.00 °

Axial distance

Distance between the axis of rotation of the sample rack and the swing axis of the robotic arm.

for 814 USB Sample Processor:

Input range	100.0 to 300.0 mm
Default value	166.0 mm

for 815 Robotic USB Sample Processor XL:

Input range	100.0 to 300.0 mm
Default value	196.0 mm

for 810 Sample Processor:

Input range	100.0 to 300.0 mm
Default value	166.0 mm

11.5.3 Properties – Swing Head

Instrument list: **Sample Processor ▶ Edit ▶ Swing Head ▶ Properties**

Edit device / Properties		
Sample Processor / Tower 1 / Swing Head 786		
Serial number 01108		
Ext. pos.	Angle [°]	Work [mm]
1	60.00	100
2	60.00	0
3	60.00	0
4	60.00	0
Ext. pos. 1-4	Robotic arm	Adjustment data
Edit		

The dialog **Edit device / Properties** shows a list of all external positions with the assigned swing angle and the specific work position for each.

[Ext. pos. 1-4]

Edit the settings that apply for all external positions (see "Properties – External positions 1-4", page 111).

[Robotic arm]

Edit the robotic arm settings (see "Properties – Robotic arm", page 112).

[Adjustment data]

Display the overview of the EEPROM data of the Swing Head. This dialog cannot be edited.

[Edit]

Edit specific settings of the selected external position (see "Properties – External position", page 113).

Properties – External positions 1-4

Swing Head: **Properties** ► **Ext. pos. 1-4**

Properties / Edit external positions 1-4

Sample Processor / Tower 1 / Swing Head 786

External positions 1-4

Shift position	<input type="text" value="0"/>	mm
Rinse position	<input type="text" value="0"/>	mm
Swing increment	<input type="text" value="10.00"/>	°

Properties which apply to all four external positions can be defined in the dialog **Properties / Edit external positions 1-4**.

Shift position

Lift position at which the robotic arm rotates to the external positions.

Input range	0 to 'Maximum stroke path' mm
Default value	0 mm
	A lift position of 0 mm corresponds to the "home position", i.e. the lift is located at the upper stop position. The maximum stroke path is defined in the properties of the tower (see "Maximum stroke path", page 109).

Rinse position

Lift position used for rinsing.



Input range	0 to 'Maximum stroke path' mm
Default value	0 mm

A lift position of **0 mm** corresponds to the "home position", i.e. the lift is located at the upper stop position. The maximum stroke path is defined in the properties of the tower (see "*Maximum stroke path*", page 109).


Swing increment

Amount by which the robotic arm can be swung relative to its current position. This parameter is used in the command **MOVE** for the settings **Destination = swing +** and **swing -**.

Input range	0.00 to 180.00 °
Default value	10.00 °

Properties – Robotic arm

Swing Head: **Properties ► Robotic arm**

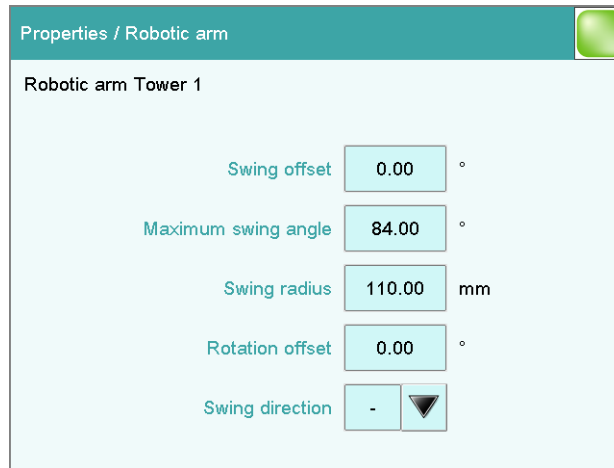

Properties of the robotic arm can be defined in the dialog **Properties / Robotic arm**.

Swing offset

The swing offset is a physical angle offset of a specific robotic arm model. The required values can be found in the leaflet for the robotic arm.

Input range	-270.00 to 270.00 °
Default value	0.00 °

Maximum swing angle

Useable swing range of the robotic arm. Each robotic arm model displays a different value on the basis of its construction. The range can also be reduced if necessary. You will find the required values in the leaflet for the robotic arm.

Input range	0.00 to 330.00 °
Default value	60.00 °

Swing radius

Maximum swing radius of the robotic arm. The swing radius is dependent on the length of the robotic arm and is, together with the axial distance (see "Axial distance," page 110), the most important variable for precise movement to a rack position. You will find the required values in the leaflet for the robotic arm.

Input range	30.00 to 300.00 mm
Default value	110.00 mm

Rotation offset

The rotation offset is the offset from the center of the tower to the center of the robotic arm. This value does not usually need to be changed. If a Swing Head is to be mounted on the tower with a lateral offset, then this value can be determined by the service technician when the rack is adjusted.

Input range	-270.00 to 270.00 °
Default value	0.00 °

Swing direction


The swing direction of the robotic arm depends on its type. For a 2-tower model, the robotic arm must be defined as right-swinging on Tower 1 and as left-swinging on Tower 2.

Selection	+ -
Default value	-

- +**
Left-swinging.
- Right-swinging.

Properties – External position

Swing Head: **Properties ► Edit**

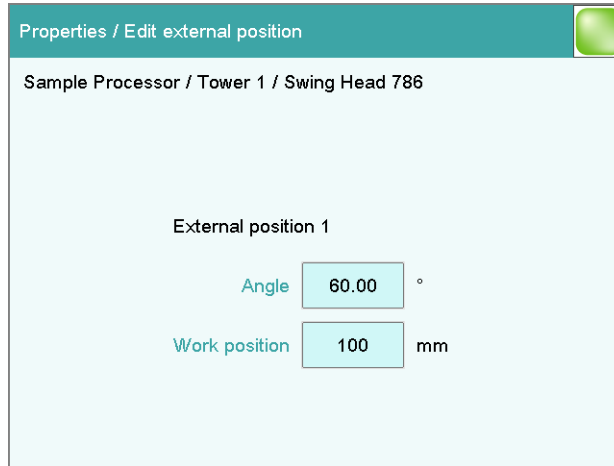

Properties which apply only to the selected external position can be defined in the dialog **Properties / Edit external position**.

Angle

Swing angle for the selected external position.

Input range	(Offset) to (Offset + max. swing range) °
Default value	60.00 °
	The offset is made up of a design-dependent angle (approx. 8...9°) together with the robotic arm offset from the robotic arm properties. The maximum swing range is also defined under the robotic arm properties (see "Properties – Robotic arm", page 112).

Work position

Work position for the selected external position.

Input range	0 to 'maximum stroke path' mm
	A lift position of 0 mm corresponds to the "home position", i.e. the lift is located at the upper stop position. The maximum stroke path is defined in the properties of the tower (see "Maximum stroke path", page 109).

11.6 Sample racks

Instrument list: **Sample Processor** ► **Edit** ► **Racks**

Edit device / Sample racks		
Rack	Positions	Code
6.2041.310	12	000001
6.2041.320	16	000010
6.2041.340	24	001000
6.2041.350	48	010000
6.2041.360	12	100000
6.2041.370	14	000011
6.2041.380	14	000101
6.2041.390	16	100001

Buttons: Load, Copy, Delete, Edit

In the dialog **Edit device / Sample racks**, you will find the list of configured racks. The attached rack is displayed in green. New racks can be loaded or created and existing ones can be edited or deleted in this dialog window.

The following data is displayed in the list:

- **Rack**
Name of the sample rack. Metrohm standard racks are designated by their order numbers.
- **Positions**
Number of positions on the rack.
- **Code**
The rack code corresponds to the arrangement of magnets on the base of the rack and is read in by the Sample Processor in order to recognize the rack.

[Load]

Load a new sample rack (see "Loading the sample rack", page 116).

[Copy]

Create a new sample rack by copying an existing rack (see "Creating a new sample rack", page 117).

[Delete]

Delete the selected sample rack from the list.

[Edit]


Edit the data of the selected sample rack (see chapter 11.6.1, page 117).

Loading the sample rack

When you receive the file for a new sample rack from your Metrohm representative, you can easily import this file into your existing system. Proceed as follows:

1 Copying the file to an external storage medium

- Copy the file to a subdirectory of "Files".

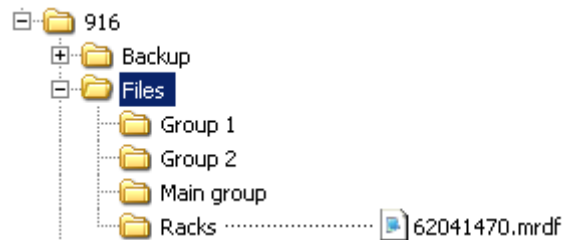

If this structure is not maintained, the new rack will not be found because the software directly accesses subdirectories of "Files".

- Plug in the external storage medium at the Ti-Touch.

2 Display the list of saved sample racks

- In the **Edit device / Sample racks** dialog, tap on the **[Load]** button.

The selection of file groups on the external storage medium is displayed. If only one group is available, then the list of the saved sample rack files will be displayed directly.

- Select the group with the desired sample rack.
- Tap on **[Show files]**.

The list with the saved sample rack files is opened.

3 Load the sample rack file

- Select the desired file.
- Tap on **[Load]**.

The new sample rack is now loaded and appears in the list of available racks.

Creating a new sample rack

Own sample racks can be simply and conveniently created by copying an existing sample rack. Proceed as follows:

1 Copy the existing rack

- In the dialog **Edit device / Sample racks**, select a sample rack which is to be used as a template.
- Tap on **[Copy]**.

The **Sample rack / Copy** dialog is displayed.

2 Enter the rack name and rack code

- In the field **New rack name**, enter a name for the new rack.
- In the field **New rack code**, **110000** is suggested. As a rule, this rack code is used for special racks. Apply this code or enter a new rack code and confirm with **[OK]**.

The new sample rack appears in the list of available racks.

11.6.1 Editing rack data

Sample rack list: **Rack ► Edit**

You can edit the data of the selected rack in the dialog **Sample rack / Edit rack data**.

Beaker radius samples

Actual radius of the sample vessels at the general sample positions of the rack.

This beaker radius may not be less than the minimum beaker radius defined in the tower properties (see "Min. beaker radius," page 109). If



the lift is to be moved to the work position, then these two values will be compared with one another.

Input range	1.0 to 100.0 mm
Selection	off
Default value	off

off

No check takes place.

Beaker sensor

Each time a sample position is to be moved to with the **MOVE** command, the beaker sensor checks whether a vessel is present. In the **MOVE** command, you define the action that takes place if the beaker sensor does not detect a vessel at the position being moved to.

Selection	Tower Robotic arm off
Default value	off

Robotic arm

A Swing Head with beaker sensor must be mounted. In addition, a suitable work position must be defined for the lift, so that the robotic arm touches the sample vessel. The work position is moved to for the purpose of beaker detection.

off

No check takes place.

1. Calibration pos.

Position of the first calibration solution for automatic calibration with a Sample Processor(see "*Definition of the calibration positions*", page 451) .

Input range	1 to "highest rack position" It is imperative that the remaining buffers/standards be placed on the rack positions directly following.
Selection	Special beaker 1...n off
Default value	off

Special beaker 1...n

If a special beaker is selected as the first calibration position, then the number of buffers used for calibration will determine the number of special beakers to be defined. Any rack position can be defined as a special beaker. It is, however, preferable to set them at high rack positions in order to be able to begin sample series at rack position 1. The special beakers are moved to in ascending order.

Rack offset

The rack offset is a production-related tolerance value between the upper and lower sections of the rack. This value is determined when the rack is adjusted and is displayed here. It can be changed if necessary.

Input range	-10.00 to 10.00 °
Default value	0.00 °

[Adjust rack]

Adjust the rack (*see chapter 11.6.2, page 123*).

[Lift pos. tower 1]/[Lift pos. tower 2]

Define rack specific lift positions (*see "Lift positions Tower 1/2", page 119*).

[Special beakers]

Define rack positions as special beakers (*see "Editing special beakers", page 120*).

Lift positions Tower 1/2

Sample rack list: **Rack ▶ Edit ▶ Lift pos. tower 1 / Lift pos. tower 2**

Edit rack data / Lift positions tower 1

Rack 6.2041.410, Code 001010, 142 Positions

Work position	<input type="text" value="120"/>	mm
Rinse position	<input type="text" value="0"/>	mm
Shift position	<input type="text" value="0"/>	mm
Special position	<input type="text" value="0"/>	mm

You can define rack-specific lift positions in the dialog **Edit rack data / Lift positions tower 1/2**. These then apply for all rack positions except those that have been defined as special beakers.

These positions can be moved to directly under manual control and with the LIFT command. Only lift positions within the maximum stroke path can be entered. This is defined in the device properties of the tower.

**NOTICE**

These lift positions can also be assigned directly in manual control after moving to the desired lift height (*see chapter 28.6.1, page 283*).

Work position

At this lift position the electrodes, stirrer and buret tips are optimally positioned for work.



Input range	0 to 'maximum stroke path' mm A lift position of 0 mm corresponds to the "home position", i.e. the lift is located at the upper stop position. The maximum stroke path is defined in the properties of the tower (<i>see "Maximum stroke path", page 109</i>).
-------------	---

Rinse position

Lift position used for rinsing.

Input range	0 to 'Maximum stroke path' mm
Default value	0 mm A lift position of 0 mm corresponds to the "home position", i.e. the lift is located at the upper stop position. The maximum stroke path is defined in the properties of the tower (<i>see "Maximum stroke path", page 109</i>).

Shift position

The lift is raised to this position with each rotational movement of the rack whenever it is located at a lower lift position.

Input range	0 to 'maximum stroke path' mm A lift position of 0 mm corresponds to the "home position", i.e. the lift is located at the upper stop position. The maximum stroke path is defined in the properties of the tower (<i>see "Maximum stroke path", page 109</i>).
-------------	---

Special position

For sample positions, you can define an additional lift position for special applications.

Input range	0 to 'maximum stroke path' mm A lift position of 0 mm corresponds to the "home position", i.e. the lift is located at the upper stop position. The maximum stroke path is defined in the properties of the tower (<i>see "Maximum stroke path", page 109</i>).
-------------	---

Editing special beakers

Sample rack list: **Rack ► Edit ► Special beakers**

Edit rack data / Special beakers	
Rack 6.2041.410, Code 001010, 142 Positions	
Special beaker	Rack position
Special beaker 1	142
Special beaker 2	0
Special beaker 3	0
Special beaker 4	0
Special beaker 5	0
Special beaker 6	0
Special beaker 7	0

A maximum of 16 rack positions can be defined as special beakers for each sample rack. The dialog **Edit rack data / Special beakers** shows a list of all special beakers with their assigned rack positions.

[Edit]

Edit the data of the selected special beaker, see the following.

Special beaker / Edit	
Special beaker 1	
Rack position	142
Work pos. tower 1	100 mm
Work pos. tower 2	0 mm
Beaker radius	off mm
Beaker sensor	Tower

You can edit the data of the selected special beaker in the dialog **Special beaker / Edit**.

Rack position

Number of rack position for selected special beaker. Any rack position can be defined as a special beaker. It is, however, preferable to set them at high rack positions in order to be able to begin sample series at rack position 1. Rack positions defined as special beakers can no longer be used as sample positions.

Input range **0 to 'maximum number of rack positions'**



Work pos. tower 1/2

Work position for the selected special beaker. One specific work position can be defined for tower 1 and tower 2.

Input range	0 to 'maximum stroke path' mm A lift position of 0 mm corresponds to the "home position", i.e. the lift is located at the upper stop position. The maximum stroke path is defined in the properties of the tower (see " <i>Maximum stroke path</i> ", page 109).
-------------	---

Beaker radius

Actual radius of the selected special beaker.

This beaker radius may not be less than the minimum beaker radius defined in the tower properties (see "*Min. beaker radius*," page 109). If the lift is to be moved to the work position, then these two values will be compared with one another.

Input range	1.0 to 100.0 mm
Selection	off
Default value	off

off

No check takes place.

Beaker sensor

Each time this special beaker is to be moved to with the **MOVE** command, the beaker sensor checks whether a vessel is present. In the **MOVE** command, you define the action that takes place if the beaker sensor does not detect a vessel at the position being moved to.

Selection	Tower Robotic arm off
Default value	off

Robotic arm

A Swing Head with beaker sensor must be mounted. In addition, a suitable work position must be defined for the lift, so that the robotic arm touches the sample vessel. The work position is moved to for the purpose of beaker detection.

off

No check takes place.

11.6.2 Rack adjustment

Instrument list: **Sample Processor** ► **Edit** ► **Racks** ► **Edit** ► **Adjust rack**


If necessary, each sample rack can be finely adjusted, i.e. the rack offset is determined in the direction of rotation. The adjustment of a rack is usually not necessary. However, if the exact positioning of a robotic arm is necessary for a high degree of accuracy (e.g. for very small sample beakers), then a fine adjustment can be carried out.

Proceed as follows:

1 Open the dialog for rack adjustment

- In the device properties of the Sample Processor, open the list of the configured sample racks.
- Select the attached rack and tap on **[Edit]**.
The **Sample rack / Edit rack data** dialog is displayed.
- Tap on the **[Adjust rack]** button.

The **Edit rack data / Adjust rack** dialog is displayed:

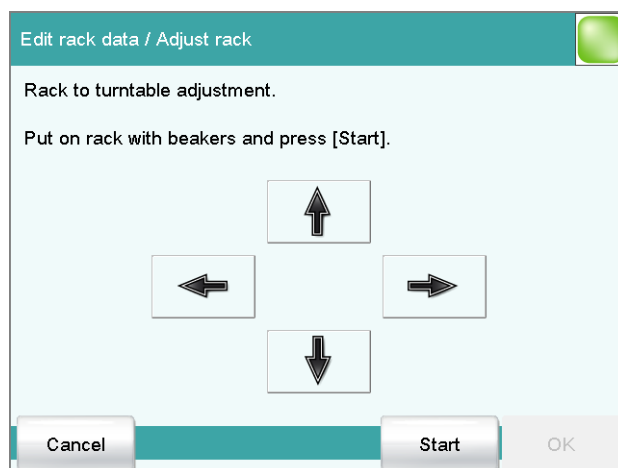

2 Carry out the adjustment


- Tap on the **[Start]** button (NOT on the **[▶]** fixed key).
The rack is being initialized. The rack then moves to position 1 and the lift is lowered to the work position.
- If necessary, the lift position can be corrected with the arrow keys **[↑]** and **[↓]**.
- Now use the arrow keys **[←]** and **[→]** to rotate the rack so that the robotic arm or the center of the titration head is positioned exactly above the center of rack position 1.
- Finish the adjustment with **[OK]**.



The lift is moved to the uppermost stop.

11.7 Printer

Instrument list: **Printer ▶ Edit**

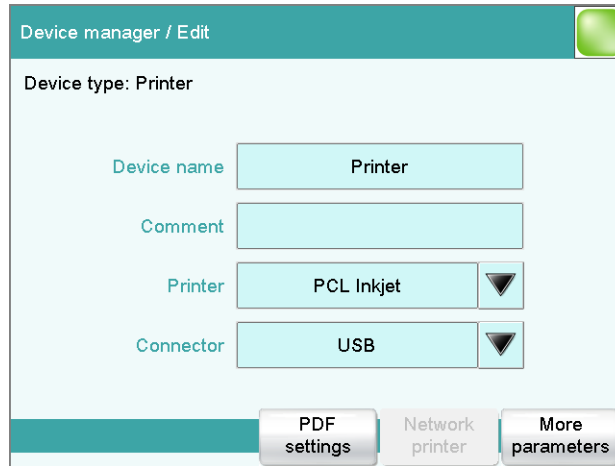

The list of devices always includes a printer, even if no corresponding device is connected. In this dialog, you also define when you would like to have a report generated as a PDF file.

Device name

Here you can enter a device name of your choice.

Entry **24 characters maximum**

Comment

Entry **24 characters maximum**

Printer

Selection of the printer type.

Selection **PCL Inkjet | PCL Laser | HP | Epson | Epson new | Canon | CUSTOM NEOS | off**

PCL Inkjet
for HP DeskJet printers.

PCL Laser
for HP LaserJet printers.

CUSTOM NEOS
Due to the narrow paper size, only reports marked with # in the "List of all printable reports" (*see chapter 27.3, page 261*) can be printed.

off
The reports will not be printed out on paper.

Connector

Selection of the connection type for the printer.

Selection	USB Ethernet
Default value	USB

USB

Local printer at a USB interface.

Ethernet

Network printer.



NOTICE

The **Ti-Touch** prints the reports with a fixed resolution of 300 dpi. If you are using a printer with a resolution of 360 dpi (or a multiple thereof, e.g. an Epson), then the text will be printed out somewhat smaller than with printers with a resolution of 300 dpi (or a multiple thereof, e.g. a Canon or HP).

11.7.1 PDF settings

Printer: **Edit ► PDF settings**

The settings for saving a report as a PDF file are defined in this dialog.

Memory

Memory location where the PDF file will be saved. The report will be saved in the directory *PDF_Report*. This directory is created when a PDF file is generated. The file name is appended with the date and time when the determination was started and the date and time when the PDF was generated (e.g.: **Identification 1-20130320-141044-20130320-141220**). The first date and time set indicates the start of the determination, the second date and time set indicates the generation of the PDF file. This prevents previously created PDF files from being overwritten each time a new file is created, e.g. when a result is recalculated and the report is generated again.

All three memory locations are listed as possible selections, even if they cannot be accessed at the moment.

Selection	off External memory 1 External memory 2 Shared memory
Default value	off

off

The report will not be saved as a PDF file.

Shared memory

The report will be saved in a shared directory on the network. The shared directory is selected in the **Edit device / Shared memory** dialog (see chapter 11.3.3, page 100).

Copy or extract content allowed

on | off (Default value: **on**)

If this option is activated, then content can be copied or deleted from the PDF file.

Printing allowed

on | off (Default value: **on**)

If this option is activated, then the PDF file can be printed.

Change the document allowed

on | off (Default value: **off**)

If this option is activated, then the PDF file can be edited.

Add or change comments allowed

on | off (Default value: **on**)

If this option is activated, then comments can be added to the PDF file.

11.7.2 Network printer

Printer: **Edit ► Network printer**

If you have connected your Ti-Touch to your network, then you can specify a network printer for your reports in this dialog.



NOTICE

If you have shared a memory location on a PC within your network and if you have configured the Ti-Touch accordingly (see chapter 11.3.3, page 100), then the settings for the parameters **Domain**, **User** and **Password** will be applied and can then no longer be modified in this dialog.

**CAUTION**

If the computer on which the network printer is configured does not have the same subnet as the Ti-Touch, then a WINS server must be present. The computer must be entered in this server. In order for the communication to work, the SMB protocol version 1 (SMBv1) must be activated. Since this function is not standard anymore in Windows 10, it has to be activated manually. From firmware version 5.916.0044, the SMB protocol version 2 and 3 is also supported.

Print server

Host name of the print server or the computer on which the network printer is configured. If you are working with Windows, then you will find the host name as follows: In the Windows **command prompt** window, enter the command **ipconfig -all**. The host name is listed together with other parameters of the print server or the computer.

Do not under any circumstances enter an IP address in this input field.

Entry	max. 32 characters
Default value	empty

Share name

Share name of the network printer on the above-specified computer. Please note that the Share name of a network printer often does not match the name of the printer. The share name can be found on the Release tab in the Properties dialog of the released network printer.

Entry	max. 32 characters
Default value	empty

Domain

Network domain in which the above-specified computer is located. If you are working with Windows, then you will find the domain name as follows: In the Windows **command window**, enter the command **ipconfig -all**. The domain is listed together with other parameters of the computer. Leave the field empty if the computer is not located within a domain.

Entry	max. 32 characters
Default value	empty

User

User name of the user authorized to access the shared network printer.

Entry	max. 32 characters
Default value	empty

Password

Password of the user configured on the computer.

Entry	max. 32 characters
Default value	empty

11.7.3 More options

Paper size

Selection of the paper size.



NOTICE

The paper size selected will be ignored when the report is printed on a **CUSTOM NEOS** printer.

Selection	A4 (210 mm x 297 mm) Letter (216 mm x 279 mm)
Default value	A4 (210 mm x 297 mm)

Color

on | off (Default value: **off**)

If this option is activated, then the report will be printed out in color.

Spooler

on | off (Default value: **on**)

If this option is activated, then you can continue to work while the printer is printing. If this option is deactivated, then the Ti-Touch will be blocked while the printout is being produced.

11.8 Balance

List of devices: **Balance ► Edit**

You define the type of balance and its connector in the dialog **Device manager / Edit**.

Device name

Here you can enter a device name of your choice.

Entry **24 characters maximum**

Comment

Entry **24 characters maximum**

Balance type

If you have connected a balance, then you have to specify the balance type here.

Selection	AND Mettler Mettler AT Mettler AX Ohaus Precisa Sartorius Shimadzu
Default value	Sartorius

RS-232

RS-232 interface to which the balance is connected. The interface parameters are adjusted in the **Edit device / Port parameters** dialog (see chapter 11.9, page 130).

Selection	COM 1 COM 2
Default value	COM 1

COM 2

The interface is disabled.

The following table indicates the balance type that needs to be selected for the balance model:



Balance	Balance type
AND	AND
Mettler AB, AE, AG, AM, AJ, PE, PM, PJ, PR, XP, XS	Mettler
Mettler AT	Mettler AT
Mettler AX, MX, UMX, PG, AB-S, PB-S	Mettler AX
Ohaus Voyager, Explorer, Analytical Plus	Ohaus
Precisa	Precisa
Sartorius	Sartorius
Shimadzu BX, BW	Shimadzu

11.9 USB/RS-232 adapter

Instrument list: **USB/RS-232 adapter** ► **Edit** ► **COM interface** ► **Edit**

As a rule, balances are equipped with a serial RS-232 interface. To connect a balance, you will require the 6.2148.050 cable. You can adjust the interface parameters in the dialog **Edit device / Port parameters**. These settings must match the settings on the connected device.



NOTICE

If you make changes to parameter settings in this dialog, then you must switch the Ti-Touch off and back on in order for the changes to take effect.

Edit device / Port parameters

Baud rate ▼

Data bits ▼

Parity ▼

Stop bits ▼

Handshake ▼

Baud rate

Transfer rate in characters per second.

Selection	1200 2400 4800 9600 19200 38400 57600 115200
Default value	9600

Data bits

Number of data bits.

Selection	7 8
Default value	8

Parity

Type of parity testing.

Selection	even odd none
Default value	none

Stop bits

Number of stop bits.

Selection	1 2
Default value	1

Handshake

Type of data transfer protocol.

**NOTICE**

In case of communication problems, try the software handshake (**Software (XON/XOFF)**).

Selection	none Software (XON/XOFF) Hardware (DTR/CTS)
Default value	Hardware (DTR/CTS)


Software (XON/XOFF)

Use the software handshake when you send a PC/LIMS report via an RS-232 interface.

11.10 PC keyboard

Instrument list: **PC keyboard ▶ Edit**

A commercially available USB keyboard can be connected to make text and numerical input easier. It will be recognized automatically and entered in the list of devices with default settings.

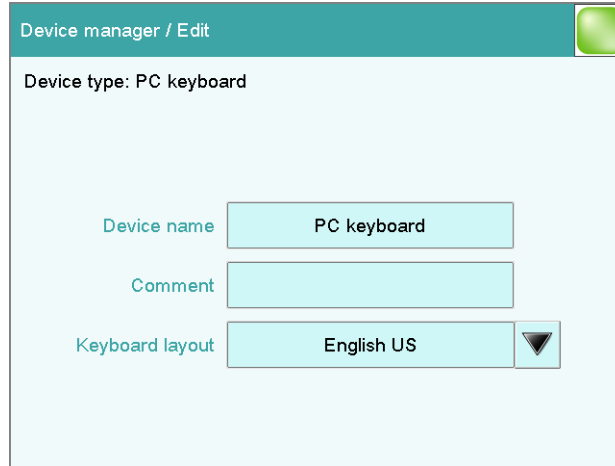

Device name

Here you can enter a device name of your choice.

Entry **24 characters maximum**

Comment

Entry **24 characters maximum**

Keyboard layout








Define the country-specific keyboard layout here.

Selection	English US German DE French FR Spanish ES German CH
Default value	English US

To enter texts and numbers with the PC keyboard, the appropriate text and number input dialog must be opened on the Ti-Touch. Only the following keys on the PC keyboard have a function:

Table 6 Key assignment

Function in the editing dialog	PC keyboard key
[Cancel]	[Esc]

Function in the editing dialog	PC keyboard key
Enter respective character	Numbers, letters and special characters + Shift
[OK]	[Enter] / [↵]
[□]	[←] (backspace)
Line break (for multi-line text input)	[Enter] / [↵]
The cursor is moved up or down by one character.	[↑] [↓]
The cursor is moved to the left or to the right by one character.	[←] [→]
	[F1]
	[F2]
	[F3]
	[F4]
	[F5]
	[F6]
	[F7]



NOTICE

The lettering of the USB keyboard may differ from above lettering, depending on the country-specific keyboard used.

11.11 Barcode reader

List of devices: **Barcode reader** ► **Edit**



Device manager / Edit

Device type: Barcode reader

Device name

Comment

Barcode input target ▼

Keyboard layout ▼

A barcode reader can be connected to read in sample data or other texts. A connected barcode reader will be recognized automatically and entered in the list of devices with default settings.

You will hear an acoustic signal as confirmation that a character string has been transmitted by the barcode reader and accepted.

NOTICE

The data will be read only if the system is in normal status, i.e. when no determination is running.

Device name

Here you can enter a device name of your choice.

Entry	24 characters maximum
-------	------------------------------

Comment

Entry	24 characters maximum
-------	------------------------------

Barcode input target

Selection of the input field for the character string read in by the barcode reader.

Selection	Active input field Method Identification 1 Identification 2 Sample size
-----------	--

Active input field

The character string is entered in the input field of the opened text- or number-input dialog.

Method

The character string is entered in the input field **Method**.

Identification 1

The character string is entered in the input field **Identification 1**.

Identification 2

The character string is entered in the input field **Identification 2**.

Sample size

The character string is entered in the input field **Sample size**. Character strings containing characters other than numbers and decimal separators will be ignored.

Keyboard layout

Specify the country-specific keyboard layout for the emulation of the PC keyboard. This setting must match the setting on the barcode reader (see documentation for the barcode reader).

Selection	English US German DE French FR Spanish ES German CH
Default value	English US

12 File manager

Main dialog: **System ► File manager**

The saved methods, determinations, sample tables, etc. are managed in the file manager. You can also create a backup of your system (all data and settings). Also, an existing backup can be reloaded.

The file memory is organized as follows:

- **Internal memory**

The following files can be stored in the internal memory:

- Methods

- **External memory**

You can use a USB flash drive as an auxiliary storage medium, for instance. The following files can be stored on an external memory:

- Backup
- Methods
- Determinations
- Sample tables
- Result tables
- Reports as *.pdf
- PC/LIMS report

12.1 Managing files

File manager: **Internal memory / External memory 1 / External memory 2**

The saved files can be organized into groups. These groups are comparable to file directories on your PC, although unlike your PC, only one level is possible.




NOTICE

The file names must be unique for each memory location, i.e. you cannot save two files with the same name, even if they are in different groups.

**NOTICE**


If you use an external storage medium with the FAT or FAT32 file system, then you can save a maximum of 999 files per group. If you find it necessary to store more than 999 files in a single group, then you must reformat the storage medium with the file system **ExFAT** (see chapter 32.11.2, page 473).

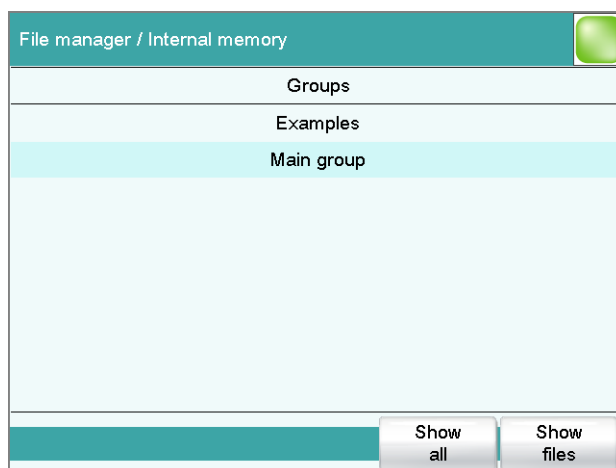**[Show all]**

Display the list of all the files in the selected memory location.

[Show files]

Display the list of all the files in the selected group.

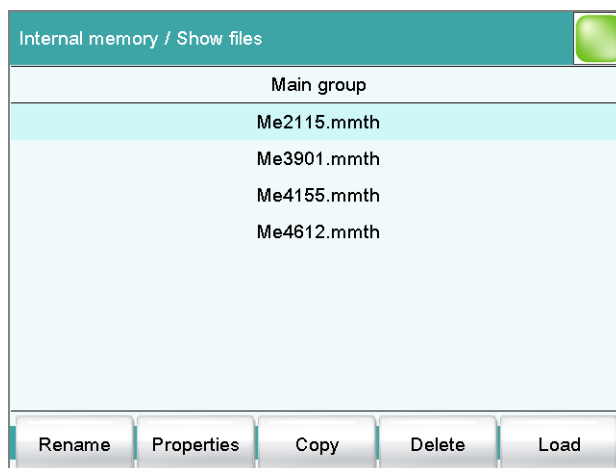**[Rename]**

Rename the selected file (see chapter 12.1.2, page 138).

[Properties]

Display the file properties (see chapter 12.1.3, page 139).

12.1.3 File properties

File manager: **Internal memory / External memory 1 / External memory 2 ► Show all / Show files ► Properties**

Detailed information concerning the file is displayed in the **Show files / Properties** dialog. They can be edited to a certain extent.

- **File name**
Name of the file.
- **File group**
Entry or selection of the group in which the file will be saved. If you enter a new name, then the file will be moved into the new group.
- **Write protection**
If this option is activated, then the file cannot be saved, deleted, moved or renamed. This is only an internal write protection feature and is independent of the write protection of the operating system on your computer.
- **Created by**
User who saved the file initially.
- **Created on**
Date and time at which the file was saved for the first time.
- **Last saved by**
User who saved the file most recently.
- **Last saved on**
Date and time at which the file has been saved most recently. Additionally, the version of the file is displayed. The version number will be increased by one each time the file is saved.
- **Size**
File size.
- **Program version**
Program version of the software with which the file was last saved.

Directory structure

The directory structure appears as follows on the external storage medium:

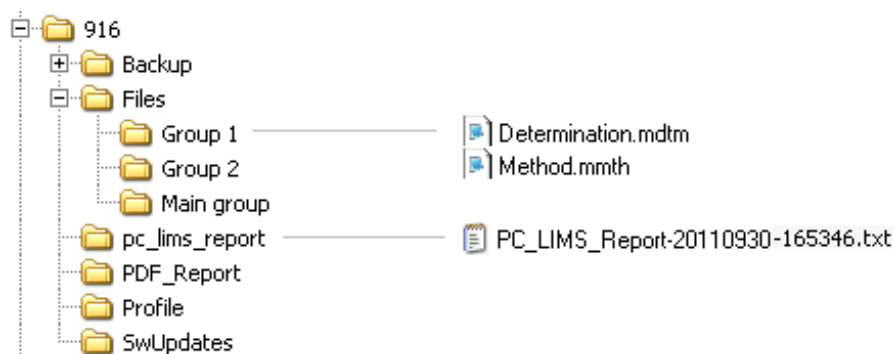


Figure 24 Directory structure on the external storage medium

Backup	All of the files of the backup are stored in this directory. The directory will be created the first time a backup is created.
Files	This directory contains all of the groups as well as the files contained in them (methods, determinations, etc.).
pc_lims_report	PC/LIMS reports are stored in this directory as TXT files. The directory will be created the first time a PC/LIMS report is printed out.
PDF_Report	PDF reports are stored in this directory. The directory will be created the first time a report is saved as a PDF file.
Profiles	The identification profile is stored in this directory. The directory will be created the first time an identification profile is generated.
SwUpdates	Files for software updates must be saved in this directory.

12.3 Creating backups / Restoring data

File manager: **External memory 1 / External memory 2**

You can use the **[Backup]** function to easily create a backup containing all the data and settings of your system. You should make a backup at regular intervals in order to avoid data loss.



NOTICE

Only **one** backup can be created on a storage medium.

If a backup has already been stored on the medium, then it will be overwritten when this function is carried out again.

12.3.1 Restoring data

You can use the function **[Restore]** to restore either a complete backup or only certain data.



CAUTION

All of the methods in the internal memory will be deleted.

Backups are not backward compatible, i.e. backups of the latest version cannot be restored using earlier versions.

Proceed as follows to restore backed-up data:

1 Select the storage medium

- Connect the storage medium on which the backup is located.
- Select the storage medium under **System ► File manager**.

2 Select data

- Tap on **[Backup]**.
The selection of the data that can be restored separately is displayed (see the following).
- Deselect the data sets which are not to be restored.
- Tap on **[Load data]**.

The selected data blocks are restored.

3 Complete the restoration

- Switch the Ti-Touch off and on again.

Restorable data blocks

The following data blocks can be loaded individually:

- **Methods**
All of the methods stored in the internal memory.
- **Sample table**
Current sample table.
- **Result table**
Current result table.
- **Current determination data**
All the data for the current determination (including the method with which the determination was carried out).
- **Control**
Settings in the control dialog.
- **User list**
Settings for each user under **System settings / User administration**.
- **System settings / User admin.**
All of the system settings, including device-specific dialog configuration and dialog options for the command list and fixed keys, device-specific settings for the user administration (login options and password options).
- **Titration**
All titration with all their data.
- **Sensors**
All sensors with all their data.
- **Device data**
All of the devices configured in the device manager, with all of the data.
- **GLP data**
All data from the GLP manager.
Deactivate if the backup is to be loaded onto a different system.
- **Common variables**
All common variables.
- **Templates**
All templates for sample data, result calculations, calibration buffers, etc.
- **Routine dialog settings**
Current routine dialog settings (see **Dialog options / Routine dialog**).
- **Sample racks**
The sample racks present in the properties of the USB sample processor.
- **Own subsequences**
All subsequences created and stored by the user.

13 GLP manager

Main dialog: **System ► GLP manager**

In the GLP manager you can document data for various GLP tests. The results of the automatic system tests carried out after switching on are also documented.


The following tests can be documented:

- GLP test "Measurement" (see chapter 13.3, page 146)
- GLP test "Titration" (see chapter 13.3, page 146)
- System validation (see chapter 13.4, page 150)

In addition, you can also:

- Create a list with your test tools (see chapter 13.2, page 145).
- Define a service interval for having routine maintenance carried out by a Metrohm service technician (see chapter 13.5, page 154).
- Define an interval for the regular performance of backups (see chapter 13.5, page 154).

Further information about quality management and validation is also given in the documentation series **Quality Management with Metrohm** and **Application Bulletin AB 252** (Validation of Metrohm titrators (potentiometric) in accordance with GLP/ISO 9001).

System / GLP manager 		
Test	Last test	Next test
Automatic system test	2010-08-30	Power on
System validation	2010-08-02	2011-01-04
GLP test "Measurement"	2010-07-01	2011-01-03
GLP test "Titration"	2010-07-02	2011-01-04

Test tools
GLP tests HW/SW
System validation
Monitoring
View test data

The table shows the last time that each test was carried out and when the next test is to be carried out. A test is entered in the list the first time that it is documented.

[Test tools]

Configuring the test tools for GLP tests (see chapter 13.2, page 145).

[GLP tests HW/SW]

Document the GLP tests "Measurement" and "Titration" (*see chapter 13.3, page 146*).

[System validation]

Document the system validation (*see chapter 13.4, page 150*).

[Monitoring]

Defining the time intervals for system maintenance and backups (*see chapter 13.5, page 154*).

[View test data]

Display the data of the selected test.

13.1 Automatic system test

The system test is carried out automatically when the Ti-Touch is switched on. The result of each individual test is shown in green if no error has occurred. If a result is shown in red, then an error occurred when the respective test was carried out. Switch the Ti-Touch off and back on again. If the error still occurs please notify Metrohm Service.

Print report at system start

on | off (Default value: **off**)

If this parameter is activated, then the result of the system test will be printed out automatically when the Ti-Touch is switched on.

13.2 Test tools

Main dialog: **System ► GLP manager ► Test tools**

You can create a list of test tools to be used in the tests in the GLP manager.

The following test tools have already been defined:

- **767 Calibrated Reference:** Device for checking measuring inputs and electrode cable.
- **822 Titr.Curve Simulator:** Curve simulator for checking the hardware and software.
- **773 pH/mV Simulator:** Device for checking measuring inputs and electrode cable.
- **868 UR Generator:** Device for checking measuring inputs and electrode cable.

[New]

Add a new test tool to the list.

[Delete]

Delete the selected test tool from the list.

[Edit]

Change the designation of the selected test tool.

13.3 GLP tests for measurement and titration

Main dialog: **System ► GLP manager ► GLP tests HW/SW ► GLP test "Measurement" / GLP test "Titration"**

You can document the GLP tests for measurements and titrations in the dialog **GLP manager / GLP tests Hardware/Software** and its subdialogs. The following procedure describes the GLP test "Measurement," but also applies for the GLP test "Titration".

Proceed as follows:

1 Open the properties dialog

- In the **System / GLP manager** dialog, tap on the **[GLP tests HW/SW]** button and then on the **[GLP test "Measurement"]** button.

2 Edit data

- Define the test method, the test results, etc. (see "Dialogs "GLP tests HW/SW / GLP test "Measurement"" and "GLP tests HW/SW / GLP test "Titration""", page 147).

3 Define the hardware used

- Tap on **[Hardware]**.

- Select the hardware used for the test (see "Dialogs "GLP test "Measurement" / Hardware" and "GLP test "Titration" / Hardware"", page 149).
- Tap on the fixed key [↩].

The dialog **GLP tests HW/SW / GLP test "Measurement"** appears again.

4 Define the test interval

- Tap on **[GLP test interval]**.
- Enter the time interval or the date for the next GLP test (see "Dialogs "GLP test "Measurement" / Test interval" and "GLP test "Titration" / Test interval"", page 148).
- Tap on the fixed key [↩].

13.3.1 Parameter description

Dialogs "GLP tests HW/SW / GLP test "Measurement"" and "GLP tests HW/SW / GLP test "Titration""

Method

Method the GLP test has been carried out with.

Entry	32 characters maximum
Selection	Selection of methods stored in the internal memory

User

User who carried out the GLP test.

Entry	24 characters maximum
Selection	Selection of configured users

Test date

Date on which the GLP was carried out.

Format: YYYY:MM:DD

Test result

Result of the test.

Selection	Test OK Test not OK
Default value	Test OK

Comment

Entry	24 characters maximum
-------	------------------------------

**[GLP test interval]**

Defining the time interval for the GLP test (see "Dialogs "GLP test "Measurement" / Test interval" and "GLP test "Titration" / Test interval", page 148).

[Hardware]

Documenting the hardware with which the GLP test was carried out (see "Dialogs "GLP test "Measurement" / Hardware" and "GLP test "Titration" / Hardware", page 149).

Dialogs "GLP test "Measurement" / Test interval" and "GLP test "Titration" / Test interval"

Test date

Date on which the GLP was carried out.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which a GLP test has to be carried out again will be monitored.

GLP test interval

If you define a time interval for the GLP test, then the date in **Next GLP test** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next GLP test

If you define a date for the next GLP test, then the **GLP test interval** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message

For all three options it is documented in the determination data (see dialog **More determination data / Messages**), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

Dialogs "GLP test "Measurement" / Hardware" and "GLP test "Titration" / Hardware"**Test tool**

Selection of the test tool. Test tools are defined at **GLP manager ► Test tools**.

Selection	Selection of configured test tools
-----------	---

Control device

Selection of the control device with which the GLP test has been carried out.

Selection	Selection of configured control devices
-----------	--

Measuring input

Selection of the measuring input used for the GLP test. The selection is not dependent on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

13.4 System validation

Main dialog: **System ► GLP manager ► System validation**

In the dialog **GLP manager / System validation** and its subdialogs, you can document the system validation results and define the time interval after which validation must be carried out again.

Proceed as follows:

1 Open the properties dialog

- In the **System / GLP manager** dialog, tap on the **[System validation]** button.

2 Edit data

- Define the method, the result, etc. (see *"GLP manager / System validation" dialog*, page 151).

3 Define the validation interval

- Tap on **[Validation interval]**.
- Enter the time interval or the date for the next system validation (see *"Dialog "System validation / Validation interval"*, page 152).
- Tap on the fixed key [**↩**].

The dialog **GLP manager / System validation** appears again.

4 Enter statistical data

- Tap on **[Test statistics]**.
- Enter the statistical data for the last system validation (see *"System validation / Test statistics" dialog*, page 153).
- Tap on the fixed key [**↩**].

The dialog **GLP manager / System validation** appears again.

5 Insert a note

- Tap on **[Note (SOP)]**.
- Enter a brief description, e.g. a summary of the SOP (standard operating procedure) according to which the system validation was carried out (see "*System validation / Note (SOP) dialog*", page 152).
- Tap on the fixed key [**↩**].

13.4.1 Parameter description

"GLP manager / System validation" dialog

Method

Method with which the system validation has been carried out.

Entry	32 characters maximum
Selection	Selection of methods stored in the internal memory

User

User who carried out the system validation.

Entry	24 characters maximum
Selection	Selection of configured users

Test date

Date on which the system validation was carried out.

Format: YYYY:MM:DD

Test result

Result of the test.

Selection	Test OK Test not OK
Default value	Test OK

Comment

Entry	24 characters maximum
-------	------------------------------

[Note (SOP)]

Enter a brief description, e.g. a summary of the SOP (standard operating procedure) according to which the system validation has been carried out.

[Validation interval]

Defining the time interval for the system validation (see "*Dialog "System validation / Validation interval"*", page 152).

[Test statistics]

Documenting the statistical data of the system validation (see *"System validation / Test statistics" dialog*, page 153).

"System validation / Note (SOP)" dialog

In this dialog, you can enter a brief text, e.g. a summary of the SOP (standard operating procedure) according to which the system validation was carried out.

Dialog "System validation / Validation interval"

Last validation

Date on which the last system validation was carried out.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which a system validation has to be carried out again will be monitored.

Validation interval

If you define a time interval for the system validation, then the date in **Next validation** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next validation

If you define a time interval for the next system validation, then the **Validation interval** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message

For all three options it is documented in the determination data (see dialog **More determination data / Messages**), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

"System validation / Test statistics" dialog

In the dialog **System validation / Test statistics**, you can document the statistical data for the last system validation.

Number (n)

Number of determinations carried out.

Input range	1 to 999999999
Default value	empty

Mean value

Mean value out of single results.

Input range	-999999999.00000 to 999999999.00000
Default value	empty

s abs

Absolute standard deviation of the results.

Input range	0.00000 to 999999999.00000
Default value	empty

s rel

Relative standard deviation of the results.

Input range	0.00000 to 100.00000 %
Default value	empty %

d rel

Systematic deviation of the results.

Input range	0.00000 to 100.00000 %
Default value	empty %

a sys

Systematic error.

Input range	0.00000 to 999999999.00000
Default value	empty

13.5 System monitoring

13.5.1 Service interval

Main dialog: **System ► GLP manager ► Monitoring ► Service interval**

In the dialog **GLP manager / Service interval**, you can define the time interval for the maintenance of the system by the Metrohm Service department. The service interval is checked each time the system is started.

Last service

Date on which the last servicing was carried out.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which system maintenance must be carried out again will be monitored.

Service interval

If you define a time interval for the system maintenance, then the date in **Next service** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next service

If you specify a date for the next system maintenance, then the **Service interval** will be tracked automatically.

Format: YYYY:MM:DD

13.5.2 Backup interval

Main dialog: **System ► GLP manager ► Monitoring ► Backup interval**

In the dialog **GLP manager / Backup interval**, you can define the time interval for backups. The backup interval is checked each time the system is started.

Last backup

Date on which the last backup was created.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which a backup has to be created again will be monitored.

Backup interval

If you define a time interval for the creation of backups, then the date in **Next backup** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next backup

If you specify a date for the next backup, then the **Backup interval** will be tracked automatically.


Format: YYYY:MM:DD

14 Common variables

Main dialog: **System ► Common variables**

You can save 25 **method-independent variables**, or common variables. These variables can be used in future calculations (as variables **CV01... CV25**). Common variables are useful, e.g. for the following applications:

- Determination of a blank value which will be taken into account during the content determination of the sample.
- Determination of the content of a standard solution, which will be taken into account during the content determination of the sample.

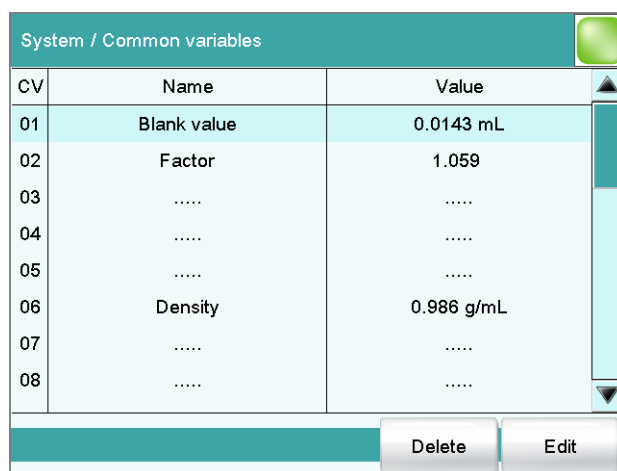

CV	Name	Value
01	Blank value	0.0143 mL
02	Factor	1.059
03
04
05
06	Density	0.986 g/mL
07
08

The designation and the value (including the unit) are displayed for every common variable in the list.

[Delete]

Delete the selected common variable from the list.

[Edit]

Edit the data of the selected common variable, see following chapter.

14.1 Editing common variables

List of common variables: **Common Variable ► Edit**

The common variables can be modified as follows:

- Edit manually, see the following.
- Automatic assignment from the determination run. A calculation result must be configured accordingly for this purpose (see chapter 14.4, page 160).

Name

Designation of the common variable.

Entry	24 characters maximum
Default value	empty

Value

Value of the common variable.

Input range	-999999999 to 999999999
Default value	empty

Unit of the common variable.

Entry	10 characters maximum
Default value	empty
Selection	% mol/L mmol/L g/L mg/L mg/mL mg/100 g ppm g mg µg mL µL mg/ piece °C µg/min mL/min µL/min
Default value	%



Next assignment

This parameter is displayed only if a validity has been defined for the common variable.

Date on which the validity of the common variable expires.

Format: YYYY:MM:DD

[Validity]

Defining the time interval for the validity of the common variable (*see chapter 14.3, page 159*).

[Properties]

Display the properties of the common variable, see following chapter.

14.2 Properties of common variables

Common variable: **Edit ► Properties**

Detailed information concerning the common variable is displayed in the **Edit common variables / Properties** dialog:

- **Status**
Status of the common variable. If the time interval for the validity has expired, then **invalid** will be displayed.
- **Method**
Method with which the value has been assigned to the common variable. If the value was entered manually, then **manual** will be displayed.
- **Method status** (only for automatic assignment from the determination run)
- **Determination status** (only for automatic assignment from the determination run)
- **Last assignment**
Date and time of the last assignment.
- **User**
User who assigned the value to the common variable.

14.3 Monitoring validity

Common variable: **Edit ► Validity**

In the dialog **Edit common variables / Validity**, you can define the time interval after which a new value must be assigned to the common variable.

Last assignment

Date on which the common variable was last assigned a value.

Format: YYYY:MM:DD

Monitoring

on | off (Default value: **off**)

If this parameter is activated, then the time interval after which the common variable must be assigned a new value will be monitored.

Validity

If you define a time interval for the validity of the common variable, then the date in **Next assignment** will be tracked automatically.

Input range	1 to 999 days
Default value	999 days

Next assignment

If you specify a date for the next assignment, then the time interval for the **Validity** will be tracked automatically.

Format: YYYY:MM:DD

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message

For all three options it is documented in the determination data (see dialog **More determination data / Messages**), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

14.4 Assigning a result automatically to a common variable

**NOTICE**

This instruction is based on the assumption that the method contains a calculation command with a calculation.

Proceed as follows to assign a result to a common variable:

1 Opening the editing dialog of the result

- In the command list, select the command **CALC**.
- Tap on the **[Edit command]** button.
- Select the result whose value is to be assigned to a common variable and tap on **[Edit]**.

The editing dialog of the result is displayed.



2 Define result options

- Tap on the **[Result options]** button.

Edit calculation / Result options
 R1 Blank value
 Variable for mean value SMN1
 Save as titer off
 Save as common variable
 Variable CV01
 More options

- Activate the parameter **Save as common variable**.
- Define **Variable = CV01...CV25**.

3 Saving the settings

Tap on the fixed keys [] or [].

In the future, the calculated result will be assigned to the selected common variable (result name, value and unit).

15 Templates

Main dialog: **System ► Templates**


You have the option of defining **system-specific templates**. You can use these templates when editing the respective data.

You can create the following templates:

- **Sample data** (see chapter 15.1, page 162)
Create sample identifications and sample assignments.
- **Custom result templates** (see chapter 15.2, page 166)
Define formulas for result calculations.
- **Input lines** (see chapter 15.3, page 170)
Define input signals on the remote interface.
- **Output lines** (see chapter 15.4, page 172)
Defining output signals on the remote interface.
- **Custom calib. buffers** (see chapter 15.5, page 175)
Define a buffer series for the calibration of pH electrodes.
- **Report header** (see chapter 15.6, page 177)
Create a system-specific report header that contains information concerning the laboratory, for example.
- **Custom electrode type** (see chapter 15.7, page 178)
Define customized limit values for the electrode test of pH electrodes.

15.1 Sample data

Main dialog: **System ► Templates ► Sample data**

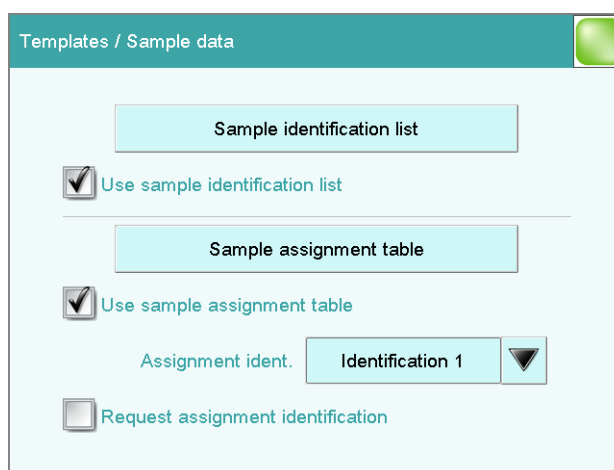

You can create the following templates in the dialog **Templates / Sample data**:

- Sample identifications (*see chapter 15.1.1, page 163*)
List with sample identifications. If you must enter the sample identification before starting a determination, then you can select the entries contained in this list.
- Sample assignments (*see chapter 15.1.2, page 164*)
You can assign a particular method to a sample identification.

Use sample identification list

on | off (Default value: **off**)

Activate this parameter so that the sample identifications defined in the list will be displayed as selections.

Use sample assignment table

on | off (Default value: **off**)

Activating/deactivating the use of the sample assignment table.



NOTICE

If this parameter is activated, the following will no longer be possible:

- Carrying out determinations without their sample identification being defined in the sample assignment table.
- Defining a method in the sample table. Previously defined methods will be ignored.

Assignment ident.

Identification which is used as assignment identification to load the correct method.

Selection	Identification 1 Identification 2
Default value	Identification 1


Request assignment identification

on | off (Default value: **off**)

If this parameter is activated, then the sample identification will be requested automatically at the start of a determination.

15.1.1 Sample identification list

Main dialog: **System ► Templates ► Sample data ► Sample identification list**

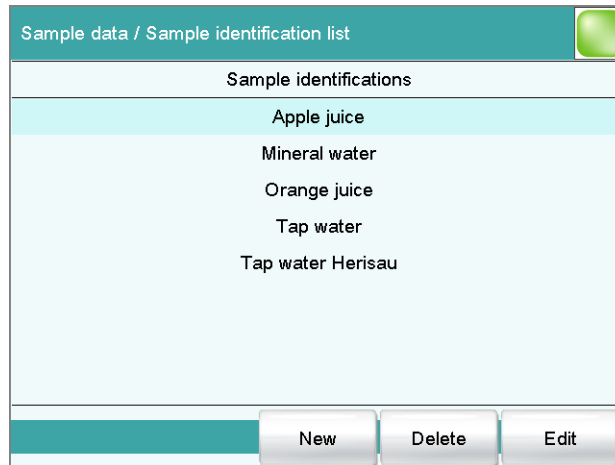

You can create a system-specific list with sample identifications in the dialog **Sample data / Sample identification list**. This list makes it easier for you to enter frequently used sample identifications, i.e. this list is available to you for entering them (main dialog, command **REQUEST**, etc.). It may be a good idea to define the unchanging part of the identification as a template and to add the variable part during sample data input.

[New]

Add a new sample identification to the list.

[Delete]

Delete the selected sample identification from the list.

[Edit]

Change the designation of the selected sample identification.

15.1.2 Sample assignment table

Main dialog: **System ► Templates ► Sample data ► Sample assignment table**

Sample ident.	Method
Apple juice	Me2115
Mineral water	Me3901
Orange juice	Me4612
Tap water *	Me4155

In the sample assignment table, a particular method is assigned to a sample identification. In this way, you ensure that your samples will be processed with the correct method; mix-ups are not longer possible. When you start a determination, you need only enter the sample identification; the method is loaded automatically.

[New]

Add a new sample assignment to the list, see following chapter.

[Delete]

Delete the selected sample assignment.

[Edit]

Edit the selected sample assignment, see following chapter.

Editing the sample assignment

Identification

Identification of the sample.



NOTICE

You can place an * as a wildcard at the beginning or end of the character string. Doing so allows you to prefix or suffix a sequential number, for example, to the identification, which will be ignored when methods are being assigned.

Entry	24 characters maximum
Selection	Selection of defined identifications in the sample identification list

Memory

Memory location the method is loaded from. All memory locations are selectable, even if they are currently not accessible.

Selection	Internal memory External memory 1 External memory 2 Shared memory
Default value	Internal memory

Shared memory

Shared directory in the network.


Method

Method that is loaded when a sample with the above-specified identification is processed. You can also enter a method that does not yet exist in the selected memory. When a determination is started there is a check whether the method is available.

Entry	32 characters maximum
Selection	Selection of methods stored in the selected memory

15.2 Custom result templates

Main dialog: **System ► Templates ► Custom result templates**

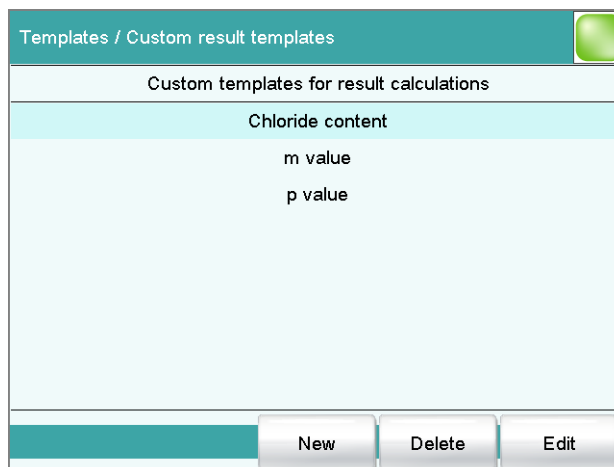

In the dialog **Templates / Custom result templates**, you can define the result calculations which can be loaded in the calculation command **CALC**. With the exception of the following points, creating a result template is identical to editing a calculation in the command **CALC** (*see chapter 29.10.1.2, page 391*):

- No result variable can be assigned to the template.
- The definition of limit values is not possible.



NOTICE

Up to nine **wildcards** can be inserted in the calculation formula, the **Variables F1...F9**. Use these variables, e.g. for the molar mass of your analyte. If you have loaded a result template with these wildcards in the calculation command, then you will automatically be prompted to enter the numerical values.

[New]

Add a new result template to the list, see following chapter.

[Delete]

Delete the selected result template from the list.

[Edit]

Edit the data of the selected result template, see following chapter.

15.2.1 Editing result templates

List of result templates: **Result template ► New / Edit**

Result name

The result name is the text which will be shown in the results display and in the report.

Entry	max. 24 characters
Default value	R

Calculation formula

Shows the calculation formula. A special editor is opened for the definition (see chapter 29.10.3, page 397).

Entry	max. 100 characters
Default value	empty

Decimal places

Number of decimal places used to display the result.

Input range	0 to 5
Default value	2

Result unit

The result unit is displayed and saved together with the result.

Entry	max. 10 characters
Selection	% mol/L mmol/L g/L mg/L mg/mL mg/100 g ppm g mg µg mL µL mg/piece °C µg/min mL/min µL/min
Default value	%

[Note]

Entering a note on the calculation.

[Note for wizard]

Enter a note on the calculation. This note is displayed when loading the result template.

[Result options]

Defining additional settings for the calculation.

"Edit result template / Result options" dialog

In the dialog **Edit result template / Result options**, settings for how to process the calculated result are defined.

Display result

on | off (Default value: **on**)

If you deactivate this parameter, the result is neither displayed in the result dialog nor printed in the result report. This can be advisable for intermediate results.

Save result in result table

on | off (Default value: **off**)

The calculated result can be saved in the result table. This may be advisable if e.g. the results of all determinations carried out on a particular day are to be displayed clearly. A maximum of nine results from a determination can be saved in the result table.

Precision

Setting, with which accuracy the result is used in additional calculations.

Selection	Round Truncate Full precision
Default value	Round

Round

The result is rounded to the defined number of decimal places (commercial rounding, in accordance with the US Pharmacopeia USP). If the digit at the first dropped decimal place is **1, 2, 3 or 4**, then it will be rounded off; if this digit is **5, 6, 7, 8 or 9**, then it will be rounded up. Negative digits will be rounded in accordance with their amount, i.e. away from zero.

Truncate

The result is cut to the number of decimal places defined.

Full precision

The result is used with full accuracy (floating point number either in "single precision" (32 bit) or in "double precision" (64 bit), according to the standard IEEE 754).

15.3 Input lines

Main dialog: **System ▶ Templates ▶ Input lines**

Signal name	Input signal
Cond OK	*****1*
End1	****1***
End2	*1*****
EndMeter	***11***
Ready*	**1****1
Ready1	*****1
Ready2	**1*****
Sample ready	***1****

In the dialog **Templates / Input lines**, you can define the input signals at the remote interface as a template. You can select these templates in the command **SCAN**. The list can contain a maximum of 20 templates.

Table 7 List of predefined input signals

Signal name	Input signal	Function
Cond OK	*****1*	queries "Cond OK" condition of device.
End1	****1***	waits for the EOD pulse from Device 1 (Titrino, Titrande).
End2	*1*****	waits for the EOD pulse from Device 2.
EndMeter	***11***	waits for the EOD pulse from 780/781 pH/Ion meter (stirrer 1 will be switched on during the waiting period).
Ready*	**1****1	queries the "Ready" condition of Devices 1 and 2 (Titrino, Titrande). The status of devices working in parallel can be scanned with this signal. In this case the "Ready" line of both devices must be statically set (i.e. constantly) at the end of a determination. Devices which only transmit a brief impulse cannot be controlled in parallel.
Ready1	*****1	queries the "Ready" condition of Device 1.
Ready2	**1*****	queries the "Ready" condition of Device 2.

Signal name	Input signal	Function
Sample ready	***1***	waits for a stepping pulse, of, for instance, a connected Sample Processor, as soon as it is ready.

[New]

Add a new input signal to the list, see following chapter.

[Delete]

Delete the selected input signal from the list.

[Edit]

Edit the data of the selected input signal, see following chapter.

15.3.1 Editing the input signal

List of input signals: **Input signal ► New / Edit**

Signal name

Designation of the template.

Entry	24 characters maximum
-------	------------------------------

Input signal

Entering the desired bit pattern:

Entering the bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status

The input lines are always numbered from right to left, i.e. with the signal *******1** line 0 is expected to be active.




NOTICE

We recommend masking lines that are of no interest or for which no defined condition can be predicted with an asterisk (*).

Entry	Bit pattern of exactly 8 characters
Default value	*****

15.4 Output lines

Main dialog: **System ► Templates ► Output lines**

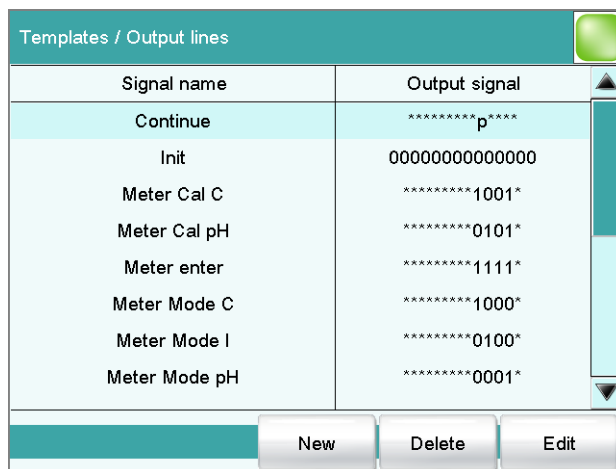

In the dialog **Templates / Output lines**, you can define the output signals at the remote interface as a template. You can select these templates in the command **CTRL**. The list can contain a maximum of 20 templates.

Table 8 List of predefined output signals

Signal name	Output signal	Function
Continue	*****p****	sends a stepping pulse to the connected Sample Processor.
Init	00000000000000	initializes the remote interface.
Meter Cal C	*****1001*	switches the 781 pH/Ion meter to concentration calibration.
Meter Cal pH	*****0101*	switches the 780/781 pH/Ion meter to pH calibration and starts the calibration.
Meter enter	*****1111*	simulates the [Enter] key of the 780/781 pH/Ion meter; mandatory for pH calibration in order to start the measurement of the second buffer.
Meter Mode C	*****1000*	switches the 781 pH/Ion meter to concentration measurement.
Meter Mode I	*****0100*	switches the 780/781 pH/Ion meter to voltametric measurement with polarization current and starts the measurement.
Meter Mode pH	*****0001*	switches the 780/781 pH/Ion meter to pH measurement and starts the measurement.

Signal name	Output signal	Function
Meter Mode T	*****0010*	switches the 780/781 pH/Ion meter to temperature measurement and starts the measurement.
Meter Mode U	*****0011*	switches the 780/781 pH/Ion meter to voltage measurement and starts the measurement.
Start device*	*****p****p	starts Devices 1 and 2 (e.g. Titrino, Titrand, etc.) *).
		The signal is transmitted as a short pulse of 200 ms.
Start device1	*****p	starts Device 1 (e.g. Titrino, Titrand, etc.) *).
		The signal is transmitted as a short pulse of 200 ms.
Start device2	*****p*****	starts Device 2 (e.g. Titrino, Titrand, etc.) *).
		The signal is transmitted as a short pulse of 200 ms.
Start Dos*	****p*p*****	starts Dosimat at Devices 1 and 2 (Titrino via "activate").
		The signal is transmitted as a short pulse of 200 ms.
Start Dos1	*****p*****	starts Dosimat at Device 1 (Titrino via "activate").
		The signal is transmitted as a short pulse of 200 ms.
Start Dos2	****p*****	starts Dosimat at Device 2 (Titrino via "activate").
		The signal is transmitted as a short pulse of 200 ms.

*) In the case of the 780/781 pH/Ion meter, a result report is triggered.

[New]

Add a new output signal to the list, see following chapter.

[Delete]

Delete the selected output signal from the list.

[Edit]

Edit the data of the selected output signal, see following chapter.

15.4.1 Editing the output signal

List of input signals: **Output signal** ► **New / Edit**

Signal name

Designation of the template.

Entry	24 characters maximum
-------	------------------------------

Output signal

Entering the desired bit pattern:

Entering the bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status
- p = set pulse

The output lines are always numbered from right to left, i.e. with the signal *******1** line 0 is set.



NOTICE

We recommend masking lines that are of no interest or for which no defined condition can be predicted with an asterisk (*).

Entry	Bit pattern of exactly 14 characters
Default value	*****


Pulse length

Duration of the sent pulses.

Input range	100 to 1000 ms
Default value	200 ms

15.5 Custom calibration buffers

Main dialog: **System ► Templates ► Custom calib. buffers**

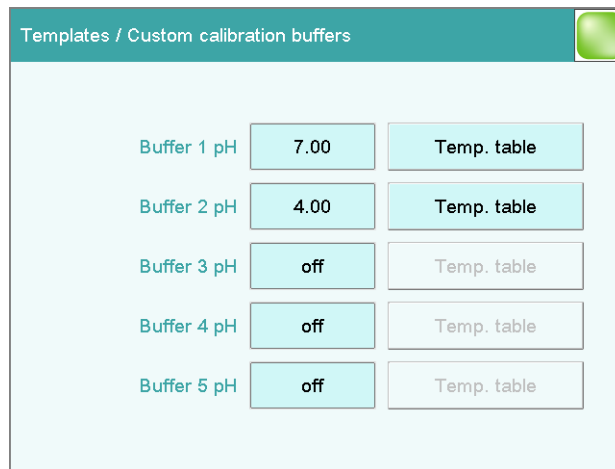

In the **Templates / Custom calibration buffers** dialog, you can define a buffer series of your own for pH calibrations with a maximum of five calibration buffers. Automatic buffer recognition is active for this buffer.

Buffer 1 pH

pH value of the first calibration buffer.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the .

Input range	-20.000 to 20.000
Selection	off
Default value	off

Buffer 2...5 pH

See **Buffer 1 pH**.

[Temp. table]

Display the temperature table for the selected buffer, see following chapter.

15.5.1 Defining calibration buffers

The **Custom calibration buffers / Temperature table** dialog lists the pH values of the buffer in the temperature range 0 - 95 °C.



Custom calibration buffers / Temperature table	
Temperature in °C	pH value
0.0	off
5.0	off
10.0	7.06
15.0	7.04
20.0	7.02
25.0	7.00
30.0	6.99
35.0	off

[Edit pH value]

Enter the pH value for the selected temperature.

"Temperature table / Edit pH value" dialog

In this dialog, the pH values of the buffer are defined at various temperatures. Enter the pH values for the temperature range in which you will carry out your pH calibration and pH measurement. If you do not know the pH values at individual temperatures, they will be calculated automatically by means of linear interpolation.

Temperature in °C

Input range	0.0 to 95.0 °C (Increment: 5.0)
-------------	--

pH value

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the .

Input range	-20.000 to 20.000
Selection	off
Default value	off

15.6 Report header

Main dialog: **System ► Templates ► Report header**

In the dialog **Templates / Report header**, you can create a system-specific report header that contains information concerning the laboratory, for example. The report header can contain a maximum of four lines of 46 characters each and will always be printed out in front of the standard report header. The standard report header contains the instrument type, serial number and program version of the software, instrument name and printing date.

Print logo

on | off (Default value: **off**)

If this parameter is activated, then the logo defined below will be printed out in the right margin.

Logo

Selection	Metrohm logo Custom logo
Default value	Metrohm logo

Custom logo

Select this setting if your logo is to be printed out in the report header. How to create a logo is described below.

Applying the custom logo

You can create a graphics file of your own and use it as a logo in the report header in place of the Metrohm logo.



NOTICE

Observe the size of the graphics object. The maximum size of your graphics file is 64 kB. We recommend a maximum of 500 × 200 pixels.

Proceed as follows to create your own logo:

1 Creating the logo

- Create a graphics file with the desired logo in an image processing program.
- Save the graph as a JPG file under the file name "CustomImage.jpg".



2 Copying the file to an external storage medium

- Copy the file "CustomImage.jpg" to the directory "916".
If this structure is not maintained, the file will not be found because the software directly accesses the "916" directory.
- Plug in the external storage medium at the Ti-Touch.

3 Loading the graphics file

- In the **Templates / Report header** dialog, tap on the **[Load]** button.
The file is copied to the internal memory.
- Switch the instrument off and back on again.

4 Applying the custom logo

- In the **Templates / Report header** dialog, activate the **Print logo** option.
- Define **Logo = Custom logo**.

Your logo will now be printed in the report header of all reports in the future.

15.7 Custom electrode type

Main dialog: **System ► Templates ► Custom electrode type**

For the electrode test of pH electrodes, you can define your own values instead of using the limit values that are saved. This can be helpful for special applications, e.g. when other quality requirements are demanded of the pH electrode or when working with a reference electrode that pushes the offset potential U_{off} beyond the required -15 to $+15$ mV (standard electrode).

Limit values can be defined for the following electrode ratings:

- Excellent electrode
- Good electrode
- Usable electrode

Lower limit U_{off}

Lower limit value for the offset voltage, i.e. the voltage at $\text{pH} = 7.0$. The value applies for all ratings.

Input range	-999 to 999 mV
Default value	-15 mV

Upper limit Uoff

Upper limit value for the offset voltage, i.e. the voltage at pH = 7.0. The value applies for all ratings.

Input range	-999 to 999 mV
Default value	15 mV

15.7.1 Limit values for the electrode rating

You can define limit values for the three electrode ratings **Excellent electrode**, **Good electrode** and **Usable electrode**. These values represent maximum values that may not be exceeded (exception: **Min. slope** = minimum value).

Streaming potential

The streaming potential is the difference between the measured voltages in stirred and unstirred solution.

for very good electrodes:

Input range	-999.9 to 999.9 mV
Default value	2.5 mV

for good electrodes:

Input range	-999.9 to 999.9 mV
Default value	3.0 mV

for usable electrodes:

Input range	-999.9 to 999.9 mV
Default value	4.0 mV

Drift

The drift in stirred solution is determined for each buffer. The sum of these values is compared with this value here.

for very good electrodes:

Input range	0.1 to 9.9 mV/min
Default value	2.0 mV/min

for good electrodes:

Input range	0.1 to 9.9 mV/min
Default value	2.5 mV/min

for usable electrodes:

Input range	0.1 to 9.9 mV/min
Default value	3.0 mV/min



Min. slope

Minimum slope of the pH electrode.

for very good electrodes:

Input range	0.1 to 999.9 %
Default value	96.5 %

for good electrodes:

Input range	0.1 to 999.9 %
Default value	96.0 %

for usable electrodes:

Input range	0.1 to 999.9 %
Default value	95.0 %

Max. slope

Maximum slope of the pH electrode.

for very good electrodes:

Input range	0.1 to 999.9 %
Default value	101.0 %

for good electrodes:

Input range	0.1 to 999.9 %
Default value	102.0 %

for usable electrodes:

Input range	0.1 to 999.9 %
Default value	103.0 %

Response time

The voltage measured after three minutes in stirred solution serves as a comparison value for determining the response time. The response time is the time after which the measured voltage is within ± 1 mV of this comparison value.

for very good electrodes:

Input range	0 to 999 s
Default value	45 s

for good electrodes:

Input range	0 to 999 s
Default value	50 s



for usable electrodes:

Input range	0 to 999 s
Default value	60 s

16 Methods


16.1 Creating a new method

Proceed as follows to create a new method:

1 Opening the method table

- In the main dialog, tap on **[Load method]**.
- Tap on **[New method]**.

The method table with the stored templates opens:

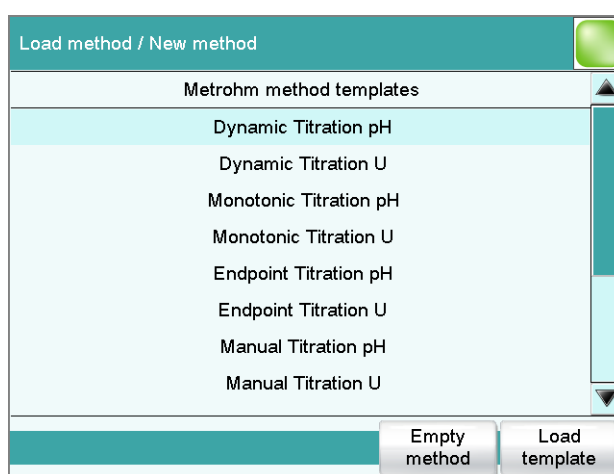

2 Loading the method

- Tap on **[Empty method]**.

or

- Select the desired template and tap on **[Load template]**.

The method is now loaded. **New method** is displayed in the main dialog in the title line.

If a new method has been created, then the individual parameters can be modified with **[Edit parameters]**.

16.2 Saving a method

If you modify method parameters, then you can save these as your own method.

To save a method, proceed as follows:

1 Open the command sequence

- In the main dialog, tap on **[Edit parameters]**.

The command sequence of the method is displayed:

Parameters / Sequence		
Current method: New method		
01	MET pH	Monotonic pH titration
02	CALC	Calculation
03	REPORT	Report
04	...	

Save method Method options Insert command Delete command Edit command

2 Modifying/applying the method name

- Tap on **[Save method]**.
A method name will be suggested for new methods. If the method has already been saved once, then the method name will be displayed.

Sequence / Save method	
Memory	Internal memory ▼
Group	Main group ▼
File name	Me2115 ▼

Cancel Save

Applying the name:

- Tap on **[Save]**.

The method will be saved and the command sequence is displayed.

Entering a new name:

- Tap on the input field **File name**.
The text editor opens.
- Enter a new file name (max. 32 characters) and confirm with **[OK]**.
- Tap on **[Save]**.

The method will be saved and the command sequence is displayed.


16.3 Loading a method

Proceed as follows to load a method:

1 Opening the method table

- In the main dialog, tap on **[Load method]**.
The selection of memory locations is displayed. This selection will be skipped if only the internal memory is accessible.
- Select the memory location where the desired method is saved.
The selection of file groups is displayed (*see chapter 12.1, page 136*). If only one group is available, then the method table will be displayed directly.
- Select the group with the desired method.
- Tap on **[Show files]**.

The method table with the stored methods opens:

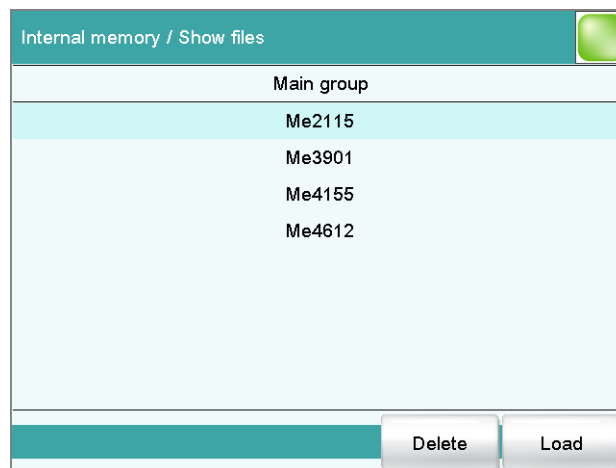

2 Selecting the method

- Select the desired method.

3 Loading the method

- Tap on **[Load]**.

The method is now loaded. The method name is displayed in the main dialog in the title line.



NOTICE

The data of the current determination will be deleted when you load a method.

16.4 Editing a method

Main dialog: **Edit parameters**

Parameters / Sequence		
Current method: New method		
01	MET pH	Monotonic pH titration
02	CALC	Calculation
03	REPORT	Report
04	...	

Save method Method options Insert command Delete command Edit command

In the dialog **Parameters / Sequence**, you can define and edit the method commands which are executed successively during a determination.

The following data is displayed in the command list for each command:

- Line number
- Name
- Comment
You can adjust the comment in accordance with your requirements.

[Save method]

Save the current method (*see chapter 16.2, page 183*).

[Method options]

Define various settings which refer to the entire method, e.g. statistics, automatic saving of determinations, settings for sample data, etc. (*see chapter 16.5, page 187*).

[Insert command]

Insert a new method command. It is inserted before the selected command.

**NOTICE**

Not all commands are available for subsequences. Commands which cannot be inserted into subsequences are disabled.

[Delete command]

Delete the selected method command.

[Edit command]

Edit the selected method command.

16.4.1 Inserting a command

A method is comprised of individual commands. When you start a determination, the commands are executed one after the other.

The following constraints apply:


- Maximum of 99 commands
- Maximum of nine commands for titrations, measurements and calibrations

Inserting a command

To insert a command in a method, proceed as follows:

1 Open the command overview

- In the main dialog, tap on **[Edit parameters]**.
- Tap on the **[Insert command]** button.

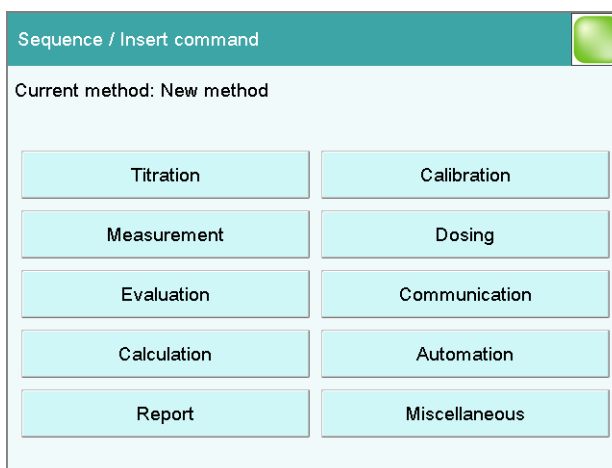

This dialog contains all of the commands which can be inserted into a method run, organized into thematic groups.

2 Select a command group

- Tap on the desired command group.

In the case of **[Report]**, the command is inserted directly into the command list. For all other command groups, the available commands will be displayed (titration / measuring commands, etc.).


3 Inserting a command

Tap on the desired command.

The dialog is closed and the command is displayed in the command list.

16.5 Method options

Main dialog: **Edit parameters ► Method options**

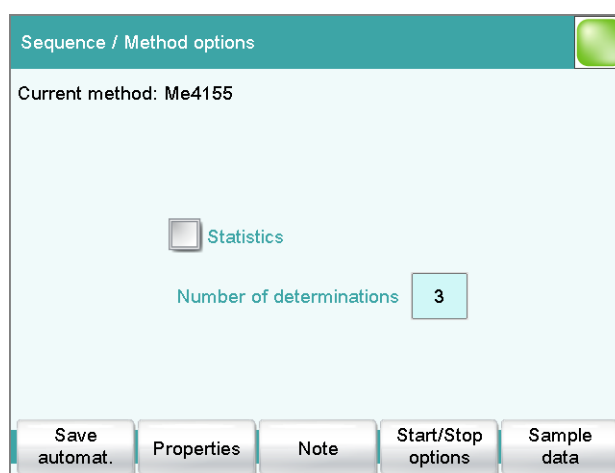

Method options are settings which apply to the method as a whole and not to an individual command.

The following settings are available:

- Activate/deactivate statistics calculations
- Save determination automatically
- Sign a method electronically
- Check the method when the determination starts
- Define a fixed sample size
- etc.



Statistics

on | off (Default value: **off**)

If this parameter is activated, then statistics calculations will be carried out for all of the defined results. The prerequisite here, however, is that a variable for the mean value is defined in the result options (see "Dialog "Edit calculation / Result options"", page 393).

Number of determinations

The number of determinations that are carried out for the statistics calculations.

If an additional determination has to be added to the determination series, because one determination has been incorrect, for example, then this can be accomplished in the statistical overview (see chapter 25, page 246).

Input range	2 to 20
Default value	3

[Save automat.]

Define the settings for automatically saving the determination (see chapter 16.5.6, page 197).

[Properties]

Display the properties of the current method and sign electronically (see chapter 16.5.4, page 195).

[Note]

Enter a note on the method (see chapter 16.5.5, page 196).

[Start/Stop options]

Edit functions which are carried out immediately after the method is started or stopped, see chapter Start options, page 188 and chapter Stop options, page 190.

[Sample data]

Specify method-specific settings for the sample data, e.g. designation of the sample identifications, limits for the sample size (see chapter 16.5.3, page 191).

16.5.1 Start options

Main dialog: **Edit parameters ► Method options ► Start/Stop options ► Start options**

In the dialog **Method options / Start options**, you can configure the functions which are carried out when a determination is started.

Method check at start

on | off (Default value: **on**)

If this option is activated, then the following tests are carried out at the method start.

- Are the required control and peripheral devices connected?
- Are the titrant and the sensor on hand?
- Have the monitoring intervals expired?
- etc.



NOTICE

We recommend that the method check only be deactivated if the titration must begin immediately after the start of the determination (e.g. with very rapid reactions such as enzyme-catalyzed reactions). If the method check is deactivated, it could happen that the determination run will be interrupted, e.g. because devices, titrants or sensors are not on hand.

The following settings can be implemented for a connected Sample Processor:

Rack reset

on | off (Default value: **off**)

If this option is activated, then the sample rack is initialized at the time of the method start. The following actions are carried out:

- The lift is moved upwards.
- The sample rack is rotated to the starting position.
- The rack code of the rack attached is being read out and the corresponding rack data is transferred to the Sample Processor.
- The sample variable is reset to the value 1.

Check rack

Definition of the sample rack which must be attached at the time of the method start. This ensures that the method will be carried out only with this sample rack.



NOTICE

The sample rack can, however, only be checked if the **Rack reset** option is activated.

Selection	no Selection of the configured sample racks
Default value	no

Increase sample variable automatically

on | **off** (Default value: **on**)

The sample variable describes the current position of the sample on the rack of the Sample Processor. If this option is activated, then the sample variable is automatically increased by 1 at the end of each determination.

This option must be deactivated if the sample variable is to be altered specifically with a **SAMPLE** command.

Request rack position at start

on | **off** (Default value: **off**)

If this option is activated, then the rack position of the first sample to be processed is requested.

16.5.2 Stop options

Main dialog: **Edit parameters** ► **Method options** ► **Start/Stop options** ► **Stop options**

In the dialog **Method options / Stop options**, you can define the actions which are carried out when a method is canceled.

The method can be canceled as follows:

- Manual stop with the fixed key []
- Stop because of an error
- Stop by remote signal via the Control Remote Box

Switch off pumps

on | **off** (Default value: **on**)

If this option is activated, then all of the connected pumps will be switched off.

Switch off stirrers

on | **off** (Default value: **on**)

If this option is activated, then all of the connected stirrers will be switched off.

Set remote lines

Selection of the signal out of the templates or entering the required bit pattern. Templates are defined under **System** ► **Templates** ► **Output lines**.

**NOTICE**

A line set active is not being reset automatically, not even at the end of the determination.

Entering the bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status
- p = set pulse

The output lines are always numbered from right to left, i.e. with the signal *****1 line 0 is set. With a pulse, the length is set to 200 ms. If you wish to set pulses with other lengths, you have to define a corresponding template.

Entry	Bit pattern containing exactly 14 characters or a max. of 24 characters for the name of the template
Default value	*****
Selection	Selection of the templates defined

16.5.3 Sample data

Main dialog: **Edit parameters ► Method options ► Sample data**

You can define the following method-specific settings in the dialog

Method options / Sample data:

- Modify the designation for the sample identifications
- Define a fixed sample size
- Define limits for the sample size
 - If you define limits, these will be monitored at the start and at the end of the determination.

Below you will find an example demonstrating the effects on the main dialog of the settings in this dialog (see "*Modifying sample data*", page 193).

Identification 1, Identification 2

on | off (Default value: **on**)

If this option is activated, then the input field for the sample identification in the main dialog will be displayed.

Designation

A designation of your own choosing can be defined for the input field for each method.

Entry	16 characters maximum
-------	------------------------------

Fixed sample size

on | off (Default value: **off**)

Activate this parameter if the same sample amount is always to be used for all determinations. If this is case, then you can define the sample size here. It will be displayed in the main dialog, but it can no longer be changed there.

Sample size

This parameter can only be edited when **Fixed sample size** is activated.

Input range	-999999999 to 999999999
Default value	1.0

Unit of sample size.

Selection	g mg µg mL µL pieces
Default value	g



NOTICE

You can also use the parameter **Fixed sample size** if you would like to establish a default for the sample size.

Example: For a particular determination, you require 10 mL of sample. Now and again, however, you have too little sample available and you must perform the determination with 5 mL, for instance. In this case, proceed as follows:

1. Activate the parameter **Fixed sample size**.
2. Define the value and unit for the sample size, in the above example, for instance, **[10 mL]**.
3. Deactivate the parameter **Fixed sample size** again.
4. Save method.

⇒ Each time you load this method, 10 mL is entered as the sample size, but in the example mentioned you can enter the lesser quantity of 5 mL.

[Sample size limits]

This button is accessible only if **Fixed sample size** is deactivated.

Define the limit values for the sample size (see "Defining limit values for the sample size", page 194).

Modifying sample data

This instruction is for the purpose of clarifying the settings which are possible in this dialog. The following settings should be made:

- Change the designation for the input field of the first sample identification.
- Hide the input field for the second sample identification.
- Define a fixed sample size.

Proceed as follows:

1 Change the designation for the input field

- Tap on the input field **Identification 1**.
The text editor opens.
- Change the designation in **Batch** and confirm with **[OK]**.
- Deactivate the parameter **Identification 2**.


2 Define a fixed sample size

- Activate the parameter **Fixed sample size**.
- Enter the value **10 mL** as **Sample size**.

3 Displaying the main dialog

Tap on the fixed key [].

The data is saved and the adjusted main dialog appears.

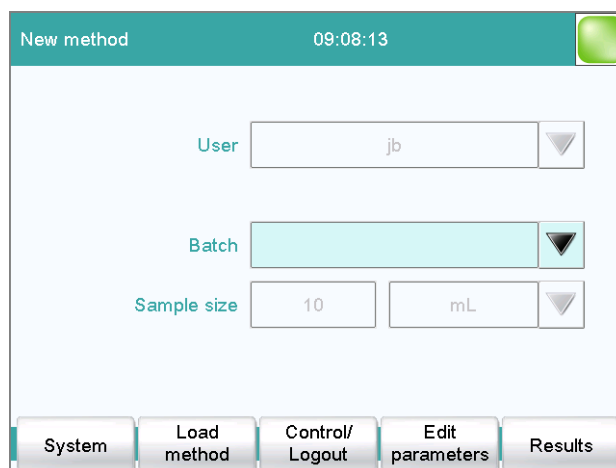

Only Identification 1 is shown with the title "Batch." The fixed sample size "10 mL" is shown, but cannot be edited.

Defining limit values for the sample size

In the dialog **Sample data / Sample size limits**, you can define the limit values for the sample size. These values are not monitored during sample data input, but rather:

- when the determination is started.
- when the determination is finished.
- upon automatic sample data query via the command **REQUEST**.
- when the determination is being recalculated.

An entry is made in the message list automatically when these limits are infringed (see chapter 24.2, page 233).

Monitoring sample size limits

on | off (Default value: **off**)

If this parameter is activated, then the sample size will be monitored.

Lower limit

Input range	-999999999 to 999999999
Default value	0

Upper limit

Input range	-999999999 to 999999999
Default value	999999999

Display message

on | off (Default value: **off**)

If this parameter is activated, a message is displayed in case the limits are infringed. You can select whether you want to continue with the determination or cancel the run.

If this parameter is deactivated, then the message will be entered only in the message list of the determination.

Timeout

The message is displayed during this time. Afterwards, the determination is continued automatically.

Input range	0 to 999999 s
Default value	30 s
Selection	off

off

The determination will not be continued until after the message has been confirmed.

16.5.4 Method properties

Main dialog: **Edit parameters** ► **Method options** ► **Properties**

Detailed information concerning the method is displayed in the **Method options / Properties** dialog and the methods can be signed electronically here:

- **Method status**

Current status of the method.

- **new**
The method has been newly created and not yet saved.
- **saved**
The current method has been saved.
- **modified**
The current method has been modified.
- **reviewed**
The current method has been signed at level 1.
- **released**
The current method has been signed at level 2.

- **Created by**

User who created the method.

- **Created on**

Date and time at which the method was created.

- **Last saved by**

User who saved the method most recently.

- **Last saved on**

Date and time at which the method was most recently saved. The version of the method is displayed as well. The version number will be increased by one each time the file is saved.

The following data is displayed only if the option **Saving modified method** under **Login options / Modification options** is activated:

- **Modified by**

User who modified and saved the method most recently.

- **Modified on**

Date and time at which the method was most recently modified and saved.

- **Reason**

Reason for the modification of the method.

- **Comment**

Comment for the modification of the method.

The following information is only shown if the method has been signed at level 1:

- **Reviewed by**

User who signed the method at level 1.



- **Reviewed on**
Date and time at which the method has been signed at level 1.
- **Reason**
Reason for the signing of the method.
- **Comment**
Comment for the signing of the method.

The following information is only shown if the method has been signed at level 2:

- **Released by**
User who signed the method at level 2.
- **Released on**
Date and time at which the method has been signed at level 2.
- **Reason**
Reason for the signing of the method.
- **Comment**
Comment for the signing of the method.

[Delete signature]

Delete all of the signatures for the method or the determination. This button is inactive if level 2 has not yet been signed (*see chapter 17.2, page 200*).

[Sign]

Sign the method or determination. This button is active only when working with activated login function and password (*see chapter 17.1, page 199*).

16.5.5 Note

Main dialog: **Edit parameters ► Method options ► Note**

In this dialog, you can enter a short text, e.g. important information about carrying out the determination.

[Display options]

Definition when the note is being displayed.

Automatically after loading the method

on | off (Default value: **off**)

If this parameter is activated, the note is displayed when loading the method. It can otherwise only be read in this dialog.

16.5.6 Saving a determination automatically

Main dialog: **Edit parameters** ► **Method options** ► **Save automat.**

In the dialog **Method options / Save automatically**, you can specify whether the determination is to be saved automatically and/or whether a PC/LIMS report is to be created.

Save determination automatically

on | off (Default value: **off**)

If this parameter is activated, then the determination will be saved automatically.

Memory

Memory location where the determination will be saved.

Selection	External memory 1 External memory 2 Shared memory
Default value	External memory 1

Shared memory

The determination will be saved in a shared directory on the network. The shared directory is selected in the **Edit device / Shared memory** dialog (see chapter 11.3.3, page 100).

Group

Directory where the determination will be saved.

Entry	32 characters maximum
Selection	Selection of available directories empty
Default value	empty

File name

File name of the determination. The file name is always extended to include the date and the time of day (YYYYMMDD-hhmmss) in order to ensure that the file names are unique for all determinations.

Entry	max. 16 characters
Selection	Identification 1 Identification 2 Method
Default value	Identification 1

Identification 1

The first 16 characters of the text which was entered in the main dialog for Identification 1 + YYYYMMDD-hhmmss.

Identification 2

The first 16 characters of the text which was entered in the main dialog for Identification 2 + YYYYMMDD-hhmmss.

**Method**

The first 16 characters of the text, of the method name +
YYYYMMDD-hhmmss.

Write protection

on | off (Default value: **on**)

If this parameter is activated, then the file cannot be saved, deleted, or renamed. This is only an internal write protection feature and is independent of the write protection of the operating system on your computer. This write protection feature protects the saved determination data against accidental modifications or modifications by unauthorized persons.

Create PC/LIMS report

on | off (Default value: **off**)

If this parameter is activated, then a machine-readable report with all of the important data for the determination is created, a so-called PC/LIMS report.

The settings are implemented in the device manager (*see chapter 11.3.2, page 99*).

When you save the PC/LIMS report as a file, the file name will be generated automatically as follows: **PC_LIMS_Report**-first 16 characters of the value of the parameter **File name-Date-Time.txt**. If the value of the parameter **File name** contains one or more of the following characters, then these will be automatically converted in the file name into the character "_": / \ : * ? " < > |.

17 Electronic signatures

17.1 Signing methods/determinations electronically

Methods: **Main dialog ▶ Edit parameters ▶ Method options ▶ Properties ▶ Sign**

Determinations: **Results dialog ▶ More data ▶ Properties ▶ Sign**

You can sign methods or determinations in the dialog **Electronic signature**. The following information is saved along with each electronic signature:

- User (complete name)
- Date and time
- Reason
- Comment

Methods can only be signed if they have the status **saved**. The status of a method can be found in the dialog **Method options / Properties**.

Signature

Selection of the level at which the determination should be signed.

Selection	Review (signature level 1) Release (signature level 2)
	A maximum of three signatures from up to three different users can be given at each level. However, it is not possible for a user to sign at both level 1 and level 2.

Review (signature level 1)

If the method has never been signed before, then only level 1 can be selected.

Release (signature level 2)

If the method has been signed three times at level 1 or already once at level 2 then only level 2 can be selected.

User

Identification of the user who signs.

Password

Password of the user.



Reason

Selection of the reasons for the signature. The list with reasons is compiled in the user administration (*see chapter 7.3.6, page 56*).

Selection	Selection of configured reasons
-----------	--

Comment

Entry	24 characters maximum
-------	------------------------------

[Change password]

Change the current password.

[Sign]

Sign the method or determination with the data entered above.

17.2 Deleting electronic signatures

If a released method must be revised, e.g. in order to adapt to altered conditions, then the signatures must be deleted. The method can otherwise not be modified. In the case of released determinations, the signatures can also be deleted. Signatures, however, cannot be deleted until after the method/determination has been signed at Level 2.

Signatures are deleted automatically in the following cases:

- A method signed at Level 1 is being altered.
- A determination signed at Level 1 is being recalculated.

User

Identification of the user who deletes the signatures.

Password

Password of the user.

Reason

Selection of the reasons for the deletion of the signature. The list with reasons is compiled in the user administration (*see chapter 7.3.6, page 56*).

Selection	Selection of configured reasons
-----------	--

Comment


Entry	24 characters maximum
-------	------------------------------

[Delete signature]

Delete all the signatures for the method or the determination. The method is assigned the status **saved**.

18 Control

Main dialog: **Control**

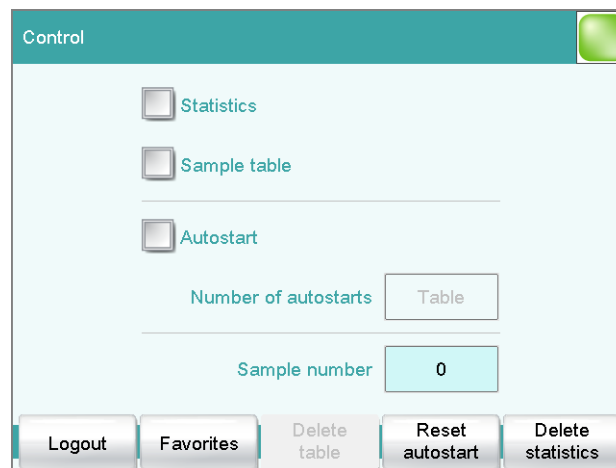

In the dialog **Control**, the settings for the execution of a single determination or of one sample series are defined.

If you work with the login function activated, you can log off the system in this dialog. The login dialog will then be shown immediately.

Statistics

on | off (Default value: **off**)

With this option, you can enable or disable the statistics calculations for individual determinations (see "Statistics", page 203).



NOTICE

The parameter **Statistics** in the method options will not be changed as a result. This setting here will be adjusted automatically to match the one in the method options.

Sample table

on | off (Default value: **off**)

If this option is activated, the sample data for a sample series can be entered in a table (see chapter 21, page 212).

Autostart

on | off (Default value: **off**)

If this option is activated, a new determination is started automatically at the end of a determination. This continues until the number specified has been reached (siehe **Number of autostarts**).



NOTICE

If you deactivate this option while a determination is running, then that determination will be ended, although the next one will however also not be started. This can be used to cancel a sample series.

Number of autostarts

Number of automatic starts.

Input range	1 to 9999
Selection	Table
Default value	Table

Table

The number of automatic starts corresponds to the number of samples in the sample table.

Sample number

The sample number is primarily used for documentation. Each time a determination is started, it is increased by one. The sample number can, however, also be entered manually. Each time the Ti-Touch is switched on, the value is reset to zero.

Input range	0 to 9999
Default value	0

[Logout]

User logout (visible only when work is done with activated login function).

[Favorites]

Save methods/sample tables as a favorite (see chapter 19, page 205). Depending on whether you are working with or without login function, user-specific or common favorites can be created. On the main dialog, a button is created for each favorite. This way, a method or a sample table can be started with one single mouse click.

[Delete table]

Delete the sample table completely.

[Reset autostart]

Reset the value of the autostart counter to zero.

[Delete statistics]

Delete all statistics data.

The statistics data should be deleted manually if a sample series has been canceled and a new sample series is to be started.

The statistics data is deleted automatically in following cases:

- when all of the determinations of the determination series have been carried out and a new determination has been started afterwards.
- when a new method is loaded.
- when a determination is loaded (the method with which the determination was carried out is loaded simultaneously with the determination).

Statistics

In the dialog **Control**, you can deactivate the statistics calculations for individual determinations. This option is required primarily in order to insert the processing of an "**urgent sample**" during the processing of a sample series when the urgent sample is not to be included in the statistics calculations.

Interrupting a determination series for an "urgent sample"

If the same method is required for the "urgent sample" as is being used for the sample series, then you need only deactivate the option **Statistics** and then reactivate it after the determination is completed. Proceed as follows if you need to process the "urgent sample" with a different method:

1 Deactivate statistics calculation

- Deactivate the option **Statistics**.

2 Saving the determination

- Save the current determination of the sample series (*see chapter 24.6, page 242*).

3 Carrying out the determination

- Load the method for the "urgent sample".
- Carry out the determination.

4 Load the last determination of the sample series

- Reload the previously saved last determination.

The determination, the method used for it and the current statistical data are loaded. The option **Statistics** is activated again and the

value of the statistics counter corresponds to that in effect before the interruption.

5 Continuing the sample series


- Carry out the remaining determinations in the sample series.

19 Favorites

Main dialog: **Control ► Favorites**

Main dialog: **System ► System settings ► User admin. ► Edit ► Favorites**

This chapter describes how you can create favorites. Favorites are comparable to the favorites/bookmarks in your Internet browser. In the main dialog, a button is created for each favorite. By pressing a single key, you can trigger an action without having to navigate through different dialogs.

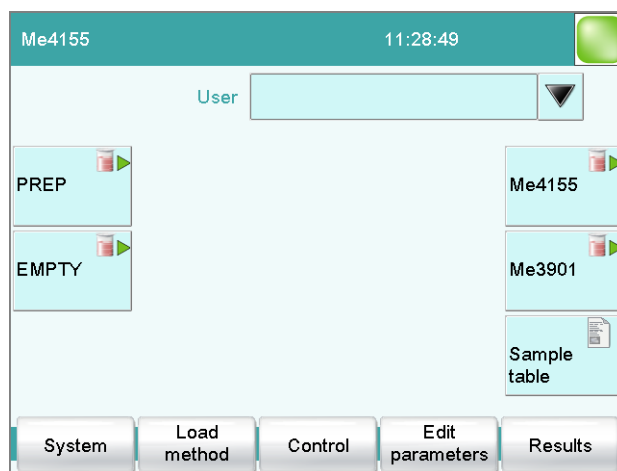

The following objects can be saved as favorites:

- Methods
- Sample tables

User-specific or common favorites can be created, depending on whether you are working with or without the login function.

- Operation with login function:
If you are working with the login function activated, **user-specific favorites** can be created. These can only be used by a certain user. User-specific favorites are created as follows:
 - in the user administration by a user with administrator rights (**System ► System settings ► User admin. ► Edit**).
 - in the dialog **Control** for the user who is logged in.
- Operation without login function:
If you are working without the login function, **common favorites** can be created. These favorites are available for all users. Common favorites are created in the **Control** dialog.



19.1 Creating favorites

Control / Favorites	
Name	Position
PREP	1
EMPTY	6
Me4155	5
Me3901	10

The list of favorites gives the designation and position of the button on the main dialog of each favorite configured.

[New]

Add a new favorite to the list, see following chapter.

[Delete]

Delete the selected favorite from the list.

[Edit]

Edit the properties of the selected favorite, see following chapter.

19.1.1 Editing favorites

Favorites / Edit	
Position	5 ▼
Name	Me4155
Type	Method ▼
Memory	Internal memory ▼
Method / Sample table	Me4155 ▼



Position

On the main dialog, a button is created for each favorite. These buttons are ordered in three rows at fixed positions. The position 1 can be found top left.

Input range	1 to 14
Default value	1

Name

The designation of the favorite is used for unambiguous identification.

Entry	24 characters maximum
-------	------------------------------

Type

Definition as to whether the favorite represents a single method or a complete sample table.

Selection	Method Sample table
Default value	Method

Memory

Memory location the method or the sample table is loaded from. Only the currently accessible memory locations are selectable.

Selection	Internal memory External memory 1 External memory 2 Shared memory
Default value	Internal memory

Shared memory

The favorite will be saved in a shared directory on the network. The shared directory is selected in the **Edit device / Shared memory** dialog (see chapter 11.3.3, page 100).

Method / Sample table

Method or sample table of the favorite.

Selection	Selection of stored methods/sample tables
-----------	--

[More options]

Parameterize the autostart function, see following chapter.

More options

The autostart function is parameterized under **[More options]**.



Autostart

on | off (Default value: **off**)

If this option is activated, a new determination is started automatically at the end of a determination. This continues until the number specified has been reached (siehe **Number of autostarts**).

Number of autostarts

This option can only be edited when **Autostart** is activated.

Number of automatic starts.

Input range	1 to 9999
Default value	1
Selection	Table

Table

The number of automatic starts corresponds to the number of samples in the sample table.

20 Sample data

You can enter the sample data (identification, sample size, etc.) in a variety of ways:

- Directly in the main dialog.
- Using the sample table. This is particularly useful with sample series. The sample table is a table in which the sample data for up to 999 samples can be entered (*see chapter 21, page 212*).
- Automatic request immediately after the start of the determination (*see chapter 20.2, page 210*).

You can also send the sample size and the unit from a connected balance in any case. With some balances, the sample identification and method can be also sent (*see chapter 32.5, page 445*).

20.1 Entering sample data in the main dialog

For a sample, you can enter the sample data directly in the main dialog, even while the determination is running (*see chapter 23, page 223*).

Identification 1

Sample identification. The sample identification can be used in calculations as the variable **CI1**. If you have defined a sample identification list (*see chapter 15.1.1, page 163*) then the entries can be selected here.

Entry	24 characters maximum
Default value	empty



Identification 2

Sample identification. The sample identification can be used in calculations as the variable **CI2**. If you have defined a sample identification list (see chapter 15.1.1, page 163) then the entries can be selected here.

Entry	24 characters maximum
Default value	empty

Sample size

Sample size. The value of the sample size can be used in calculations as the variable **C00**.

Input range	-99999999 to 9999999999
Default value	1.0

Unit of sample size.

Entry	6 characters maximum
Selection	g mg µg mL µL pieces
Default value	g

Unit defined in method:

The unit defined in the method is ignored. The unit entered in the sample table is used.

Balance connected and configured:

Value from the connected balance.

Fixed sample size defined in method:

The value will be entered in the corresponding line when the method is loaded. The existing entry will be overwritten.

20.2 Requesting sample data at the start of the determination

The sample data can be automatically requested immediately after the start of the determination in order to ensure that the sample data entry is not forgotten. This automatic inquiry is requisite for reweighing.

Me2115

01 REQUEST Data request

Identification 1

Sample size mL

Continue

To accomplish this, you must insert and configure the command **REQUEST** in the method (see chapter 29.15.3, page 433).

If the **Hold sequence** parameter is activated, then the run will be paused and must be continued with **[Continue]** after the sample data has been entered. If the **Hold sequence** parameter is deactivated, then the determination will continue in the background until the measured data is required for further processing. This dialog will be displayed until the entering of the sample data is confirmed with **[Continue]**, even if the titration/measurement is already completed. This ensures that the sample data is available for calculations.

If a fixed sample size has been defined in the method (see chapter 16.5.3, page 191), then it will be displayed at the time of the entry but it will not be editable.

21 Sample table

21.1 General

Main dialog: **Sample table**

The sample table is a table in which the sample data for up to 999 samples can be entered. If you are processing larger-sized sample series, it makes sense to enter the sample data in this table.


Activating the sample table

Proceed as follows to enable the sample table:

1 Opening the control dialog

- In the main dialog, tap on **[Control]**.


The control dialog is displayed:

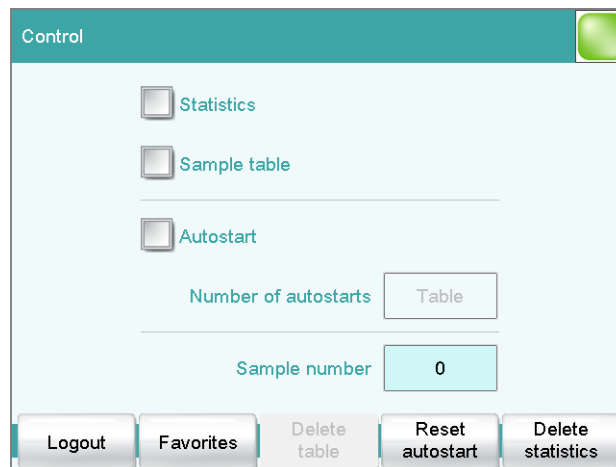

2 Activating the sample table

- Activate the **Sample table** option.
- Tap on the fixed key **[🏠]**.

The main dialog is displayed:

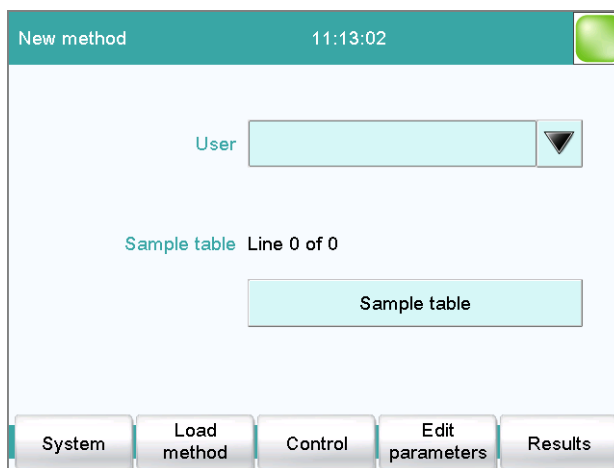

The display will show the number of determinations already carried out and the total number of sample lines containing data. The sample table is still empty in this example.

The sample table contains numbered lines; each line represents one determination. In the standard settings, the first sample identification and the sample size of each sample are displayed. In the properties, you define the columns to be displayed (*see chapter 21.3, page 217*).

Sample table		
No.	Identification 1	Sample size
1	#2370015	1.0 g
2	#2370015	1.0 g
3	#2370016	1.0 g
4	#2370016	1.0 g
5	...	

[Load/ Save]

Load a saved sample table (*see "Loading a sample table", page 215*) or save the current sample table (*see "Saving a sample table", page 214*).

[Properties]

Edit the properties of the sample table (*see chapter 21.3, page 217*).

[Insert line]

Insert a new line above the line selected.

[Delete]

Delete the selected line from the sample table.

[Edit]

Edit the data of the selected line (see chapter 21.2, page 215).

Saving a sample table**NOTICE**

Sample tables can only be saved on external storage media or on a shared file system.

Proceed as follows to save a sample table:

1 Opening the dialog

- In the **Sample table** dialog, tap on the **[Load/ Save]** button.

The **Sample table / Load/Save** dialog is displayed.

If no external storage medium is plugged in or if no file system is shared, then **[Save]** is disabled.

2 Modify/apply a name

- Tap on **[Save]**.
A name is suggested for new sample tables. If the sample table has already been saved, the previous name will be displayed.

Applying the name:

- Tap on **[Save]**.

The sample table is saved.

Entering a new name:

- Tap on the input field **File name**.
The text editor opens.
- Enter a new file name (max. 32 characters) and confirm with **[OK]**.
- Tap on **[Save]**.

The sample table is saved.

Loading a sample table

Proceed as follows to load a sample table:

1 Display the list of saved sample tables

- In the **Sample table** dialog, tap on the **[Load/ Save]** button.
The **Sample table / Load/Save** dialog is displayed.
If no external storage medium is plugged in or if no file system is shared, then **[Load]** is disabled.
- Tap on **[Load]**.
The selection of memory locations is displayed. This selection will be skipped if only one memory can be accessed.
- Select the memory location where the desired sample table is stored.
The selection of file groups is displayed (*see chapter 12.1, page 136*). If only one group is available, then the list of the saved sample tables will be displayed directly.
- Select the group with the desired sample table.
- Tap on **[Show files]**.

The list with the saved sample table is opened.

2 Select the sample table

- Select the desired sample table.

3 Loading a sample table

- Tap on **[Load]**.

The sample table is now loaded.

21.2 Editing sample data

Sample table / Edit

Line number - 1 +

Method Me2115 ▼

Identification 1 #2370015

Identification 2

Sample size 1.0 g ▼



At the very top you will see the line number of the selected line. In this example, the sample data of the first line is displayed. You can scroll between individual data sets with the keys **[-]** and **[+]**.

Method

Method used for processing the sample.

If you use the sample assignment table (*see chapter 15.1.2, page 164*), then the method can no longer be defined here.

Selection	Selection of stored methods empty
Default value	empty

empty

The currently loaded method is used.

Identification 1

Sample identification. The sample identification can be used in calculations as the variable **CI1**. If you have defined a sample identification list (*see chapter 15.1.1, page 163*) then the entries can be selected here.

Entry	24 characters maximum
Default value	empty

Identification 2

Sample identification. The sample identification can be used in calculations as the variable **CI2**. If you have defined a sample identification list (*see chapter 15.1.1, page 163*) then the entries can be selected here.

Entry	24 characters maximum
Default value	empty

Sample size

Sample size. The value of the sample size can be used in calculations as the variable **C00**.

Input range	-99999999 to 9999999999
Default value	1.0

Unit of sample size.

Entry	6 characters maximum
Selection	g mg µg mL µL pieces
Default value	g

Unit defined in method:

The unit defined in the method is ignored. The unit entered in the sample table is used.

Balance connected and configured:

Value from the connected balance.

Fixed sample size defined in method:


The value will be entered in the corresponding line when the method is loaded. The existing entry will be overwritten.



NOTICE

The sample size limits defined in the method are not monitored when the sample data is being entered in the sample table.

21.3 Properties

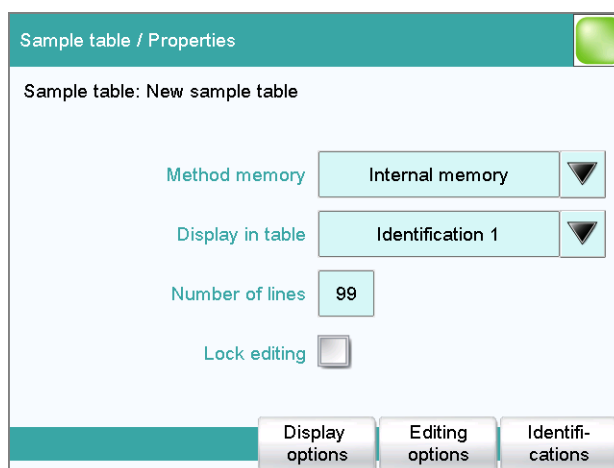

Method memory

Memory location the method is loaded from. All memory locations are selectable, even if they are currently not accessible.

Selection	Internal memory External memory 1 External memory 2 Shared memory
Default value	Internal memory

Shared memory

Shared directory in the network.

Display in table

Selection of the columns to be displayed in the sample table.

Selection	Method Identification 1 Identification 2
Default value	Identification 1

**Method**

Only the assigned method is displayed for each sample.

Identification 1

The first sample identification and the sample size are displayed for each sample.

Identification 2

The second sample identification and the sample size are displayed for each sample.

Number of lines

Maximum number of lines which can be present in the sample table.

Input range	2 to 999
Default value	99

Lock editing

on | off (Default value: **off**)

If this option is activated, then the sample data can no longer be altered.

Display options

Configure the sample data to be displayed in the dialog **Sample table / Edit** (see *"Properties / Display options" dialog*, page 218).

Editing options

Specify the settings for the processing of the sample data (see *"Properties / Editing options" dialog*, page 218).

Identifications

Define the designations for the input fields of Identification 1 and Identification 2 (see *"Dialog Properties / Identifications"*, page 219).

"Properties / Display options" dialog

Sample table: **Properties ► Display options**


In this dialog, you can configure which of the following sample data can be edited in the sample table:

- **Method**
- **Identification 1**
- **Identification 2**
- **Sample size**

"Properties / Editing options" dialog

Sample table: **Properties ► Editing options**

In this dialog, you can define the settings so that the input of the sample data becomes simpler and faster.

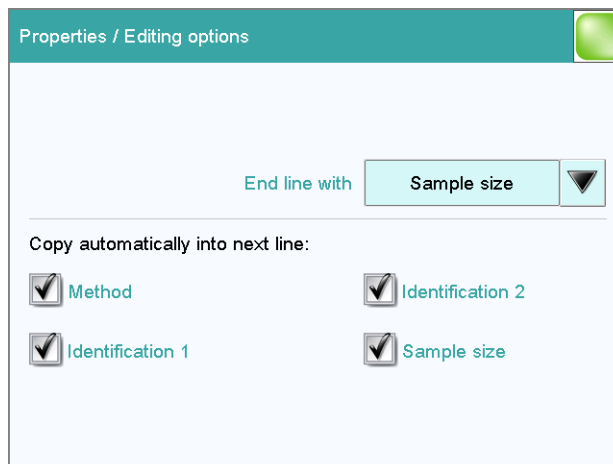

End line with

Selection as to with which input the editing dialog for the next sample is automatically displayed.

Selection	Sample size Identification 1 Identification 2 manual
Default value	Sample size

manual

The editing dialog for the next sample can be displayed with the **[New sample]** button in the editing dialog.

If, for example, you have to enter similar data for each sample, then you can automatically copy the data which is identical for each sample into the next line:

- **Method**
- **Identification 1**
- **Identification 2**
- **Sample size**

Dialog "Properties / Identifications"

Sample table: **Properties ► Identifications**

In the dialog **Properties / Identifications**, you can alter the designations for the sample identifications in accordance with specific methods.

22 Determination run

This chapter provides a description of how to carry out a determination, i.e. how to process a single sample or a sample series. The current sample data is used for calculations.

22.1 Carrying out a single determination

When you carry out a determination, you can enter the sample data in various ways (see chapter 20, page 209). The following instructions describe how to enter the sample data in the main dialog.

To carry out a single determination, proceed as follows:

1 Loading the method

- See Chapter "Loading a method", Page 184.

2 Preparing the sample



NOTICE

Calculate the amount of the sample in such a way that a titrant consumption of 10...90% of the cylinder volume will result.

- Weigh in or measure the sample in a sample vessel.
- Add solvent.
- Place the sample vessel on the stirrer.
- Immerse the electrode and buret tip in the solution.

3 Entering the sample size

- Tap on the input field **Sample size** in the dialog. The number editor opens.
- Enter the sample size and confirm with **[OK]**.

4 Starting a determination

- Tap on the fixed key [▶].

Sequence of the determination


- After a determination is started, a check is made as to
 - whether all of the devices required are connected and ready for operation.
 - whether all titrants and sensors are present.
 - whether the monitoring intervals for titrants, sensors and common variables are still valid.
 - if a sample assignment table is activated (*see chapter 15.1.2, page 164*):
 - whether the entered sample identification is defined in the sample assignment table. If this is the case, then the assigned method will be loaded automatically and the actual determination will be started.



NOTICE

Manual titrations (MAT command):

Start of titration: Because of the technical implementation involved, a few microliters are already dosed at the start of a manual titration. The possibility of deviation from the actual endpoint volume is excluded, however, because the already dosed volume is taken into account.

End of titration: A manual titration must always be ended with the **[Cancel command]** button (NOT with the [] fixed key) or with the **[FILL]** key on the Manual Dosing Controller. This is the only way to ensure that subsequent method commands will be carried out.

22.2 Performing a sample series

You can use the following functions with sample series:

- **Statistics for multiple determinations**
The statistics calculations are enabled in the method options (*see chapter 16.5, page 187*).
- **Sample table**
You can enter the sample data for an entire sample series in the sample table (*see chapter 21, page 212*).


23 Live modifications

23.1 Editing the sample data of the running determination

The sample data can be entered or modified in the main dialog while a determination is running. The sample data entered at the end of the determination is always used in calculations.

Proceed as follows to edit the sample data:

1 Display the main dialog

- Tap on the fixed key [].

The main dialog is displayed. The determination continues to run in the background.

2 Edit the sample data

- Edit the sample data and apply with **[OK]**.

3 Display the live dialog

- Tap on **[Live display]**.

The live dialog is displayed once again.



NOTICE

Make sure that the input dialogs are closed before the determination is ended. Otherwise the determination will have to be recalculated.

23.2 Editing the sample table while a determination is running

You can insert new lines or delete existing ones or edit sample data while a determination is running.



NOTICE

To ensure that no problems arise during the run and that the current data is always available for the calculation, we recommend that the input dialogs always be closed.

Editing the sample table

Proceed as follows to edit the sample table:

1 Display the main dialog

- Tap on the fixed key [].

The main dialog is displayed. The determination continues to run in the background.

2 Select the sample data

- Tap on **Sample table**.
The sample table is displayed. The first line always contains the sample data of the running determination.
- Select the desired line.
- Tap on the button **[Edit]**.

3 Edit the sample data

- Edit the sample data and apply with **[OK]**.



NOTICE

In addition to the sample data, the method can also be modified, except in cases where the determination is running.


4 Display the live dialog

- Tap on the fixed key [].
The main dialog is displayed.
- Tap on **[Live display]**.

The live dialog is displayed once again.

23.3 Live display

Main dialog: **Live display**

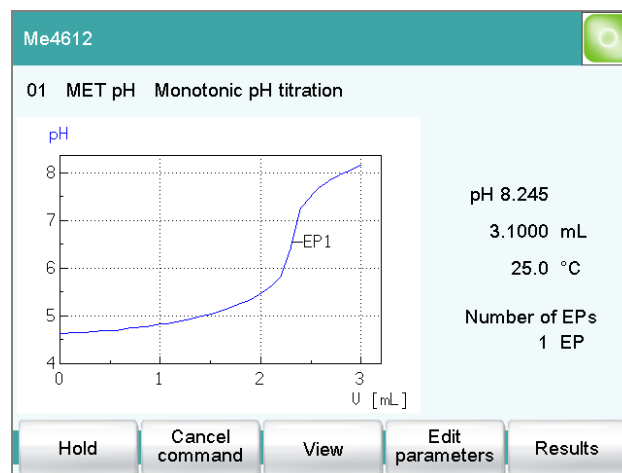


The current method and the system status are displayed in the title line. The currently running command is displayed directly underneath the title line.

The following functions are available to you while a determination is running:

- **[Hold]**
Pause the determination. The label changes to **[Continue]**. This can be used to continue the run.
- **[Cancel command]**
Cancel the current method command. The next command is started.
- **[View]** (only for titrations/measurements)
To modify the settings for the curve and measured value display, see below.
- **[Sample data]** (visible only if a titration or a measurement is not currently running)
Switch to the main dialog in order to edit the sample data.
- **[Edit parameters]**
Edit the method options and live parameters (see chapter 23.4, page 227).
- **[Results]**
Display the results dialog (see chapter 24, page 229).

Dialog "View"

Main dialog: **Live display** ► **View**

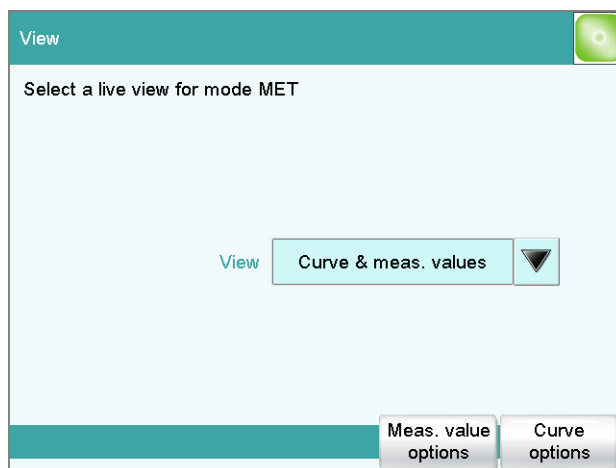

View

Type of curve display.

Selection	Curve Curve & meas. values Measured values Curve & sample data
Default value	Curve & meas. values

Curve

Only the curve is displayed.

Curve & meas. values

The measured values are displayed to the right next to the curve.

Measured values

Only the measured values are displayed.

Curve & sample data

The sample data is displayed to the right next to the curve.

[Meas. value options]

Define settings for the display of the measured values.

A maximum of three different values can be displayed at the same time. The selection depends on the titration/measuring mode. The settings are saved separately for each mode.

[Curve options]

Define the settings for the curve display (*see "Curve options", page 243*). The settings are saved separately for each titration/measuring mode and apply not only for the live display but also for the curve display in the results dialog.

23.4 Live parameters

Live display: **Edit parameters**

Certain method parameters can be edited while a determination is being carried out. The modified parameters are taken into account at once. If you modify, for instance, the start conditions after the start volume has been dosed, then these modifications will not be taken into account until the next determination.



NOTICE

If the live modifications are also to be used in the future, then the method must be saved after the determination has been completed. The live modifications are discarded as soon as you load a new method.

If you are working with the sample table and a different method is used for the next sample, then the live modifications will be discarded.

Proceed as follows to modify method parameters during an ongoing determination:

1 Select the method command

- Tap on **Edit parameters**.
The command sequence is displayed.

Parameters / Sequence		
Current method: Me4612		
01	REQUEST	Data request
02	MET pH	Monotonic pH titration
03	CALC	Calculation
04	REPORT	Report
05	...	

Save method Method options Insert command Delete command Edit command

[Edit command] is enabled only for those commands for which the live parameters can be edited.


All of the functions except for the method properties can be edited with **[Method options]**.

- Select the desired method command.
- Tap on **[Edit command]**.

**2 Edit the method parameters**

- Change the desired parameters accordingly.

3 Display the live dialog


- Tap on the fixed key [].
The main dialog is displayed.
- Tap on **[Live display]**.

The live dialog is displayed once again.

24 Results and more determination data

Main dialog: **Results**

Detailed specifications concerning the current determination are displayed in the **Results** dialog:

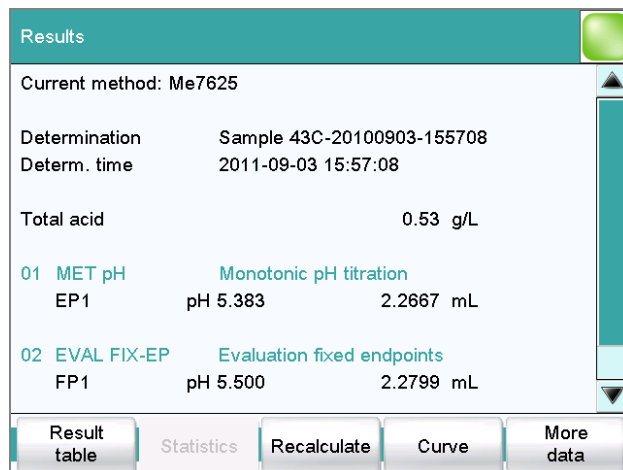

- **Current method**
Method with which the determination was carried out.
- **Determination**
Name of the determination. Displayed only if the determination has been saved.
- **Determination time**
Date and time at which the determination was carried out.
- **Results**
Calculated results. If monitoring of the result limits is activated, then the result will be highlighted in color:
 - Green lettering: The result is within the limits
 - Red lettering: The result is outside the limits
- **Commands**
List of the data-generating commands (titrations, measurements, etc.). The most important data is displayed for each command.

[Result table]

Open result table (*see chapter 26, page 251*). The result table is a table in which the results for several determinations can be saved.

[Statistics]

Display the statistical overview of a determination series (*see chapter 25, page 246*).

[Local com. var.]

(available only with loaded determinations)

Display the list of the common variables defined at the time of the determination (see chapter 24.3, page 234).

[Messages]

Display list with all of the messages which occurred during the determination (see chapter 24.2, page 233).

[View data]

Display all of the details for the selected command.

24.1.1 Details

Results dialog: **More data ► View data**

More determination data / View data		
01 MET pH Monotonic pH titration		1M
Stop criterion	Stop volume reached	MTS
Titration duration	48.9 s	MCD
Start volume	1.0000 mL	MSV
End volume	3.0000 mL	MCV
Temp. measurement	Pt1000	MTM
Initial measured value	3.507	MIM
Initial temperature	27.8 °C	MIT
Start duration	6.5 s	MSD
Start measured value	3.682	MSM
Start temperature	27.8 °C	MST
End measured value	6.812	MCM

All of the variables generated by the determination are listed in the **More determination data / View data** dialog. These can be used in calculations.

[Calibration data]

(available only with loaded determinations)

Display calibration data of the sensor used which was valid at the time of the determination (see "Calibration and titrant data", page 233).

[Titrant data]

(available only with loaded determinations)

Display titrant data which was valid at the time of the determination (see "Calibration and titrant data", page 233).

[Endpoint details]

Display details for each endpoint found (see "Endpoint details", page 232).

[Measuring point list]

Display measuring point list of titrations and measurements (see "Measuring point list", page 232).

Measuring point list

Results dialog: **More data ▶ View data ▶ Measuring point list**

Results dialog: **Curve ▶ Display curve ▶ Measuring point list**

View data / Measuring point list		
01 MET pH Monotonic pH titration		1M
Volume [mL]	Meas.value [pH]	Delta MV [pH]
1.00000	3.682	0.000
1.10000	3.713	0.032
1.20000	3.749	0.036
1.30000	3.790	0.040
1.40000	3.836	0.046
1.50000	3.889	0.053
1.60000	3.950	0.061

View

The individual measured values are listed in the **View data / Measuring point list** dialog. You can configure which columns are displayed (**[View]** button). These settings are saved separately for each mode (DET, MET, etc.).

Column 1 / Column 2 / Column 3

Quantity displayed in the columns 1, 2 and 3. Both the selection of the quantities and the default setting depend on the mode.

Selection	Volume Measured value ERC Delta meas.value Time Temperature Volume drift Signal drift Index
-----------	--

Endpoint details

Results dialog: **More data ▶ View data ▶ Endpoint details**

Results dialog: **Curve ▶ Display curve ▶ Endpoint details**

Detailed information concerning each equivalence point or endpoint determined is displayed in the **View data / Endpoint details** dialog.

The volume, the measured value, the equivalence point criterion ERC (only DET, MET), the time and the temperature are specified for each equivalence point or endpoint determined.

The equivalence point criterion determined (ERC = Equivalence point Recognition Criterion) is a measure for the size of the jump in the titration

curve. The ERC determined is compared with the EP criterion defined in the method for the evaluation of the titration curve *see "Evaluation and equivalence point criterion with DET", page 303 and "Evaluation and equivalence point criterion with MET", page 319*). The ERC is thus an important quantity when you need to adjust the parameters for the evaluation of the equivalence points.

Calibration and titrant data

Results dialog: **More data ▶ View data ▶ Calibration data**

Results dialog: **More data ▶ View data ▶ Titrant data**

The calibration data or titrant data which was current at the time the determination was performed is displayed in the dialogs **View data / Calibration data** and **View data / Titrant data**.



NOTICE

Recalculation:

The concentration and titer of the titrant can be modified retroactively. The values displayed here are used for the recalculation of a loaded determination. The values saved in the dialog **System / Titrants** are not changed as a result.

24.2 Messages

Results dialog: **More data ▶ Messages**

If messages occurred during a determination, they will be displayed in the **More determination data / Messages** dialog. The following data for the message is indicated under **[Details]**:

- **Message number**
Unique identification number.
- **Time**
Moment that the message occurred.
- **Message**
Message text.

- **Method status**
Current status of the method.
 - **new**
The method has been newly created and not yet saved.
 - **saved**
The current method has been saved.
 - **modified**
The current method has been modified but not saved yet (e.g. live modification).
 - **reviewed**
The current method has been signed at level 1.
 - **released**
The current method has been signed at level 2.
- **Determination**
Name of the determination. Displayed only with saved determinations.
- **End of determin.**
Way in which the determination was ended.
 - **Regular without errors**
The determination was automatically ended at the end of the method.
 - **Manual stop**
The determination has been canceled with the fixed key [].
 - **Error**
The determination has been canceled due to an error.
- **Status of deter.**
 - **original**
The determination has not been recalculated.
 - **recalculated**
The determination has been recalculated.
 - **loaded**
The determination has been loaded.
 - **saved**
The determination has been saved. The version number indicates how many times the file has been saved.
 - **reviewed**
The determination has been signed at level 1.
 - **released**
The determination has been signed at level 2.

The following data is displayed only if the option **Recalculating determination** under **Login options / Modification options** is activated:

- **Recalculated by**
User who recalculated the determination.
- **Recalculated on**
Date and time at which the determination was recalculated.



- **Reason**
Reason for the recalculation of the determination.
- **Comment**
Comment for the recalculation of the determination.

The following information is displayed only if the determination has been signed at level 1:

- **Reviewed by**
User who signed the determination at level 1.
- **Reviewed on**
Date and time at which the determination was signed at level 1.
- **Reason**
Reason for the signing of the determination.
- **Comment**
Comment for the signing of the determination.

The following information is displayed only if the determination has been signed at level 2:

- **Released by**
User who signed the determination at level 2.
- **Released on**
Date and time at which the determination was signed at level 2.
- **Reason**
Reason for the signing of the determination.
- **Comment**
Comment for the signing of the determination.
- **Used devices**
Hardware that was used for the determination.
 - Ti-Touch with program version of the software and the serial number.
 - Measuring inputs with ADC type und serial number.
 - MSB connector with peripheral devices connected to it (dosing device, stirrer).

[Delete signature]

Delete all of the signatures for the method or the determination. This button is inactive if level 2 has not yet been signed (*see chapter 17.2, page 200*).

[Sign]

Sign the method or determination. This button is active only when working with activated login function and password (*see chapter 17.1, page 199*).


24.5 Loading a determination

Proceed as follows to load a determination:

1 Opening the determination list

- In the results dialog, tap on **[More data]**.
- Tap on **[Load/ Save]**.
- Tap on **[Load]**.
The selection of memory locations is displayed. This selection will be skipped if only one external storage medium is available.
- Select the memory location where the desired determination is saved.
The selection of file groups is displayed (*see chapter 12.1, page 136*). If only one group is available, then the determination list will be displayed directly.
- Select the group with the desired determination.
- Tap on **[Show files]**.

The determination list with the stored determination opens:

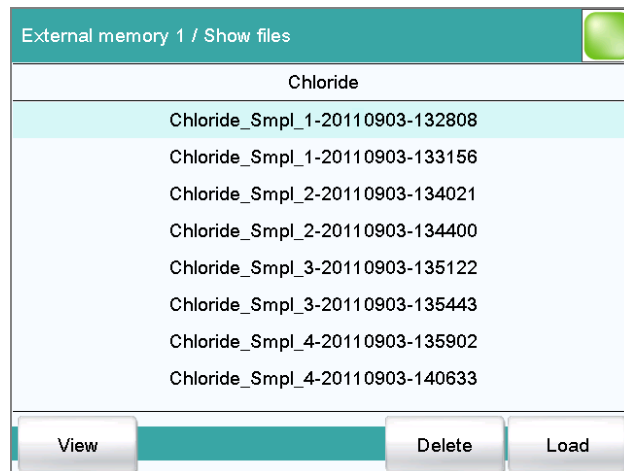

2 Selecting a determination

- Select the desired determination.

3 Loading the determination

- Tap on **[Load]**.

The determination is now loaded. The properties are displayed in the results dialog.

**NOTICE**

When you load a determination, the method with which the determination was carried out will be loaded automatically.

For this reason, save any modifications which may have been made to the method currently loaded before you load a determination.

24.5.1 Determination list

Results dialog: **More data ► Load/ Save ► Load ► Show files ► View**

Show files / View		
Sorted by Determ. time		10
No.	Identification 1	Result 1
1	Blank methanol	0.01 %
2	Blank methanol	0.02 %
3	Calcium hardness	2.015 mmol/L
4	Calcium hardness	1.981 mmol/L
5	Titer of NaOH	0.99615
6	Titer of NaOH	0.99417
7	Titer of NaOH	0.98729

In the **Show files / View** dialog you can depict the individual determinations in greater detail. The sorting criterion and the number of saved determinations are displayed at the very top. The list contains numbered lines; each line represents one determination. In the default settings, the first sample identification and the first calculated result of each determination are displayed. In the properties, you define which data is displayed in the columns and the criteria according to which the determinations are sorted (see "Properties", page 239). In addition, you can define search filters to aid you in finding the desired determinations (see "Filter criteria", page 240).

[Previous 100]

Display the previous 100 determinations.

[Next 100]

Display the next 100 determinations.

[Properties]

Configure the display of the columns and define filter criteria.

[Delete]


Delete the selected file.

[Load]

Load the selected file.

Properties

Determination list: **Properties**

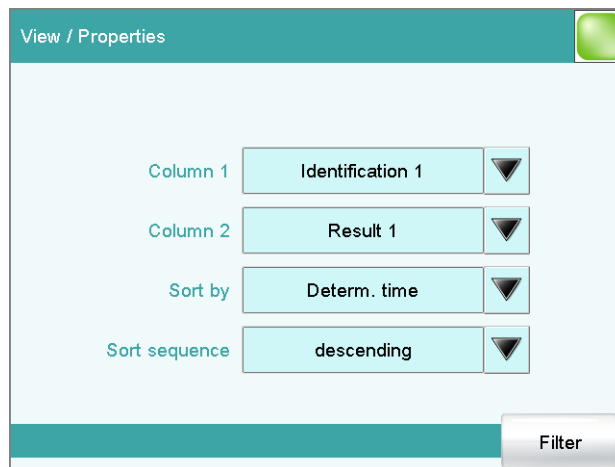

In the **View / Properties** dialog, you can define which data is displayed in the columns and the criteria according to which the determinations are to be sorted.

Column 1

Selection of what is to be displayed in the first column.

Selection	Identification 1 Identification 2 Determ. time Result name Method Sample size User
Default value	Identification 1

Column 2

Selection of what is to be displayed in the second column.

Selection	Result 1...9 The numbering of the results corresponds to the order in which the results were calculated in the determination run.
-----------	---

Sort by

Selection according to which criterion the list is to be sorted.



Selection	Identification 1 Identification 2 Determ. time Result Method Sample size User
Default value	Determ. time

Identification 1

Sort alphabetically according to the first sample identification.

Identification 2

Sort alphabetically according to the second sample identification.

Determ. time

Sort chronologically according to the determination time.

Result

Sort according to the numerical value of the result.

Method

Sort alphabetically according to the method name.

Sample size

Sort according to the numerical value of the sample size.

User

Sort alphabetically according to the user who carried out the determination.

Sort sequence


Selection	ascending descending
Default value	descending

[Filter]

Specify the filter criteria for the list.

Filter criteria

Determination list: **Properties ▶ Filter**

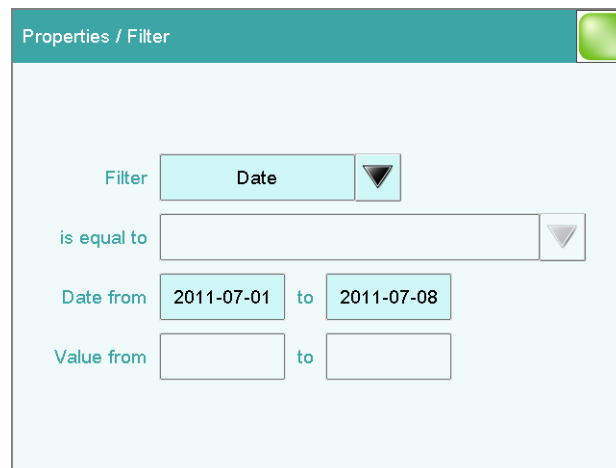

You can set filters in the **Properties / Filter** dialog. A filter is a rule defining what is to be displayed in the determination list. If a filter has been set,

this is indicated with a corresponding symbol in the **Show files / View** dialog.

Filter

Define the criterion according to which the list is to be filtered. The only lines which will be displayed are those which fulfil the filter criterion.

Selection	no filter Identification 1 Identification 2 Date Result Method Sample size User
Default value	no filter

no filter

The list itself will not be filtered.

Identification 1

In the input field **is equal to**, select or enter the first sample identification according to which filtering is to be carried out.

Identification 2

In the input field **is equal to**, select or enter the second sample identification according to which filtering is to be carried out.

Date

In the input fields **Date from ... to**, specify the period of time according to which filtering is to be carried out.

Result

In the input fields **Value from ... to**, specify the value range according to which filtering is to be carried out.

Method

In the field **is equal to**, select or enter the method name according to which filtering is to be carried out.

Sample size

In the input fields **Value from ... to**, specify the value range according to which filtering is to be carried out.

User

In the field **is equal to**, select or enter the user according to whom filtering is to be carried out.



NOTICE

The filter can be applied to empty fields as well, for example, if no sample identification has been entered. In this case, select only the desired filter criterion and leave the fields below empty.

24.6 Saving a determination

Results dialog: **More data ▶ Load/ Save ▶ Save**

Determinations can be saved as follows:


- Manually in this dialog.
The suggested default setting for file names is the first 16 characters of Identification 1, followed by the determination time in the format YYYYMMDD-hhmmss. Extending the file name with the determination time has the advantage that the file names of all determinations are unique.
- Automatically at the end of the determination (*see chapter 16.5.6, page 197*).

The following data is stored for each determination:

- Calculated results
- Measuring point lists, equivalence points, curves and variables which were generated during the determination
- Determination properties
- Statistics (if carried out)
- Method with which the determination was carried out
- Common variables (values and properties of all common variables at the time of the determination)
- Titrant data of the titrant used
- Calibration data of the sensor used

24.7 Curves

Results dialog: **Curve**

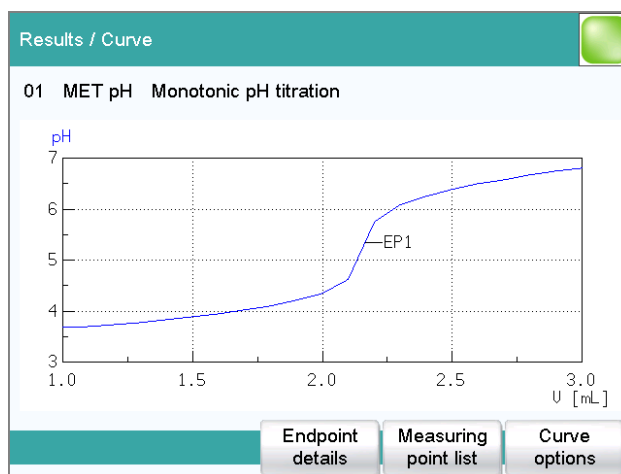

The curve of the current determination is displayed in the **Results / Curve** dialog. You can configure the curve display.

[Endpoint details]

Display details for each endpoint found (see "Endpoint details", page 232).

[Measuring point list]


Display measuring point list of titrations and measurements (see "Measuring point list", page 232).

[Curve options]

Define the settings for the curve display (see "Curve options", page 243). The settings are saved separately for each titration/measuring mode and apply not only for the live display but also for the curve display in the results dialog.

Curve options

Results dialog: **Curve ▶ Curve options**

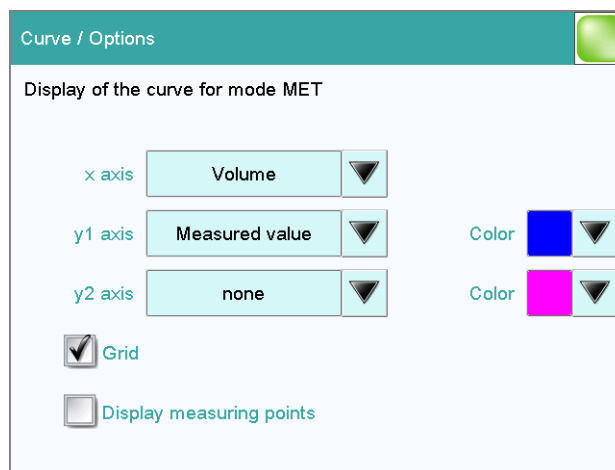

You can make the settings for the curve display in the **Curve / Options** dialog. These settings are saved separately for each mode (DET, MET, etc.) and also apply for the live display. Both the selection of the quantities and the default setting depend on the mode.

x axis

Quantity which is shown on the x axis of the graph.

Selection	Volume Measured value ERC Delta meas.value Time Temperature Volume drift Signal drift
-----------	--

y1 axis

Quantity which is shown on the y1 axis of the graph.

Selection	Volume Measured value ERC Delta meas.value Time Temperature Volume drift Signal drift
-----------	--

y2 axis

Quantity which is shown on the y2 axis of the graph.

Selection	none Volume Measured value ERC Delta meas.value Time Temperature Volume drift Signal drift
Default value	none

Color

Selection of a color for the curves.

Grid

on | off (Default value: **on**)

If this parameter is activated, grid lines are shown.

Display measuring points

on | off (Default value: **off**)

If this parameter is activated, the curve is shown with the single measured values.

24.8 Recalculation and reevaluation

Results dialog: **Recalculate**

You can use the **[Recalculate]** function to recalculate and reevaluate determinations. This is necessary in the following cases:

- If you have modified the evaluation parameters in a way that changes the recognized endpoints or equivalence points.
- If you have modified the calculations.
- If you have modified the variables used in calculations, e.g. sample size, titer or common variables.



NOTICE

Neither the recalculation nor the reevaluation can be undone.

We therefore recommend that the original determination be saved beforehand (*see chapter 24.6, page 242*). The best way of doing this is to use the **Save determination automatically** function (*see chapter 16.5.6, page 197*).

The following method commands are executed once again with the **[Recalculate]** function:

- DET and MET titrations:
The potentiometric evaluation will be carried out again.
- EVAL commands:
All evaluations are carried out again. You can also insert new evaluations retroactively.
- CALC command:
All calculations are carried out again. You can also define new calculations retroactively.
If you subsequently modify the titrant in a method, then the variables TITER and CONC of the "new" titrant will be used. If a result is assigned to the variable TITER in a calculation, then the titer of the "new" titrant will also be overwritten.



NOTICE

If you modify the calibration data of the sensor used in the **System / Sensors** dialog, these modifications will not be taken into account. Changing the sensor in the method also has no influence on the measured data.

If a determination has been recalculated, this will be documented in the determination properties (see chapter 24.4, page 234).

Recalculation/reevaluation of loaded determinations

With loaded determinations, it is assumed that they were carried out a long time ago and that the **titrant data** and **common variables** of the system no longer correspond to those of the determination. This is why the most important titrant data of the titrants used and the list of **common variables** are stored with each determination. This data will be used if you recalculate a loaded determination. If you wish to recalculate the result with a corrected **common variable** or with a corrected titer, you must therefore modify the respective values in the **More determination data / Local common variables** or **More determination data / View data** dialog.

If a result is assigned to the variable TITER or to a **common variable**, then the titer or **common variable** which is currently valid in the system will not be overwritten until after a query has been confirmed.

If you modify the titrant retroactively in the method, this will not be taken into account.

The recalculated results of loaded determinations are not saved in the result table.

25 Statistics

Main dialog: **Results ► Statistics**


In the dialog **Results**, you can use **[Statistics]** to display the statistical overview of a determination overview. You can statistically evaluate a maximum of nine results calculated in a determination. A statistical series can contain a maximum of 20 determinations.

Activating the statistics function

Proceed as follows to activate the statistics function:


1 Editing the method options

- In the main dialog, tap on **[Edit parameters]**.
- Tap on **[Method options]**.

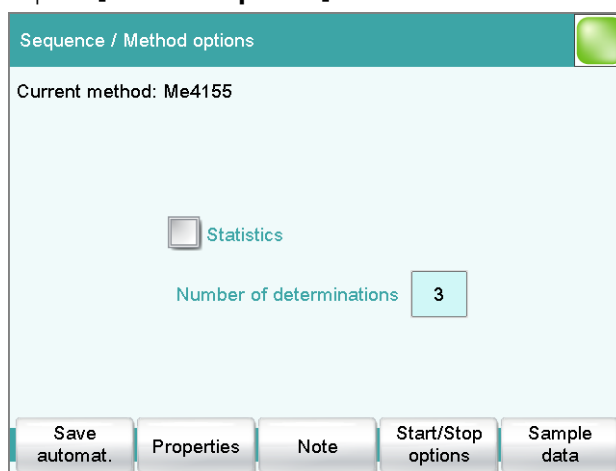

- Activate the **Statistics** option.
- Under **Number of determinations**, enter the desired number of determinations.

The results for which statistical calculations are performed are defined in the calculation command **CALC** (see "Variable for mean value", page 168).

Results / Statistics 	
Current method: Me3901 3 of 3	
Result name	Mean value
Content (mmol/L)	100.31 mmol/L
Volume of HCl	5.0 mL
<div style="display: flex; justify-content: space-between;"> Add determ. Details </div>	

In this dialog, the method with which the determination was carried out is displayed in the first line. The second line shows how many determinations have already been carried out and the total number of determinations to be carried out. The name of the result and the mean value are shown for each result.

[Add determ.]

Add a further determination to the determination series (*see chapter 25.3, page 249*).

[Details]

Display the statistics results for the selected result and the single results from all determinations, see next chapter.



NOTICE

If you recalculate a determination, the statistics for this determination are also recalculated.

25.1 Displaying details for a result

Results dialog: **Statistics ▶ Details**

Statistics / Details		
Result name: Content (mmol/L)		SMN1
Mean value	100.31 mmol/L	n=03
s abs	0.236 mmol/L	
s rel	0.24 %	
No.	Sample size	Result
1	5.0 mL	100.56 mmol/L
2	4.5 mL	100.09 mmol/L
3	5.5 mL	100.28 mmol/L
Sample data		Determ. on/off
		Result on/off

In the dialog **Statistics / Details**, the following details are displayed:

- Result name
- Mean value
On the right-hand side of the dialog the number of single results from which the mean value was calculated is displayed. In this example, it is 3.
- Absolute standard deviation (**s abs**)
- Relative standard deviation (**s rel**)
- Result and sample size of each single determination
Results that could not be calculated are indicated with **invalid** and ignored.

[Sample data]

Display the sample data of the selected determination.

[Determ. on/off]

Remove all of the results of the selected determination from the statistics. All of the entries of this line are marked with an asterisk (*), the statistics will be recalculated automatically.

[Result on/off]

Remove the selected result from the statistics. The result is marked with an asterisk (*), the statistics will be recalculated automatically.

**NOTICE**

If the mean value has been assigned to a common variable or to the variable TITER (see command **CALC**), then the determination must be recalculated manually (see chapter 24.8, page 244).

25.2 Deleting statistical data

In the dialog **Control**, you can delete all of the statistics data manually with the function **[Delete statistics]**. The statistics data is deleted automatically in the following cases:

- when all of the determinations of the determination series have been carried out and a new determination has been started afterwards.
- when a new method is loaded (even if it is the same method)
If the **sample table** or the **sample assignment table** is used, the statistics data is only deleted when another method is loaded.

25.3 Adding a determination to a determination series

You can use the function **[Add determ.]** to add an additional sample to a determination series, e.g. because a determination was faulty and had to be removed from the statistics.

Results / Statistics	
Current method: Me3901 3 of 3 (+2)	
Result name	Mean value
Content (mmol/L)	100.31 mmol/L
Volume of HCl	5.0 mL
<div style="display: flex; justify-content: space-between; align-items: center;"> Add determ. Details </div>	

The number of determinations by which the set statistics counter has been increased is shown above the list in brackets. This modification is also displayed in the main dialog and in the statistics report.



NOTICE

The number of samples defined in the method options does not, however, change as a result.

26 Result table

Main dialog: **Results ► Result table**

Results / Result table		
Sorted by Determ. time		6 of 100
No.	Identification 1	Result 1
1	Content of Ca #32	0.101 mol/L
2	Content of Ca #32	0.092 mol/L
3	Content of Ca #32	0.098 mol/L
4	Titer NaOH	0.9946
5	Titer NaOH	0.9942
6	Titer NaOH	0.9967

Buttons: Load/ Save, Properties, Delete all, Delete, Details

The result table is suitable for displaying in detail the results of all of the determinations performed on a single day, for instance. You can save the results of up to 200 determinations. A maximum of 9 results from a single determination can be saved in the result table. You define which results from this table are to be saved using the **CALC** command. Because the size of the display is not sufficient to display all of the data at the same time, it is practical to print out the result table.

The sorting criterion and the number of determinations from which results are saved in the result table are displayed at the top. The list contains numbered lines; each line represents one determination. In the default settings, the first sample identification and the first calculated result of each determination are displayed. In the properties, you define which data is displayed in the columns and the criteria according to which the determinations are sorted (*see chapter 26.1, page 252*). In addition, you can define search filters to aid you in finding the desired determinations (*see "Filter criteria", page 254*).

[Load/ Save]

Load a saved result table (*see chapter 26.3, page 256*) or save the current result table (*see chapter 26.2, page 256*).

[Properties]

Configure the display of the columns and define filter criteria.

[Delete all]

Delete the result table completely.



[Delete]

Delete the selected line from the result table.

[Details]

Display details concerning the selected determination.


"Result table / Details" dialog

The following determination data is displayed in this dialog:

- **User**
User who carried out the determination.
- **Method**
Method with which the determination was carried out.
- **Determ. time**
Date and time at which the determination was carried out.
- Sample data (Identification 1 and 2 and sample size).
- All of the results of this determination saved in the result table.

26.1 Properties

Result table: **Properties**

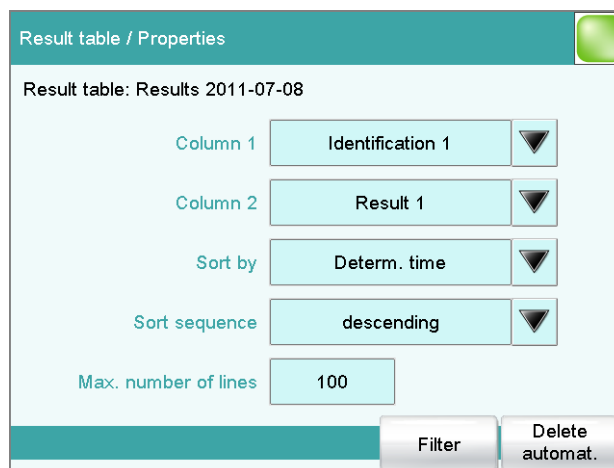

In the **Result table / Properties** dialog, you can select the data to be displayed with each determination as well as the sorting criterion for the list.

Column 1

Selection of what is to be displayed in the first column.

Selection	Identification 1 Identification 2 Determ. time Result name Method Sample size User
Default value	Identification 1

Column 2

Selection of what is to be displayed in the second column.

Selection	Result 1...9 The numbering of the results corresponds to the order in which the results were calculated in the determination run.
-----------	---

Sort by

Selection according to which criterion the list is to be sorted.

Selection	Identification 1 Identification 2 Determ. time Result Method Sample size User Determ. time
Default value	Determ. time

Identification 1

Sort alphabetically according to the first sample identification.

Identification 2

Sort alphabetically according to the second sample identification.

Determ. time

Sort chronologically according to the determination time.

Result

Sort according to the numerical value of the result.

Method

Sort alphabetically according to the method name.

Sample size

Sort according to the numerical value of the sample size.

User

Sort alphabetically according to the user who carried out the determination.

Sort sequence

Selection	ascending descending
Default value	descending

Max. number of lines

If the result table contains the quantity of lines defined here, then no further results can be saved. If this is the case, save the result table and create a new one.

Input range	10 to 200
Default value	100

[Filter]

Specify the filter criteria for the list.

[Delete automat.]

Define the settings for the automatic deletion of the result table (see *"Deleting the result table automatically"*, page 255).

Filter criteria

Result table: **Properties ► Filter**

You can set filters in the **Properties / Filter** dialog. A filter is a rule defining what is to be displayed in the result table. If a filter has been set, this is indicated with a corresponding symbol in the **Results / Result table** dialog.

Filter

Define the criterion according to which the list is to be filtered. The only lines which will be displayed are those which fulfil the filter criterion.

Selection	no filter Identification 1 Identification 2 Date Result Method Sample size User
Default value	no filter

no filter

The list itself will not be filtered.

Identification 1

In the input field **is equal to**, select or enter the first sample identification according to which filtering is to be carried out.

Identification 2

In the input field **is equal to**, select or enter the second sample identification according to which filtering is to be carried out.

Date

In the input fields **Date from ... to**, specify the period of time according to which filtering is to be carried out.

Result

In the input fields **Value from ... to**, specify the value range according to which filtering is to be carried out.

Method

In the field **is equal to**, select or enter the method name according to which filtering is to be carried out.

Sample size

In the input fields **Value from ... to**, specify the value range according to which filtering is to be carried out.

User

In the field **is equal to**, select or enter the user according to whom filtering is to be carried out.

**NOTICE**

The filter can be applied to empty fields as well, for example, if no sample identification has been entered. In this case, select only the desired filter criterion and leave the fields below empty.

Deleting the result table automatically

Result table: **Properties ► Delete automat.**

In the **Properties / Delete automatically** dialog, you define when the contents of the result table is to be automatically deleted.

Delete oldest line only

on | off (Default value: **off**)

If this option is activated at the same time that the result table is full, the oldest entry will be deleted as soon as the results of a new determination are saved.

Time when result table is deleted:**Switch on**

on | off (Default value: **off**)

If this option is activated, then the result table will be deleted each time you switch on the Ti-Touch.

Start a new sample series

on | off (Default value: **off**)

If this option is activated, then the result table will be deleted when you start a new sample series (with activated autostart function!).

Save result table

on | off (Default value: **on**)

If this option is activated, the result table will be deleted as soon as you have saved it.

26.2 Saving the result table



NOTICE

When you save the result table, the complete determination will not be saved with each entry. If you wish to save the determination data for each determination, use the **Save determination automatically** function (see chapter 16.5.6, page 197).

26.3 Loading the result table

Proceed as follows to load a result table:

1 Opening the current result table

- In the results dialog, tap on **[Result table]**.
- Tap on **[Load/ Save]**.
- Tap on **[Load]**.

The selection of memory locations is displayed. This selection will be skipped if only one external storage medium is available.

- Select the memory location where the desired result table is saved.

The selection of file groups is displayed (see chapter 12.1, page 136). If only one group is available, then the list of the saved result tables will be displayed directly.

- Select the group with the desired result table.
- Tap on **[Show files]**.

The list with the saved result tables opens.

2 Selecting the result table

- Select the desired result table.


3 Loading the result table

- Tap on **[Load]**.


The result table is now loaded and will be displayed immediately.

27 Printing

Reports can be printed out as follows:

- Manually with the [] fixed key in this dialog. The fixed key is disabled while a determination is running.
- Automatically in the determination run with the command **REPORT** (see chapter 29.11, page 403).

How to configure a printer is described in *Chapter 11.7, page 124*.

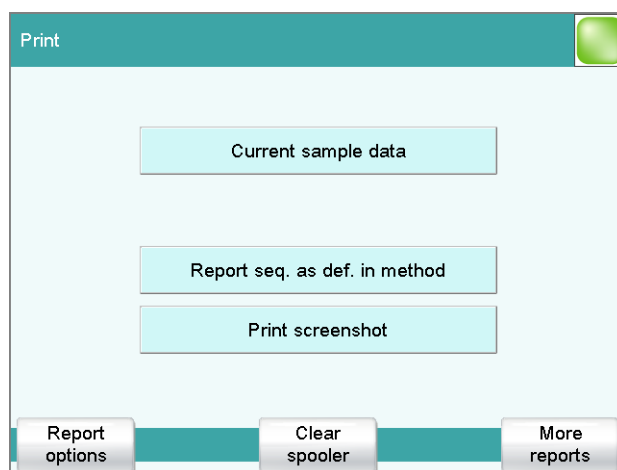


This screenshot shows the printing dialog associated with the main dialog. The sample data is entered in the main dialog, which is why you can print out the sample data from there in a context-sensitive manner with [**Current sample data**].

The following buttons are also available in the other print dialogs:

[Report seq. as def. in method]

Print out all of the reports which are defined in the loaded method.

[Print screenshot]

Activate the "Print Screen" function. The current screen contents will be printed out the next time the fixed key [] is tapped on.

[Report options]

Defining the general settings for the report output.


[Clear spooler]

Delete all outstanding printing orders.


[More reports]

Open the list of available reports.

More reports

Fixed key []: **More reports**

The dialog **Print / More reports** displays a list of all of the available reports.

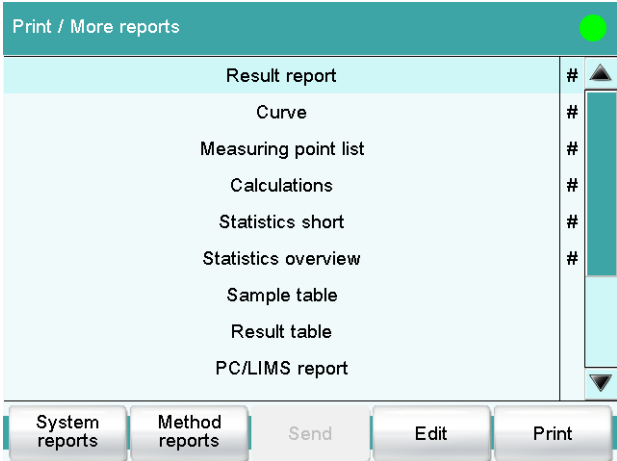

NOTICE

If you are using a **CUSTOM Neos** printer, you can print only the reports marked with a #. The designation on the reports is only visible once the **CUSTOM NEOS** printer is selected (see screenshot above).

[System reports]

Opening the list of reports for the system settings.



NOTICE

If you are using a **CUSTOM NEOS** printer, you cannot print system reports.

[Method reports]

Opening the list of reports for the current method.



NOTICE

If you are using a **CUSTOM NEOS** printer, you cannot print method reports.

[Send]

Create a PC/LIMS report.

[Edit]

Edit the settings of the selected report.

[Print]

Print the selected report.

27.1 General report options

Fixed key []: **Report options**

The settings for the report printout can be adjusted in the dialog **Print / Report options** to indicate, for example, whether a report header, a signature line or a frame is to be printed.

Report header

Defining the output of the report header. The report header contains general information such as device type, serial number, program version and the printing date.



NOTICE

You can also define your own report header, which is printed in addition to this report header (**System ► Templates ► Report header**).

Selection	off once on each page
Default value	on each page

off

The report header will not be printed.

once

The report header will only be printed on the first page.

on each page

The report header will be printed on every page.

Signature line

Output of a special line for date and signature. This line will be printed at the very bottom of every page.

Selection	off once on each page
Default value	off

off

No signature line will be printed.

**once**

The signature line will only be printed on the last page.

on each page

The signature line will be printed on every page.

Frame

on | off (Default value: **on**)

If this parameter is activated, a frame is printed as lateral marking.

27.2 Settings of the individual reports

Fixed key []: **More reports ► Report ► Edit**

Command REPORT: **Report ► Edit**

Settings can be edited for the following reports (for details, see online help):

- **Result report**
- **Curve**
Definition of the curve size, display of the individual measuring points and grid lines, etc.
- **Measuring point list**
Definition of the method command to which the measuring point list is to be printed.
- **Calculations**
Definition of the accuracy with which the variables used as well as the results calculated are printed.
- **Statistics short**
Definition of the time at which the report is to be printed (with each determination or only at the end of a sample series).
- **Statistics overview**
Definition of the time at which the report is to be printed (with each determination or only at the end of a sample series).
- **Sample table**
- **Result table**
Definition of the time at which the report is to be printed (with each determination, at the end of a sample series or at the end of the sample table).
- **Used devices**
- **Variables**
- **Monitoring**

27.3 List of all printable reports

The following reports can be printed out with the fixed key :



NOTICE

If you have selected a **CUSTOM NEOS** printer, you can print only the reports that are marked with a # on the display.

Report	Contents
Result report	# Report with determination properties, sample data, calculated results, etc. If there are multiple determinations, the statistics will also be printed out.
Curve	# Curve report. The settings for the curve print-out are adopted from the curve options of the live display.
Measuring point list	# Measuring point list report.
Calculations	# Details concerning the calculations carried out (parameter settings in accordance with the CALC command and the calculated results).
Used devices	The devices used for the determination, as displayed in the More determination data / Properties dialog.
Variables	All determination variables, as displayed in the results dialog.
Monitoring	Details for the monitored quantities (STAT only).
Statistics	
Statistics short	# Summary of the statistics calculations. The number of determinations, the mean value, and the absolute and relative standard deviations are printed out for each result.
Statistics overview	# Detailed statistical overview. The sample data and all individual results are printed out for each determination. The number of determinations, the mean value, and the absolute and relative standard deviations are printed out for each result.
PC/LIMS	




Report	Contents
PC/LIMS report	Machine-readable report with all of the important data for a determination. This report can be saved as a TXT or UTF-8 file or sent to a terminal program or a LIMS via an RS-232 interface. The definition is carried out in the device manager (<i>see chapter 11.3.2, page 99</i>).
Sample data	
Current sample data	Sample data of the current determination, <i>context-sensitive only</i> from the main dialog.
Sample table	List of all determinations in the sample table with the respective sample data, as entered in the sample table.
Result table	
Result table	List of all determinations in the result table with results and with the determination data, as saved in the result table.
More reports ► Method reports	
Method sequence	Method properties and the list of all method commands.
Parameters full	Method properties and options, all method commands with all parameters. All parameters which no longer have their default settings will be printed in bold . All parameters which have been modified in comparison to the stored version of the method are indicated by an *.
Titration & measuring param.	Method properties; titration, measurement and calibration commands with all parameters. All parameters which no longer have their default settings will be printed in bold . All parameters which have been modified in comparison to the stored version of the method are indicated by an *.
Modified parameters	Method properties, all method parameters which have been modified in comparison to the stored version of the method.
Non default parameters	Method properties, all method parameters which no longer have any default settings.
More reports ► System reports	
System settings	
System settings	Settings for acoustic signals and accuracy of the measured value display.

Report	Contents
Dialog options	Settings for routine dialog and expert dialog.
User list	List with all of the users configured in the system, together with their data, <i>context-sensitive only</i> from the user list.
Login options	Login options, <i>context-sensitive only</i> from the respective dialog.
Titrants	
Titrant list	List of all the titrants configured in the system.
All titrant data short	The most important titrant data of all titrants (name, concentration, titer, last titer determination).
All titrant data full	All of the titrant data of all of the titrants (working life, titer options, data concerning the exchange unit/dosing unit).
Titrant data short	The most important titrant data of the selected titrant, <i>context-sensitive only</i> from the editing dialog.
Titrant data full	All titrant data of the selected titrant, <i>context-sensitive only</i> from the editing dialog.
Sensors	
Sensor list	List of all the sensors configured in the system.
All sensor data short	The most important sensor data for all sensors (name, calibration data).
All sensor data full	All of the sensor data for all of the sensors (working life, complete calibration data, calibration interval).
Sensor data short	The most important sensor data of the selected sensor, <i>context-sensitive only</i> from the editing dialog.
Sensor data full	All of the sensor data of the selected sensor, <i>context-sensitive only</i> from the editing dialog.
Device manager	
Device list	List of all devices configured in the system.
All device properties	Properties of all the devices configured in the system.
Device properties	Properties of the selected device, <i>context-sensitive only</i> from the editing dialog.
GLP manager	
GLP data	All data stored in the GLP manager.
Common variables	

Report	Contents
Common variable list	List of all the common variables defined in the system, together with their most important data (name, value, status).
All common variable properties	Properties of all common variables (name, value, validity, status).
Common variable properties	Properties of the selected common variable (name, value, validity, status), <i>context-sensitive only</i> from the editing dialog.
Local Common variable list	List with the common variables available at the time of the determination, together with their most important data (name, value, status), <i>context-sensitive only</i> from the dialog "More determination data / Local common variables".
Local common var. properties	Properties of the selected common variable (name, value, validity, status), <i>context-sensitive only</i> from the dialog "More determination data / Local common variables".
Templates	
Templates sample data	Sample identification list and sample assignment table.
Result template list	List with all of the user-generated result templates.
All result templates details	Details of the all of the user-generated result templates (calculation formula, result options, note).
Result template details	Details of the selected result template, <i>context-sensitive only</i> from the editing dialog.
Input/Output lines	List with all of the defined input and output lines at the remote interface (name, bit pattern).
Custom calibration buffers	Temperature tables for all defined custom calibration buffers.
Rack tables	
Sample rack list	List with all of the sample racks configured in the system, together with designation, number of positions and rack code.

28 Manual control

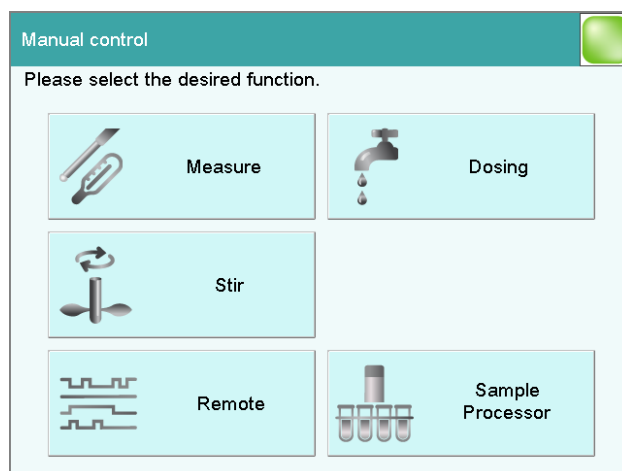

The following functions are available in the manual control:

[Measure]

Carrying out manual measurements (*see chapter 28.2, page 267*).

Measuring modes:

- **pH** (potentiometric pH measurement)
- **U** (potentiometric voltage measurement)
- **Ipol** (voltametric measurement with selectable polarization current)
- **Upol** (amperometric measurement with selectable polarization voltage)
- **T** (temperature measurement)

[Dosing]



NOTICE

Only active if a Dosino or dosing unit is connected.

Manual dosing (*see chapter 28.3, page 269*).

The following dosing functions are available:

- Preparing the exchange unit or dosing unit
- Emptying the dosing unit
- Filling the dosing cylinder of the exchange unit/dosing unit
- Dosing a specified volume
- Dosing continuously

[Stir]

Controlling the stirrer manually (*see chapter 28.4, page 278*).

[Remote]

Scanning the input lines on the remote interface and setting the output lines (see chapter 28.5, page 280).


[Sample Processor]

Controlling the Sample Processor manually (see chapter 28.6, page 281).


28.1 Opening and closing the manual control

Opening the manual control

1 Opening the dialog window

- Tap on the fixed key [].

Manual control opens:

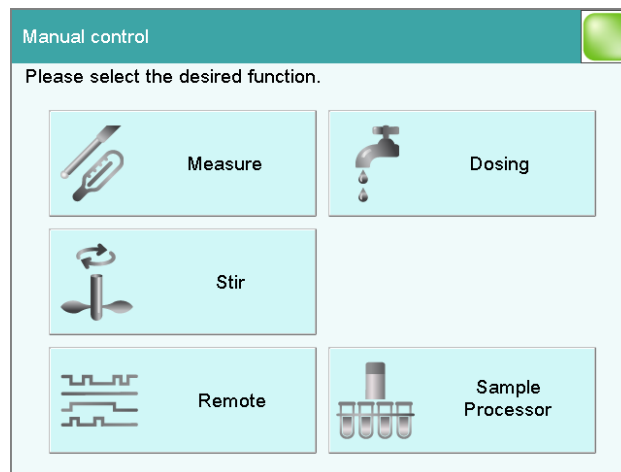

2 Select the function

- Tap on the button for the desired function.

If the function can be carried out by several devices, a dialog to select the device will appear first; otherwise the dialog for the selected function will appear immediately.

Closing the manual control

- #### 1
- Tap on the fixed key [].



NOTICE

Manual control can also be exited when a manual sequence has been started but has not yet ended. The fact that a sequence has been started in manual control can be recognized by the corresponding symbol in the title line (see chapter 6.2.3, page 38).

28.2 Measuring

Manual control ► Measure

Manual measurements can be carried out with the function **[Measure]**.

Proceed as follows:

1 Selecting a sensor

- Select the desired sensor from the sensor list.
All of the sensors available in the sensor list are displayed. Sensors are defined under **System ► Sensors**.
- Select the measuring input at which the sensor is connected.

2 Select a measuring mode

- Select the desired measuring mode.
Only those measuring modes are displayed which are advisable for the selected sensor.


3 Enter a measuring temperature

- Enter the measuring temperature if no temperature sensor is connected. If a temperature sensor is connected, then the temperature will be measured automatically. This temperature is used for automatic temperature compensation with pH measurements.

4 Starting the measurement

- Tap on **[Start]**.

The measured value display is opened:

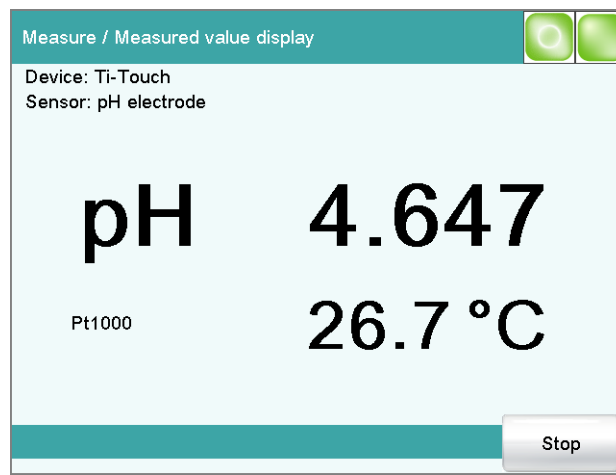

The ongoing measurement is canceled with **[Stop]**. The label switches once again to **[Start]**.

28.2.1 Parameter description

"Manual control / Measure" dialog

Sensor

Selection of the sensor from the sensor list. Sensors are defined under **System ► Sensors**.

Selection	Selection of configured sensors
-----------	---------------------------------

Measuring input

Selection of the measuring input the sensor is connected to. The selection depends on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Measuring mode

Selection of the measuring mode. Only those measuring modes are displayed which are advisable for the selected sensor.

Selection	pH U Ipol Upol T
pH	potentiometric pH measurement
U	potentiometric voltage measurement
Ipol	voltametric measurement with selectable polarization current
Upol	amperometric measurement with selectable polarization voltage
T	temperature measurement

Temperature

Temperature entered manually. If a temperature sensor is connected, then the temperature will be measured continuously. This value is used for temperature correction in pH measurements.

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

[Info sensor]

Display information on the sensor.

28.3 Dosing


Manual control ► Dosing

The following dosing functions are available:

- Preparing the exchange or dosing unit (*see chapter 28.3.3, page 276*)
- Emptying the dosing unit (*see chapter 28.3.4, page 277*)
- Filling the dosing cylinder of the exchange unit/dosing unit (*see chapter 28.3.5, page 277*)
- Dosing a specified volume (*see chapter 28.3.2, page 273*)
- Dosing continuously (*see chapter 28.3.1, page 272*)

Selecting the dosing device

If several dosing devices are connected to a control instrument, then the dialog for selecting a dosing device will be opened.

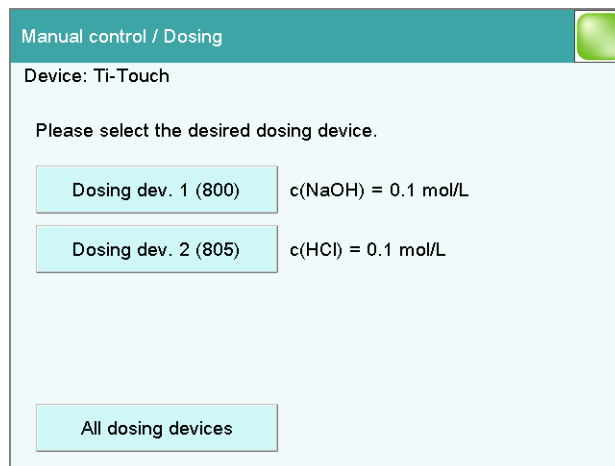

[Dosing device X]

Selection of the dosing device. The MSB connector and the type of the connected dosing device are displayed.

Various information is shown next to the button:

- Titrant
For dosing devices of the type 8XX with an intelligent exchange unit or dosing unit the titrant name is read off the data chip.
- Cylinder volume
For dosing devices of the type 8XX with an exchange unit without data chip only the cylinder volume is displayed.
- ---
For exchange or dosing units which are not or not correctly attached.
- Manual busy
When a dosing function is being carried out.


[All dosing devices]

Carrying out the functions **[Prepare]**, **[Empty]** and **[Fill]** at the same time with several dosing devices of the control device.


Selecting the dosing function

If only one dosing device is connected, then the properties of this dosing device will be shown directly. This information varies according to the type of dosing device:

- Type 8XX with integrated data chip:
If an exchange or dosing unit with integrated data chip is attached, the data stored on the data chip is displayed.

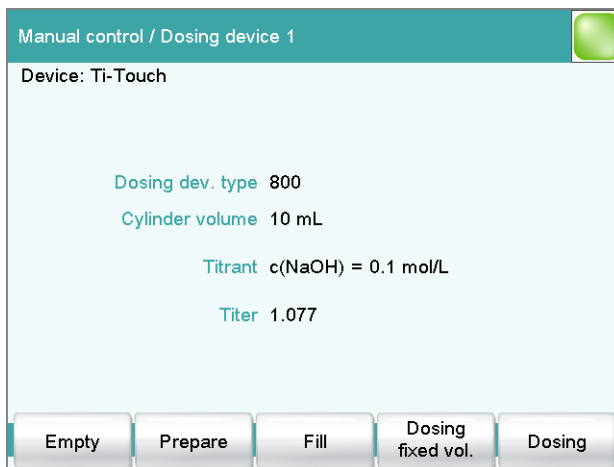

- Type 8XX without integrated data chip:
If an exchange unit without integrated data chip is attached, then the titrant can be selected from the titrant list. We recommend always selecting the titrant. This ensures that the parameters defined for the titrant for the preparing are used. Titrants are defined under **System ► Titrants**.

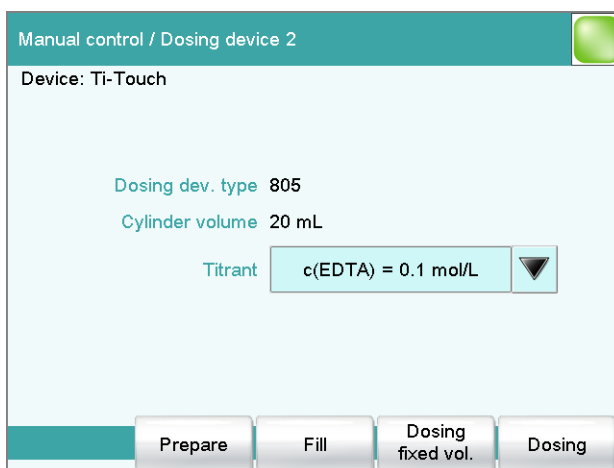**[Empty]**

Empty the cylinder and the tubings of the dosing unit (see chapter 28.3.4, page 277).

[Prepare]

Rinse the cylinder and the tubings of the exchange / dosing unit (see chapter 28.3.3, page 276).

[Fill]

Fill the cylinder of the exchange / dosing unit (see chapter 28.3.5, page 277).

[Dosing fixed vol.]


Dose a specified volume (see chapter 28.3.2, page 273).

[Dosing]

Dose continuously (see chapter 28.3.1, page 272).

28.3.1 Continuous dosing

Manual control: **Dosing ► Dosing**

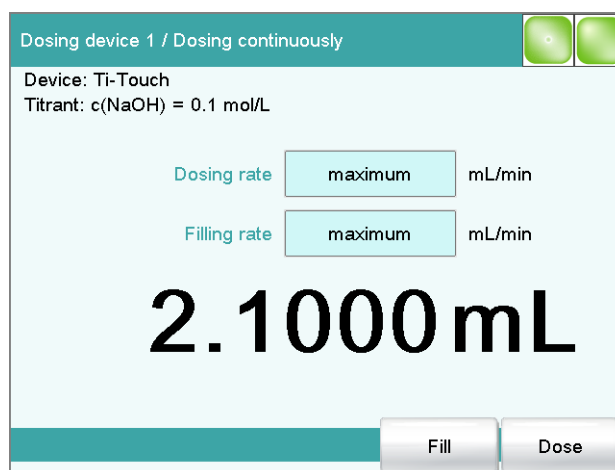

With the function **[Dosing]**, dosing is continuous for as long as you hold the button **[Dose]** pressed down.

Proceed as follows:

1 Configure the dosing function**NOTICE**

- The dosing and filling rates should be decreased for viscous liquids.
- The maximum dosing and filling rates depend on the cylinder volume (see chapter 32, page 442).

- Enter the dosing rate.
- Enter the filling rate.

2 Start dosing

- Tap on **[Dose]**.

The volume dosed is displayed. When the volume of one cylinder has been dosed, the dosing cylinder will be refilled automatically.


3 Fill the cylinder

- Tap on **[Fill]**.

The dosing cylinder is filled. The displayed volume value will be reset to 0.0000 mL.



NOTICE

If you exit the dialog with the fixed key [>], the dosing cylinder is filled automatically.

Parameter description

Dosing rate

Rate at which it is dosed.

Input range	0.01 to 166.00 mL/min
Selection	maximum dynamic
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

dynamic

This setting can only be selected when the dosing cylinder is filled. The dosing is being carried out faster and faster until the maximum dosing rate is reached (starting with 1 mL/min and then doubling the dosing rate every 1.5 s).

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 32, page 442*).

28.3.2 Dosing fixed volumes

Manual control: **Dosing ► Dosing fixed vol.**

Dosing device 1 / Dosing fixed volume

Device: Ti-Touch
Titrant: c(NaOH) = 0.1 mol/L

Volume 2.500 mL

Dosing rate 50 mL/min

Filling rate 50 mL/min

Fill automatic.

Fill Start

You can dose a particular volume with the **[Dosing fixed vol.]** function.
Proceed as follows:

1 Configure the dosing function




NOTICE

- The dosing and filling rates should be decreased for viscous liquids.
 - The maximum dosing and filling rates depend on the cylinder volume (*see chapter 32, page 442*).
- Enter the desired volume.
 - Enter the dosing rate.
 - Enter the filling rate.

2 Start dosing

- Tap on **[Start]**.

The volume display is shown. When the volume of one cylinder has been dosed, the dosing cylinder will be refilled automatically.

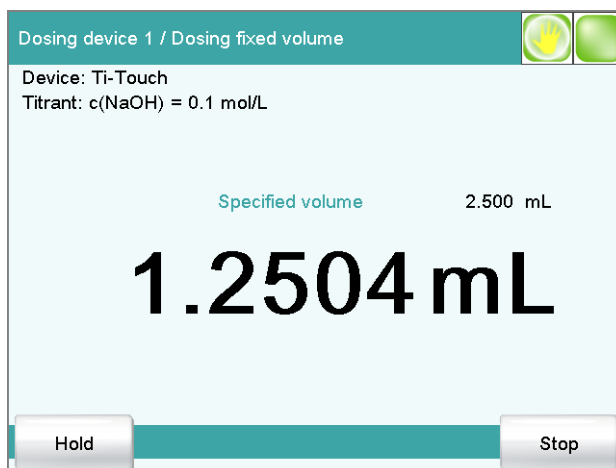

Continuous dosing is paused with **[Hold]**. The label changes to **[Continue]**. It can be used to resume the run.

3 Fill the cylinder

With the default settings (see "Fill automatic.", page 276) the dosing cylinder is filled automatically.

Otherwise:

- Tap on **[Fill]**.

The dosing cylinder is filled. The displayed volume value will be reset to 0.0000 mL.

Parameter description

Volume

Volume which is dosed.

Input range	0.00000 to 99999.9 mL
Default value	0.10000 mL

Dosing rate

The rate at which dosing takes place.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (see chapter 32, page 442).

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (see chapter 32, page 442).

Fill automatic.

on | off (Default value: **on**)

If this parameter is activated, then the dosing cylinder is being automatically filled at the end of the dosing.

28.3.3 Preparing

Manual control: **Dosing ► Prepare**

The **Prepare** function is used to rinse and fill the cylinder and tubings of the exchange or dosing unit air bubble-free. You should carry out this function before the first determination or once a day.

If the titrant is selected (see "Selecting the dosing function", page 270), the parameters defined for the titrant for preparing/emptying and the tubing parameters will be used (dosing unit only). If the titrant is not selected, default parameters will be used (see Chapter 32.2.2, page 443 and Chapter 32.3.2, page 444).

The following two figures show the live display of an exchange unit and a dosing unit:

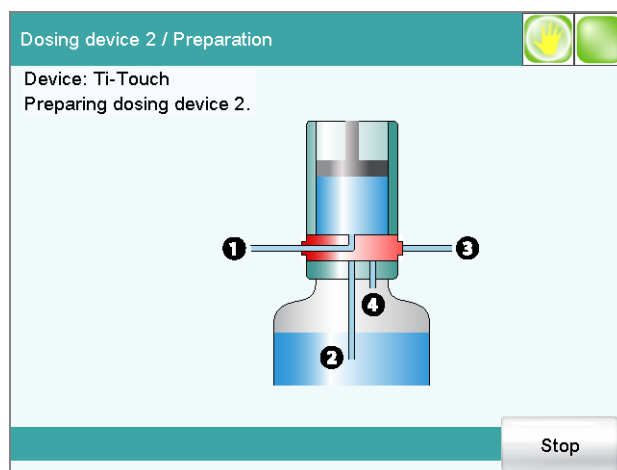


Figure 25 Live display "Preparing the dosing unit"

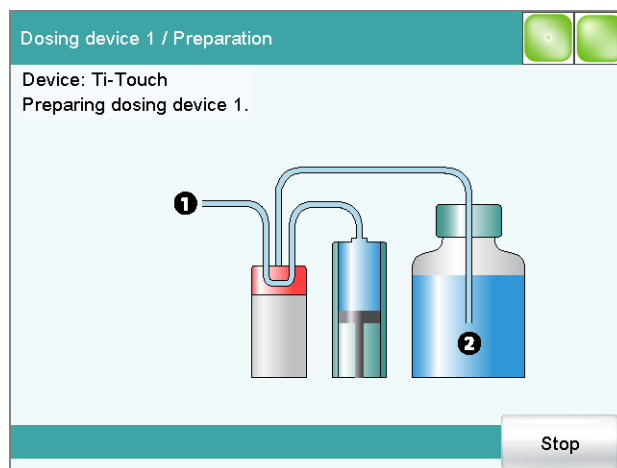


Figure 26 Live display "Preparing the exchange unit"

28.3.4 Emptying

Manual control: **Dosing ► Empty**



NOTICE

The function **Empty** is possible only with dosing units.

The cylinders and the tubings of the dosing unit are emptied with this function.

If the titrant is selected (see "Selecting the dosing function", page 270), the parameters defined for the titrant for preparing/emptying and the tubing parameters will be used. If the titrant is not selected, default parameters will be used (see Chapter 32.3.2, page 444 and Chapter 32.2.2, page 443).

28.3.5 Filling

Manual control: **Dosing ► Fill**

You can use the function **[Fill]** to fill the dosing cylinder of the exchange unit/dosing unit manually. The filling rate cannot be configured; the maximum filling rate will be used.

28.3.6 Replacing reagent



NOTICE


You can easily change the reagent in a dosing unit without no contact with the chemicals using the functions **Prepare** and **Empty**.

Proceed as follows:

- 1 Execute the function **[Empty]**.
- 2 Attach the dosing unit to the bottle with a suitable solvent.
- 3 Execute the function **[Prepare]**.
The last traces of the reagent in the cylinder and tubings are removed with the solvent.
- 4 Execute the function **[Empty]** again.
- 5 Attach the dosing unit to the bottle with the new reagent.
- 6 Execute the function **[Prepare]**.
The dosing unit can now be used.

28.4 Stirring

Manual control ► Stir

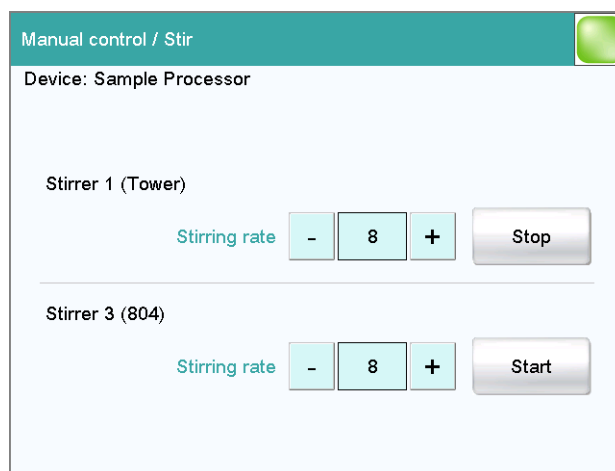

The MSB connector and type of stirrer is indicated for each connected stirrer.

You can control a connected stirrer manually with the function **[Stir]**.

Proceed as follows:

1 Setting the stirring rate

- Tap on the [-] or [+] button.
Each time one of the buttons is tapped, the stirring rate is reduced or increased by one level.

2 Switching on the stirrer

- Tap on [Start].

The stirrer is started.

3 Switching off the stirrer

- Tap on [Stop].

The stirrer is stopped.

Parameter description

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 32.4, page 444*.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8



NOTICE

If a stirrer is used in the determination run, it can still be manually controlled at the same time. For example, you can use the manual control to reduce the stirring rate of a stirrer which has been started in the determination run.

28.5 Remote

Manual control ► Remote

With the function **[Remote]**, you can define manual output signals to the remote interface of a connected Remote Box. It is not possible to define input signals; the current status of the input lines is, however, displayed.

Inputs

Current status of the input lines.

Outputs

Selection of the signal out of the templates or entering the required bit pattern. Templates are defined under **System ► Templates ► Output lines**.

Entering the bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status
- p = set pulse

The output lines are always numbered from right to left, i.e. with the signal **00000000000001** line 0 is set. With a pulse, the length is set to 200 ms. If you wish to set pulses with other lengths, you have to define a corresponding template.


Entry	Bit pattern containing exactly 14 characters or a max. of 24 characters for the name of the template
Default value	00000000000001
Selection	Selection of the templates defined

[Set]

Set a defined output signal.

28.6 Sample Processor

Manual control ► Sample Processor

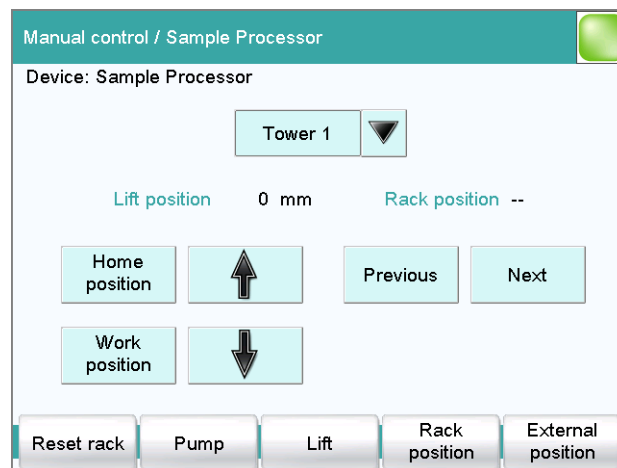

With the function **[Sample Processor]**, you can control a connected Sample Processor manually. The following functions are available:

- Rotating the sample rack
- Moving the lift
- Defining specific lift positions (work position, rinse position, etc.)
- Switching pumps on and off
- Initializing the rack
- Defining external positions outside of the sample rack (only with Swing Head)

Tower for the manual control.

Selection	Tower 1 Tower 2
Tower 1	View from the front, the right tower.
Tower 2	View from the front, the left tower. This selection is only available for Sample Processors with two towers.

Lift position

Current lift position.

Rack position

Current rack position. If the rack is not set to a valid position, e.g. after a rack reset, -- is displayed to indicate this.



NOTICE

Note that the rack has to be set to a valid position for all lift movements! Otherwise a corresponding error message will be displayed.

Move to a valid rack position with **[Previous]** or **[Next]**. Alternatively, you can carry out a MOVE command.

[Home position]

Move the lift to the home position (corresponds to lift position 0 mm).

[Work position]

Move the lift to the work position of the current rack position or the external position.

The work position can directly be defined here (see *"Assigning lift positions"*, page 284) or in the device manager (see *"Lift positions Tower 1/2"*, page 119).

[↑]

Continuously move the lift upwards as long as the button is pressed down. The lift rate used is defined in the dialog **Sample Processor / Lift**.

[↓]

Continuously move the lift downwards as long as the button is pressed down. The lift rate used is defined in the dialog **Sample Processor / Lift**.

[Previous]

Move to the previous rack position.

[Next]

Move to the next rack position.

[Reset rack]

Initialize the rack. This function is identical with the command **RACK**. The following actions are carried out:

- The lift is moved upwards.
- The sample rack is rotated to the starting position.
- The rack code of the rack attached is being read out and the corresponding rack data is transferred to the Sample Processor.
- The sample variable is being reset to the value 1.

[Pump]

Switch on and off the pumps mounted or connected to the tower.

[Lift]

Move the lift to any or a specific lift position and assigning lift positions (see chapter 28.6.1, page 283).

[Rack position]

Move the rack to any sample position or to defined positions which are defined as special beakers (see chapter 28.6.2, page 285).

[External position]

This button is only enabled when a Swing Head is mounted to the Sample Processor.

Swing the robotic arm to an external position outside the sample rack and assign the swing angle and the lift position (see chapter 28.6.3, page 286).

28.6.1 Moving the lift

Manual control: **Sample Processor ▶ Lift**

The screenshot shows a software interface for controlling the lift. The title bar reads 'Sample Processor / Lift'. Below the title bar, it says 'Tower 1'. The main area contains three rows of controls: 'Current lift pos.' with a value of '0 mm'; 'Move to lift pos.' with a dropdown menu showing 'Work position' and a unit indicator 'mm'; and 'Lift rate' with a value of '25' and a unit indicator 'mm/s'. At the bottom of the dialog, there are two buttons: 'Assign lift pos.' on the left and 'Start' on the right.

In the dialog **Sample Processor / Lift**, you can move the lift to any lift position. You can also define frequently used positions (work position, rinse position, etc.) as specific lift positions and then to move to them with ease (see "Assigning lift positions", page 284). This definition is possible in this dialog for the attached rack, and in the device manager for all of the racks in the list.

Current lift pos.

Current lift position.

Move to lift pos.

Selecting a predefined lift position or enter any lift position.

Input range	0 to 'maximum stroke path' mm A lift position of 0 mm corresponds to the "home position", i.e. the lift is located at the upper stop position. The maximum stroke path is defined in the properties of the tower (<i>see "Maximum stroke path", page 109</i>). If a higher value is entered, an error message will be displayed.
Selection	Work position Shift position Rinse position Special position
Default value	Work position

Lift rate

Rate at which the lift is moved in the manual control.

Input range	5 to 25 mm/s
Default value	25 mm/s

[Assign lift pos.]

Assign the current lift position to a preset lift position.

Assigning lift positions

Specific lift positions are stored separately for every sample rack used as well as for tower 1 and tower 2. The following specific lift positions are available:

- **General rack positions**
For general rack positions you can define one work position, one shift position and one rinse position as well as one special position.
- **Special beaker positions**
A specific work position can be defined for each special beaker. The definitions of the general rack positions are used for the shift position, rinse position and special position of the affected tower.
- **External positions** (only with robotic arm and mounted Swing Head, *see Chapter 28.6.3, page 286*)
For each of the four possible external positions, a specific work position can be defined. Shift position and rinse position can only be defined for all four external positions commonly. A special position is not possible.

Proceed as follows to assign the current lift position to a specific lift position:

1 Move to lift position

- Enter the desired lift position in mm and tap on [**Start**].

The lift moves to the desired lift position.

2 Assign lift position

- Tap on **[Assign lift pos.]**.
The **Lift / Assign lift position** dialog is displayed:

Lift / Assign lift position	
Current lift pos. 100 mm	
Position name	Position height
Work position	130 mm
Shift position	50 mm
Rinse position	100 mm
Special position	0 mm
Assign	

- Select the desired specific lift position and tap on **[Assign]**.
The current lift position is assigned to the specific lift position.

28.6.2 Moving to a rack position

Manual control: **Sample Processor ► Rack position**

Sample Processor / Rack position	
Tower 1	
Current rack pos. 1	
Rack position	1 ▼
Shift rate	Special beaker 1 %/s
Shift direction	auto ▼
Swing rate	55 %/s
Start	

In the dialog **Sample Processor / Rack position**, you can move to any position on the attached rack.

Current rack pos.

Current rack position. If the rack is not set to a valid position, e.g. after a rack reset, -- is displayed to indicate this.

Rack position

Desired rack position.



Input range	1 to Number of positions on the rack attached.
Default value	1
Selection	Special beaker 1...16

Special beaker 1...16

The selection depends on how many rack positions are defined as special beakers (see "Editing special beakers", page 120).

Shift rate

Rate at which the sample rack is moved.

Input range	3 to 20 °/s
Default value	20 °/s

Shift direction

Direction in which the rack is moved.

Selection	auto + -
Default value	auto

auto

A shift direction, with which the shorter way has to be passed, is automatically selected.

+

Counterclockwise rotation.

-

Clockwise rotation.


Swing rate

Rate at which the robotic arm is swung in the manual control.

Input range	10 to 55 °/s
Default value	55 °/s

28.6.3 External positions

Manual control: **Sample Processor ► External position**

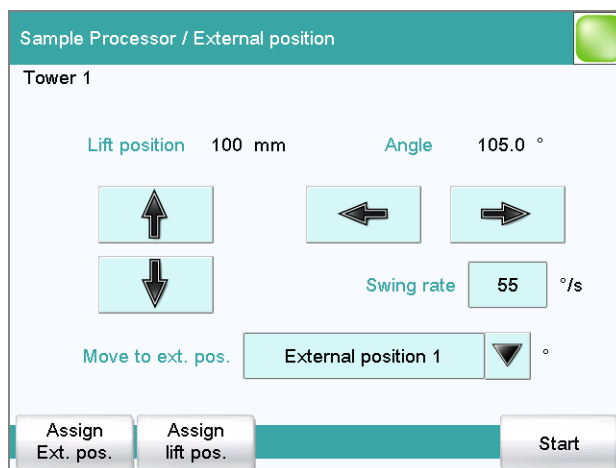

NOTICE

This dialog is accessible only if you have a Swing Head with robotic arm attached to the Sample Processor.

In the dialog **Sample Processor / External position**, you can swivel the robotic arm to an external position outside of the sample rack. As with the rack positions, here you can also define up to four frequently used external positions as specific positions. You can assign specific lift positions to these positions, analogous to the rack positions.

External positions are required if you have, for example, a measuring vessel or a titration cell mounted in addition to the rack.

Lift position

Current lift position.

Angle

Current angle position of the robotic arm.

[↑]

Only enabled, when the robotic arm has been moved to a specific external position with **[Start]**.

Continuously move the lift upwards as long as the button is pressed down. The lift rate used is defined in the dialog **Sample Processor / Lift**.

[↓]

Only enabled, when the robotic arm has been moved to a specific external position with **[Start]**.

Continuously move the lift downwards as long as the button is pressed down. The lift rate used is defined in the dialog **Sample Processor / Lift**.



[←]

Swing the robotic arm in clockwise direction as long as the button is pressed down.

[⇒]

Swing the robotic arm in counterclockwise direction as long as the button is pressed down.

Swing rate

Rate at which the robotic arm is swung in the manual control.

Input range	10 to 55 °/s
Default value	55 °/s

Move to ext. pos.

Selecting a predefined position or entering any swing angle.

Input range	(Offset) to (Offset + max. swing range) ° The offset is made up of a design-dependent angle (approx. 8...9°) together with the robotic arm offset from the robotic arm properties. The maximum swing range is also defined under the robotic arm properties (see " <i>Properties – Robotic arm</i> ", page 112).
Selection	External position 1...4

[Assign Ext. pos.]

Assign the current angle position of the robotic arm to an external position as swing angle.

[Assign lift pos.]

Assign the current lift position to a preset lift position.

Assigning swing angles and lift positions

For each of the four possible external positions, a specific work position can be defined. Shift position and rinse position can only be defined for all four external positions commonly. A special position is not possible.

Proceed as follows to assign a swing angle to an external position and to assign the current lift position to a specific lift position:

1 Move to the external position

- Swing the robotic arm to the desired position with the arrow keys [←] or [⇒].

2 Assign the external position

- Tap on **[Assign Ext. pos.]**.
The **External position / Assign swing angle** dialog is displayed:

External position / Assign swing angle	
Current robotic arm angle 98.0 °	
External position	Angle
1	105.0 °
2	60.0 °
3	60.0 °
4	60.0 °
Assign	

- Select the desired external position and tap on **[Assign]**.

The current angle position of the robotic arm is assigned to the external position.

3 Move to the external position

- Tap on [**↩**].
The next higher-level dialog is displayed.
- Select **Move to ext. pos. = External position X** (X = 1...4) and tap on **[Start]**.

The robotic arm swings to the desired position.


4 Move to lift position

- Move the lift to the desired height with the arrow keys [**↑**] or [**↓**].

5 Assign lift position

- Tap on **[Assign lift pos.]**.

The **Lift / Assign lift position** dialog is displayed:

Lift / Assign lift position 	
Current lift pos. 100 mm	
Position name	Position height
Work position	130 mm
Shift position	50 mm
Rinse position	100 mm
Special position	0 mm
Assign	

- Select the desired specific lift position.
- Tap on **[Assign]**.

The current lift position is assigned to the specific lift position.

29 Parameters



NOTICE

For most of the numerical parameters, the result of a calculation can also be utilized in place of a number. A description of this can be found in the appendix (*see chapter 32.7, page 449*).

29.1 Dynamic equivalence point titrations (DET)

29.1.1 Start conditions

The parameters that are carried out before the start of titration are defined under **[Start conditions]**. The start conditions are processed in the following sequence:

1. Start volume
2. Start measured value
3. Start slope
4. Pause

Start volume

Volume that is dosed prior to the start of the titration.

Input range	0.00000 to 9999.99 mL
Default value	0.00000 mL

Dosing rate

Rate at which the start volume is dosed.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

Pause

Waiting time, e.g. for the electrode to settle down after the start or a reaction time after the dosing of a start volume. The time entered here is only running when all start conditions have been fulfilled.

Input range	0 to 999999 s
Default value	0 s



More start conditions

Start meas. value

Before starting the titration, it will be dosed until this measured value is reached. If the start measured value is reached by the dosing of a start volume, then the titration starts directly.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measuring mode pH:

Input range	-20.000 to 20.000
Selection	off
Default value	off

Measuring mode U, Ipol:

Input range	-2000.0 to 2000.0 mV
Selection	off
Default value	off

Measuring mode Upol:

Input range	-200.0 to 200.0 μA
Selection	off
Default value	off

Start slope

Before starting the titration, it will be dosed until this slope (measured value per volume) is reached. If the start slope is reached by the dosing of a start volume, then the titration starts directly.

Measuring mode pH:

Input range	00.000 to 9,999 pH/mL
Selection	off
Default value	off

Measuring mode U, Ipol:

Input range	0 to 999 mV/mL
Selection	off
Default value	off

Measuring mode Upol:

Input range	0 to 99 μA/mL
Selection	off
Default value	off

Dosing rate

Rate to be dosed with until the start measured value or the start slope is reached.

Input range	0.01 to 166.00 mL/min
Default value	5.00 mL/min
Selection	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

Initial measured value

The [**Initial meas. value**] is measured before carrying out the **start conditions**. This value is used for the **titration direction determination**, which is necessary for the **start measured value** and **stop measured value** parameters.

Signal drift

Maximum permissible drift for the measured value acceptance, i.e. maximum change of the measured value per minute. For this value, a suitable maximum waiting time is automatically calculated except you already have modified this waiting time.

Measuring mode pH, U and Ipol:

Input range	0.1 to 999.0 mV/min
Selection	off
Default value	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed.

Measuring mode Upol:

Input range	0.01 to 99.90 µA/min
Selection	off
Default value	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed.

Min. waiting time

The measured value is not accepted until the minimum waiting time has elapsed, even if the signal drift has already been reached. The minimum waiting time is only important for drift-controlled measurements.

measurement = automatic or **continuous** is defined (see sensor dialog of the command). This value is used for temperature correction in pH measurements.

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

User-defined parameters

These parameters are only accessible when **Titration rate = user**.

Meas. point density

A small value means small volume increments, i.e. a high measuring point density. The curve then shows all the finest details which also include noise; this could cause unwanted equivalence points to be found. A larger value, i.e. a smaller measuring point density, permits quicker titrations. If you are using a dosing device with a small cylinder volume then a smaller measuring point density value may be beneficial. However, you should also set a smaller signal drift and a higher EP criterion at the same time.

Input range	0 to 9
Default value	4

Min. increment

This smallest permitted volume increment is added at the start of the titration and with steep curves in the region of the equivalence point. Very small values should only be used if a low titrant consumption is expected; otherwise unwanted equivalence points could be evaluated.

Input range	0.05 to 999.90 µL
Default value	10.00 µL

Max. increment

A maximum volume increment should be selected in the following cases:

- when titration consumption is very low up until the equivalence point is reached.
- when a start volume is dosed up until shortly before the equivalence point is reached.
- when the change of direction in the jumping range is very abrupt, because otherwise it could easily happen that an excessively large volume will be dosed in the region of the equivalence point.

The value should not be less than 1/100 cylinder volume.

Input range	0.1 to 9999.9 µL
Selection	off
Default value	off

**NOTICE**

It is not advisable to select similar volumes for the minimum and the maximum increment. Monotonic equivalence point titration (MET) is appropriate for these applications.

Dosing rate

Rate at which the volume increments are dosed.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

Signal drift

Maximum permissible drift for the measured value acceptance, i.e. maximum change of the measured value per minute. This type of titration is often referred to as equilibrium titration. For this value, a suitable maximum waiting time is automatically calculated except you already have modified this waiting time.

**NOTICE**

A constant measured value is often only reached after a certain time, as mixing and the reaction itself require a certain time. The response time of an electrode can also increase with time, i.e. reaching a constant measured value takes longer and longer. Drift-controlled measured value acceptance is particularly advisable in such cases, as the measured values are only accepted when equilibrium has almost been reached.

Measuring mode pH, U and Ipol:

Input range	0.1 to 999.0 mV/min
Default value	50.0 mV/min
Selection	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed. This can be useful when the titration reaction proceeds slowly or the electrode is slow to respond.

Measuring mode Upol:

Input range	0.01 to 99.90 $\mu\text{A}/\text{min}$
Default value	50.00 $\mu\text{A}/\text{min}$
Selection	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed. This can be useful when the titration reaction proceeds slowly or the electrode is slow to respond.

Min. waiting time

The measured value is not accepted until the minimum waiting time has elapsed, even if the signal drift has already been reached. The minimum waiting time is only important for drift-controlled measurements.

Input range	0 to 999999 s
Default value	0 s

Max. waiting time

If the signal drift has been switched off or has not yet been reached, then the measured value will be accepted when the maximum waiting time has elapsed. As default value, a suitable waiting time for the signal drift is automatically calculated.

Input range	0 to 999999 s
Default value	26 s

Table 9 Default values of the predefined titration rates for DET

	Titration rate		
	slow	optimal	fast
Meas. point density	2	4	6
Min. increment	10.00 μL	10.00 μL	30.00 μL
Max. increment	off	off	off
Dosing rate	maximum	maximum	maximum
Signal drift			
– pH, U and Ipol	20.0 mV/min	50.0 mV/min	80.0 mV/min
– Upol	20.0 $\mu\text{A}/\text{min}$	50.0 $\mu\text{A}/\text{min}$	80.0 $\mu\text{A}/\text{min}$
Min. waiting time	0 s	0 s	0 s
Max. waiting time	38 s	26 s	21 s



29.1.3 Stop conditions

The conditions for canceling the titration are defined under [Stop conditions].

Stop volume

The titration is stopped when the specified volume has been dosed since the start of the titration. This volume should be adjusted to the size of the titration vessel in order to prevent the contents from running over.

Input range	0.0000 to 9999.99 mL
Default value	100,000 mL
Selection	off

Stop meas. value

The titration is canceled when the specified measured value has been reached since the start of the titration.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measuring mode pH:

Input range	-20.000 to 20.000
Selection	off
Default value	off

Measuring mode U, Ipol:

Input range	-2000.0 to 2000.0 mV
Selection	off
Default value	off

Measuring mode Upol:

Input range	-200.0 to 200.0 µA
Selection	off
Default value	off

Stop EP

The titration is stopped when the specified number of equivalence points has been found.

Input range	1 to 9
Default value	9
Selection	off

Volume after EP

This volume will be added when the number of equivalence points defined under **Stop EP** has been found. The curve shape after the equivalence point can also be seen this way.

Input range	0.01000 to 9999.99 mL
Selection	off
Default value	off

Stop time

The titration is stopped when the specified time has elapsed since the start of the titration.

Input range	0 to 999999 s
Selection	off
Default value	off

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 32, page 442*).

29.1.4 Potentiometric evaluation

The parameters for the evaluation of the titration curve are defined under **[Potentiometr. evaluation]**.

EP criterion

The Equivalence point Recognition Criterion (ERC) found is compared with this value. Equivalence points whose ERC is less than the value defined here will be ignored.

Input range	0 to 200
Default value	5

EP recognition

This parameter allows you to filter out only the equivalence points that are being sought.

Selection	all greatest last ascending descending off
Default value	all

**all**

All equivalence points will be recognized.

greatest

Only the equivalence point with the greatest ERC value, i.e. the steepest jump, will be recognized.

last

Only the last equivalence point will be recognized.

ascending

Only equivalence points with a positive slope of the titration curve will be recognized.

descending

Only equivalence points with a negative slope of the titration curve will be recognized.

off

No evaluation takes place.

Set windows

This parameter allows to recognize equivalence points only in a particular range (so-called window) of the curve. Equivalence points outside this window will not be recognized. A maximum of 9 windows can be defined. Setting windows is advisable when interference and unnecessary equivalence points are to be suppressed.

Selection	Measured value Volume off
Default value	off

Measured value

The windows are defined on the measured value axis.

Volume

The windows are defined on the volume axis.

off

The entire titration curve is being evaluated.

Setting windows

The list of defined windows is opened with **[Set window]**.

When the list is opened for the first time, a window over the entire measured value range or volume range is already defined. The windows must not overlap, they only may lie next to one another. For each window, own criteria can be defined for the equivalence point recognition.

**NOTICE**

When you define a result variable instead of a numerical value as limit value (see chapter 32.7, page 449), then just one window can be defined.

[New]

Define a new window. This is only possible when not the entire range is covered yet.

[Delete]

Delete the selected window.

[Edit]

Edit the settings of the selected window.

Lower limit

Measured value or volume for the lower limit.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measured value (measuring mode pH):

Input range	-20.000 to 20.000
Default value	-20.000

Measured value (measuring mode U, I_{pol}):

Input range	-2000.0 to 2000.0 mV
Default value	-2000.0 mV

Measured value (measuring mode U_{pol}):

Input range	-200.00 to 200.00 µA
Default value	-200.00 µA

Volume:

Input range	0.00000 to 9999.99 mL
Default value	0.00000 mL

Upper limit

Measured value or volume for the upper limit.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.



Measured value (measuring mode pH):

Input range	-20.000 to 20.000
Default value	20.000

Measured value (measuring mode U, I_{pol}):

Input range	-2000.0 to 2000.0 mV
Default value	2000.0 mV

Measured value (measuring mode U_{pol}):

Input range	-200.00 to 200.00 µA
Default value	200.00 µA

Volume:

Input range	0.00000 to 9999.99 mL
Default value	9999.99 mL

EP criterion

The Equivalence point Recognition Criterion (ERC) found is compared with this value. Equivalence points whose ERC is less than the value defined here will be ignored.

Input range	0 to 200
Default value	5

EP recognition

This parameter allows you to filter out only the equivalence points that are being sought.

Selection	first greatest last ascending descending
Default value	first

first

Only the first equivalence point will be recognized.

greatest

Only the equivalence point with the greatest ERC value, i.e. the steepest jump, will be recognized.

last

Only the last equivalence point will be recognized.

ascending

Only equivalence points with a positive slope of the titration curve will be recognized.

descending

Only equivalence points with a negative slope of the titration curve will be recognized.

Only one equivalence point will be recognized in a window. The numbering of the equivalence points (EP) is defined by the numbering of the win-

dows (e.g. EP2 in window 2), so that even if EPs are missing, the calculations will still be carried out with the correctly assigned EP volumes.

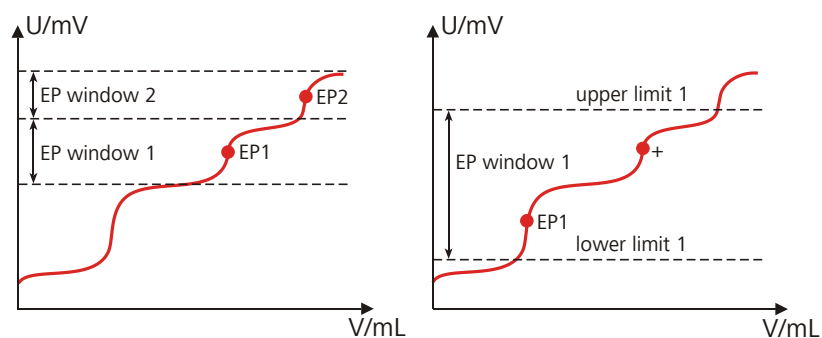


Figure 27 Equivalence point recognition and equivalence point numbering within windows

In the first example in the above figure, two equivalence points are recognized in two different windows (EP1 in window 1 and EP2 in window 2). In the second example, two equivalence points are found in one window, but only the first one is recognized. In order to ensure that the user recognizes that more than one equivalence point was found in the set window, EP1 is marked with a "+" in the result view. In addition, a corresponding message is entered in the message list.

Evaluation and equivalence point criterion with DET

The equivalence points (EP) are localized in a way similar to the Tubbs method [1][2]. The volume value of the equivalence point (V_E) is shifted from the inflection point (see arrow) towards the smaller circle of curvature for real asymmetric titration curves.

[1] C. F. Tubbs, *Anal. Chem.* **1954**, 26, 1670–1671.

[2] E. Bartholomé, E. Biekert, H. Hellmann, H. Ley, M. Weigert, E. Weise, *Ullmanns Encyklopädie der technischen Chemie*, Vol. 5, Verlag Chemie, Weinheim, 1980, p. 659.

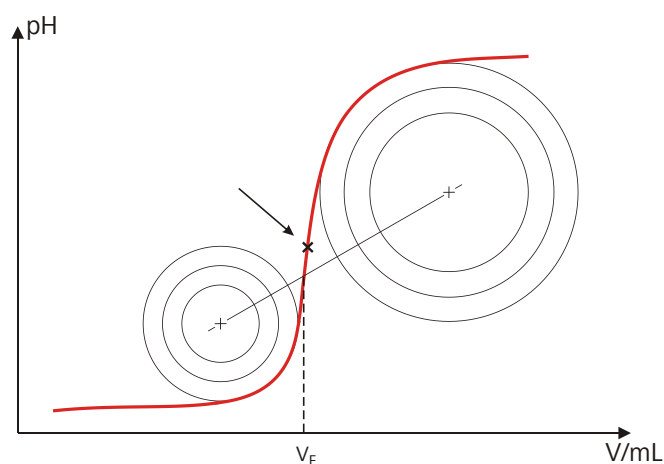


Figure 28 Tubbs method for determining the equivalence point



The figure shows that the evaluation still requires measured values from the measuring point list even after the equivalence point.

For the recognition of the EPs found, the set EP criterion is compared to the ERC (Equivalence point Recognition Criterion) found. The ERC is the first derivative of the titration curve combined with a mathematical function that is more sensitive for flat jumps than for steeper ones. EPs whose ERC is smaller than the defined EP criterion will not be recognized. The ERC is displayed in the results dialog for each discovered and recognized EP. If you adjust the EP criterion later in order to recognize more or fewer EPs, then you can reevaluate the determination (**[Recalculate]** function in the results dialog).

29.1.5 Control instrument

The control instrument the determination is carried out with is selected under **[Control device]**. Control instruments are defined under **System ► Device manager**.



NOTICE

This button will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control instrument.

Control device

Selection of the control instrument from the list of devices. Only those devices are displayed which are able to carry out the command.

Selection	Selection of configured control instruments
-----------	--

29.1.6 Sensor

The parameters for the sensor are edited under **[Sensor]**.

Measuring input

Selection of the measuring input the sensor is connected to. The selection is not dependent on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Sensor

Selection of the sensor from the sensor list. The selection depends on the measuring mode. Sensors are defined under **System ► Sensors**. You can also enter a sensor name which is not contained in the sensor list. When a determination is started there is a check whether the sensor is contained in the sensor list.

Selection	Selection of configured sensors
-----------	--

I(pol)

The polarization current is the current that is applied to a polarizable electrode during voltametric measurement. This parameter is available only with I(pol) determinations.

The values that are actually adjustable may vary from the values specified here. The binding values can be found in the manual, in the *Technical specifications* chapter.

Input range	-125.0 to 125.0 μA (Increment: 2.5)
Default value	5.0 μA
Selection	-1.0 μA 1.0 μA

U(pol)

The polarization voltage is the voltage applied to the polarizable electrode during an amperometric measurement. This parameter is available only with U(pol) determinations.

The values that are actually adjustable may vary from the values specified here. The binding values can be found in the manual, in the *Technical specifications* chapter.

Input range	-1,250 to 1,250 mV (Increment: 25)
Default value	400 mV

Electrode check

on | off (Default value: **off**)

For the following electrodes, an electrode check can be carried out:

- pH electrodes
- Metal electrodes
- Ion-selective electrodes

A check is made that the electrode is properly connected and that no short-circuit is present. The electrode check is carried out when this command is started. Note that this electrode check has nothing to do with the electrode test (command **ELT**).

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

continuous

A temperature sensor must be connected. The temperature is measured continuously.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 32.4, page 444*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

Switch off automatically

on | off (Default value: **on**)

If this parameter is activated, the stirrer will be switched off automatically at the end of the titration, measurement, etc.

29.2 Monotonic equivalence point titrations (MET)

29.2.1 Start conditions

The parameters that are carried out before the start of titration are defined under **[Start conditions]**. The start conditions are processed in the following sequence:

1. Start volume
2. Start measured value
3. Start slope
4. Pause

Start volume

Volume that is dosed prior to the start of the titration.

Input range	0.00000 to 9999.99 mL
Default value	0.00000 mL



Dosing rate

Rate at which the start volume is dosed.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

Pause

Waiting time, e.g. for the electrode to settle down after the start or a reaction time after the dosing of a start volume. The time entered here is only running when all start conditions have been fulfilled.

Input range	0 to 999999 s
Default value	0 s

More start conditions

Start meas. value

Before starting the titration, it will be dosed until this measured value is reached. If the start measured value is reached by the dosing of a start volume, then the titration starts directly.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measuring mode pH:

Input range	-20.000 to 20.000
Selection	off
Default value	off

Measuring mode U, I_{pol}:

Input range	-2000.0 to 2000.0 mV
Selection	off
Default value	off

Measuring mode U_{pol}:

Input range	-200.0 to 200.0 μA
Selection	off
Default value	off

Start slope

Before starting the titration, it will be dosed until this slope (measured value per volume) is reached. If the start slope is reached by the dosing of a start volume, then the titration starts directly.

Measuring mode pH:

Input range	00.000 to 9,999 pH/mL
Selection	off
Default value	off

Measuring mode U, Ipol:

Input range	0 to 999 mV/mL
Selection	off
Default value	off

Measuring mode Upol:

Input range	0 to 99 µA/mL
Selection	off
Default value	off

Dosing rate

Rate to be dosed with until the start measured value or the start slope is reached.

Input range	0.01 to 166.00 mL/min
Default value	5.00 mL/min
Selection	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

Initial measured value

The **[Initial meas. value]** is measured before carrying out the **start conditions**. This value is used for the **titration direction determination**, which is necessary for the **start measured value** and **stop measured value** parameters.

Signal drift

Maximum permissible drift for the measured value acceptance, i.e. maximum change of the measured value per minute. For this value, a suitable maximum waiting time is automatically calculated except you already have modified this waiting time.

Measuring mode pH, U and Ipol:

Input range	0.1 to 999.0 mV/min
-------------	----------------------------

optimal

For all standard titrations. The parameters have been optimized for the most frequent applications.

fast

For fast and less critical titrations.

user

The individual titration parameters can be modified.

**NOTICE**

Select **optimal** as titration rate when you are developing a new titration method. This parameter is suitable for a lot of titrations and only needs adaptation in special cases.

The settings of the individual titration rates are listed in *table 10, page 313*.

Temperature

Temperature entered manually. The temperature is being continuously measured when a temperature sensor is connected and when **Temp. measurement = automatic** or **continuous** is defined (see sensor dialog of the command). This value is used for temperature correction in pH measurements.

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

User-defined parameters

These parameters are only accessible when **Titration rate = user**.

Volume increment

The volume that is dosed at each dosing step. High accuracy requires using the correct volume increment.

A good guideline for the volume increment is 0.1 mL. A total of 20 dosing steps is recommended until the equivalence point is reached.

With greater consumption, the titration time can be optimized using the start conditions. Smaller volume increments (smaller than 0.1 mL) are required to determine blank values or to ensure accuracy with highly asymmetric curves. Larger volume increments are advisable for flatter curves.



NOTICE

The accuracy of the evaluation cannot be increased by using smaller increments as the measured value changes between two measuring points are then of the same order of magnitude as the noise.

Input range	0.00005 to 999.900 mL
Default value	0.10000 mL

Dosing rate

Rate at which the volume increments are dosed.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

Signal drift

Maximum permissible drift for the measured value acceptance, i.e. maximum change of the measured value per minute. This type of titration is often referred to as equilibrium titration. For this value, a suitable maximum waiting time is automatically calculated except you already have modified this waiting time.



NOTICE

A constant measured value is often only reached after a certain time, as mixing and the reaction itself require a certain time. The response time of an electrode can also increase with time, i.e. reaching a constant measured value takes longer and longer. Drift-controlled measured value acceptance is particularly advisable in such cases, as the measured values are only accepted when equilibrium has almost been reached.

Measuring mode pH, U and I_{pol}:

Input range	0.1 to 999.0 mV/min
Default value	50.0 mV/min
Selection	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed. This can be useful when the titration reaction proceeds slowly or the electrode is slow to respond.

Measuring mode Upol:

Input range	0.01 to 99.90 $\mu\text{A}/\text{min}$
Default value	50.00 $\mu\text{A}/\text{min}$
Selection	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed. This can be useful when the titration reaction proceeds slowly or the electrode is slow to respond.

Min. waiting time

The measured value is not accepted until the minimum waiting time has elapsed, even if the signal drift has already been reached. The minimum waiting time is only important for drift-controlled measurements.

Input range	0 to 999999 s
Default value	0 s

Max. waiting time

If the signal drift has been switched off or has not yet been reached, then the measured value will be accepted when the maximum waiting time has elapsed. As default value, a suitable waiting time for the signal drift is automatically calculated.

Input range	0 to 999999 s
Default value	26 s

Table 10 Default values of the predefined titration rates for MET

	Titration rate		
	slow	optimal	fast
Volume increment	0.05000 mL	0.10000 mL	0.20000 mL
Dosing rate	maximum	maximum	maximum
Signal drift			
– pH, U and I _{pol}	20.0 mV/min	50.0 mV/min	80.0 mV/min
– U _{pol}	20.0 $\mu\text{A}/\text{min}$	50.0 $\mu\text{A}/\text{min}$	80.0 $\mu\text{A}/\text{min}$
Min. waiting time	0 s	0 s	0 s
Max. waiting time	38 s	26 s	21 s

Volume after EP

This volume will be added when the number of equivalence points defined under **Stop EP** has been found. The curve shape after the equivalence point can also be seen this way.

Input range	0.01000 to 9999.99 mL
Selection	off
Default value	off

Stop time

The titration is stopped when the specified time has elapsed since the start of the titration.

Input range	0 to 999999 s
Selection	off
Default value	off

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 32, page 442*).

29.2.4 Potentiometric evaluation

The parameters for the evaluation of the titration curve are defined under **[Potentiometr. evaluation]**.

EP criterion

The Equivalence point Recognition Criterion (ERC) found is compared with this value. Equivalence points whose ERC is less than the value defined here will be ignored.

Measuring mode pH:

Input range	0.10 to 9.99
Default value	0.50

Measuring mode U, I_{pol}:

Input range	1 to 999 mV
Default value	30 mV

Setting windows

The list of defined windows is opened with **[Set window]**.

When the list is opened for the first time, a window over the entire measured value range or volume range is already defined. The windows must not overlap, they only may lie next to one another. For each window, own criteria can be defined for the equivalence point recognition.



NOTICE

When you define a result variable instead of a numerical value as limit value (see chapter 32.7, page 449), then just one window can be defined.

[New]

Define a new window. This is only possible when not the entire range is covered yet.

[Delete]

Delete the selected window.

[Edit]

Edit the settings of the selected window.

Lower limit

Measured value or volume for the lower limit.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measured value (measuring mode pH):

Input range	-20.000 to 20.000
Default value	-20.000

Measured value (measuring mode U, I_{pol}):

Input range	-2000.0 to 2000.0 mV
Default value	-2000.0 mV

Measured value (measuring mode U_{pol}):

Input range	-200.00 to 200.00 µA
Default value	-200.00 µA

Volume:

Input range	0.00000 to 9999.99 mL
Default value	0.00000 mL



Upper limit

Measured value or volume for the upper limit.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measured value (measuring mode pH):

Input range	-20.000 to 20.000
Default value	20.000

Measured value (measuring mode U, I_{pol}):

Input range	-2000.0 to 2000.0 mV
Default value	2000.0 mV

Measured value (measuring mode U_{pol}):

Input range	-200.00 to 200.00 µA
Default value	200.00 µA

Volume:

Input range	0.00000 to 9999.99 mL
Default value	9999.99 mL

EP criterion

The Equivalence point Recognition Criterion (ERC) found is compared with this value. Equivalence points whose ERC is less than the value defined here will be ignored.

Measuring mode pH:

Input range	0.10 to 9.99
Default value	0.50

Measuring mode U, I_{pol}:

Input range	1 to 999 mV
Default value	30 mV

Measuring mode U_{pol}:

Input range	0.1 to 99.9 µA
Default value	2.0 µA

EP recognition

This parameter allows you to filter out only the equivalence points that are being sought.

Selection	first greatest last ascending descending
Default value	first

first

Only the first equivalence point will be recognized.

greatest

Only the equivalence point with the greatest ERC value, i.e. the steepest jump, will be recognized.

last

Only the last equivalence point will be recognized.

ascending

Only equivalence points with a positive slope of the titration curve will be recognized.

descending

Only equivalence points with a negative slope of the titration curve will be recognized.

Only one equivalence point will be recognized in a window. The numbering of the equivalence points (EP) is defined by the numbering of the windows (e.g. EP2 in window 2), so that even if EPs are missing, the calculations will still be carried out with the correctly assigned EP volumes.

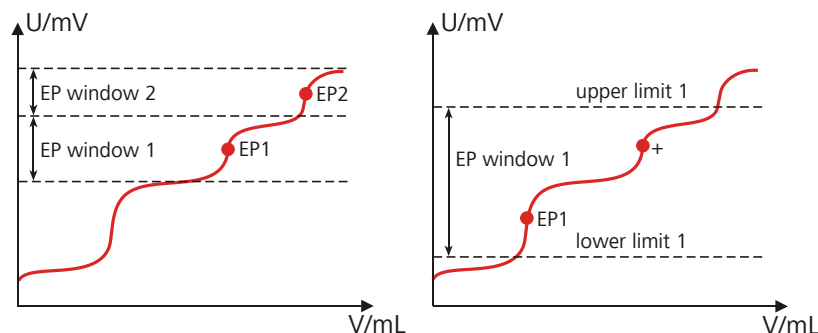


Figure 29 Equivalence point recognition and equivalence point numbering within windows

In the first example in the above figure, two equivalence points are recognized in two different windows (EP1 in window 1 and EP2 in window 2). In the second example, two equivalence points are found in one window, but only the first one is recognized. In order to ensure that the user recognizes that more than one equivalence point was found in the set window, EP1 is marked with a "+" in the result view. In addition, a corresponding message is entered in the message list.

Evaluation and equivalence point criterion with MET

The equivalence points (EPs) are localized by a method based on the Fortuin method which has been adapted by Metrohm for numerical methods. A search is made for the largest measured value change (Δ_n). The exact EP is determined by using an interpolation factor ρ which depends on the Δ values before and after Δ_n .

$$V_{EP} = V_0 + \rho \cdot \Delta V$$



- V_{EP} : EP volume
- V_0 : Dosed total volume before Δ_n
- ΔV : Volume increment
- ρ : Interpolation factor according to Fortuin

For the recognition of the EPs found, the set EP criterion is compared to the ERC (Equivalence point Recognition Criterion) found. The ERC is the sum of the measured value changes before and after the jump:

$$|\Delta_{n-2}| + |\Delta_{n-1}| + |\Delta_n| + |\Delta_{n+1}| + |\Delta_{n+2}|$$

In certain cases only three or only one summand is taken into account.

EPs whose ERC is smaller than the defined EP criterion will not be recognized. The ERC is displayed in the results dialog for each discovered and recognized EP. If you adjust the EP criterion later in order to recognize more or fewer EPs, then you can reevaluate the determination (**[Recalculate]** function in the results dialog).

29.2.5 Control instrument

The control instrument the determination is carried out with is selected under **[Control device]**. Control instruments are defined under **System ► Device manager**.



NOTICE

This button will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control instrument.

Control device

Selection of the control instrument from the list of devices. Only those devices are displayed which are able to carry out the command.

Selection	Selection of configured control instruments
-----------	--

29.2.6 Sensor

The parameters for the sensor are edited under **[Sensor]**.

Measuring input

Selection of the measuring input the sensor is connected to. The selection is not dependent on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Sensor

Selection of the sensor from the sensor list. The selection depends on the measuring mode. Sensors are defined under **System ► Sensors**. You can also enter a sensor name which is not contained in the sensor list. When a determination is started there is a check whether the sensor is contained in the sensor list.

Selection	Selection of configured sensors
-----------	---------------------------------

I(pol)

The polarization current is the current that is applied to a polarizable electrode during voltametric measurement. This parameter is available only with I(pol) determinations.

The values that are actually adjustable may vary from the values specified here. The binding values can be found in the manual, in the *Technical specifications* chapter.

Input range	-125.0 to 125.0 μA (Increment: 2.5)
Default value	5.0 μA
Selection	-1.0 μA 1.0 μA

U(pol)

The polarization voltage is the voltage applied to the polarizable electrode during an amperometric measurement. This parameter is available only with U(pol) determinations.

The values that are actually adjustable may vary from the values specified here. The binding values can be found in the manual, in the *Technical specifications* chapter.

Input range	-1,250 to 1,250 mV (Increment: 25)
Default value	400 mV

Electrode check

on | off (Default value: **off**)

For the following electrodes, an electrode check can be carried out:

- pH electrodes
- Metal electrodes
- Ion-selective electrodes

A check is made that the electrode is properly connected and that no short-circuit is present. The electrode check is carried out when this command is started. Note that this electrode check has nothing to do with the electrode test (command **ELT**).

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

continuous

A temperature sensor must be connected. The temperature is measured continuously.

automatic

If a temperature sensor is connected then the temperature will be measured continuously. Otherwise, the temperature entered manually will be used (see dialog of the titration and measuring parameters).

off

The temperature will not be measured. The temperature entered manually is used (see dialog of the titration and measuring parameters).

29.2.7 Dosing device

The parameters for the dosing device are edited under **[Dosing device]**.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the correct data (titer, concentration, etc.) is always used for the calculation. Titrants are defined under **Sys-tem ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without an integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place.

29.2.8 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 32.4, page 444*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

Switch off automatically

on | off (Default value: **on**)

If this parameter is activated, the stirrer will be switched off automatically at the end of the titration, measurement, etc.

29.3 Endpoint titrations (SET)

29.3.1 Start conditions

The parameters that are carried out before the start of titration are defined under **[Start conditions]**.

Pause 1

Waiting time, e.g. for the electrode to settle down before a start volume is added.

Input range	0 to 999999 s
Default value	0 s

Start volume

Volume that is dosed prior to the start of the titration.

Input range	0.00000 to 9999.99 mL
Default value	0.00000 mL

Dosing rate

Rate at which the start volume is dosed.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

Pause 2

Waiting time, e.g. for the electrode to settle down after the start or a reaction time after the dosing of a start volume.

Input range	0 to 999999 s
Default value	0 s

Initial measured value

The **[Initial meas. value]** is measured before carrying out the **start conditions**. This value is used for the **titration direction determination** when the titration direction is set to **auto** and only 1 EP is defined.

Signal drift

Maximum permissible drift for the measured value acceptance, i.e. maximum change of the measured value per minute. For this value, a suitable maximum waiting time is automatically calculated except you already have modified this waiting time.

Measuring mode pH, U and I_{pol}:

Input range	0.1 to 999.0 mV/min
Selection	off
Default value	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed.

Measuring mode U_{pol}:

Input range	0.01 to 99.90 µA/min
Selection	off
Default value	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed.

Min. waiting time

The measured value is not accepted until the minimum waiting time has elapsed, even if the signal drift has already been reached. The minimum waiting time is only important for drift-controlled measurements.

Input range	0 to 999999 s
Default value	0 s

Max. waiting time

If the signal drift has been switched off or has not yet been reached, then the measured value will be accepted when the maximum waiting time has elapsed. As default value, a suitable waiting time for the signal drift is automatically calculated.

Input range	0 to 999999 s
Default value	1 s

29.3.2 Control parameters

The control parameters for endpoint 1 and endpoint 2 are defined under **[Control parameters]**.

Endpoint 1 at

Measured value for the first endpoint.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.



Measuring mode pH:

Input range	-20.000 to 20.000
Selection	off
Default value	off

Measuring mode U and I_{pol}:

Input range	-2000.0 to 2000.0 mV
Selection	off
Default value	off

Measuring mode U_{pol}:

Input range	-200.00 to 200.00 µA
Selection	off
Default value	off

Titration rate

Three predefined sets of parameters can be selected for the titration rate.

Selection	slow optimal fast user
Default value	optimal

slow

For steep titration curves for which dosing must be carried out in small steps at the endpoint.

optimal

For all standard titrations. The parameters have been optimized for the most frequent applications.

fast

For flat titration curves for which the endpoint is reached only slowly.

user

The individual titration parameters can be modified.

The settings of the individual titration rates are listed in *table 11, page 328*.

User-defined parameters

These parameters are only accessible when **Titration rate = user**.

Dynamics

This parameter defines the control range before the specified endpoint. Individual volume steps are dosed in the control range; the dosing is finely controlled. The closer the endpoint, the slower the dosing until the dosing rate defined under **Min. rate** has been reached. The larger the control range, the slower the titration. Outside of the control range, dosing is carried out continuously; the dosing rate is defined under **Max. rate**.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measuring mode pH:

Input range	0.001 to 20.000
Default value	2.000
Selection	off

Measuring mode U and I_{pol}:

Input range	0.1 to 2000.0 mV
Default value	100.0 mV
Selection	off

Measuring mode U_{pol}:

Input range	0.01 to 200.00 µA
Default value	10.00 µA
Selection	off

Max. rate

Rate at which dosing is carried out outside of the control range.

Input range	0.01 to 166.00 mL/min
Default value	10.00 mL/min
Selection	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

Min. rate

Rate at which dosing is carried out at the very beginning of the titration and in the control range at the end of the titration. This parameter has a decisive influence on the titration rate and thus also on the accuracy. The smaller the selected minimum rate, the slower the titration.

Input range	0.01 to 9999.00 µL/min
Default value	25.00 µL/min

Table 11 Default values of the predefined titration rates for SET

	Titration rate		
	slow	optimal	fast
Dynamics			
– pH	5.000	2.000	0.500
– U und I _{pol}	300.0 mV	100.0 mV	30.0 mV
– U _{pol}	40.00 µA	10.00 µA	5.00 µA
Max. rate	1.00 mL/min	10.00 mL/min	maximum
Min. rate	5.00 µL/min	25.00 µL/min	50.00 µL/min

Stop criterion

Stop criterion

The titration is stopped when the endpoint has been reached and this stop criterion has been fulfilled. If no stop criterion has been selected then the titration will not be stopped. The stop conditions (*see chapter 29.3.4, page 330*) always lead to a stop, even if the stop criterion has not been reached.



NOTICE

In older instructions the delay time was usually defined as the stop criterion. The same delay time may however result in different stopping time points – because of different smallest possible volume increments (depending on the cylinder volume). In contrast, if the stop drift is used, the titration will always be stopped at the same curve slope dV/dt .

Selection	drift time off
Default value	drift

drift

The titration is stopped when the stop drift has been reached.

time

The titration is stopped when the endpoint has been exceeded during a certain period of time (**Delay time**).

off

The titration will not be stopped until the stop conditions have been fulfilled.

Stop drift

This parameter can only be edited with **Stop criterion = drift**.

The titration is stopped when the endpoint and the stop drift have been reached.

Input range	1 to 999 $\mu\text{L}/\text{min}$
Default value	20 $\mu\text{L}/\text{min}$

Delay time

This parameter can only be edited with **Stop criterion = time**.

When the endpoint has been reached, the specified time is allowed to elapse after the last dosing and the titration is then stopped.

Input range	0 to 999 s
Default value	10 s

Endpoint 2

The control parameters for the second endpoint are defined under **[Endpoint 2]**. The parameters and input ranges are identical with those for the first endpoint.

29.3.3 Titration parameters

Under **[Titration parameters]**, the parameters influencing the run of the entire titration are defined.

Titration direction

With the setting "auto", the titration direction is determined automatically from the initial measured value and the set endpoint. It is recommended that you specify whether the titration direction is positive or negative. If two endpoints have been set, then the titration direction will be defined automatically. In this case, the setting will be ignored.

Selection	+ - auto
Default value	auto

+

Positive measured value change by the addition of titrant, i.e. in the direction of a higher pH value, greater voltage or greater current.

-

Negative measured value change by the addition of titrant, i.e. in the direction of a lower pH value, lesser voltage or lesser current.

auto

The titration direction is determined automatically from the initial measured value and the set endpoint.

Extraction time

Minimum duration of the titration. The titration will not be stopped during the extraction time, even if the endpoint has already been reached. The titration is however stopped if a stop condition is fulfilled during this time



(see chapter 29.3.4, page 330). The entry of an extraction time may be advisable, e.g. for the titration of sparingly soluble samples.

Input range	0 to 999999 s
Default value	0 s

Temperature

Temperature entered manually. The temperature is being continuously measured when a temperature sensor is connected and when **Temp. measurement = automatic** or **continuous** is defined (see sensor dialog of the command). This value is used for temperature correction in pH measurements.

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

Time interval MP

Time interval for entering a measuring point in the measuring point list. The measuring point list is limited to 1000 measuring points.

Input range	0.1 to 999999.0 s
Default value	2.0 s

29.3.4 Stop conditions

The conditions for canceling the titration are defined under **[Stop conditions]**, if this does not occur automatically. This could be the case when the endpoint set is not reached or if the stop criterion (see "Stop criterion", page 328) is not fulfilled.

Stop volume

The titration is stopped when the specified volume has been dosed since the start of the titration. This volume should be adjusted to the size of the titration vessel in order to prevent the contents from running over.

Input range	0.00000 to 9999.99 mL
Default value	100,000 mL
Selection	off

Stop time

The titration is stopped when the specified time has elapsed following the termination of the start conditions.

Input range	1 to 999999 s
Selection	off
Default value	off

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum


The maximum filling rate depends on the cylinder volume (*see chapter 32, page 442*).

29.3.5 Conditioning

The conditions required for conditioning are defined under **[Conditioning]**.

Conditioning

on | off (Default value: **off**)

If this parameter is activated, then the first time the titration is started the working medium will be titrated to the endpoint with the specified control parameters. The status is kept stable. The actual method run does not begin until [] has been pressed once more.

Start drift

Conditioning OK will be displayed as soon as this drift has been reached and the titration can be started.

Input range	1 to 999 µL/min
Default value	20 µL/min

Drift correction

The endpoint volume can be corrected by taking the drift value into account. For this, the drift is multiplied with the drift correction time and this value is then subtracted from the endpoint volume. The drift correction time is the time interval between the end of conditioning and the end of the determination.

Selection	auto manual off
Default value	off

auto

The value of the current drift is automatically applied at the start of the titration.

manual

If the drift is known throughout a longer period of time, this can be entered manually.

off

No drift correction takes place.

Drift value

This parameter can only be edited with **Drift correction = manual**.

Drift for manual drift correction.

Input range	0.0 to 99.9 µL/min
Default value	0.0 µL/min


Measured value display

on | off (Default value: **off**)

If this parameter is activated, the currently measured value is displayed during the conditioning.

Conditioning options [Cond. options]

Cond. stop volume

Maximum permissible volume that can be dosed during conditioning. Conditioning is stopped when the specified volume is dosed. If conditioning is continued by pressing [] once again, then the titrant volume that has already been dosed will not be taken into account; i.e. the dosing starts again at zero. The stop volume should be adjusted to the size of the titration cell in order to prevent any overflow.

Input range	0.00000 to 9999.99 mL
Default value	20.0000 mL
Selection	off

Cond. stop time

Maximum permissible time over which conditioning may take place. Conditioning is stopped when the specified time has elapsed.

Input range	1 to 999999 s
Selection	off
Default value	off

Delay reconditioning

on | off (Default value: **off**)

If this parameter is activated, the conditioning will not start before all the method commands have been processed. Otherwise, the conditioning will immediately start after the titration command.



NOTICE

This parameter must be activated if commands follow which must not run simultaneously with the conditioning.

Example: Emptying the titration cell followed by adding new working medium.

29.3.6 Cell

The monitoring of the reagent used is activated or deactivated under **[Cell]**.

Reagent monitoring

Selection of the reagent from the list of reagents. The selection depends on the titration mode. Reagents are defined under **System ► Reagents**.

Selection	Selection of configured reagents off
Default value	off

off

The reagent monitoring is not active.

29.3.7 Control instrument

The control instrument the determination is carried out with is selected under **[Control device]**. Control instruments are defined under **System ► Device manager**.



NOTICE

This button will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control instrument.

Control device

Selection of the control instrument from the list of devices. Only those devices are displayed which are able to carry out the command.

Selection	Selection of configured control instruments
-----------	--

29.3.8 Sensor

The parameters for the sensor are edited under **[Sensor]**.

Measuring input

Selection of the measuring input the sensor is connected to. The selection is not dependent on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Sensor

Selection of the sensor from the sensor list. The selection depends on the measuring mode. Sensors are defined under **System ► Sensors**. You can also enter a sensor name which is not contained in the sensor list. When a



determination is started there is a check whether the sensor is contained in the sensor list.

Selection	Selection of configured sensors
-----------	--

I(pol)

The polarization current is the current that is applied to a polarizable electrode during voltametric measurement. This parameter is available only with I(pol) determinations.

The values that are actually adjustable may vary from the values specified here. The binding values can be found in the manual, in the *Technical specifications* chapter.

Input range	-125.0 to 125.0 μA (Increment: 2.5)
Default value	5.0 μA
Selection	-1.0 μA 1.0 μA

U(pol)

The polarization voltage is the voltage applied to the polarizable electrode during an amperometric measurement. This parameter is available only with U(pol) determinations.

The values that are actually adjustable may vary from the values specified here. The binding values can be found in the manual, in the *Technical specifications* chapter.

Input range	-1,250 to 1,250 mV (Increment: 25)
Default value	400 mV

Electrode check

on | off (Default value: **off**)

For the following electrodes, an electrode check can be carried out:

- pH electrodes
- Metal electrodes
- Ion-selective electrodes

A check is made that the electrode is properly connected and that no short-circuit is present. The electrode check is carried out when this command is started. Note that this electrode check has nothing to do with the electrode test (command **ELT**).

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

continuous

A temperature sensor must be connected. The temperature is measured continuously.

automatic

If a temperature sensor is connected then the temperature will be measured continuously. Otherwise, the temperature entered manually will be used (see dialog of the titration and measuring parameters).

off

The temperature will not be measured. The temperature entered manually is used (see dialog of the titration and measuring parameters).

29.3.9 Dosing device

The parameters for the dosing device are edited under **[Dosing device]**.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the correct data (titer, concentration, etc.) is always used for the calculation. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without an integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place.



29.3.10 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 32.4, page 444*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

Switch off automatically

on | off (Default value: **on**)

If this parameter is activated, the stirrer will be switched off automatically at the end of the titration, measurement, etc.

29.4 STAT titrations (STAT)

29.4.1 Start conditions

The parameters that are carried out before the start of titration are defined under **[Start conditions]**.

Pause 1

Waiting time, e.g. for the electrode to settle down before a start volume is added.

Input range	0 to 999999 s
Default value	0 s

Start volume

Volume that is dosed prior to the start of the titration.

Input range	0.00000 to 9999.99 mL
Default value	0.00000 mL

Dosing rate

Rate at which the start volume is dosed.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

Pause 2

Waiting time, e.g. for the electrode to settle down after the start or a reaction time after the dosing of a start volume.

Input range	0 to 999999 s
Default value	0 s

Initial measured value

The **[Initial meas. value]** is measured before carrying out the **start conditions**. This value is used for the **titration direction determination** when the titration direction is set to **auto**.

Signal drift

Maximum permissible drift for the measured value acceptance, i.e. maximum change of the measured value per minute. For this value, a suitable maximum waiting time is automatically calculated except you already have modified this waiting time.



Input range	0.1 to 999.0 mV/min
Selection	off
Default value	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed.

Min. waiting time

The measured value is not accepted until the minimum waiting time has elapsed, even if the signal drift has already been reached. The minimum waiting time is only important for drift-controlled measurements.

Input range	0 to 999999 s
Default value	0 s

Max. waiting time

If the signal drift has been switched off or has not yet been reached, then the measured value will be accepted when the maximum waiting time has elapsed. As default value, a suitable waiting time for the signal drift is automatically calculated.

Input range	0 to 999999 s
Default value	1 s

29.4.2 Control parameters

The parameters for the control point are defined under **[Control parameters]**.

Control point at

Measured value for the control point.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measuring mode pH:

Input range	-20.000 to 20.000
Selection	off
Default value	off

Measuring mode U:

Input range	-2000.0 to 2000.0 mV
Selection	off
Default value	off

Titration rate

Three predefined sets of parameters can be selected for the titration rate.

Selection	50 µL/min 100 µL/min 500 µL/min user
Default value	100 µL/min

50 µL/min

For reactions with an expected reaction rate in the range of 50 µL/min.

100 µL/min

For reactions with an expected reaction rate in the range of 100 µL/min.

500 µL/min

For reactions with an expected reaction rate in the range of 500 µL/min.

user

The individual titration parameters can be modified.

The settings of the individual titration rates are listed in *table 12, page 340*.

User-defined parameters

These parameters are only accessible when **Titration rate = user**.

Dynamics

The control range defines the measured value range before the specified control point. This control point has a vital influence on the titration rate and thus also on the measuring accuracy.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measuring mode pH:

Input range	0.001 to 20.000
Default value	1.000
Selection	off

Measuring mode U:

Input range	0.1 to 2,000.0 mV
Default value	60.0 mV
Selection	off

Max. rate

The maximum permissible rate at which dosing takes place.

Rule of thumb for the maximum rate in mL/min:

0.005 · expected reaction rate in µL/min

Input range	0.01 to 166.00 mL/min
Default value	0.75 mL/min
Selection	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

Min. rate

The rate at which dosing is carried out at the very beginning of the titration and in the control range at the end of the titration. This parameter has a decisive influence on the behavior at the control point. The minimum rate also has a decisive influence on how exactly the control point can be kept constant. The smaller the selected minimum rate, the longer it takes until the control point is reached for the first time.

The following rule of thumb should be considered in order to obtain a dosing at the control point as constant (regular) as possible:

Min. rate in $\mu\text{L}/\text{min} = 0.1 \cdot \text{expected reaction rate in } \mu\text{L}/\text{min}$

Input range	0.01 to 9999.00 $\mu\text{L}/\text{min}$
Default value	10.00 $\mu\text{L}/\text{min}$

Table 12 Default values of the predefined titration rates for STAT

	Reaction rate		
	50 $\mu\text{L}/\text{min}$	100 $\mu\text{L}/\text{min}$	500 $\mu\text{L}/\text{min}$
Dynamics			
– pH	1.000	1.000	1.000
– U	60.0 mV	60.0 mV	60.0 mV
Max. rate	0.25 mL/min	0.75 mL/min	2.00 mL/min
Min. rate	5.00 $\mu\text{L}/\text{min}$	10.00 $\mu\text{L}/\text{min}$	40.00 $\mu\text{L}/\text{min}$

Dosing and control range

Outside the control range the dosing rate is mainly characterized by the maximum rate. The parameters **Max. rate** and **Dynamics** should be optimized together in such a way that the titration does not overshoot too much when the control point is reached. The Dynamics should be selected in such a way that the measured value remains within the control range while the control point is maintained. A slightly larger control range should be defined for slow reactions (e.g. pH = 3, U = 180 mV). The set-point value is often set with a SET pretitration before the substrate is added. This means that the STAT titration can be started already with a small deviation from the control point.

29.4.3 Titration parameters

Under **[Titration parameters]**, the parameters influencing the run of the entire titration are defined.

Titration direction

With the setting "auto", the titration direction is determined automatically from the initial measured value and the set control point. It is recommended that you specify whether the change of the measured value is positive or negative.

Selection	+ - auto
Default value	auto

+

Positive measured value change by the addition of titrant, i.e. in the direction of a higher pH value or greater voltage.

-

Negative measured value change by the addition of titrant, i.e. in the direction of a lower pH value or lesser voltage.

auto

The titration direction is determined automatically from the initial measured value and the set control point.

Temperature

Temperature entered manually. The temperature is being continuously measured when a temperature sensor is connected and when **Temp. measurement = automatic** or **continuous** is defined (see sensor dialog of the command). This value is used for temperature correction in pH measurements.

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

Time interval MP

Time interval for entering a measuring point in the measuring point list. The measuring point list is limited to 1000 measuring points.

Input range	0.1 to 999999.0 s
Default value	2.0 s

Start time

The measured values are not entered in the measuring point list until this time has elapsed since the start of the titration.

Input range	0 to 999999 s
Default value	0 s

Start meas. value

The measured values are not entered in the measuring point list until this start measured value has been reached.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measuring mode pH:

Input range	-20.000 to 20.000
Selection	off
Default value	off

Measuring mode U:

Input range	-2000.0 to 2000.0 mV
Selection	off
Default value	off

Start rate

The measured values are not entered in the measuring point list until the dosing rate has fallen below this value. Checking this parameter is not started until 10 s after the start of the command.

Input range	0.01 to 166.00 mL/min
Selection	off
Default value	off



NOTICE

If several of the parameters **Start time**, **Start meas. value** or **Start rate** have been defined, then all of them must be fulfilled for measured values to be entered in the measuring point list.

29.4.4 Stop conditions

The conditions for canceling the titration are defined under **[Stop conditions]**.

Stop volume

The titration is stopped when the specified volume has been dosed since the start of the titration. This volume should be adjusted to the size of the titration vessel in order to prevent the contents from running over.

Input range	0.00000 to 9999.99 mL
Default value	100,000 mL
Selection	off

Stop time

The titration is stopped when the specified time has elapsed.

Input range	0 to 999999 s
Selection	off
Default value	off

Since

This parameter can only be edited when a stop time has been defined.

This parameter defines the start of the stop time.

Selection	Start Control point Last dosing
-----------	--

Start

The stop time starts after the start volume has been dosed and the following pause time has elapsed.

Control point

The stop time starts as soon as the control point has been reached for the first time.

Last dosing

The stop time starts after the last increment has been dosed, i. e. the stop time is reset to zero again after each dosing step.

Stop rate

The titration is stopped when the current titration rate is lower than the value entered. Checking this parameter is not started until 10 s after the start of the command.

Input range	0.01 to 166.00 µL/min
Selection	off
Default value	off

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 32, page 442*).

29.4.5 Monitoring

Under **[Monitoring]**, the parameters for monitoring the following quantities are defined:

- Measured value
- Dosing rate (an average rate)
- Temperature

Monitoring is not used for parameterization of the control behavior. Remote signals or RS-232 commands can be assigned to the monitored quantities (see "Communication", page 348). These remote signals or RS-232 commands are sent if a limit has been infringed. This can be used, for example, to switch a thermostat on/off if necessary.

The parameters are identical for all monitored quantities.

Monitoring

on | off (Default value: **off**)

Enabling and disabling the monitoring function. The monitoring of the dosing rate is not started until 10 s after the start of the titration.

Lower limit

When the monitored quantity falls below this value, the action defined in the following is triggered.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measuring mode pH:

Input range	-20.000 to 20.000
Default value	-20.000

Measuring mode U:

Input range	-2000.0 to 2000.0 mV
Default value	-2000.0 mV

Dosing rate:

Input range	0.00 to 166.00 mL/min
Default value	0.00 mL/min

Temperature:

Input range	-20.0 to 150.0 °C
Default value	-20.0 °C

Lower hysteresis

The lower hysteresis represents the tolerance range for the lower limit (see figure 32, page 348). When **Action = Wait** the dosing does not continue until the monitored quantity has once again exceeded the lower limit incl. the value entered here.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measuring mode pH:

Input range	0.000 to 20.000
Default value	0.020

Measuring mode U:

Input range	0.0 to 2000.0 mV
Default value	2.0 mV

Dosing rate:

Input range	0.00 to 166.00 mL/min
Default value	0.20 mL/min

Temperature:

Input range	0.0 to 150.0 °C
Default value	0.2 °C

Upper limit

When this value is exceeded, the action defined in the following is triggered.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measuring mode pH:

Input range	-20.000 to 20.000
Default value	20.000

Measuring mode U:

Input range	-2000.0 to 2000.0 mV
Default value	2000.0 mV

Dosing rate:

Input range	0.00 to 166.00 mL/min
Default value	166.00 mL/min

*Temperature:*

Input range	-20.0 to 150.0 °C
Default value	150.0 °C

Upper hysteresis

The upper hysteresis represents the tolerance range for the upper limit (see figure 32, page 348). When **Action = Wait**, the dosing does not continue until the monitored quantity has once again fallen below the upper limit incl. the value entered here.

The values that are actually measurable may vary from the values specified here. The binding measurement value ranges can be found in the manual, in the "Technical specifications" chapter.

Measuring mode pH:

Input range	0.000 to 20.000
Default value	0.020

Measuring mode U:

Input range	0.0 to 2000.0 mV
Default value	2.0 mV

Dosing rate:

Input range	0.00 to 166.00 mL/min
Default value	0.20 mL/min

Temperature:

Input range	0.0 to 150.0 °C
Default value	0.2 °C

Action

Selection of the action which is carried out when the limits of the monitored quantity are infringed.

Selection	None Exit method Cancel command Pause Wait
-----------	---

None

No action will be taken.

Exit method

The method is stopped immediately.

Cancel command

The current command is stopped und the next command in the method will be started.

Pause

Adding reagent is stopped until the dosing is continued manually with **[Continue]**. If several quantities are monitored, the other quantities will still be monitored in the background.

Wait

Adding reagent is stopped until the value of the monitored quantity is within the limits again (incl. hysteresis). The dosing is continued automatically. If several quantities are monitored, the other quantities will still be monitored in the background.

**NOTICE**

If the lower limit is infringed while monitoring the dosing rate, the settings **Action = Pause** and **Action = Wait** do not make sense because the average dosing rate becomes smaller and smaller during the waiting time. The valid range can never be reached again.

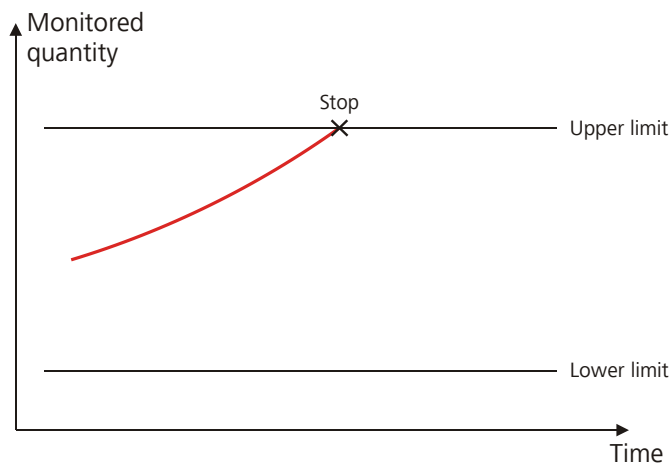


Figure 30 Action "Exit method" or "Cancel command"

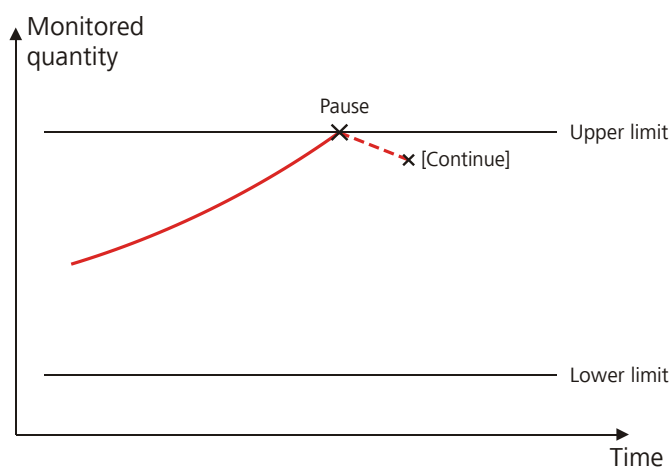


Figure 31 Action "Pause"

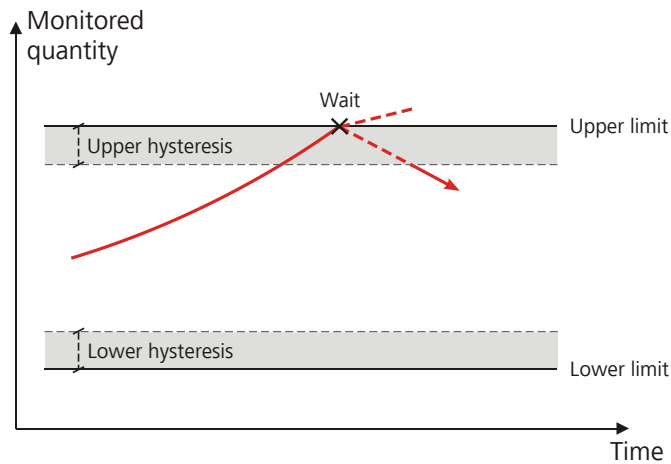


Figure 32 Action "Wait"

Communication

Under **[Communication]**, remote signals or RS-232 commands can be assigned to the monitored quantities. These remote signals or RS-232 commands are always sent, independent of the action defined for a limit infringement (see "Action", page 346).

[New]

Define a new assignment.

[Delete]

Delete the selected assignment.

[Edit]

Edit the settings for the selected assignment.

Monitored quantity

Selection of the monitored quantity to which a remote signal or an RS-232 command is to be assigned.

Selection	Measured value Temperature Dosing rate any
-----------	---

any

If several quantities are monitored, the remote signal or RS-232 command set is sent as soon as a limit of one of the monitored quantities is exceeded.

Infringed limit

Indicating the limit whose infringement causes the remote signal or RS-232 command defined in the following to be sent.

Selection	upper lower any back in range
-----------	--

upper

The remote signal or the RS-232 command is sent when the upper limit is exceeded.

lower

The remote signal or the RS-232 command is sent when the lower limit is exceeded.

any

The remote signal or the RS-232 command is sent when the upper or the lower limit is infringed.

back in range

The remote signal or the RS-232 command is sent when the value of the monitored quantity is within the limits again (incl. hysteresis).

Interface

Selection of the Remote Box or the serial interface via which the remote signal or the RS-232 command is sent.

Selection	Remote Box 1...4 COM1...8
-----------	------------------------------------

Output signal

This parameter can only be edited when a Remote Box is selected under **Interface**.

Selection of the signal out of the templates or entering the required bit pattern. Templates are defined under **System ► Templates ► Output lines**.



NOTICE

A line set active is not being reset automatically, nor at the end of the determination nor when the monitored quantity is in the range again.

Entering a bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status
- p = set pulse

The output lines are always numbered from right to left, i.e. with the signal *******1** line 0 is set. With a pulse, the length is set to 200 ms. If it is required to set pulses with other lengths, you have to define a corresponding template.

Entry	Bit patterns containing exactly 14 characters or a max. of 24 characters for the name of the template
Default value	*****

Selection **Selection of the templates defined**

Character string

This parameter can only be edited when a serial interface (COM) is selected under **Interface**.

Input of the RS-232 command as character string. All characters of the ASCII code page can be used. Control characters (Esc, FF etc.) have to be entered as a three-place, decimal ASCII code, starting with a slash. Each command is automatically terminated with the ASCII characters **CR** and **LF**.

Entry	ASCII string with a max. of 24 characters
Default value	empty

29.4.6 Control instrument

The control instrument the determination is carried out with is selected under **[Control device]**. Control instruments are defined under **System ► Device manager**.



NOTICE

This button will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control instrument.

Control device

Selection of the control instrument from the list of devices. Only those devices are displayed which are able to carry out the command.

Selection	Selection of configured control instruments
-----------	--

29.4.7 Sensor

The parameters for the sensor are edited under **[Sensor]**.

Measuring input

Selection of the measuring input the sensor is connected to. The selection is not dependent on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Sensor

Selection of the sensor from the sensor list. The selection depends on the measuring mode. Sensors are defined under **System ► Sensors**. You can also enter a sensor name which is not contained in the sensor list. When a

determination is started there is a check whether the sensor is contained in the sensor list.

Selection	Selection of configured sensors
-----------	--

Electrode check

on | off (Default value: **off**)

For the following electrodes, an electrode check can be carried out:

- pH electrodes
- Metal electrodes
- Ion-selective electrodes

A check is made that the electrode is properly connected and that no short-circuit is present. The electrode check is carried out when this command is started. Note that this electrode check has nothing to do with the electrode test (command **ELT**).

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

continuous

A temperature sensor must be connected. The temperature is measured continuously.

automatic

If a temperature sensor is connected then the temperature will be measured continuously. Otherwise, the temperature entered manually will be used (see dialog of the titration and measuring parameters).

off

The temperature will not be measured. The temperature entered manually is used (see dialog of the titration and measuring parameters).

29.4.8 Dosing device

The parameters for the dosing device are edited under **[Dosing device]**.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1



Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the correct data (titer, concentration, etc.) is always used for the calculation. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without an integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place.

Tandem dosing

A second dosing device can be defined under **[Tandem dosing]** in order to enable uninterrupted dosing. Dosing is carried out with a combination of two dosing devices so that the second dosing device is dosing while the first one is being filled and vice versa.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	off

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without integrated data chip, the cylinder volume and the type of dosing drive are checked. For

the selected titrant, only the time interval for the GLP test of the buret unit is checked at the start of the determination.



NOTICE

The concentration, validity of the titer and the working life of the titrant are ignored.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place.

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 32, page 442*).

However, in order to ensure uninterrupted dosing, the following points must be taken into account:

- Use the highest possible filling rate in order to keep the filling times as short as possible. The filling rate should be decreased for viscous liquids.
- When you use two buret units with different dosing cylinders, the filling rate for the larger sized cylinder must be at the minimum:

$$v_{2,Fill} \geq v_{1,Fill} \cdot \frac{V_{Cyl.2}}{V_{Cyl.1}}$$

$v_{2,Fill}$ = filling rate in mL/min for the larger sized cylinder

$v_{1,Fill}$ = filling rate in mL/min for the smaller cylinder

$V_{Cyl.2}$ = cylinder volume in mL of the buret unit of the second dosing device

$V_{Cyl.1}$ = cylinder volume in mL of the buret unit of the first dosing device

Example:

Dosing device 1: volume = 20 mL, filling rate = 50 mL/min

Dosing device 2: volume = 50 mL



Filling rate $2 \geq 50 \text{ mL/min} \cdot 50 \text{ mL} / 20 \text{ mL} \geq 125 \text{ mL/min}$

- The dosing rate must not exceed 75% of the filling rate of the smaller cylinder. These values are listed in the following table, valid at maximum filling rate:

Table 13 Maximum dosing rate for different dosing cylinders

Cylinder volume	maximum dosing rate	
	Exchange unit	Dosing unit
1 mL	2.25 mL/min	–
2 mL	–	5.00 mL/min
5 mL	11.25 mL/min	12.50 mL/min
10 mL	22.50 mL/min	25.00 mL/min
20 mL	45.00 mL/min	50.00 mL/min
50 mL	112.50 mL/min	124.50 mL/min

29.4.9 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 32.4, page 444*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

Switch off automatically

on | off (Default value: **on**)

If this parameter is activated, the stirrer will be switched off automatically at the end of the titration, measurement, etc.

29.5 Manual titrations (MAT)


A manual titration can be carried out with a Manual Dosing Controller or with the buttons **[Dose]** and **[Cancel command]**.

In order for the manual titration to be carried out with the 6.2107.120 Manual Dosing Controller, the latter must be connected to the MSB connector of the 916 Ti-Touch via the 6.2148.010 Remote Box MSB.

Manual titration using the Manual Dosing Controller

- 1 Press the **[GO]** key to start the method.
- 2 Press the **[GO]** key during the **MAT** command for dosing.
- 3 End the MAT command with the **[FILL]** key.

Manual titration using the buttons

- 1 Press the  key to start the method.
- 2 Press the **[Dose]** button during the MAT command for dosing.
- 3 End the MAT command with the **[Cancel command]** button.



NOTICE

Start of titration: Because of the technical implementation involved, a few microliters are already dosed at the start of a manual titration. The possibility of deviation from the actual endpoint volume is excluded, however. The already dosed volume is taken into account.

End of titration: A manual titration must always be ended with the **[Cancel command]** button (NOT with the **[]** fixed key) or with the **[FILL]** key on the Manual Dosing Controller. This is the only way to ensure that subsequent method commands will be carried out.

29.5.1 Dosing parameters

Under **[Dosing parameters]**, the parameters influencing the run of the manual titration are defined.

Dosing ramp

The dosing ramp is a delay in the dosing rate at the time of the start of a dosing step. It is advantageous primarily for manual titrations with indicator, when small volume steps are to be dosed before the equivalence point. The selectable delay time determines after which amount of time the defined dosing rate is to be reached.

Input range	0 to 10 s
Default value	2 s

Dosing rate

The rate at which dosing takes place.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

Time interval MP

Time interval for entering a measuring point in the measuring point list. The measuring point list is limited to 1000 measuring points.

Input range	0.1 to 999999.0 s
Default value	5.0 s

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 32, page 442*).

Temperature

Temperature entered manually. The temperature is being continuously measured when a temperature sensor is connected and when **Temp. measurement = automatic** or **continuous** is defined (see sensor dialog

of the command). This value is used for temperature correction in pH measurements.

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

29.5.2 Control instrument

The control instrument the determination is carried out with is selected under **[Control device]**. Control instruments are defined under **System ► Device manager**.



NOTICE

This button will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control instrument.

Control device

Selection of the control instrument from the list of devices. Only those devices are displayed which are able to carry out the command.

Selection	Selection of configured control instruments
-----------	--

29.5.3 Sensor

The parameters for the sensor are edited under **[Sensor]**.

Measuring input

Selection of the measuring input the sensor is connected to. The selection is not dependent on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Sensor

Selection of the sensor from the sensor list. The selection depends on the measuring mode. Sensors are defined under **System ► Sensors**. You can also enter a sensor name which is not contained in the sensor list. When a determination is started there is a check whether the sensor is contained in the sensor list.

Selection	Selection of configured sensors
-----------	--

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

continuous

A temperature sensor must be connected. The temperature is measured continuously.

automatic

If a temperature sensor is connected then the temperature will be measured continuously. Otherwise, the temperature entered manually will be used (see dialog of the titration and measuring parameters).

off

The temperature will not be measured. The temperature entered manually is used (see dialog of the titration and measuring parameters).

29.5.4 Dosing device

The parameters for the dosing device are edited under **[Dosing device]**.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the correct data (titer, concentration, etc.) is always used for the calculation. Titrants are defined under **Sys-tem ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without an integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place.

Tandem dosing

A second dosing device can be defined under **[Tandem dosing]** in order to enable uninterrupted dosing. Dosing is carried out with a combination of two dosing devices so that the second dosing device is dosing while the first one is being filled and vice versa.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	off

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, only the time interval for the GLP test of the buret unit is checked at the start of the determination.



NOTICE

The concentration, validity of the titer and the working life of the titrant are ignored.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place.

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

**maximum**

The maximum filling rate depends on the cylinder volume (*see chapter 32, page 442*).

However, in order to ensure uninterrupted dosing, the following points must be taken into account:

- Use the highest possible filling rate in order to keep the filling times as short as possible. The filling rate should be decreased for viscous liquids.
- When you use two buret units with different dosing cylinders, the filling rate for the larger sized cylinder must be at the minimum:

$$v_{2,Fill} \geq v_{1,Fill} \cdot \frac{V_{Cyl.2}}{V_{Cyl.1}}$$

$v_{2,Fill}$ = filling rate in mL/min for the larger sized cylinder

$v_{1,Fill}$ = filling rate in mL/min for the smaller cylinder

$V_{Cyl.2}$ = cylinder volume in mL of the buret unit of the second dosing device

$V_{Cyl.1}$ = cylinder volume in mL of the buret unit of the first dosing device

Example:

Dosing device 1: volume = 20 mL, filling rate = 50 mL/min

Dosing device 2: volume = 50 mL

Filling rate 2 \geq 50 mL/min \cdot 50 mL / 20 mL \geq 125 mL/min

- The dosing rate must not exceed 75% of the filling rate of the smaller cylinder. These values are listed in the following table, valid at maximum filling rate:

Table 14 Maximum dosing rate for different dosing cylinders

Cylinder volume	maximum dosing rate	
	Exchange unit	Dosing unit
1 mL	2.25 mL/min	–
2 mL	–	5.00 mL/min
5 mL	11.25 mL/min	12.50 mL/min
10 mL	22.50 mL/min	25.00 mL/min
20 mL	45.00 mL/min	50.00 mL/min
50 mL	112.50 mL/min	124.50 mL/min

29.5.5 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 32.4, page 444*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

Switch off automatically

on | off (Default value: **on**)

If this parameter is activated, the stirrer will be switched off automatically at the end of the titration, measurement, etc.

29.6 Measurements (MEAS)

29.6.1 Measuring parameters

Under **[Measuring parameters]**, the parameters influencing the run of the entire measurement are defined.

Signal drift

Maximum permissible drift for the measured value acceptance, i.e. maximum change of the measured value per minute. For this value, a suitable maximum waiting time is automatically calculated except you already have modified this waiting time.



NOTICE

A constant measured value is often only reached after a certain time, as mixing and the reaction itself require a certain time. The response time of an electrode can also increase with time, i.e., reaching a constant measured value takes longer and longer. Drift-controlled measured value acceptance is particularly advisable in such cases, as the measured values are only accepted when equilibrium has almost been reached.

Measuring mode pH, U and I_{pol}:

Input range	0.1 to 999.0 mV/min
Default value	10.0 mV/min
Selection	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed.

Measuring mode U_{pol}:

Input range	0.01 to 99.90 µA/min
Default value	10.00 µA/min
Selection	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed.

Measuring mode T:

Input range	0.1 to 999.0 °C/min
Default value	0.5 °C/min
Selection	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed.

Min. waiting time

The measured value is not accepted until the minimum waiting time has elapsed, even if the signal drift has already been reached. The minimum waiting time is only important for drift-controlled measurements.

Input range	0 to 999999 s
Default value	0 s

Max. waiting time

If the signal drift has been switched off or has not yet been reached, then the measured value will be accepted when the maximum waiting time has elapsed. As default value, a suitable waiting time for the signal drift is automatically calculated.

Input range	0 to 999999 s The default value depends on the measuring mode.
Selection	off

off

The measurement is continued endlessly.

Stop meas. value

The measurement is canceled when the specified measured value has been reached since the start of the measurement.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the .

Measuring mode pH:

Input range	-20.000 to 20.000
Selection	off
Default value	off

Measuring mode U, Ipol:

Input range	-2000.0 to 2000.0 mV
Selection	off
Default value	off

Measuring mode Upol:

Input range	-200.00 to 200.00 µA
Selection	off
Default value	off



Measuring mode T:

Input range	-20.0 to 150.0 °C
Selection	off
Default value	off

Temperature

This parameter is not available with the **MEAS T** command (temperature measurement).

Temperature entered manually. The temperature is being continuously measured when a temperature sensor is connected and when **Temp. measurement = automatic** or **continuous** is defined (see sensor dialog of the command). This value is used for temperature correction in pH measurements.

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

Time interval MP

Time interval for entering a measuring point in the measuring point list. The measuring point list is limited to 1000 measuring points.

Input range	0.1 to 999999.0 s
Default value	2.0 s

29.6.2 Control instrument

The control instrument the determination is carried out with is selected under **[Control device]**. Control instruments are defined under **System ► Device manager**.



NOTICE

This button will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control instrument.

Control device

Selection of the control instrument from the list of devices. Only those devices are displayed which are able to carry out the command.

Selection	Selection of configured control instruments
-----------	--

29.6.3 Sensor

The parameters for the sensor are edited under **[Sensor]**.

The parameters are valid for all measuring modes except for **Temp. measurement**. This parameter is not available with the **MEAS T** command (temperature measurement).

Measuring input

Selection of the measuring input the sensor is connected to. The selection is not dependent on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Sensor

Selection of the sensor from the sensor list. The selection depends on the measuring mode. Sensors are defined under **System ► Sensors**. You can also enter a sensor name which is not contained in the sensor list. When a determination is started there is a check whether the sensor is contained in the sensor list.

Selection	Selection of configured sensors
-----------	--

I(pol)

The polarization current is the current that is applied to a polarizable electrode during voltametric measurement. This parameter is available only with I(pol) determinations.

The values that are actually adjustable may vary from the values specified here. The binding values can be found in the manual, in the *Technical specifications* chapter.

Input range	-125.0 to 125.0 μA (Increment: 2.5)
Default value	5.0 μA
Selection	-1.0 μA 1.0 μA

U(pol)

The polarization voltage is the voltage applied to the polarizable electrode during an amperometric measurement. This parameter is available only with U(pol) determinations.

The values that are actually adjustable may vary from the values specified here. The binding values can be found in the manual, in the *Technical specifications* chapter.

Input range	-1,250 to 1,250 mV (Increment: 25)
Default value	400 mV



Electrode check

on | off (Default value: **off**)

For the following electrodes, an electrode check can be carried out:

- pH electrodes
- Metal electrodes
- Ion-selective electrodes

A check is made that the electrode is properly connected and that no short-circuit is present. The electrode check is carried out when this command is started. Note that this electrode check has nothing to do with the electrode test (command **ELT**).

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

continuous

A temperature sensor must be connected. The temperature is measured continuously.

automatic

If a temperature sensor is connected then the temperature will be measured continuously. Otherwise, the temperature entered manually will be used (see dialog of the titration and measuring parameters).

off

The temperature will not be measured. The temperature entered manually is used (see dialog of the titration and measuring parameters).

29.6.4 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rota-

tional speed is specified in *chapter 32.4, page 444*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

Switch off automatically

on | off (Default value: **on**)

If this parameter is activated, the stirrer will be switched off automatically at the end of the titration, measurement, etc.

29.7 Calibration of sensors (CAL)

29.7.1 Calibration parameters (CAL pH)

Under **[Calibration parameters]**, the parameters influencing the run of the entire calibration are defined.

Signal drift

Maximum permissible drift for the measured value acceptance, i.e. maximum change of the measured value per minute. For this value, a suitable maximum waiting time is automatically calculated except you already have modified this waiting time.

Input range	0.1 to 999.0 mV/min
Default value	2.0 mV/min
Selection	off

off

Measured value acceptance will take place after the maximum waiting time has elapsed.

Min. waiting time

The measured value is not accepted until the minimum waiting time has elapsed, even if the signal drift has already been reached. The minimum waiting time is only important for drift-controlled measurements.

Input range	0 to 999999 s
Default value	10 s

Max. waiting time

If the signal drift has been switched off or has not yet been reached, then the measured value will be accepted when the maximum waiting time has



elapsed. As default value, a suitable waiting time for the signal drift is automatically calculated.

Input range	0 to 999999 s
Default value	110 s
Selection	off

Temperature

Temperature entered manually. The temperature is being continuously measured when a temperature sensor is connected and when **Temp. measurement = automatic** or **continuous** is defined (see sensor dialog of the command).

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

Sample Processor

If the calibration is carried out with a Sample Processor, the buffers or standards are automatically changed. This is why no request for the calibration temperature appears when starting the calibration. The value entered above is used.

Selection	off Remote USB
Default value	off

off

For calibrations without Sample Processor. During the calibration, a request for changing the buffer/standard is displayed.

Remote

For automatic calibrations with Sample Processors, which are connected via the Remote Box.

USB

For automatic calibrations with a Sample Processor that is connected via USB(see "Definition of the calibration positions", page 451).

Buffers

The buffer type and the number of buffers is defined under **[Buffers]**.

Buffer type

Selection of a predefined buffer series or definition of special buffers. In the case of predefined buffer series, the instrument automatically recognizes which buffer is involved.

Selection	Metrohm NIST DIN Fisher Fluka Basel Mettler Merck Tit. Beckman Radiometer Custom Special Merck Cer. Baker Hamilton Precisa
Default value	Metrohm

Custom

Definition of a custom buffer series. The buffer series is defined under **System ► Templates ► Custom calib. buffers.**

Special

Up to five calibration buffers can be defined in the method. The automatic buffer detection is not activated in this case. The buffers must be measured precisely in the specified sequence.

Merck Cer.

Reference temperature = 25 °C. When using Merck CertiPUR buffers (20 °C) the buffer type **Merck Tit.** must be selected.

Number of buffers

This parameter is not visible when **Buffer type = Special** is defined.

Number of buffers that are used for calibration. If calibration is accomplished with more than two buffers, then they can be used repeatedly in order to give them more statistical weight. The first two buffers must however always be different from one another.

Selection	1 2 3 4 5
Default value	2

Buffer 1 pH

This parameter is visible only when **Buffer type = Special**.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the .

Input range	-20.000 to 20.000
Default value	7.000

Buffer 2 pH

This parameter is visible only when **Buffer type = Special**.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the .

Input range	-20.000 to 20.000
Default value	4.000
Selection	off

Buffer 3 pH

This parameter is visible only when **Buffer type = Special**.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the .

Input range	-20.000 to 20.000
Selection	off
Default value	off

Buffer 4 pH

See **Buffer 3 pH**.

Buffer 5 pH

See **Buffer 3 pH**.

Stirrer control

The parameters for the stirrer control during the calibration are edited under **[Stirrer control]**. Ensure that a stirrer has been selected in the **Edit command / Stirrer** dialog.

**NOTICE**

Deactivate the **Stir solution during measurement** parameter when you carry out the calibration with a Sample Processor. The stirrer control for the tower stirrer is ineffective with these instruments.

Stir solution during measurement

on | off (Default value: **on**)

If this parameter is activated, the stirrer will be switched on and off automatically at the start or end of the measurement.

Stir before meas.

This parameter can only be edited when the parameter **Stir solution during measurement** is deactivated.

If the stirrer is switched off during the measurement, it will be stirred for the time entered before measuring.

Input range	0 to 999999 s
Default value	0 s

Pause before meas.

This parameter can only be edited when the parameter **Stir solution during measurement** is deactivated.

If it is stirred before the measurement, a waiting time can be defined here before starting the measurement. No stirring or measuring occurs during this time.

Input range	0 to 999999 s
Default value	0 s

29.7.2 Control instrument

The control instrument the determination is carried out with is selected under **[Control device]**. Control instruments are defined under **System ► Device manager**.



NOTICE

This button will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control instrument.

Control device

Selection of the control instrument from the list of devices. Only those devices are displayed which are able to carry out the command.

Selection	Selection of configured control instruments
-----------	--

29.7.3 Sensor

The parameters for the sensor are edited under **[Sensor]**.

Measuring input

Selection of the measuring input the sensor is connected to. The selection is not dependent on whether the control device has one or two measuring interfaces.

Selection	1 2
Default value	1

Sensor

Selection of the sensor from the sensor list. The selection depends on the measuring mode. Sensors are defined under **System ► Sensors**. You can also enter a sensor name which is not contained in the sensor list. When a determination is started there is a check whether the sensor is contained in the sensor list.

Selection	Selection of configured sensors
-----------	--

Electrode check

on | off (Default value: **off**)

For pH electrodes, an electrode check can be carried out. A check is made that the electrode is properly connected and that no short-circuit is present. The electrode check is carried out when this command is started. Note that this electrode check has nothing to do with the electrode test (command **ELT**).

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

continuous

A temperature sensor must be connected. The temperature is measured continuously.

automatic

If a temperature sensor is connected then the temperature will be measured continuously. Otherwise, the calibration temperature entered manually will be used.

off

The temperature will not be measured. The calibration temperature entered manually is used. The temperature is requested after the start, except the determination is carried out with a Sample Processor.

29.7.4 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to $+15$. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 32.4, page 444*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

29.8 Electrode test for pH electrodes (ELT)

An automatic electrode test can be carried out for pH electrodes which will permit a qualification of the electrode used. Differentiated specifications regarding characteristic measurement properties of your pH electrode (slope, response time, etc.) will be obtained as the result. The concluding evaluation of the electrode is carried out by testing these results against the specified limit values. Once the electrode test has been successfully completed, the calibration data is stored with the sensor data of the sensor used.

These limit values are saved in the software for the following types of pH electrodes:

- Standard electrodes
- Gel electrodes
- Electrodes with a nonaqueous reference electrolyte

You can, however, also define limits of your own choosing (*see chapter 15.7, page 178*). The precise classification of the measuring chain is determined by the reference system used. You will find an overview of these electrode types and the corresponding limit values in *table 15, page 377*.

Faulty electrodes often exhibit certain combinations of measurement results of the electrode test which fall outside the intended limits. This makes it possible to have more precise data concerning possible causes of errors and recommended remedial measures.

Buffers

You will require flawless buffers in the pH values 4, 7 and 9 for carrying out the electrode test. We recommend for this purpose the appropriate Metrohm buffers. The electrode test can, however, in principle also be carried out with any of the buffer sets contained in the memory.

The buffers must be measured in the following sequence:

1. pH 9
2. pH 4
3. pH 7



NOTICE

Ensure in particular the flawlessness of the pH 9 buffer. It can readily deviate from its specified pH value as the result of absorption of CO₂ from the ambient atmosphere, thus leading to incorrect test results.



Stirrer

A stirrer has to be connected to the titrator for the electrode test.

Temperature

Take care to ensure that the entire electrode test is carried out at a consistently uniform temperature. Work should be accomplished at room temperature wherever possible, in light of the fact that temperature exercises a great influence on response time. The preset limit values are calibrated to 25 °C.

29.8.1 Electrode test parameters

Under **[Electrode test param.]**, the parameters influencing the sequence of the electrode test are defined.

Buffer type

Selecting a predefined buffer series.

Selection	Metrohm NIST DIN Fluka Basel Mettler Merck Tit. Radiometer Merck Cer. Baker Hamilton Precisa
Default value	Metrohm

Merck Cer.

Reference temperature = 25 °C. When using Merck CertiPUR buffers (20 °C) the buffer type **Merck Tit.** must be selected.

Sample Processor

If the calibration is carried out with a Sample Processor, the buffers or standards are automatically changed. This is why no request for the calibration temperature appears when starting the calibration. The value entered above is used.

Selection	off Remote USB
Default value	off

off

For calibrations without Sample Processor. During the calibration, a request for changing the buffer/standard is displayed.

Remote

For automatic calibrations with Sample Processors, which are connected via the Remote Box.

USB

For automatic calibrations with a Sample Processor that is connected via USB(see *"Definition of the calibration positions", page 451*).

Electrode type

Selection of the electrode type.

Selection	Standard Gel Non-aqueous Custom
Default value	Standard

Standard

Electrodes which contain e.g. KCl solution as reference electrolyte.

Gel

Electrodes which contain Idrolyte as reference electrolyte.

Non-aqueous

Electrodes, which contain a non-aqueous reference electrolyte, e.g. TEABr in ethylene glycol or LiCl in ethanol.

Custom

Under **System ► Templates** a custom electrode type can be defined.

Temperature

Temperature entered manually. The temperature is being continuously measured when a temperature sensor is connected and when **Temp. measurement = automatic** or **continuous** is defined (see sensor dialog of the command).

Input range	-20.0 to 150.0 °C
Default value	25.0 °C

29.8.2 Control instrument

The control instrument the determination is carried out with is selected under **[Control device]**. Control instruments are defined under **System ► Device manager**.

**NOTICE**

This button will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control instrument.

Control device

Selection of the control instrument from the list of devices. Only those devices are displayed which are able to carry out the command.

Selection	Selection of configured control instruments
-----------	--

29.8.3 Sensor

The parameters for the sensor are edited under **[Sensor]**.

Measuring input

Selection of the measuring input the sensor is connected to. The selection is not dependent on whether the control device has one or two measuring interfaces.



Selection	1 2
Default value	1

Sensor

Selection of the sensor from the sensor list. The selection depends on the measuring mode. Sensors are defined under **System ► Sensors**. You can also enter a sensor name which is not contained in the sensor list. When a determination is started there is a check whether the sensor is contained in the sensor list.

Selection	Selection of configured sensors
-----------	--

Temp. measurement

Type of temperature measurement.

Selection	continuous automatic off
Default value	automatic

continuous

A temperature sensor must be connected. The temperature is measured continuously.

automatic

If a temperature sensor is connected then the temperature will be measured continuously. Otherwise, the calibration temperature entered manually will be used.

off

The temperature will not be measured. The calibration temperature entered manually is used. The temperature is requested after the start, except the determination is carried out with a Sample Processor.

29.8.4 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Stirring rate

Setting the stirring rate. It can be set in steps of –15 to +15. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 32.4, page 444*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

29.8.5 Additional information

Table 15 Limit values of the three electrode types

	Rating		
	Excellent electrode	Good electrode	Usable electrode
Electrode type "Standard"			
Streaming potential	2.5 mV	3.0 mV	4.0 mV
Drift	2.0 mV/s	2.5 mV/s	3.0 mV/s
Min. slope	96.5%	96.0%	95.0%
Max. slope	101.0%	102.0%	103.0%
Response time	45 s	50 s	60 s
Lower limit Uoff	-15 mV		
Upper limit Uoff	15 mV		
Electrode type "Gel"			
Streaming potential	3.0 mV	3.5 mV	4.5 mV
Drift	2.5 mV/s	3.0 mV/s	4.0 mV/s
Min. slope	96.5%	96.0%	95.0%
Max. slope	101.0%	102.0%	103.0%
Response time	60 s	75 s	90 s
Lower limit Uoff	-20 mV		
Upper limit Uoff	20 mV		
Electrode type "Non-aqueous"			
Streaming potential	3.0 mV	4.5 mV	6.0 mV
Drift	5.0 mV/s	7.0 mV/s	9.0 mV/s
Min. slope	88.0%	80.0%	70.0%
Max. slope	120.0%	130.0%	140.0%
Response time	60 s	75 s	90 s
Lower limit Uoff	-50 mV		
Upper limit Uoff	100 mV		



Table 16 Possible errors during the electrode test

Test criterion	Message	Remedy
<p>Buffer pH 9:</p> <p>The drift in stirred solution is > 1 mV.</p>	General problem	<ul style="list-style-type: none"> ▪ Connect an electrode. ▪ Replace the defective electrode cable. ▪ Clean the diaphragm (see leaflet for the electrode). ▪ Replace the electrode.
<p>Buffer pH 9:</p> <p>-10 mV < U(not stirred) < 10 mV</p> <p>and</p> <p>The sum of the drift values after 1, 2, 3 and 4 min is < 12 mV/min.</p>	Short circuit	Replace the electrode.
<p>A response time does not fulfill the limit value for the rating Usable electrode.</p>	Glass membrane / Diaphragm	Clean the diaphragm (see leaflet for the electrode).
<p>The pH values are not defined for all the buffers at the temperature measured.</p>	No buffer data	Repeat the electrode test at a temperature at which the pH values of all the buffers are defined.
<p>All the slopes fulfill the limit values for the rating Usable electrode.</p> <p>and</p> <p>U_{off} is outside the preset limits.</p>	Unsuitable reference electrode	<ul style="list-style-type: none"> ▪ Repeat the electrode test with a suitable reference system. ▪ Replace the contaminated reference electrolyte. ▪ Select Electrode type = Custom and adjust the limits for U_{off}.
<p>The streaming potential is too high.</p>	Contaminated diaphragm	Clean the diaphragm (see leaflet for the electrode).
<p>Two slopes do not fulfill the limit values for the rating Usable electrode.</p>	Wrong buffer	Repeat the electrode test with buffers pH 4, 7 and 9.
<p>All the slopes do not fulfill the limit values for the rating Usable electrode.</p>	Partial short circuit	<ul style="list-style-type: none"> ▪ Check the temperature sensor. ▪ Enter the correct calibration temperature.

29.9 Evaluations (EVAL)

Various **additional evaluations** (EVAL commands) can be carried out for titrations and measurements. The evaluation always refers to the last titration or measurement prior to the EVAL command. Only those evaluation commands which are available for the last titration or measurement prior to the EVAL command can be inserted in the list of commands. If a titration or measuring command is deleted before the EVAL command, then it will be shown in red in the list of commands, as the reference is missing.

EVAL commands can also be inserted in the method run at a later time and the determination can be reevaluated (**[Recalculate]** function in the results dialog).

Additional evaluations for DET titrations

The following additional evaluations are possible:

- **EVAL FIX-EP** (fixed endpoints)
Measured quantities:
 - Measured value
 - Time
 - Volume
- **EVAL pK/HNP** (pK value / half neutralization potential HNP)
This evaluation is only possible for the measuring modes **pH** and **U**.
- **EVAL MIN/MAX** (minimum/maximum)
- **EVAL BREAK** (break points)

Additional evaluations for MET titrations

The following additional evaluations are possible:

- **EVAL FIX-EP** (fixed endpoints)
Measured quantities:
 - Measured value
 - Time
 - Volume
- **EVAL pK/HNP** (pK value / half neutralization potential HNP)
This evaluation is only possible for the measuring modes **pH** and **U**.
- **EVAL MIN/MAX** (minimum/maximum)
- **EVAL BREAK** (break points)

Additional evaluations for SET titrations

The following additional evaluations are possible:



- **EVAL FIX-EP** (fixed endpoints)
Measured quantities:
 - Measured value
 - Time
 - Volume
- **EVAL MIN/MAX** (minimum/maximum)
- **EVAL RATE** (average dosing rate)

Additional evaluations for STAT titrations

The following additional evaluations are possible:

- **EVAL FIX-EP** (fixed endpoints)
Measured quantities:
 - Measured value
 - Time
 - Volume
- **EVAL MIN/MAX** (minimum/maximum)
- **EVAL RATE** (average dosing rate)

Additional evaluations for measurements

The following additional evaluations are possible:

- **EVAL FIX-EP** (fixed endpoints)
Measured quantities:
 - Measured value
 - Time
- **EVAL MIN/MAX** (minimum/maximum)
- **EVAL BREAK** (break points)

29.9.1 Fixed endpoint evaluation (EVAL FIX-EP)

For the fixed endpoint evaluation, the associated values are interpolated from the measuring point list for one quantity (measured value, volume, etc.). Up to nine fixed endpoints can be evaluated with one command.

Fixed quantity

Selection of the quantity to which the associated value is interpolated from the measuring point list.

Selection	Measured value Time Volume
Default value	Measured value

Fixed EP1 at

The fixed endpoint must lie between the first and the final entry in the measuring point list.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the .

Measured value (measuring mode pH):

Input range	-20.000 to 20.000
Selection	off
Default value	off

Measured value (measuring mode U, Ipol):

Input range	-2,000.0 to 2,000.0 mV
Selection	off
Default value	off

Measured value (measuring mode Upol):

Input range	-200.00 to 200.00 µA
Selection	off
Default value	off

Measured value (measuring mode T):

Input range	-20.0 to 150.0 °C
Selection	off
Default value	off

Time:

Input range	0.0 to 999999.9 s
Selection	off
Default value	off

Volume:

Input range	0.00000 to 9,999.99 mL
Selection	off
Default value	off

Fixed EP2 at...Fixed EP9 at

See **Fixed EP1 at**.

29.9.2 **pK value and half neutralization potential evaluation (EVAL pK/HNP)**

The pK value (pH measuring mode) or the half neutralization potential (U measuring mode) can be determined for DET and MET titrations.

The activities of conjugated acid-base pairs are linked by the so-called Henderson-Hasselbalch equation:

$$\text{pH} = \text{pK}_a + \log (a_B/a_A)$$

If the activities of the acid and the conjugated base are equal ($a_A = a_B$), then $\text{pH}=\text{pK}_a$. This is the value at the half neutralization point and can be extrapolated from the titration curve. A careful pH calibration is necessary for pK evaluations. Nonetheless, the determined pK value is only an

approximation, as the ionic strength is not taken into account. In order to obtain more accurate values, titrations must be carried out with decreasing ionic strength and the results extrapolated to the ionic strength zero. pK evaluation in aqueous solution is limited to the range $3.5 < pK < 10.5$ because of the leveling effect of strong acids and the lack of jumps with very weak acids. pK values of mixtures of acids and polyvalent acids can also be determined.

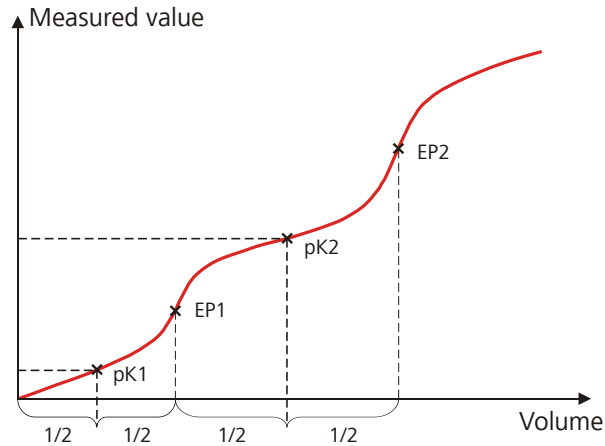


Figure 33 Evaluation of pK value / half neutralization potential

In non-aqueous solutions the half neutralization potential (HNP) is frequently used instead of the pK value. The HNP is evaluated in the same way as the pK value.

No parameters can be edited for the command **EVAL pK /HNP**.



NOTICE

If a start volume is to be added then it must be smaller than the half endpoint volume.

29.9.3 Minimum and maximum evaluation (EVAL MIN/MAX)

For the minimum or maximum measured value the associated volume, time and temperature are interpolated from the measuring point list. The evaluation begins as soon as the slope of the curve exceeds a particular threshold value.

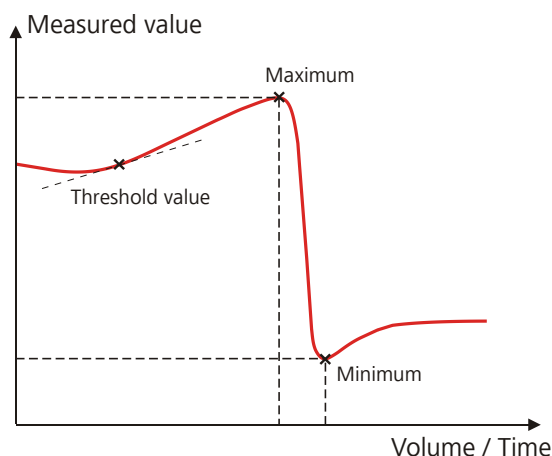


Figure 34 Evaluation of minimum and maximum

Evaluation

Selection whether the minimum or the maximum of the titration curve is being evaluated. With a command either the minimum or the maximum can be evaluated. If both values are needed, a second command has to be defined in the method.

Selection	Minimum Maximum
Default value	Maximum

Threshold value

The evaluation of the minimum or maximum begins as soon as the slope of the curve exceeds the set threshold value. Use a lower threshold value if the minimum or maximum is not found.

Measuring mode pH:

Input range	0.1 to 20
Default value	1.0
	With titrations, the unit is pH/mL, with measurements pH/s.

Measuring mode U, Ipol:

Input range	1.0 to 2,000.0
Default value	25.0
	With titrations the unit is mV/mL, with measurements mV/s.

Measuring mode Upol:

Input range	0.5 to 10.0
Default value	5.0
	With titrations the unit is μ A/mL, with measurements μ A/s.

Set windows

This parameter allows to evaluate break points only in a certain range (so-called window) of the curve. Break points outside this window will not be evaluated. Only one window can be defined. Within this window however, all break points are evaluated.

Selection	off Measured value Time Volume
Default value	off

off

The entire titration curve is being evaluated.

Measured value

The window is defined on the measured value axis.

Time

The window is defined on the time axis.

Volume

The window is defined on the volume axis.

Setting windows

The lower and upper limit of the window is defined under **[Set window]**.

Lower limit

Measured value, time or volume for the lower limit.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the .

Measured value (measuring mode pH):

Input range	-20.000 to 20.000
Default value	-20.000

Measured value (measuring mode U, Ipol):

Input range	-2,000.0 to 2,000.0 mV
Default value	-2,000.0 mV

Measured value (measuring mode Upol):

Input range	-200.00 to 200.00 µA
Default value	-200.00 µA

Measured value (measuring mode T):

Input range	-20.0 to 150.0 °C
Default value	-20.0 °C

Time:

Input range	0.0 to 999999.9 s
Default value	0.0 s

Volume:

Input range	0.00000 to 9,999.99 mL
Default value	0.00000 mL

Upper limit

Measured value, time or volume for the upper limit.

The values that are actually measurable may vary from the values specified here, depending on the hardware. The binding measurement value ranges can be found in the .

Measured value (measuring mode pH):

Input range	-20.000 to 20.000
Default value	20.000

Measured value (measuring mode U, Ipol):

Input range	-2,000.0 to 2,000.0 mV
Default value	2,000.0 mV

Measured value (measuring mode U_{pol}):

Input range	-200.00 to 200.00 µA
Default value	200.00 µA

Measured value (measuring mode T):

Input range	-20.0 to 150.0 °C
Default value	150.0 °C

Time:

Input range	0.0 to 999,999.9 s
Default value	999,999.9 s

Volume:

Input range	0.00000 to 9,999.99 mL
Default value	9,999.99 mL

29.9.5 Rate evaluation (EVAL RATE)

The rate evaluation provides the median dosing rate in a defined range (so-called window). The median dosing rate is determined by means of linear regression through a minimum of three measuring points. If the dosing is carried out with only one dosing device and if more than one cylinder volume is dosed, then the median dosing rate will be reduced as a result of the filling time.

When the command is opened for the first time, a window over the entire time range is already defined. In contrary to the potentiometric evaluation

of a DET/MET titration the individual windows may overlap. A maximum of 9 windows can be defined.

[New]

Define a new window.

[Delete]

Delete the selected window.

[Edit]

Edit the settings of the selected window.

Lower limit

Time for the lower limit.

Input range	0 to 999999 s
Default value	0 s

Upper limit

Time for the upper limit.

Input range	0 to 999999 s
Default value	999999 s

29.10 Calculations

The following calculation commands are available:

CALC	Defining calculations of a determination.
CALC LIVE	Defining the calculation whose result will be displayed in the live display (so-called live result) during a titration or measurement.

29.10.1 Calculations (CALC)

Calculations are defined with the command **CALC**. A method may contain a maximum of nine calculation commands. A maximum of nine calculations can be defined per command. A series of variables (raw data from the determination, previously calculated results, etc.) is available for the calculations.

The command contains the following options, among others:

- Save result as titer
- Save result as common variable
- Save result in result table
- Define result limits



Sequence / Edit command	
02 CALC	Calculation
Result	Result name
R1	Content
New Delete Edit	

The result name is specified in the list for each calculation.

[New]

Define a new calculation or load an existing template, see following chapter.

[Delete]

Delete the selected calculation from the list.

[Edit]

Edit the data of the selected calculation (see chapter 29.10.1.2, page 391).

29.10.1.1 Loading a result template


The most common calculations are already available, the so-called result templates, in order to facilitate the definition of a new calculation. You also have the option, however, of defining templates of your own (see dialog **System / Templates**).

Proceed as follows to load an existing template:

1 Select a result template


- In the **Sequence / Edit command** dialog, tap on the **[New]** button.

The list of Metrohm result templates is displayed:

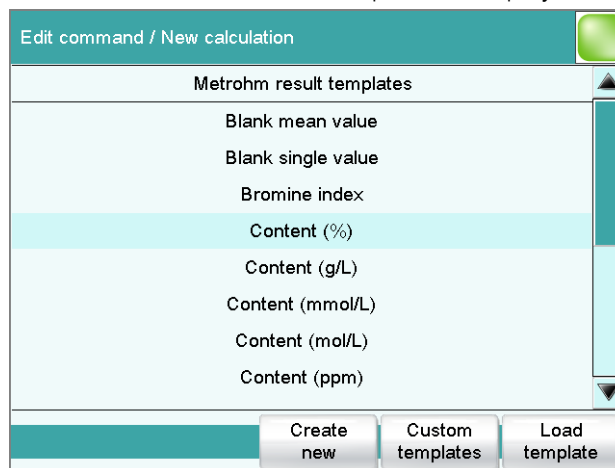

- Select the desired Metrohm result template and tap on **[Load template]**.
or
Tap on **[Custom templates]** and select a template you generated yourself.

If a statistic variable SMN# (# = 1..9) is defined in the template which is already used in another calculation, then the following message will be displayed:

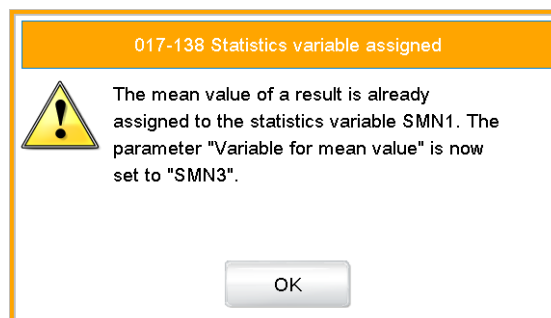

2 Confirm the message

- Tap on **[OK]**.

The Note dialog is displayed. This note is defined in the dialog **Edit result template / Note for wizard**.

New calculation / Load template

Enter the following variable:

F1= Molar mass for one equivalent of the substance that is determined.

Please consider the information in the note under edit result.

Cancel Back Next

3 Enter the numerical values of the wildcards

- Tap on **[Next]**.
The list of wildcards is displayed:

New calculation / Load template

F1= 35.45 F6=

F2= F7=

F3= F3=

F4= F9=

F5=

Cancel Back Next

- Enter the numerical values of the wildcards.
- Tap on **[Next]**.

The editing dialog of the new calculation is displayed:

The calculation can now be edited further, e.g. define result limits, see following chapter.

29.10.1.2 Editing a calculation

Result name

The result name is the text which will be shown in the result display and in the report. The default name corresponds to the result variable.

Entry	max. 24 characters
Default value	R1...R9

Calc. formula R1...R9

Shows the calculation formula. A special editor is opened for the definition (see chapter 29.10.3, page 397). The designation **R1...R9** corresponds to the result variable. Under **[Result variable]** this can later be modified.

Entry	max. 100 characters
Default value	empty

Decimal places

Number of decimal places used to display the result.

Input range	0 to 5
Default value	2

Result unit

The result unit is displayed and saved together with the result.

Entry	max. 10 characters
Selection	% mol/L mmol/L g/L mg/L mg/mL mg/100 g ppm g mg µg mL µL mg/piece °C µg/min mL/min µL/min
Default value	%

**[Note]**

Entering a note on the calculation.

[Result variable]

Modifying the result variable.

[Result limits]

Defining the limits for the monitoring of the result.

[Result options]

Defining additional settings for the calculation.

Dialog "Edit calculation / Note"

In this dialog you can enter a short text, e.g. to describe the variables used.

Dialog "Edit calculation / Result variable"

When a new calculation is created, the result variable will be automatically issued. In this dialog, however, it can be modified at any time.

Result variable

To each calculation belongs an unambiguous result variable. With this result variable you can use this result in additional calculations.

Selection	R1 R2 R3 R4 R5 R6 R7 R8 R9
	The selection contains only those result variables which are not issued in this calculation command yet.

Dialog "Edit calculation / Result limits"

For each result, limit values can be defined. These result limits are monitored when the calculation is carried out. If monitoring is activated, the results are shown as follows in the result display:

- Green, if the result falls within the limit values
- Red, if the result lies outside the limit values

Monitoring result limits

on | off (Default value: **off**)

Enabling and disabling the monitoring function.

Lower limit

When the monitored quantity falls below this value, the action defined in the following is triggered.

Input range	0 to 9999999999
Default value	0

Upper limit

When this value is exceeded, the action defined in the following is triggered.

Input range	0 to 9999999999
Default value	9999999999

Action

Selection of the action, which is carried out when the result limits are exceeded.

Selection	Display message Document message Cancel determination
Default value	Display message

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the result limits have been exceeded.

Cancel determination

The determination is stopped.

Dialog "Edit calculation / Result options"

In the dialog **Edit calculation / Result options**, settings for how to process the calculated result are defined.

Variable for mean value

If the statistic calculations have been activated (see method options), the mean value of the single results will be saved as variable SMN1 to SMN9. As default value, always the first free variable is being displayed.

Selection	SMN1 ... SMN9 off
-----------	----------------------------

off

For the result not statistic calculations are carried out.

Save as titer

The result can be saved as titer for the selected titrant.



NOTICE

The titer is assigned to the titrant defined in the last titration command before the calculation. Please note that the CALC command containing the titer assignment is inserted after the titration command with which the titer is determined.

Selection	Single value Mean value off
Default value	off

Single value

The result of the current determination is saved as titer.

Mean value

If statistic calculations are carried out for the result, the current mean value of the determination series is saved.

off

The result is not saved as titer.

Save as common variable

on | off (Default value: **off**)

The calculated result can be saved as a method-independent variable, a so-called common variable. The result is then also available in other methods for calculations. In all cases, only the single value will be saved, even if the statistics function is enabled.

Variable

Selection of the common variable to which the result is being assigned.

Selection	CV01 ... CV25
-----------	----------------------

Display result

on | off (Default value: **on**)

If you deactivate this parameter, the result is neither displayed in the result dialog nor printed in the result report. This can be advisable for intermediate results.

Save result in result table

on | off (Default value: **off**)

The calculated result can be saved in the result table. This may be advisable if e.g. the results of all determinations carried out on a particular day are to be displayed clearly. A maximum of nine results from a determination can be saved in the result table.

Precision

Setting, with which accuracy the result is used in additional calculations.

Selection	Round Truncate Full precision
Default value	Round

Round

The result is rounded to the defined number of decimal places (commercial rounding, in accordance with the US Pharmacopeia USP). If the digit at the first dropped decimal place is **1, 2, 3 or 4**, then it will be rounded off; if this digit is **5, 6, 7, 8 or 9**, then it will be rounded up. Negative digits will be rounded in accordance with their amount, i.e. away from zero.

Truncate

The result is cut to the number of decimal places defined.

Full precision

The result is used with full accuracy (floating point number either in "single precision" (32 bit) or in "double precision" (64 bit), according to the standard IEEE 754).

29.10.2 Calculations (CALC LIVE)

The command **CALC LIVE** can be used to define a calculation, the result of which will be shown in the live display during a titration or measurement. The current volume or measured value, respectively, is used for this purpose. This is helpful, e.g. for Karl Fischer titrations, in order to be able to track the water content directly during the titration.



CAUTION

The command must be inserted directly before the corresponding titration or measuring command.

This command is identical with the **CALC** command except for the following differences:

- The number of variables is limited.
- There are no result options.
- The result cannot be monitored.
- The result variable cannot be modified.
- The result will only be displayed in the live display. It appears neither in the results dialog nor in a report.

Result name

With live calculations no custom result name can be defined, the designation **LR** cannot be modified.

Calculation formula

Shows the calculation formula. A special editor is opened for the definition (see chapter 29.10.3, page 397).

Entry	max. 100 characters
Default value	empty

Decimal places

Number of decimal places used to display the result.

Input range	0 to 5
Default value	2

Result unit

The result unit is displayed and saved together with the result.

Entry	max. 10 characters
Selection	% mol/L mmol/L g/L mg/L mg/mL mg/100 g ppm g mg µg mL µL mg/piece °C µg/min mL/min µL/min
Default value	%

[Note]

Entering a note on the calculation.

Dialog "Edit calculation / Note"

In this dialog you can enter a short text, e.g. to describe the variables used.

Displaying the live result

Proceed as follows to display the live result in the live display:

1 Starting a determination

- Tap on [▶].

The determination starts and the live display is shown:

2 Define the display options

- Tap on the **[View]** button.
- Tap on the **[Meas. value options]** button.
A maximum of three measured values can be shown in the live display.
- Select the option **Live result** at one of the three parameters.

3 Display the live display


- Tap twice on [↵].

The current result is displayed,

29.10.3 Formula editor

The formulas for the calculations are entered with the formula editor. The maximum formula length is 100 characters.

The formula editor contains buttons for entering numbers, mathematical operators and variables. The variables can be selected from lists. This is advisable in order to avoid spelling mistakes. The formula editor is equipped with an automatic syntax check, which is activated when applying the formula. The standard rules of priority apply for the calculation operations.

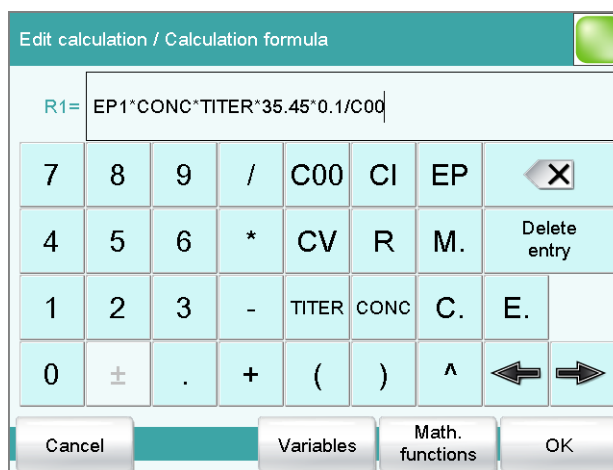

Key	Description
C00	Sample size
CI	Sample identification CI# (# = 1 - 2)
EP	Volume of endpoint EP# (# = 1...9)
CV	Common Variable CV# (# = 1...25)
R	Result R# (# = 1 - 9)
TITER	Titer of the titrant selected in the previous command
CONC	Concentration of the titrant selected in the previous command
□	The character in front of the cursor is deleted.
Delete entry	The content of the input field is deleted completely.

Key	Description
⇐	The cursor within the input field is shifted to the left by one character.
⇒	The cursor within the input field is shifted to the right by one character.

If the method contains more than one data-generating command (titration, measurement etc.) before the calculation command, then you will have to indicate the command identification in front of the variable:

Variable	Description
#M.	Titration or measuring command
#C.	Calculation command
#E.	Evaluation command

"#" stands for a sequential number that you must enter manually. The number for each command identification will be counted separately and does not necessarily correspond to the number of the command line.

Example: The variable **1M.TITER** corresponds to the titer of the titrant selected in the first titration command.



CAUTION

The command identifications are not adjusted automatically when additional commands are subsequently inserted into the method. Check the formula for correctness in such cases.

[Variables]

Opening the list of additional variables (*see chapter 29.10.3.1, page 398*).

[Math. functions]

Opening the list with mathematical functions (*see chapter 29.10.3.2, page 399*).

29.10.3.1 Variables

All of the variables which can be used for calculations are divided up according to subject in the dialog **Calculation formula / Variables**:

- **System variable**
List of the system-specific variables. These variables describe the current status of the system.
- **Result and statistics variables**
List of the variables which are required for calculating results and statistics for the current determination.

- **Common variables**

List of common variables. All of the common variables are listed, no matter whether they possess a value or not.

You will find a detailed description of all of the variables in *chapter 29.10.3.3, page 400*.

29.10.3.2 Mathematical functions

The following mathematical functions can be utilized in a calculation formula:

Function	Description
y^z	Power function Example: $4^2 = 16$
SQRT(X)	square root of X Example: $\sqrt{EP1}$
ABS(X)	absolute value of X Example: ABS(C00); in order, e.g. with reweighings, to convert the negative sample size to a positive value for later calculations
LN(X)	natural logarithm of X
LOG(X)	decimal logarithm of X
FRAC(X)	Fraction of X Example: $FRAC(2.5971) = 0.5971$
INT(X)	integer part of X Example: $INT(2.5971) = 2$
TST(X,Y)	Test function If invalid variables (e.g. missing endpoints) occur in a calculation, then these can be replaced with a valid value by using this function. In this way invalid results can be avoided. <ul style="list-style-type: none"> ▪ Syntax: <ul style="list-style-type: none"> – X = variable to be tested – Y = substitute value

29.10.3.3 Variable list

The following table contains all of the variables which can be used for calculations. For variables having an index (e.g. 'EP1'), the index must be entered manually. In the following table, this index is characterized with the character "#".

Variable	Description
C00	Sample size
CI1, CI2	Sample identifications The sample identifications can be used in calculations only if numerical values are entered.
DD	Duration of the entire determination
Titration	
TITER	Titer of the titrant selected in the titration command
CONC	Concentration of the titrant selected in the titration command
Titration, measurements, calibrations	
EP#	Volume/quantity of endpoint EP# (# = 1 - 9)
EC#	Charge at endpoint EP# (# = 1 - 9)
EM#	Measured value of endpoint EP# (# = 1...9)
EF#	ERC of endpoint EP# (# = 1 - 9)
ET#	Temperature at endpoint EP# (# = 1...9)
ED#	Time at endpoint EP# (# = 1...9)
ESI#	Recognition of endpoint EP# (# = 1 - 9) Endpoint found = 1, no endpoint = 0
FMN	Mean gas flow during MEAS T/F ON to MEAS T/F OFF
TMN	Mean temperature during MEAS T/F ON to MEAS T/F OFF
TMI	Minimum temperature during MEAS T/F ON to MEAS T/F OFF
TMA	Maximum temperature during MEAS T/F ON to MEAS T/F OFF
MIM	Initial measured value, i.e. measured value prior to the processing of the start conditions
MIT	Initial temperature, i.e. temperature prior to the processing of the start conditions
MSA	Volume for start volume
MSP	Volume for start measured value
MSS	Volume for start slope
MSV	Volume for all the start conditions
MSD	Duration of start conditions
MSM	Start measured value, i.e. measured value after the processing of the start conditions
MST	Start temperature, i.e. temperature after the processing of the start conditions

Variable	Description
MCD	Total duration of the titration, measurement or calibration
MTM	Type of temperature measurement (Pt1000, NTC or manually) Format = Text
MDD	Duration of effective dosing, i.e. without filling times, pauses
MDC	Drift for drift correction
DDC	Time for drift correction
MCV	End volume, i.e. total dosed volume at the end of the titration
MCQ	End quantity, i.e. total amount of removed water or of generated bromine, respectively, at the end of the titration (in µg)
MCC	Final charge, i.e. total charge in mAs at the end of the titration
MCM	End measured value
MCT	End temperature
MMP	Number of measuring points in measuring point list
MTS	Stop type Format = Text
LV	Currently dosed volume of the ongoing determination
LM	Current measured value of the ongoing determination
LD	Current duration of the ongoing titration or measuring command
LT	Current temperature of the ongoing determination
Sensor	
MEN	Electrode zero point pH(0) or E(0)
MSL	Electrode slope
MVA	Electrode variance (mathematically only defined for three standards or more);
MCL	Cell constant of conductivity measuring cell
Evaluations	
FP#	Volume of fixed endpoint FP# (# = 1 - 9)
FM#	Measured value of fixed endpoint FP# (# = 1 - 9)
FT#	Temperature at fixed endpoint FP# (# = 1 - 9)
FD#	Time at fixed endpoint FP# (# = 1 - 9)
HP#	Volume of pK value / half neutralization potential HP#
HM#	Measured value of pK value / half neutralization potential HP#
HT#	Temperature at pK value / half neutralization potential HP#
HD#	Time at pK value / half neutralization potential HP#
XIP	Volume/quantity of minimum measured value




Variable	Description
XIM	Minimum measured value
XIT	Temperature at minimum measured value
XID	Time at minimum measured value
XAP	Volume/quantity of maximum measured value
XAM	Maximum measured value
XAT	Temperature at maximum measured value
XAD	Time at maximum measured value
BP#	Volume of break point BP# (# = 1 - 9)
BM#	Measured value of break point BP# (# = 1 - 9)
BT#	Temperature at break point BP# (# = 1 - 9)
BD#	Time at break point BP# (# = 1 - 9)
RD#	Mean rate in time slot # (# = 1 - 9)
RDS#	Standard deviation for RD# (# = 1 - 9)
RDC#	Correlation coefficient for RD# (# = 1 - 9)
RM	Mean rate for whole range
RMS	Standard deviation for RM
RMC	Correlation coefficient for RM
Results, statistics	
R#	Result (# = 1–9)
SMN#	Mean value of result R# (# = 1–9)
SSA#	Absolute standard deviation of SMN# (# = 1 - 9)
SSR#	Relative standard deviation of SMN# in % (# = 1 - 9)
SNR#	Number of results from which the mean value SMN# was calculated (# = 1 - 9)
SSD	Number of determinations that are carried out for the statistics calculations
SNT	Statistics status (statistics activated = 1, statistics deactivated = 0)
Common variables	
CV#	Common variable (# = 1...25)
System variables	
%RN	Sample number, i.e. number that is increased by one at each start of a determination
%SC	Start counter, see dialog More determination data / Properties
%AS	Autostart status (autostart activated = 1, autostart deactivated = 0)
%AC	Autostart counter, i.e. number of autostarts already carried out
%AD	Autostart setpoint counter, i.e. number of autostarts to be carried out

Variable	Description
%SS	Status of the sample table (sample table activated = 1, sample table deactivated = 0)
%SL	Number of the current sample table line
%SE	Last sample in the sample table processed (yes = 1, no = 0)

29.11 Reports (REPORT)

The reports to be printed out are defined with the command **REPORT**. A maximum of ten reports can be defined for each command. The following reports can be inserted, among others:

- Result report
- Calculation report
- Parameter report
- System reports (system settings, sensor list, GLP data etc.)
- Method reports (method run, method parameters etc.)

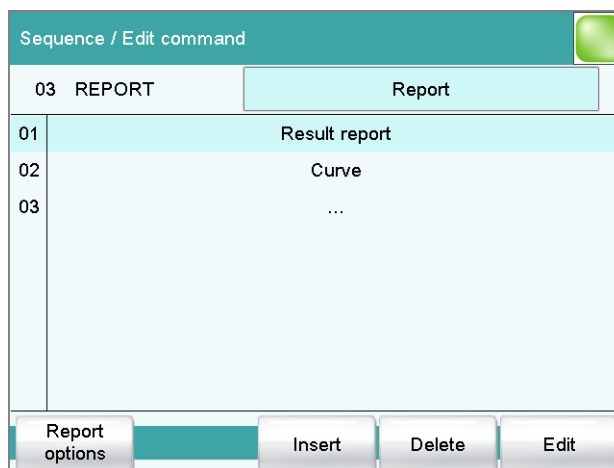

[Report options]

Defining the general settings for the report output.

[Insert]

Add a new report to the list.

[Delete]

Delete the selected report from the list.

[Edit]

Edit the settings of the selected report.

29.11.1 General report options

The settings for the report printout can be adjusted in the dialog **Edit command / Report options** to indicate, for example, whether a report header, a signature line or a frame is to be printed.

Report header

Defining the output of the report header. The report header contains general information such as device type, serial number, program version and the printing date.



NOTICE

You can also define your own report header, which is printed in addition to this report header (**System ► Templates ► Report header**).

Selection	off once on each page
Default value	on each page

off

The report header will not be printed.

once

The report header will only be printed on the first page.

on each page

The report header will be printed on every page.

Signature line

Output of a special line for date and signature. This line will be printed at the very bottom of every page.

Selection	off once on each page
Default value	off

off

No signature line will be printed.

once

The signature line will only be printed on the last page.

on each page

The signature line will be printed on every page.

Frame

on | off (Default value: **on**)

If this parameter is activated, a frame is printed as lateral marking.

29.11.2 Settings of the individual reports

Fixed key []: **More reports ► Report ► Edit**

Command REPORT: **Report ► Edit**

Settings can be edited for the following reports (for details, see online help):

- **Result report**
- **Curve**
Definition of the curve size, display of the individual measuring points and grid lines, etc.
- **Measuring point list**
Definition of the method command to which the measuring point list is to be printed.
- **Calculations**
Definition of the accuracy with which the variables used as well as the results calculated are printed.
- **Statistics short**
Definition of the time at which the report is to be printed (with each determination or only at the end of a sample series).
- **Statistics overview**
Definition of the time at which the report is to be printed (with each determination or only at the end of a sample series).
- **Sample table**
- **Result table**
Definition of the time at which the report is to be printed (with each determination, at the end of a sample series or at the end of the sample table).
- **Used devices**
- **Variables**
- **Monitoring**

29.11.3 List of reports

The following reports can be inserted in the command **REPORT**:



NOTICE

If you are using a **CUSTOM NEOS** printer, you can print only the reports that are marked with a **#** on the display.

Report	Contents
Result report	# Report with determination properties, sample data, calculated results, etc. If there are multiple determinations, the statistics will also be printed out.



Report	Contents
Curve	# Curve report.
Measuring point list	# Measuring point list report.
Calculations	# Details concerning the calculations carried out (parameter settings in accordance with the CALC command and the calculated results).
Used devices	The devices used for the determination, as displayed in the More determination data / Properties dialog.
Variables	All determination variables, as displayed in the results dialog.
Monitoring	Details for the monitored quantities (STAT only).
Statistics	
Statistics short	# Summary of the statistics calculations. The number of determinations, the mean value, and the absolute and relative standard deviations are printed out for each result.
Statistics overview	# Detailed statistical overview. The sample data and all individual results are printed out for each determination. The number of determinations, the mean value, and the absolute and relative standard deviations are printed out for each result.
Sample data	
Sample table	List of all determinations in the sample table with the respective sample data, as entered in the sample table.
Result table	
Result table	List of all determinations in the result table with results and with the determination data, as saved in the result table.
More reports ► Method reports	
Method sequence	Method properties and the list of all method commands.
Parameters full	Method properties and options, all method commands with all parameters. All parameters which no longer have their default settings will be printed in bold . All parameters which have been modified in comparison to the stored version of the method are indicated by an *.
Titration & measuring param.	Method properties; titration, measurement and calibration commands with all parameters. All parameters which no longer have their default settings will be printed in bold . All parameters which have been modified in comparison to the stored version of the method are indicated by an *.
Modified parameters	Method properties, all method parameters which have been modified in comparison to the stored version of the method.
Non default parameters	Method properties, all method parameters which no longer have any default settings.
More reports ► System reports	

Report	Contents
System settings	
System settings	Settings for acoustic signals and accuracy of the measured value display.
Dialog options	Settings for routine dialog and expert dialog.
Titrants	
Titrant list	List of all the titrants configured in the system.
All titrant data short	The most important titrant data of all titrants (name, concentration, titer, last titer determination).
All titrant data full	All of the titrant data of all of the titrants (working life, titer options, data concerning the exchange unit/dosing unit).
Sensors	
Sensor list	List of all the sensors configured in the system.
All sensor data short	The most important sensor data for all sensors (name, calibration data).
All sensor data full	All of the sensor data for all of the sensors (working life, complete calibration data, calibration interval).
Device manager	
Device list	List of all devices configured in the system.
All device properties	Properties of all the devices configured in the system.
GLP manager	
GLP data	All data stored in the GLP manager.
Common variables	
Common variable list	List of all the common variables defined in the system, together with their most important data (name, value, status).
All common variable properties	Properties of all common variables (name, value, validity, status).
Templates	
Templates sample data	Sample identification list and sample assignment table.
Result template list	List with all of the user-generated result templates.
All result templates details	Details of the all of the user-generated result templates (calculation formula, result options, note).
Input/Output lines	List with all of the defined input and output lines at the remote interface (name, bit pattern).
Custom calibration buffers	Temperature tables for all defined custom calibration buffers.
Rack tables	
Sample rack list	List with all of the sample racks configured in the system, together with designation, number of positions and rack code.
Miscellaneous	
Form feed	If this entry is inserted between two reports, then these will each be printed out on a separate page.

29.12 Dosing and Liquid Handling

The following dosing commands are available:

Preparing the exchange unit or dosing unit (PREP)	Rinsing the cylinder and the tubings of the exchange unit / dosing unit.
Emptying the dosing unit (EMPTY)	Emptying the cylinder and the tubings of the dosing unit.
Dosing a fixed volume (ADD)	Dosing a specified volume.
Liquid Handling (LQH)	Carrying out complex dosing tasks with a Dosing unit.

29.12.1 Preparing an exchange or dosing unit (PREP)

The **PREP** command is used to rinse and fill air bubble-free the cylinder and tubings of the exchange or dosing unit. You should carry out this function before the first determination or once a day.

Control device

This parameter will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control device. Control devices are defined under **System ► Device manager**.

Selection of the control device from the list of devices.

Selection	Selection of configured control devices
-----------	--

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the parameters defined for the titrant for the preparing as well as the tubing parameters (dosing unit only) are used. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether

the type of dosing drive matches. For buret units without integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place. The command is carried out with default parameters (see Chapter 32.2.2, page 443 and Chapter 32.3.2, page 444).

Cylinder volume to

This parameter can only be edited with **Titrant = not defined**.

Selection of the port via which the titrant is ejected. This setting is only relevant for dosing units, in case of preparing an exchange unit, this setting is ignored.

Selection	Port 1 Port 2 Port 3 Port 4
Default value	Port 1

29.12.2 Emptying a dosing unit (EMPTY)

The **EMPTY** command is used to empty the cylinder and tubings of the dosing unit.

Control device

This parameter will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control device. Control devices are defined under **System ► Device manager**.

Selection of the control device from the list of devices.

Selection	Selection of configured control devices
-----------	--

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the parameters defined for the titrant for the preparing as well as the tubing parameters (dosing unit only) are used.

Titration reagents are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place. The command is carried out with default parameters (see Chapter 32.2.2, page 443 and Chapter 32.3.2, page 444).

Air inlet

This parameter can only be edited with **Titrant = not defined**.

Selection of the port via which air is aspirated. This setting is only relevant for dosing units, in case of preparing an exchange unit, this setting is ignored.

Selection	Port 1 Port 2 Port 3 Port 4
Default value	Port 4

29.12.3 Dosing a specified volume (ADD)

You can dose a specified volume with the command **ADD**.

29.12.3.1 Dosing parameters

The parameters for the dosing are defined under **[Dosing parameters]**.

Volume

Volume which is dosed.

Input range	0.00000 to 99999.9 mL
Default value	10.0000 mL

Dosing rate

The rate at which dosing takes place.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing rate depends on the cylinder volume (*see chapter 32, page 442*).

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 32, page 442*).

29.12.3.2 Control instrument

The control instrument the determination is carried out with is selected under **[Control device]**. Control instruments are defined under **System ► Device manager**.

**NOTICE**

This button will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control instrument.

Control device

Selection of the control instrument from the list of devices. Only those devices are displayed which are able to carry out the command.

Selection	Selection of configured control instruments
-----------	--

29.12.3.3 Dosing device

The parameters for the dosing device are edited under **[Dosing device]**.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the correct data (titer, concentration, etc.) is always used for the calculation. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the



titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without an integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place.

Tandem dosing

A second dosing device can be defined under **[Tandem dosing]** in order to enable uninterrupted dosing. Dosing is carried out with a combination of two dosing devices so that the second dosing device is dosing while the first one is being filled and vice versa.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	off

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether the type of dosing drive matches. For buret units without integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, only the time interval for the GLP test of the buret unit is checked at the start of the determination.

**NOTICE**

The concentration, validity of the titer and the working life of the titrant are ignored.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined

No check takes place.

Filling rate

Rate at which the dosing cylinder is filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum filling rate depends on the cylinder volume (*see chapter 32, page 442*).

However, in order to ensure uninterrupted dosing, the following points must be taken into account:

- Use the highest possible filling rate in order to keep the filling times as short as possible. The filling rate should be decreased for viscous liquids.
- When you use two buret units with different dosing cylinders, the filling rate for the larger sized cylinder must be at the minimum:

$$v_{2,Fill} \geq v_{1,Fill} \cdot \frac{V_{Cyl.2}}{V_{Cyl.1}}$$

$v_{2,Fill}$ = filling rate in mL/min for the larger sized cylinder

$v_{1,Fill}$ = filling rate in mL/min for the smaller cylinder

$V_{Cyl.2}$ = cylinder volume in mL of the buret unit of the second dosing device

$V_{Cyl.1}$ = cylinder volume in mL of the buret unit of the first dosing device

Example:

Dosing device 1: volume = 20 mL, filling rate = 50 mL/min

Dosing device 2: volume = 50 mL

Filling rate 2 \geq 50 mL/min \cdot 50 mL / 20 mL \geq 125 mL/min



- The dosing rate must not exceed 75% of the filling rate of the smaller cylinder. These values are listed in the following table, valid at maximum filling rate:

Table 17 Maximum dosing rate for different dosing cylinders

Cylinder volume	maximum dosing rate	
	Exchange unit	Dosing unit
1 mL	2.25 mL/min	–
2 mL	–	5.00 mL/min
5 mL	11.25 mL/min	12.50 mL/min
10 mL	22.50 mL/min	25.00 mL/min
20 mL	45.00 mL/min	50.00 mL/min
50 mL	112.50 mL/min	124.50 mL/min

29.12.3.4 Stirrer

The parameters for the stirrer are edited under **[Stirrer]**.

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4 off
Default value	1

off

No stirrer will be used.

Stirring rate

Setting the stirring rate. It can be set in steps of –15 to +15. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 32.4, page 444*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	–15 to 15
Default value	8

Switch off automatically**on | off** (Default value: **on**)

If this parameter is activated, the stirrer will be switched off automatically at the end of the titration, measurement, etc.

29.12.4 Liquid Handling (LQH)

The command **LQH** can be used to carry out Liquid Handling tasks with a dosing device of the type **Dosino 800**. The manifold options of the Dosino can be exploited in full in combination with a USB Sample Processor. The four ports of the dosing units can be used at will as output or input ports. This means that not only simple dosing and filling procedures are possible. Complex Liquid Handling tasks such as pipetting or sample transfers can be carried out without difficulty. Several LQH and automation commands are required for this purpose which are best combined into subsequences.

29.12.4.1 Liquid Handling parameters

The parameters for the Liquid Handling function are defined under **[Liquid Handling param.]**.

Function

Type of Liquid Handling function.

Selection	Dose Fill Aspirate Eject Exchange position Change port Compensate End volume
-----------	---

Dose

The given volume is dosed. The dosing cylinder neither is automatically filled before nor afterwards.

Fill

The dosing cylinder is filled. The valve disc then remains at the selected port.

Aspirate

Liquid is aspirated. The dosing cylinder neither is automatically filled before nor afterwards. The aspirating volume has to be aspirated with a single piston stroke.

Eject

The whole cylinder content is ejected. The piston is, in contrast to the function **End volume**, lowered to the stop point, i.e. past the maximum volume mark. This function should only be carried out for preparing a dosing unit before the pipetting procedure.

Exchange position

The dosing cylinder is filled. This means that, for example, air can be aspirated via Port 4. The valve disc is then rotated to Port 2 and the dosing drive can be removed from the dosing unit.

Change port

The stopcock is moved to the given port, but no piston movement takes place.

Compensate

Because of the fact that the dosing units are interchangeable, the coupling of the Dosino piston rod (spindle) exhibits a low mechanical tolerance that can be noticed when the piston changes its direction of movement. This tolerance can be compensated with this function. A short piston movement is first made in the same direction as the previous movement, which is then followed by a piston movement in the reverse direction.

End volume

The whole cylinder content is ejected. The piston is lowered to the maximum volume mark i.e. until 10'000 pulses have been dosed. This command should be used for pipetting functions for emptying the cylinder.

In-/Outlet

Selection of the port via which the Liquid Handling command is carried out.

Selection	Port 1	Port 2	Port 3	Port 4
Port 1	Default value, when Function = Dose, Aspirate and End volume .			
Port 2	Default value, when Function = Fill, Exchange position and Change port .			
Port 4	Default value, when Function = Eject and Compensate .			

Volume

This parameter can only be edited with **Function = Dose** and **Aspirate**.

Volume which is dosed or aspirated.

Function = Dose:

Input range	0.00000 to 99999.9 mL
Default value	1.00000 mL

Function = Aspirate:

Input range	0.00000 to 50.0000 mL
Default value	1.00000 mL

Flow rate

This parameter can only be edited with **Function = Dose, Fill, Aspirate, Eject, Exchange position** and **End volume**.

Rate at which it is dosed or filled.

Input range	0.01 to 166.00 mL/min
Selection	maximum
Default value	maximum

maximum

The maximum dosing and filling rates depend on the cylinder volume (see chapter 32, page 442).

29.12.4.2 Control instrument

The control instrument the determination is carried out with is selected under **[Control device]**. Control instruments are defined under **System ► Device manager**.



NOTICE

This button will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control instrument.

Control device

Selection of the control instrument from the list of devices. Only those devices are displayed which are able to carry out the command.

Selection	Selection of configured control instruments
-----------	--

29.12.4.3 Dosing device

The parameters for the dosing device are edited under **[Dosing device]**.

Dosing device

Selection of the MSB connector to which the dosing device is connected. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Titrant

Selection of the titrant from the titrant list. We recommend always selecting the titrant. This ensures that the correct data (titer, concentration, etc.) is always used for the calculation. Titrants are defined under **System ► Titrants**. You can also enter a name which is not contained in the titrant list. When a determination is started there is a check whether the solution is contained in the list.

For buret units with integrated data chip, a check is made in the method run to verify whether the correct titrant has been attached and whether



the type of dosing drive matches. For buret units without an integrated data chip, the cylinder volume and the type of dosing drive are checked. For the selected titrant, the validity of the titer, the working life of the titrant and the time interval for the GLP test of the buret unit are checked at the start of the determination.

Selection	Selection of configured titrants not defined
Default value	not defined

not defined
No check takes place.

29.13 Communication

The following communications commands are available:

- Scanning remote lines (SCAN)** Defining remote signals which are awaited before the next method command is started.
- Setting remote lines (CTRL)** Defining remote signals.
- Receiving RS-232 commands (SCAN RS)** Defining RS-232 commands which are awaited before the next method command is started.
- Sending RS-232 commands (CTRL RS)** Defining the RS-232 commands which are sent.

29.13.1 Scanning remote lines (SCAN)

The command **SCAN** can be used to define input signals at the remote interface which are awaited before the next method command is started.

Control device

This parameter will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control device. Control devices are defined under **System ▶ Device manager**.

Selection of the control device from the list of devices.

Selection	Selection of configured control devices
-----------	--

Remote Box

Selection of the MSB connector the Remote Box is connected to. The selection does not depend on the number of MSB connectors of the control instrument. The remote signal defined is requested on this Remote Box.

Selection	1 2 3 4
Default value	1

Input signal

Selection of the signal out of the templates or entering the required bit pattern. Templates are defined under **System ► Templates ► Input lines**.

Entering a bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status

The input lines are always numbered from right to left, i.e. with the signal *******1** line 0 is expected to be active.



NOTICE

We recommend masking lines that are of no interest or for which no defined condition can be predicted with an asterisk (*).

Entry	Bit patterns containing exactly 8 characters or a max. of 24 characters for the name of the template
Default value	*****
Selection	Selection of the templates defined

Timeout

When this time interval has expired without recognizing the remote signal, the action defined in the following is triggered.

Input range	0 to 999 s
Default value	0 s

Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.



Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

29.13.2 Setting remote lines (CTRL)

The command **CTRL** can be used to define output signals at the remote interface.

Control device

This parameter will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control device. Control devices are defined under **System ► Device manager**.

Selection of the control device from the list of devices.

Selection	Selection of configured control devices
-----------	--

Remote Box

Selection of the MSB connector the Remote Box is connected to. The selection does not depend on the number of MSB connectors of the control instrument. The remote signal defined is sent via this Remote Box.

Selection	1 2 3 4
Default value	1

Output signal

Selection of the signal out of the templates or entering the required bit pattern. Templates are defined under **System ► Templates ► Output lines**.



NOTICE

A line set active is not being reset automatically, not even at the end of the determination.

Entering a bit pattern:

- 0 = line inactive
- 1 = line active
- * = retain line status
- p = set pulse

The output lines are always numbered from right to left, i.e. with the signal *******1** line 0 is set. With a pulse, the length is set to 200 ms. If it is required to set pulses with other lengths, you have to define a corresponding template.



NOTICE

We recommend masking lines that are of no interest or for which no defined condition can be predicted with an asterisk (*).

Entry	Bit patterns containing exactly 14 characters or a max. of 24 characters for the name of the template
Default value	*****
Selection	Selection of the templates defined

29.13.3 Scanning the RS-232 interface (SCAN RS)

The command **SCAN RS** can be used to define RS-232 commands which are awaited before the next method command is started.

Serial port

Selection of the serial interface the peripheral device is connected to. The RS-232 command defined is requested on this interface.

Selection	COM1 COM2 COM3 COM4 COM5 COM6 COM7 COM8
Default value	COM1

Character string

Input of the RS-232 command as character string. All characters of the ASCII code page can be used. Control characters (Esc, FF etc.) have to be entered as a three-place, decimal ASCII code, starting with a slash. Each command is automatically terminated with the ASCII characters **CR** and **LF**. The ***** can be used for one or more characters.

Entry	ASCII string with a max. of 24 characters
Default value	!.R

Timeout

When this time interval has expired without recognizing the RS-232 command, the action defined in the following is triggered.

Input range	0 to 999 s
Default value	0 s



Action

Selection of the action which is carried out when the time interval has expired.

Selection	Display message Document message Cancel determination
Default value	Display message For all three options it is documented in the determination data (see dialog More determination data / Messages), that the time interval has been expired.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Document message

In the determination data it will be documented, that the time interval has been expired.

Cancel determination

The determination is stopped.

29.13.4 Defining RS-232 commands (CONTROL RS)

The command **CONTROL RS** can be used to define the RS-232 commands which are sent.

Serial port

Selection of the serial interface the peripheral device is connected to. The RS-232 command defined is sent via this interface.

Selection	COM1 COM2 COM3 COM4 COM5 COM6 COM7 COM8
Default value	COM1

Character string

Input of the RS-232 command as character string. All characters of the ASCII code page can be used. Control characters (Esc, FF etc.) have to be entered as a three-place, decimal ASCII code, starting with a slash. Each command is automatically terminated with the ASCII characters **CR** and **LF**.

Entry	ASCII string with a max. of 24 characters
Default value	&M;\$G

29.14 Automation

29.14.1 Rotating sample rack (MOVE)

The command **MOVE** is used to move to a rack position or an external position.

Tower

Selection of the tower, with which the command is carried out. Both Tower 1 and Tower 2 can always be selected, even if your Sample Processor has only one tower.

Selection	1 2
Default value	1

1

View from the front, the right tower.

2

View from the front, the left tower.

Destination

Selection of the required target position.

Selection	Sample Ext. position Special beaker Rack position Current sample + Current sample - Next position Prev. position Calibration pos. rotate + rotate - swing + swing -
Default value	Sample

Sample

Rack position, which is defined by the sample variable. The sample variable is automatically increased by 1 after every method run or it can selectively be modified with the command **SAMPLE**.

Ext. position

One of the four external positions which can be moved to with the Swing Head. Enter the number of the external position (**1...4**) in the field to the right.

Special beaker

Special beakers, which are defined in the rack table, can be directly moved to. Enter the number of the special beaker (**1...16**) in field to the right.

Rack position

Any rack position. Enter the number of the rack position (**1...999**) in the field to the right.

**Current sample +**

Starting from the current sample (defined by the sample variable) the rack is moved forward by the number of rack positions (**1...999**) entered in the field to the right.

Current sample -

Starting from the current sample (defined by the sample variable) the rack is moved backward by the number of rack positions (**1...999**) entered in the field to the right.

Next position

Starting from the current rack position, the rack is moved forward by one position.

Prev. position

Starting from the current rack position, the rack is moved backward by one position.

Calibration pos.

For automatic calibrations with a Sample Processor (see "*Definition of the calibration positions*", page 451).

rotate +

Moving the rack forward by a certain increment. The rotation increment is defined in the properties of the tower.

rotate -

Moving the rack backward by a certain increment. The rotation increment is defined in the properties of the tower.

swing +

Swinging the robotic arm outward by a certain increment (towards higher angular degrees). The swing increment is defined in the properties of the Swing Head.

swing -

Swinging the robotic arm towards the center of the rack by a certain increment (towards lower angular degrees). The swing increment is defined in the properties of the Swing Head.

Beaker test action

This parameter can only be edited with **Destination = Sample, Next position** or **Prev. position**.

Selection of the action which is carried out when the beaker sensor does not detect any vessel at the sample position moved to. In addition, enable the beaker sensor in the rack table.

Selection	Rotate rack Display message
Default value	Display message

Rotate rack

The rack is moved to the next vessel on a sample position. **Exception:** If the sample variable is newly defined in the method before this command (command **SAMPLE**), then the option **Display message** is automatically valid.

Display message

A message is displayed. You can select whether you want to continue with the determination or cancel the run.

Options**Shift rate**

Rate at which the sample rack is moved.

Input range	3 to 20 °/s
Default value	20 °/s

Shift direction

Direction in which the rack is moved.

Selection	auto + -
Default value	auto

auto

A shift direction, with which the shorter way has to be passed, is automatically selected.

+

Counterclockwise rotation.

-

Clockwise rotation.

Swing rate

Rate at which the robotic arm is swung.

Input range	10 to 55 °/s
Default value	55 °/s

29.14.2 Moving the lift (LIFT)

The lift of a Sample Processor is moved with the command **LIFT**. The movement can, however, only be carried out if the Sample Processor is located at a valid rack position. This will not be the case, for example, after a Rack Reset.

Tower

Selection of the tower, with which the command is carried out. Both Tower 1 and Tower 2 can always be selected, even if your Sample Processor has only one tower.



Selection	1 2
Default value	1

1

View from the front, the right tower.

2

View from the front, the left tower.

Lift position

Selecting a predefined lift position or enter any lift position. The predefined lift positions are defined in the device properties of the Sample Processor. They can, however, also be assigned in the manual control.

Input range	0 to 235 mm
Selection	Work position Shift position Rinse position Special position Home position
Default value	Work position

Home position

The lift is being located at the upper stop position, i.e. at the position 0 mm.

Lift rate

Speed at which the lift is moved.

Input range	5 to 25 mm/s
Default value	20 mm/s

29.14.3 Controlling pumps (PUMP)

The **PUMP** command is used to control the pumps mounted on or connected to the tower of the sample changer.

Tower

Selection of the tower, with which the command is carried out. Both Tower 1 and Tower 2 can always be selected, even if your Sample Processor has only one tower.

Selection	1 2
Default value	1

1

View from the front, the right tower.

2

View from the front, the left tower.

Pump

Selection of the pump.

Selection	1 2 1 + 2
Default value	1

1

Pump 1 of the selected tower is switched.

2

Pump 2 of the selected tower is switched.

1 + 2

Both pumps of the selected tower are switched at the same time.

Status/Duration

Switching the pump(s) on and off.

Input range	0 to 999 s (Increment: 1) The pump is running during this time.
Selection	on off
Default value	on

on

The pump is switched on. The pump is running until it is explicitly switched off. If the determination is canceled with the key [], the pump is also switched off.

off

The pump is switched off.

29.14.4 Resetting the rack (RACK)

The following actions are carried out with the command **RACK**:

- Rack, lift and robotic arm are being reset.
- The rack code of the rack attached is being read out and the corresponding rack data is transferred to the Sample Processor.
- The sample variable is reset to the value 1.

Check rack

on | off (Default value: **off**)

Activate this parameter when the rack attached is to be checked. For this you, however, have to additionally select the rack in the dialog **Method options / Start options** at the option **Check rack**.

29.14.5 Defining the sample variable (SAMPLE)

The sample variable describes the current position of the sample on the rack of the Sample Processor. It is automatically increased by 1 after the end of a method run. In the following cases, the sample variable is automatically reset to the value 1:

- if the autostart counter is reset.
- if the sample rack is reset.

The sample variable can be altered specifically with the command **SAMPLE**.

Sample variable

Modifying the sample variable.

Selection	= + -
Default value	+

=

The sample variable corresponds to the number entered in the field **Value**.

+

The sample variable corresponds to the current value plus the number entered in the field **Value**.

-

The sample variable corresponds to the current value minus the number entered in the field **Value**.

Value

Value by which the current sample variable is to be modified.



NOTICE

The sample variable can only be set on rack positions which are not defined as special beakers.

Input range	1 to 999
Default value	1

29.14.6 Creating a subsequence (SUBSEQ)

It is useful, particularly with more complex tasks, e.g. rinsing sequences or Liquid Handling, to combine individual command sequences into a single subsequence. Subsequences are managed the same as method commands. A maximum of 99 commands can be combined to form a single subsequence. There is no limit to the number of subsequences within one method.

[Load/ Save]

Loading a saved subsequence or saving the current subsequence.

[Subseq. options]

Defining various settings which correspond to the entire subsequence.

[Insert command]

Insert a new method command. It is inserted before the selected command.

**NOTICE**

Not all commands are available for subsequences. Commands which cannot be inserted into subsequences are disabled.

[Delete command]


Delete the selected method command.

[Edit command]

Edit the selected method command.

29.14.6.1 Loading / saving a subsequence

[Load/ Save] is used to load stored subsequences or to save new subsequences. This means that they are available for all methods and only have to be created once.

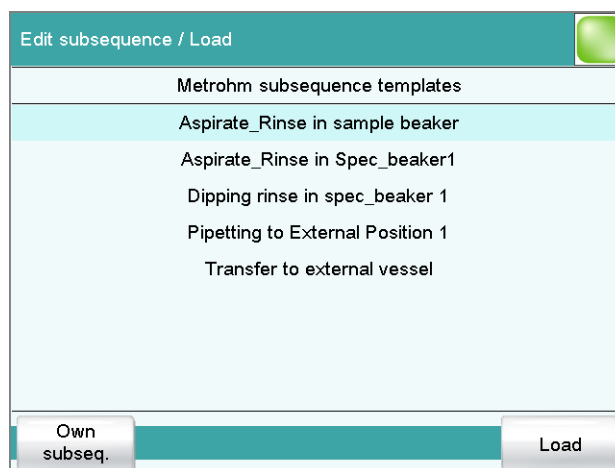Loading a subsequence**[Own subseq.]**

Opening the list of own subsequences.

[Load]

Loading the selected subsequence.

Saving a subsequence

Custom-made subsequences are saved in the internal memory.



NOTICE

In order to be able to use the subsequences also for other titration systems, you have to create a backup. You can only restore the subsequences from this backup on another system.

File name

File name of the subsequence.

Entry	max. 32 characters
-------	---------------------------

[Save]

Saving the subsequence to the internal memory.

29.14.6.2 Subsequence options

In this dialog, the settings can be defined which correspond to the entire subsequence.

Subsequence type

Selection	Sample sequence Start sequence End sequence Stop sequence Cond. sequence
Default value	Sample sequence

Sample sequence

The subsequence is carried out with every determination.

Start sequence

This subsequence is only carried out at the start of a sample series when the autostart counter = 1.

End sequence

This subsequence is only carried out at the last sample of a series when the autostart counter has reached the setpoint value (number of autostarts).

Stop sequence

The stop sequence is only carried out when canceling a method because of one of the following cases: Manual stop with [], stop because of an error, stop via remote signal.

Cond. sequence

This subsequence is carried out directly before the conditioning. This gives you the possibility to start e.g. a connected Polytron® high-frequency homogenizer before the conditioning as well as to set its stirring rate. Use the command **CONTROL RS** for this.

Cycles

Number of directly consecutive repetitions of the subsequence.

Input range	0 to 999
Default value	1
Selection	Calibration

Calibration

For automatic calibrations with a Sample Processor (*see chapter 32.8, page 450*).

[Note]

Entering a note on the subsequence.

Dialog "Subsequence options / Note"

A short text can be entered in this dialog, e.g. for the description of the subsequence commands or for the application purpose.

[Display options]

Definition when the note is being displayed.

Automatically after loading the subsequence

on | off (Default value: **off**)

If this parameter is activated, the note is displayed when loading the subsequence. It can otherwise only be read in this dialog.

29.15 Miscellaneous commands**29.15.1 Controlling a stirrer (STIR)**

A stirrer is controlled with the command **STIR**.

Control device

This parameter will only be shown when, in addition to the Ti-Touch, a Sample Processor has been configured as control device. Control devices are defined under **System ► Device manager**.

Selection of the control device from the list of devices.

Selection	Selection of configured control devices
-----------	--

Stirrer

Selection of the MSB connector the stirrer is connected to. The selection does not depend on the number of MSB connectors of the control instrument.

Selection	1 2 3 4
Default value	1

Status/Duration

Switching the stirrer on or off.

Input range	0 to 999 s (Increment: 1)
Selection	on off
Default value	on

on

The stirrer is switched on. The stirrer runs until it is explicitly switched off. If the determination is canceled with the key [], the stirrer is also switched off.

off

The stirrer is switched off.

Stirring rate

Setting the stirring rate. It can be set in steps of -15 to +15. The default setting **8** corresponds to 1000 rpm. The formula for calculating the rotational speed is specified in *chapter 32.4, page 444*. The optimum stirring rate can be tested in the manual control.

The algebraic sign of the stirring rate changes the direction in which the stirring is done. When the stirrer is viewed from above, this means:

- "+": counterclockwise rotation
- "-": clockwise rotation

Input range	-15 to 15
Default value	8

29.15.2 Pausing the method run (WAIT)

The method run can be paused with the command **WAIT**.

Hold sequence

on | off (Default value: **off**)

If this parameter is activated, then the method run is paused until it is continued manually.

Waiting time

This parameter can only be edited when **Hold sequence** is deactivated.

The method run is automatically continued after this waiting time.

Input range	0 to 99999 s
Default value	30 s

Message

on | off (Default value: **off**)

If this parameter is activated, the text message defined in the following is displayed during the waiting time.

Message text

This parameter can only be edited when **Message** is activated.

Text which is displayed during the waiting time.

Entry	max. 28 characters
Default value	empty

29.15.3 Scan data (REQUEST)

The following data can be scanned in the method run with the command **REQUEST**:

- Sample data
 - Sample identification 1
 - Sample identification 2
 - Sample size (value and unit)
- Common variable

You can select thereby whether the sequence should be paused or resumed in the background.

Sample ident.

Selection of the sample identification that is queried in the method run.

Selection	off Identification 1 Identification 2 ID1 & ID2
Default value	off

Sample size

on | off (Default value: **on**)

If this parameter is activated, then the value for the sample size will be requested.

Sample size unit

on | off (Default value: **off**)

If this parameter is activated, then the unit for the sample size will be requested.

Common variable

Selection of the common variable which is queried in the method run.

Selection	CV01...CV25 off
Default value	off

30 Operation and maintenance

30.1 System initialization

In very rare instances, a faulty file system (e.g. because of a program crash) may lead to an impairment of program functioning. The internal file system must be initialized in such cases.



CAUTION

If you carry out a system initialization, then all user data except the methods will be deleted.

We recommend creating a backup of the system at regular intervals in order to avoid data losses.

Proceed as follows for the system initialization:

1 Switching off the instrument

- Press the power switch at the rear.

2 Switching on the instrument

- Press the power switch at the rear.
An image with a frog will be displayed after a couple of seconds.
- While the image is being displayed, press the power switch again and hold it down.
- Do not release the power switch until an acoustic signal sounds.

A system message is displayed.

3 Confirm system initialization

- Confirm the message with **[Yes]**.
A second security prompt is displayed.
- Confirm this message with **[Yes]** as well.

Initialization is started. After the initialization has been completed, the 916 Ti-Touch is started automatically.

30.2 Maintenance

Maintenance

The electronic and mechanical functional groups of Metrohm instruments can and should be checked by specialist personnel from Metrohm as part of a regular preventive maintenance schedule. Please ask your local Metrohm representative regarding the precise terms and conditions involved in concluding a corresponding maintenance agreement.

For detailed information on this topic, please visit www.metrohm.com.

Cleaning

Cleaning the surfaces of the instrument

Prerequisites

- The instrument is disconnected from the power grid.

- 1 Clean the surfaces with a damp cloth.



NOTICE

Water or ethanol can be used as a cleaning medium.



NOTICE

The connectors at the rear of the instrument must only be cleaned with a dry cloth.

31 Troubleshooting

31.1 Editing methods

Problem	Cause	Remedy
The method command cannot be inserted.	No control instrument which supports the method command is configured in the device manager.	Add a control instrument which supports the desired method command to the list of devices (see chapter 11, page 95).
The additional evaluations (EVAL commands) cannot be inserted.	A titration or measuring command to which the evaluation can be applied is missing.	<ol style="list-style-type: none"> 1. Insert a titration or measuring command. 2. Insert the EVAL command directly after it.
A MEAS T/F ON command or MEAS T/F OFF command is depicted in red in the command list.	The corresponding MEAS T/F ON command or MEAS T/F OFF command is missing or not inserted in the correct order in the method run. To be able to execute a method, both MEAS T/F commands are needed in the method run.	Insert the missing MEAS T/F command and observe the logical order of the commands.
An EVAL command is depicted in red in the command list.	The associated titration or measuring command has been deleted or replaced by a different one.	<ol style="list-style-type: none"> 1. Delete the EVAL command. 2. Insert a titration or measuring command. 3. Insert the EVAL command directly after it.

31.2 Sample table


Problem	Cause	Remedy
The method cannot be entered in the sample table.	The option Use sample assignment table is activated. The method assigned to the sample identification will be loaded automatically at the start of the determination.	Deactivate the option (see chapter 15.1, page 162).




31.3 Results/Statistics

Problem	Cause	Remedy
No statistics are carried out for a result.	<i>Statistics is not activated.</i>	In the method options, activate the option Statistics (see chapter 16.5, page 187).
	<i>No statistics variable has been assigned to the calculation.</i>	In the result options (command CALC), define a variable for the mean value (see "Variable for mean value", page 168).
	<i>Statistics is not activated.</i>	In the control dialog, activate the Statistics option (see chapter 18, page 201).
A result has been removed from the statistics, the mean value has not been reassigned to the common variable or to the TITER variable.	<i>Assignment is not carried out automatically with retroactive modifications.</i>	Recalculate the determination manually (see chapter 24, page 229).
The result is not displayed in the result table.	<i>The column display in the result table is incorrectly configured.</i>	Modify the settings for the display accordingly (see chapter 26.1, page 252).
	<i>The parameter Save result in result table is not activated.</i>	In the result options (command CALC), activate the parameter (see "Save result in result table", page 169).

31.4 Printing

Problem	Cause	Remedy
Tapping on the fixed key [] has no effect.	<i>The fixed key is disabled.</i>	In the Dialog options / Fixed keys dialog, activate the Print option (see chapter 7.2, page 43).
	<i>A determination is running.</i>	Wait until the determination is finished.

31.5 Manual control

Problem	Cause	Remedy
Tapping on the fixed key [] has no effect.	<i>The fixed key is disabled.</i>	In the Dialog options / Fixed keys dialog, activate the Manual control option (see chapter 7.2, page 43).
A button in the manual control is disabled.	<i>The required hardware is either not connected or incorrectly connected.</i>	<ol style="list-style-type: none"> 1. Switch off the 916 Ti-Touch. 2. Connect the hardware correctly. 3. Switch the 916 Ti-Touch back on.
	<i>The required hardware is being used in the ongoing determination.</i>	Wait until the determination is finished.

31.6 Miscellaneous

Problem	Cause	Remedy
The 916 Ti-Touch cannot be switched on.	<i>The Ti-Touch is connected to an MSB socket on the control instrument.</i>	Plug in the Ti-Touch at the Controller connector (see manual for control instrument).

31.7 SET titration

Problem	Cause	Remedy
The titration will not be finished.	<i>The minimum dosing rate is too low.</i>	Define Titration rate = user and increase the minimum rate (Min. rate) (see chapter 29.3.2, page 325).
	<i>The stop criterion is unsuitable.</i>	Adjust the control parameters (see chapter 29.3.2, page 325): <ul style="list-style-type: none"> ▪ Increase the stop drift. ▪ Select a short delay time.
The sample is over-titrated.	<i>The control parameters are unsuitable.</i>	Adjust the control parameters (see chapter 29.3.2, page 325): <ul style="list-style-type: none"> ▪ Select Titration rate = slow. ▪ Define Titration rate = user and increase the control range. ▪ Define Titration rate = user and reduce the maximum rate (Max. rate).



Problem	Cause	Remedy
		<ul style="list-style-type: none"> Define Titration rate = user and reduce the minimum rate (Min. rate). Stir faster. Arrange the electrode and buret tip to an optimum .
	<i>The electrode responds too slowly.</i>	Replace the electrode.
The titration time is too long.	<i>The control parameters are unsuitable.</i>	Adjust the control parameters (see chapter 29.3.2, page 325): <ul style="list-style-type: none"> Select Titration rate = optimal or fast. Define Titration rate = user and reduce the control range. Define Titration rate = user and increase the maximum rate (Max. rate). Define Titration rate = user and increase the minimum rate (Min. rate).
The results are spread widely.	<i>The minimum dosing rate is too high.</i>	Define Titration rate = user and reduce the minimum rate (Min. rate) (see chapter 29.3.2, page 325).
	<i>The control range is too small.</i>	Define Titration rate = user and increase the control range.
	<i>The electrode responds too slowly.</i>	Replace the electrode.

31.8 STAT titration

Problem	Cause	Remedy
The control point is not kept stable. The measured value is at times too high and at times too low. The controller "oscillates".	<i>The control parameters are unsuitable.</i>	Adjust the control parameters (see chapter 29.4.2, page 338): <ul style="list-style-type: none"> Reduce the Max. rate and possibly also the Min. rate. Increase control range.
	<i>The stirring is not optimal.</i>	Check and optimize efficiency of the stirring.
	<i>The arrangement of electrode and buret tip is not optimum.</i>	Arrange the electrode and buret tip to an optimum (see manual of the instrument).

Problem	Cause	Remedy
	<i>The cylinder volume of the exchange/dosing unit might be too large.</i>	Use an exchange/dosing unit with a smaller cylinder volume (yields a smaller volume increment per pulse).
The control point is not reached for too long.	<i>The control parameters are unsuitable.</i>	Adjust the control parameters (see chapter 29.4.2, page 338): <ul style="list-style-type: none">▪ Reduce control range.▪ increase Min. rate.▪ increase Max. rate.

32 Appendix

32.1 Manual Dosing Controller, optional accessory for manual titration

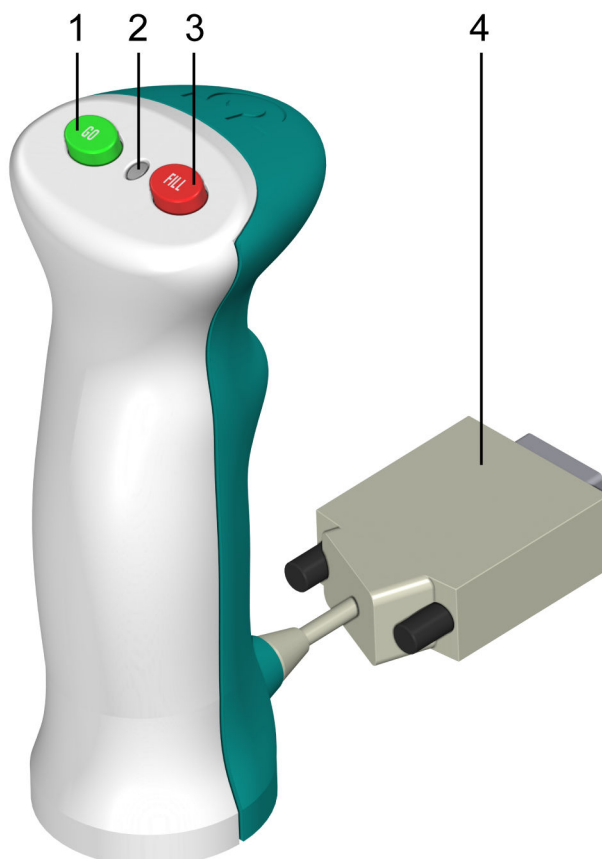


Figure 36 Manual Dosing Controller 6.2107.120 – Overview

1 Start/Dosing key [GO]

For starting the method and for dosing during the MAT command.

2 Status LED

Indicates whether the instrument is ready.

3 Cancel command button [FILL]

For canceling the running command.

4 Connector plug

With 25-pin D-sub plug.

32.2 Dosing unit

32.2.1 Maximum dosing and filling rate

The maximum dosing rate and maximum filling rate for the dosing unit depend on the cylinder volume:

Cylinder volume	Maximum rate
2 mL	6.67 mL/min
5 mL	16.67 mL/min
10 mL	33.33 mL/min
20 mL	66.67 mL/min
50 mL	166.00 mL/min

Independent of the cylinder volume, values ranging from 0.01 to 166.00 mL/min can always be entered. When the function is carried out the rate will be, if necessary, decreased automatically to the highest possible value.

32.2.2 Default parameters for preparing (PREP) and emptying (EMPTY)

The **PREP** command is used to rinse and fill the cylinder and tubings of the dosing unit air bubble-free. You should carry out this function before the first determination or once per day.

The command **EMPTY** is used to empty the cylinders and the tubings of the dosing unit.

If no titrant is selected in the command, preparing and emptying will be carried out with the following default parameters:

- The entire cylinder volume is dosed at the maximum dosing rate through Port 1.
- The cylinder is filled at the maximum filling rate through Port 2.
- The following dimensions apply for the tubings:
 - Tubing at Port 1: length = 40.0 cm, diameter = 2 mm
 - Tubing at Port 2: length = 25.0 cm, diameter = 2 mm

32.3 Exchange unit

32.3.1 Maximum dosing and filling rate

The maximum dosing rate and maximum filling rate for the exchange unit depend on the cylinder volume:

Cylinder volume	Maximum rate
1 mL	3.00 mL/min
5 mL	15.00 mL/min
10 mL	30.00 mL/min
20 mL	60.00 mL/min
50 mL	150.00 mL/min

Independent of the cylinder volume, values ranging from 0.01 to 166.00 mL/min can always be entered. When the function is carried out the rate will be, if necessary, decreased automatically to the highest possible value.

32.3.2 Default parameters for preparing (PREP)

The **PREP** command is used to rinse and fill the cylinder and tubings of the exchange unit air bubble-free. You should carry out this function before the first determination or once per day.

If no titrant is selected in the command, preparing will be carried out with the following default parameters:

- The entire cylinder volume is dosed twice at the maximum dosing rate.
- The cylinder is filled at the maximum filling rate.

32.4 Stirring rate

The stirring rate can be adjusted in steps from -15 to +15.

The approximate rotational speed for the internal magnetic stirrer (depends on the product version) can be calculated with the following formula:

$$\text{Rotational speed/min (r/min)} = 125 \cdot \text{Stirring rate}$$

Example:

Configured stirring rate: 8

Rotational speed in revolutions per minutes = $125 \cdot 8 = 1,000$

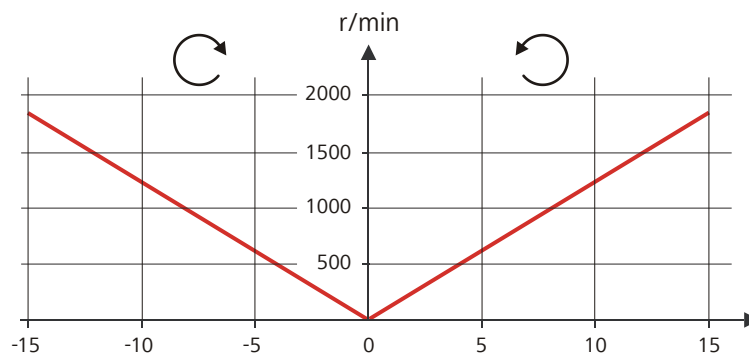


Figure 37 Rotational speed depending on the stirring rate

The information on the separately connectable 802 propeller stirrer can be found in the "802 Stirrer" manual.

32.5 Balance

The sample size and the associated unit can be sent from a connected balance. The sample size is transmitted as a number with up to ten characters (including algebraic sign and decimal point).

Sample size and unit are sent as a single character string. They are separated by a space character. The string is terminated with the ASCII characters **CR** and **LF**.

If the balance sends a negative sample size (e.g. when you are reweighing a sample), then the algebraic sign is adopted. The algebraic sign is, however, ignored for the calculations.



NOTICE

With some balances, the sample identification and the method can be sent in addition to the sample size.

Make sure that the balance does not send the sample size until the end.

Mettler AX

For the Mettler AX balance, the fields that contain the sample identification or the method must be designated as follows:

- Designation for the field with the method name: **METHOD**
- Designation for the field with sample identification 1: **ID1**
- Designation for the field with sample identification 2: **ID2**



32.6 Remote interface

The 6.2148.010 Remote Box allows devices to be controlled which cannot be connected directly to the MSB interface of the Ti-Touch.

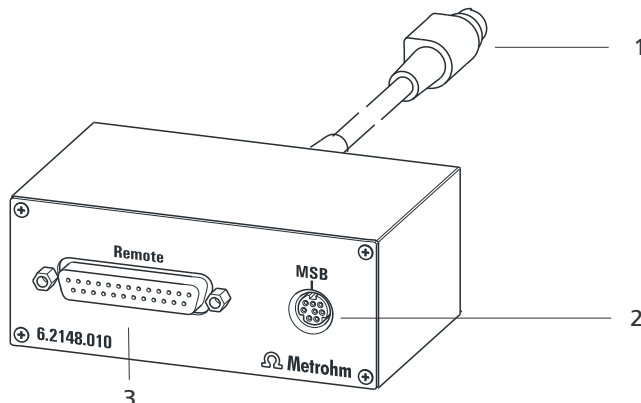


Figure 38 Connectors of the Remote Box

- | | |
|--|---|
| 1 Cable
For connecting to the Ti-Touch. | 2 MSB connector
Metrohm Serial Bus. For connecting external dosing devices or stirrers. |
| 3 Remote connector
For connecting instruments with a remote interface. | |

32.6.1 Pin assignment of the remote interface


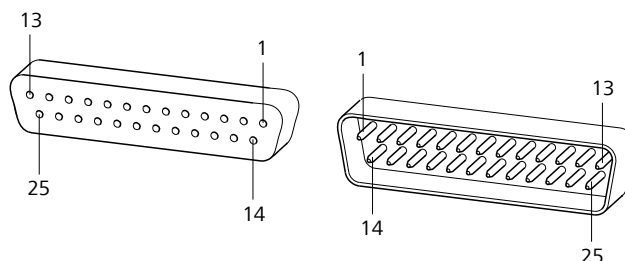



Figure 39 Pin assignment of remote socket and plug

The above figure of the pin assignment of a Metrohm remote interface applies not only for the Remote Box, but also for all Metrohm devices with 25-pin D-Sub remote connector.

Inputs

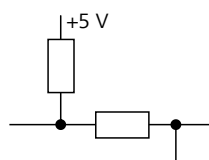


approx. 50 kΩ Pull-up

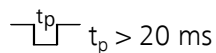

$t_p > 20 \text{ ms}$

active = low, inactive = high

Outputs

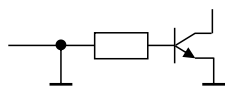

Open Collector

$t_p > 200 \text{ ms}$

active = low, inactive = high

$I_C = 20 \text{ mA}$, $V_{CE0} = 40 \text{ V}$

+5 V: maximum load = 20 mA

The following tables offer information concerning the assignment of the individual pins and their function.

Table 18 Inputs and outputs of the remote interface

Assignment	Pin No.	Function
Input 0	21	Start
Input 1	9	Stop
Input 2	22	
Input 3	10	Quit
Input 4	23	-
Input 5	11	
Input 6	24	
Input 7	12	
Output 0	5	Ready
Output 1	18	Conditioning OK
Output 2	4	Determination
Output 3	17	EOD
Output 4	3	
Output 5	16	Error
Output 6	1	
Output 7	2	Warning
Output 8	6	
Output 9	7	
Output 10	8	
Output 11	13	
Output 12	19	



Assignment	Pin No.	Function
Output 13	20	
0 volts / GND	14	
+5 volts	15	
0 volts / GND	25	

Table 19 Explanation of the individual functions

Function	Explanation
Start	The current method is started at the time of activation. $t_{pulse} > 100 \text{ ms}$
Stop	The current method is canceled (Stop) at the time of activation. $t_{pulse} > 100 \text{ ms}$
Quit	The current command in the determination run will be canceled at the time of activation. $t_{pulse} > 100 \text{ ms}$
Ready	The instrument is ready to receive a start signal.
Conditioning OK	The line is set when Conditioning with SET titration is at OK. The line remains set until the determination is started with [START] .
Determination	The instrument performs a data-generating determination.
EOD	End of Determination. Pulse ($t_{pulse} = 200 \text{ ms}$) after a determination or after a buffer during calibration using a Sample Processor.
Error	The line is set for error message display.
Warning	The line is set for warning message display.

32.7 Result variables as parameter setting

For most of the numerical parameters, the result of a calculation can also be utilized in place of a number. This result must be calculated in a preceding calculation command (command **CALC**). A typical application is the use of a relative start volume.

The following example is used to explain how this method is used to define a sample size-dependent start volume for a titration. Proceed as follows:

1 Create a new calculation

- Insert the calculation command **CALC** in front of the titration command.
- Enter the calculation formula, e.g. $R1 = C00 \cdot 3$.

01 CALC Calculation

Result name: Relative start volume

Calc. formula R1: C00*3

Decimal places: 3

Result unit: mL

Note Result variable Result limits Result options

2 Define a result variable as parameter setting

- In the titration command, enter the variable **R1** as the start volume.

02 DET pH Dynamic pH titration

Start volume mL

Dosing rate mL/min

Pause s

The selection of the result variables is displayed in the number editor.



NOTICE

If the method before the command with the result variables applied contains more than one calculation commands (possibly with result variables bearing the same name), then the result variable of the preceding calculation command will always be used.

Please note that the result must be within the input range of the parameter, as otherwise the determination run will be stopped when the parameter is accessed.

32.8 Electrode calibration with Sample Processors

Basic framework for a method for automatic calibrations

If the calibration of electrodes is carried out with a Sample Processor, then the basic framework of the method will appear as follows:

1. Subsequence with the commands required for the calibration (command **SUBSEQ**)
 - a. Command for change of buffer (command **MOVE**)
 - b. Command for moving the lift (command **LIFT**)
 - c. Calibration command (command **CAL pH**)
 - d. ...
2. ...

In the interest of general comprehensibility, other commands which are required are not listed here. A method template is available for an automatic pH calibration with subsequent measurement.

Command parameterization

The following special settings are required with the individual commands:

- Command **SUBSEQ**:
The setting **Cycles = Calibration** must be adjusted in the Subsequence options dialog.
- Command **MOVE** for changing the buffer:
The only definition made in this command is that a calibration position is to be moved to. The setting **Destination = Calibration pos.** must be adjusted for this purpose. The question of which rack positions are used as calibration positions must be defined in the rack table (see below).
- Command **CAL pH**:
The setting **Sample Processor = USB** must be adjusted in the Calibration parameters dialog.

Definition of the calibration positions

The positions of the individual buffers on the rack are defined in the rack table of the rack used. The following options are available:

- **Utilization of sample positions**
The desired rack position for the first buffer must be entered in the dialog **Sample rack / Edit rack data** for the parameter **1. Calibration pos.**. In this case, it is imperative that the remaining buffers be placed on the rack positions directly following.
- **Utilization of special beakers**
Any rack position can be defined as a special beaker. It is, however, preferable to set them at high rack positions in order to be able to begin sample series at rack position 1. The special beakers are moved to in ascending order.
The list of special beakers is displayed in the dialog **Sample rack / Edit rack data** by pressing **[Special beakers]**. The number of buffers used for calibration will determine the number of special beakers to be defined. We recommend defining special beakers 1-n. Afterwards, in the dialog **Sample rack / Edit rack data**, for the parameter **1. Calibration pos.**, the option **Special beaker 1** must be selected for the first buffer.



32.9 Stored buffer series for pH calibration

The temperature-dependent pH values of several common pH buffers are stored in the Ti-Touch for automatic buffer recognition during pH calibration. In addition to the Metrohm buffer solutions, other reference buffers and technical buffers are also included in the tables.

The following tables provide an overview of the stored pH(T) series.

pH values printed in **bold** are the values for the reference temperature of the respective buffer set.

pH values printed in *italics* are interpolated or extrapolated values, all the others correspond to the manufacturer's specifications.

32.9.1 Metrohm

Table 20 Metrohm buffer solutions

Temp. (°C)	Metrohm		
	pH 4.00	pH 7.00	pH 9.00
0	3.99	7.11	9.27
5	3.99	7.08	9.18
10	3.99	7.06	9.13
15	3.99	7.04	9.08
20	3.99	7.02	9.04
25	4.00	7.00	9.00
30	4.00	6.99	8.96
35	4.01	6.98	8.93
40	4.02	6.98	8.90
45	4.03	6.97	8.87
50	4.04	6.97	8.84
55	4.06	6.97	8.81
60	4.07	6.97	8.79
65	4.09	6.98	8.76
70	4.11	6.98	8.74
75	4.13	6.99	8.73
80	4.15	7.00	8.71
85	4.18	7.00	8.70

Temp. (°C)	Metrohm		
	pH 4.00	pH 7.00	pH 9.00
90	4.20	7.01	8.68
95	4.23	7.02	8.67



NOTICE

Update

The values of the individual buffers with the corresponding temperatures are kept up to date as far as possible.

However, they may be changed by the respective manufacturers.

32.9.2 NIST (according to DIN standard 19266, 2015)

Table 21 NIST buffer solutions

Temp. (°C)	NIST (according to DIN standard 19266, 2015-05)				
	pH 1.679	pH 4.005	pH 6.865	pH 9.180	pH 12.454
0	1.666	4.000	6.984	9.464	-
5	1.668	3.998	6.951	9.395	13.207
10	1.670	3.997	6.923	9.332	13.003
15	1.672	3.998	6.900	9.276	12.810
20	1.675	4.000	6.881	9.225	12.627
25	1.679	4.005	6.865	9.180	12.454
30	1.683	4.011	6.853	9.139	12.289
35	1.688	4.018	6.844	9.102	12.133
40	1.694	4.027	6.838	9.068	11.984
45	1.700	4.038	6.836	9.040	11.841
50	1.707	4.050	6.833	9.011	11.705
55	1.715	4.075	6.834	8.985	11.574
60	1.723	4.091	6.836	8.962	11.449
65	1.733	4.108	6.841	8.942	-
70	1.743	4.126	6.845	8.921	-
75	1.754	4.145	6.852	8.903	-
80	1.766	4.164	6.859	8.885	-



Temp. (°C)	NIST (according to DIN standard 19266, 2015-05)				
	pH 1.679	pH 4.005	pH 6.865	pH 9.180	pH 12.454
85	1.779	4.184	6.868	8.868	-
90	1.792	4.205	6.877	8.850	-
95	1.806	4.227	6.886	8.833	-



NOTICE

Update

The values of the individual buffers with the corresponding temperatures are kept as up to date as possible.

However, they may be changed by the respective manufacturers.

The NIST buffers are identical to the buffer solutions that are used in Chinese pharmacopoeia.

32.9.3 DIN (according to DIN standard 19267, 2012)

Table 22 *DIN buffer solutions*

Temp. (°C)	DIN (according to DIN standard 19267, 2012-08)					
	pH 1.09	pH 3.06	pH 4.65	pH 6.79	pH 9.23	pH 12.75
0	1.08	-	4.67	6.89	9.48	-
5	1.08	-	4.66	6.86	9.43	-
10	1.09	3.10	4.66	6.84	9.37	13.37
15	1.09	3.08	4.65	6.82	9.32	13.15
20	1.09	3.07	4.65	6.80	9.27	12.96
25	1.09	3.06	4.65	6.79	9.23	12.75
30	1.10	3.05	4.65	6.78	9.18	12.61
35	1.10	3.05	4.66	6.77	9.13	12.44
40	1.10	3.04	4.66	6.76	9.09	12.29
45	1.10	3.04	4.67	6.76	9.04	12.13
50	1.11	3.04	4.68	6.76	9.00	11.98
55	1.11	3.04	4.69	6.76	8.97	11.84
60	1.11	3.04	4.70	6.76	8.92	11.69
65	1.11	3.04	4.71	6.76	8.90	11.56

Temp. (°C)	DIN (according to DIN standard 19267, 2012-08)					
	pH 1.09	pH 3.06	pH 4.65	pH 6.79	pH 9.23	pH 12.75
70	1.11	3.04	4.72	6.76	8.88	11.43
75	1.12	3.04	4.74	6.77	8.86	11.30
80	1.12	3.05	4.75	6.78	8.85	11.19
85	1.12	3.06	4.77	6.79	8.83	11.08
90	1.13	3.07	4.79	6.80	8.82	10.99
95	-	-	-	-	-	-



NOTICE

Update

The values of the individual buffers with the corresponding temperatures are kept as up to date as possible.

However, they may be changed by the respective manufacturers.

32.9.4 Fisher

Table 23 Fisher buffer solutions (valid up to and including firmware version 5.916.0043)

Temp. (°C)	Fisher			
	pH 2.00	pH 4.00	pH 7.00	pH 10.00
0	-	4.01	7.13	10.34
5	1.98	3.99	7.10	10.26
10	1.98	4.00	7.07	10.19
15	2.02	3.99	7.05	10.12
20	2.00	4.00	7.02	10.06
25	2.00	4.00	7.00	10.00
30	2.00	4.01	6.99	9.94
35	2.02	4.02	6.98	9.90
40	2.01	4.03	6.97	9.85
45	2.01	4.04	6.97	9.81
50	2.01	4.06	6.97	9.78
55	-	4.07	6.97	9.74



Temp. (°C)	Fisher			
	pH 2.00	pH 4.00	pH 7.00	pH 10.00
60	-	4.09	6.98	9.70
65	-	4.11	6.99	9.68
70	-	4.13	7.00	9.65
75	-	4.14	7.02	9.63
80	-	4.16	7.03	9.62
85	-	4.18	7.06	9.61
90	-	4.21	7.08	9.60
95	-	4.23	7.11	9.60

Table 24 Fisher buffer solutions (valid from firmware version 5.916.0044)

Temp. (°C)	Fisher			
	pH 2.00	pH 4.00	pH 7.00	pH 10.00
0	-	4.01	7.13	10.34
5	1.98	3.99	7.10	10.26
10	1.98	4.00	7.07	10.19
15	2.02	3.99	7.05	10.12
20	2.00	4.00	7.02	10.06
25	2.00	4.00	7.00	10.00
30	2.00	4.01	6.99	9.94
35	2.02	4.02	6.98	9.90
40	2.01	4.03	6.97	9.85
45	2.01	4.04	6.97	9.81
50	2.01	4.06	6.97	9.78
55	-	4.07	6.97	9.74
60	-	4.09	6.98	9.70



NOTICE

Update

The values of the individual buffers with the corresponding temperatures are kept as up to date as possible.

However, they may be changed by the respective manufacturers.

32.9.5 Honeywell Fluka

Table 25 Buffer solutions Honeywell Fluka

Temp. (°C)	Fluka		
	pH 4.000	pH 7.000	pH 9.000
0	4.030	7.130	9.240
5	4.025	7.090	9.175
10	4.020	7.050	9.110
15	4.010	7.020	9.055
20	4.000	7.000	9.000
25	4.000	6.990	8.965
30	4.000	6.980	8.930
35	4.000	6.975	8.895
40	4.000	6.970	8.860
45	4.000	6.965	8.830
50	4.000	6.960	8.800
55	4.000	6.960	8.775
60	4.000	6.960	8.750
65	4.000	6.965	8.730
70	4.000	6.970	8.710
75	4.000	6.975	8.690
80	4.000	6.980	8.670
85	4.000	6.990	8.655
90	4.000	7.000	8.640

**NOTICE****Update**

The values of the individual buffers with the corresponding temperatures are kept as up to date as possible.

However, they may be changed by the respective manufacturers.

32.9.6 Mettler ToledoTable 26 *Mettler Toledo buffer solutions*

Temp. (°C)	Mettler Toledo				
	pH 2.00	pH 4.01	pH 7.00	pH 9.21	pH 11.00
0	2.03	4.01	7.12	9.52	11.90
5	2.02	4.01	7.09	9.45	11.72
10	2.01	4.00	7.06	9.38	11.54
15	2.00	4.00	7.04	9.32	11.36
20	2.00	4.00	7.02	9.26	11.18
25	2.00	4.01	7.00	9.21	11.00
30	1.99	4.01	6.99	9.16	10.82
35	1.99	4.02	6.98	9.11	10.64
40	1.98	4.03	6.97	9.06	10.46
45	1.98	4.04	6.97	9.03	10.28
50	1.98	4.06	6.97	8.99	10.10
55	1.98	4.08	6.98	8.96	-
60	1.98	4.10	6.98	8.93	-
65	1.98	4.13	6.99	8.90	-
70	1.99	4.16	7.00	8.88	-
75	1.99	4.19	7.02	8.85	-
80	2.00	4.22	7.04	8.83	-
85	2.00	4.26	7.06	8.81	-
90	2.00	4.30	7.09	8.79	-
95	2.00	4.35	7.12	8.77	-

**NOTICE****Update**

The values of the individual buffers with the corresponding temperatures are kept as up to date as possible.

However, they may be changed by the respective manufacturers.

32.9.7 Beckmann

Table 27 Beckmann buffer solutions (valid up to and including firmware version 5.916.0043)

Temp. (°C)	Beckmann		
	pH 4.00	pH 7.00	pH 10.01
0	4.00	7.12	10.32
5	4.00	7.09	10.25
10	4.00	7.06	10.18
15	4.00	7.04	10.12
20	4.00	7.02	10.06
25	4.00	7.00	10.01
30	4.01	6.99	9.97
35	4.02	6.99	9.93
40	4.03	6.98	9.89
45	4.05	6.98	9.86
50	4.06	6.97	9.83
55	4.08	6.98	-
60	4.09	6.98	-
65	4.11	6.99	-
70	4.12	6.99	-
75	4.14	7.00	-
80	4.16	7.00	-
85	4.18	7.01	-
90	4.19	7.02	-
95	4.21	7.03	-

Table 28 Beckmann buffer solutions (valid from firmware version 5.916.0044)

Temp. (°C)	Beckmann		
	pH 4.01	pH 7.00	pH 10.01
0	4.00	7.12	10.32
5	4.00	7.09	10.25
10	4.00	7.06	10.18
15	4.00	7.04	10.12
20	4.00	7.01	10.06
25	4.01	7.00	10.01
30	4.01	6.99	9.97
35	4.02	6.99	9.93
40	4.03	6.97	9.89
45	4.05	6.97	9.86
50	4.06	6.97	9.83



NOTICE

Update

The values of the individual buffers with the corresponding temperatures are kept as up to date as possible.

However, they may be changed by the respective manufacturers.

32.9.8 Radiometer Analytical

Table 29 Radiometer Analytical buffer solutions (valid up to and including firmware version 5.916.0043)

Temp. (°C)	Radiometer Analytical			
	pH 1.679	pH 4.005	pH 7.000	pH 9.180
0	1.666	4.000	7.118	9.464
5	1.668	3.998	7.087	9.395
10	1.670	3.997	7.059	9.332
15	1.672	3.998	7.036	9.276
20	1.675	4.001	7.016	9.225

Temp. (°C)	Radiometer Analytical			
	pH 1.679	pH 4.005	pH 7.000	pH 9.180
25	1.679	4.005	7.000	9.180
30	1.683	4.011	6.987	9.139
35	1.688	4.018	6.977	9.102
40	1.694	4.027	6.970	9.068
45	1.700	4.038	6.965	9.038
50	1.707	4.050	6.964	9.011
55	1.715	4.064	6.965	8.985
60	1.723	4.080	6.968	8.962
65	1.732	4.097	6.974	8.941
70	1.743	4.116	6.982	8.921
75	1.754	4.137	6.992	8.900
80	1.765	4.159	7.004	8.885
85	1.778	4.183	7.018	8.867
90	1.792	4.210	7.034	8.850
95	-	4.240	-	-

Table 30 Radiometer Analytical buffer solutions (valid from firmware version 5.916.0044)

Temp. (°C)	Radiometer Analytical			
	pH 1.679	pH 4.005	pH 7.000	pH 9.180
0	1.666	4.000	7.118	9.464
5	1.668	3.998	7.087	9.395
10	1.670	3.997	7.059	9.332
15	1.672	3.998	7.036	9.276
20	1.675	4.001	7.016	9.225
25	1.679	4.005	7.000	9.180
30	1.683	4.011	6.987	9.139
35	1.688	4.018	6.977	9.102
40	1.694	4.027	6.970	9.068
45	1.700	4.038	6.965	9.038
50	1.707	4.050	6.964	9.010



Temp. (°C)	Radiometer Analytical			
	pH 1.679	pH 4.005	pH 7.000	pH 9.180
55	1.715	4.064	6.965	8.985
60	1.723	4.080	6.968	8.962
65	1.732	4.097	6.974	8.941
70	1.743	4.116	6.982	8.921
75	1.754	4.137	6.992	8.900
80	1.765	4.159	7.004	8.884
85	1.778	4.183	7.018	8.867
90	1.792	4.208	7.034	8.850

**NOTICE****Update**

The values of the individual buffers with the corresponding temperatures are kept as up to date as possible.

However, they may be changed by the respective manufacturers.

32.9.9 Baker

Table 31 Baker buffer solutions

Temp. (°C)	Baker			
	pH 4.00	pH 7.00	pH 9.00	pH 10.00
0	4.00	7.13	9.23	10.30
5	4.00	7.09	9.17	10.24
10	4.00	7.05	9.10	10.17
15	4.00	7.03	9.05	10.11
20	4.00	7.00	9.00	10.05
25	4.00	6.98	8.96	10.00
30	4.01	6.98	8.91	9.96
35	4.02	6.98	8.88	9.93
40	4.03	6.97	8.84	9.89
45	4.04	6.97	8.81	9.86
50	4.05	6.96	8.78	9.82

Temp. (°C)	Baker			
	pH 4.00	pH 7.00	pH 9.00	pH 10.00
55	4.07	6.96	8.76	9.79
60	4.08	6.96	8.73	9.76
65	4.10	6.97	8.71	9.74
70	4.12	6.97	8.69	9.72
75	4.14	6.98	8.68	9.70
80	4.16	6.98	8.66	9.68
85	4.19	6.99	8.64	9.66
90	4.21	7.00	8.62	9.64



NOTICE

Update

The values of the individual buffers with the corresponding temperatures are kept as up to date as possible.

However, they may be changed by the respective manufacturers.

32.9.10 Hamilton DURACAL

Table 32 Hamilton DURACAL buffer solutions

Temp. (°C)	Hamilton DURACAL			
	pH 4.01	pH 7.00	pH 9.21	pH 10.01
0	-	-	-	-
5	4.01	7.09	9.45	10.19
10	4.00	7.06	9.38	10.15
15	4.00	7.04	9.32	10.11
20	4.00	7.02	9.26	10.06
25	4.01	7.00	9.21	10.01
30	4.01	6.99	9.16	9.97
35	4.02	6.98	9.11	9.92
40	4.03	6.97	9.06	9.86
45	4.04	6.97	9.03	9.83
50	4.06	6.97	8.99	9.79

**NOTICE****Update**

The values of the individual buffers with the corresponding temperatures are kept as up to date as possible.

However, they may be changed by the respective manufacturers.

32.9.11 Precisa

Table 33 *Precisa buffer solutions*

Temp. (°C)	Precisa		
	pH 4.00	pH 7.00	pH 9.00
0	3.99	7.11	9.27
5	3.99	7.08	9.18
10	3.99	7.06	9.13
15	3.99	7.04	9.08
20	3.99	7.02	9.04
25	4.00	7.00	9.00
30	4.00	6.99	8.96
35	4.01	6.98	8.93
40	4.02	6.98	8.90
45	4.03	6.97	8.87
50	4.04	6.97	8.84
55	4.06	6.97	8.81
60	4.07	6.97	8.79
65	4.09	6.98	8.76
70	4.11	6.98	8.74
75	4.13	6.99	8.73
80	4.15	7.00	8.71
85	4.18	7.00	8.70
90	4.20	7.01	8.68
95	4.23	7.02	8.67

**NOTICE****Update**

The values of the individual buffers with the corresponding temperatures are kept as up to date as possible.

However, they may be changed by the respective manufacturers.

32.9.12 Merck CertiPUR 20 / Titrisol

Table 34 Merck CertiPUR 20 / Titrisol buffer solutions (valid up to and including firmware version 5.916.0043)

Temp. (°C)	Merck CertiPUR 20 / Titrisol				
	pH 2.00	pH 4.00	pH 7.00	pH 9.00	pH 12.00
0	2.01	4.05	7.13	9.24	12.58
5	2.01	4.04	7.07	9.16	12.41
10	2.01	4.02	7.05	9.11	12.26
15	2.00	4.01	7.02	9.05	12.10
20	2.00	4.00	7.00	9.00	12.00
25	2.00	4.01	6.98	8.95	11.88
30	2.00	4.01	6.98	8.91	11.72
35	2.00	4.01	6.96	8.88	11.67
40	2.00	4.01	6.95	8.85	11.54
45	2.00	4.00	6.95	8.82	11.435
50	2.00	4.00	6.95	8.79	11.33
55	2.00	4.00	6.95	8.76	11.185
60	2.00	4.00	6.96	8.73	11.04
65	2.00	4.00	6.96	8.715	10.97
70	2.01	4.00	6.96	8.70	10.90
75	2.01	4.00	6.96	8.68	10.80
80	2.01	4.00	6.97	8.66	10.70
85	2.01	4.00	6.98	8.65	10.59
90	2.01	4.00	7.00	8.64	10.48
95	-	4.00	7.02	-	-

Table 35 Merck CertiPUR 20 / Titrisol buffer solutions (valid from firmware version 5.916.0044)

Merck CertiPUR 20 / Titrisol					
Article no.	109433	109435	109477	109476	109462
Temp. (°C)	pH 2.00	pH 4.00	pH 7.00	pH 9.00	pH 11.00
0	2.01	4.05	7.13	9.24	11.45
5	2.01	4.04	7.07	9.16	11.32
10	2.01	4.02	7.05	9.11	11.20
15	2.00	4.01	7.02	9.05	11.10
20	2.00	4.00	7.00	9.00	11.00
25	2.00	4.01	6.98	8.95	10.90
30	2.00	4.01	6.98	8.91	10.81
35	2.00	4.01	6.96	8.88	10.72
40	2.00	4.01	6.95	8.79	10.64
45	2.00	4.00	6.95	8.82	10.56
50	2.00	4.00	6.95	-	10.48



NOTICE

Update

The values of the individual buffers with the corresponding temperatures are kept as up to date as possible. However, they may be changed by the respective manufacturers.

32.9.13 Merck CertiPUR 25

Table 36 Merck CertiPUR 25 buffer solutions (valid up to and including firmware version 5.916.0043)

Temp. (°C)	Merck CertiPUR (25°C)			
	pH 4.00	pH 7.00	pH 9.00	pH 10.00
0	-	-	-	-
5	4.05	7.09	9.22	10.22
10	4.04	7.08	9.16	10.16
15	4.02	7.04	9.10	10.10

Temp. (°C)	Merck CertiPUR (25°C)			
	pH	pH	pH	pH
	4.00	7.00	9.00	10.00
20	4.01	7.02	9.06	10.05
25	4.00	7.00	9.00	10.00
30	3.99	6.98	8.98	9.94
35	3.98	6.98	8.93	9.90
40	3.98	6.97	8.89	9.86
45	3.98	6.97	8.86	9.80
50	3.98	6.97	8.84	9.73

Table 37 Merck CertiPUR 25 buffer solutions (valid from firmware version 5.916.0044)

Article no.	Merck CertiPUR (25°C)			
	109445	109407	109408	109409
	pH	pH	pH	pH
Temp. (°C)	4.00	7.00	9.00	10.00
0	-	-	-	-
5	4.05	7.09	9.22	10.22
10	4.04	7.08	9.16	10.16
15	4.02	7.04	9.10	10.10
20	4.01	7.02	9.05	10.05
25	4.00	7.00	9.00	10.00
30	3.99	6.98	8.96	9.94
35	3.98	6.98	8.93	9.90
40	3.98	6.97	8.89	9.86
45	3.98	6.97	8.86	9.80
50	3.98	6.97	8.84	9.73



NOTICE

Update

The values of the individual buffers with the corresponding temperatures are kept as up to date as possible. However, they may be changed by the respective manufacturers.


32.10 Using AuditTrailViewer

When you work in accordance with the *FDA Guidance 21 CFR Part 11*, logs with detailed user activities (Audit Trail) are important. The *AuditTrailViewer* supports you in your efforts to compile, filter and export logs of this type as TXT files.

First you must install the *AuditTrailViewer* on your PC. Afterwards, you must create a backup of your Ti-Touch (see chapter 12.3, page 142). The Audit Trail is generated automatically when you create the backup. The file name of the Audit Trail is **log.madt**. In the following you will find detailed information about how you must proceed.

32.10.1 Installing AuditTrailViewer

You will find the file for installing the *AuditTrailViewer* on the USB flash drive provided.

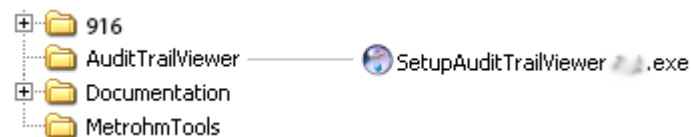

Proceed as follows:

- 1 Double-click on the file **SetupAuditTrailViewer 1.2.2.exe**.
The dialog **Welcome to the AuditTrailViewer Setup Wizard** is displayed.
- 2 Click on **[Next >]**.
The **License Agreement** dialog is displayed.
- 3 Click on **[I Agree]** to accept the contractual conditions.
The **Choose Install Location** dialog is displayed. A memory location for the *AuditTrailViewer* is suggested in this dialog.
- 4 If you accept the suggested memory location, click on **[Next >]**.
or
Select a different memory location and click on **[Next >]**.
The **Confirm Installation** dialog is displayed.
- 5 Click on **[Install]** to install the *AuditTrailViewer*.

As soon as the installation is completed, the dialog **Installation Complete** is displayed.

6 Click on **[Next >]**.

The dialog **Completing the AuditTrailViewer Setup Wizard** is displayed.

7 If the *AuditTrailViewer* should not be started immediately, deactivate the check box **Run AuditTrailViewer**.


8 Click on **[Finish]**.

If the preceding instruction step has been skipped, then the *AuditTrailViewer* will start.

32.10.2 Opening the Audit Trail

Create a backup of the data and of the settings of your Ti-Touch (see chapter 12.3, page 142).

1 Open the backup on your PC.

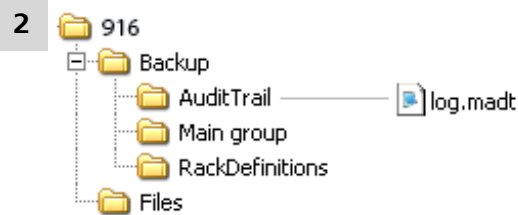

Open the file **log.madt** with the *AuditTrailViewer*.



AuditTrailViewer dialog window









File Filter Info					
No.	Date	User	Category	Action	Details
1	2011-10-31 11:30:22	Johnson	Method	New	01 Dynamic Titration pH
2	2011-10-31 11:30:34	Johnson	Method	Delete command	New method V0 03 REPORT
3	2011-10-31 11:30:40	Johnson	Method	Edit	New method V0 01 DET pH Stirrer off
4	2011-10-31 11:30:43	Johnson	Method	Edit	New method V0 01 DET pH Titrant not defined
5	2011-10-31 11:30:43	Johnson	Method	Edit	New method V0 01 DET pH Titrant
6	2011-10-31 11:30:48	Johnson	Method	Load result template	New method V0 Content (g/L)
7	2011-10-31 11:30:52	Johnson	Method	Edit	New method V0 Calc. formula EP1*CONC*TITER*3/C00
8	2011-10-31 11:31:08	Johnson	System	Message	003-908 Number of autostarts Yes/OK
9	2011-10-31 11:31:14	Johnson	Method	Start	Start key pressed
10	2011-10-31 11:31:15	Johnson	Method	Start	New method V0 1.0 g
11	2011-10-31 11:31:28	Johnson	Method	Manual stop	New method V0
12	2011-10-31 11:31:36	Johnson	System	Message	002-907 Method modified Yes/OK
13	2011-10-31 11:31:36	Johnson	Security	Logout	
14	2011-10-31 11:31:55	Meier	Security	Login message	Wrong password
15	2011-10-31 11:31:57	Johnson	System	Message	002-102 Wrong password Yes/OK
16	2011-10-31 11:31:58	Meier	Security	Log in	
17	2011-10-31 11:32:27	Meier	Method	Load	KFT 2 V1 Internal memory
18	2011-10-31 11:32:32	Meier	Security	Logout	
19	2011-10-31 11:32:45	Meier	Security	Login message	Wrong password
20	2011-10-31 11:32:47	Meier	System	Message	002-102 Wrong password Yes/OK
21	2011-10-31 11:33:01	Meier	Security	Change password	
22	2011-10-31 11:33:03	Meier	Security	Login message	Wrong password
23	2011-10-31 11:33:05	Meier	System	Message	002-102 Wrong password Yes/OK
24	2011-10-31 11:33:08	Meier	Security	Log in	
25	2011-10-31 11:33:12	Meier	Method	Start	Start key pressed
26	2011-10-31 11:33:13	Meier	Method	Start	KFT 2 V1 1.0 g
27	2011-10-31 11:33:14	Meier	System	Message	009-111 Stirrer missing Yes/OK
28	2011-10-31 11:33:15	Meier	Security	Logout	
29	2011-10-31 11:33:27	Chang	Security	Log in	
30	2011-10-31 11:33:32	Chang	Method	New	01 Dynamic Titration pH
31	2011-10-31 11:33:38	Chang	Method	Delete command	New method V0 03 REPORT
32	2011-10-31 11:33:43	Chang	Method	Edit	New method V0 01 DET pH Start volume 5 mL
33	2011-10-31 11:33:48	Chang	Method	Load result template	New method V0 Content (g/L)

32.10.3 Contents of the Audit Trail

In the following you will find a description of the dialog window and of the menu bar of the *AuditTrailViewer*.

Table 38 Dialog window

Column	Content
No.	Each entry is numbered in sequence.

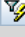Column	Content
Symbol	Classification of the entry: <ul style="list-style-type: none"> ▪  Actions that are neither relevant to safety nor alter the determination data. ▪  Actions, e.g. changes to the login options and recalculation of determinations. ▪  Errors that occur, e.g. entry of an incorrect password.
Date	Precise time of the event.
User	User who triggered the action.
Category	Category to which the entry belongs.
Action	Designation of the action.
Details	Details of the action.
Menu bar	
 Print	Print Audit Trail. An Audit Trail can be printed by means of the  symbol.
 Update	Update view. The list is updated automatically only at the time the dialog window is opened.
 Quick filter	Filter the Audit Trail.
 Show all	Display all entries once again.



32.10.4 Filtering the Audit Trail

The entries in the Audit Trail can be filtered. You can set the following filter criteria:

- **Date**
- **User**
- **Category**
- **Action**

Proceed as follows:

- 1 Click in a cell that contains the desired filter criterion (e.g. **Category** = **Method**).

- 2 Click on the  icon or the menu item **Filter ► Quick filter**.
The only entries which will be displayed are those which correspond to the desired filter criterion.
- 3 Click on the  icon or the menu item **Filter ► Show all**.
All of the entries are displayed once again.

32.10.5 Exporting the Audit Trail

You can export the Audit Trail as .txt file. This way, the Audit Trail can be archived and is available as required for inspections.

Proceed as follows:

- 1 Use the menu item **File ► Export** to open the dialog window for the export.
- 2 Enter the path and file name with the extension **.txt**.
- 3 Click on **[Open]** and then on **[Export]**.
The Audit Trail is exported.

32.11 Diagnosis

System ► Diagnosis

The electronic and mechanical functional groups of Metrohm instruments can and should be checked by specialist personnel from Metrohm as part of a regular maintenance schedule. Please ask your local Metrohm representative regarding the precise terms and conditions involved in concluding a corresponding maintenance agreement.

The following functions are available:

- **LCD test**
Check the display for faulty pixels (*see chapter 32.11.1, page 473*).
- **Format storage medium**
Format the external storage medium (*see chapter 32.11.2, page 473*).
- **Remove storage medium**
Remove the external storage medium safely (*see chapter 32.11.3, page 474*).
- **Touch adjustment**
Adjust the touch-sensitive screen (*see chapter 32.11.4, page 474*).

- **Touch screen test**
Test the function of the touch-sensitive screen (see chapter 32.11.5, page 475).
- **Software update**
Update the software of Ti-Touch and of the control instruments (see chapter 32.11.6, page 476).
- **Service**
Special functions for the service technician (see chapter 32.11.7, page 479).

32.11.1 LCD test

System ► Diagnosis ► LCD test

You can use the LCD test to check the display for faulty pixels. To do this, various test images are displayed one after the other.



NOTICE

The [↩] fixed key is used to display the preceding test picture; the test can be stopped at any time with the [🏠] fixed key.

Proceed as follows:

- 1 ▪ Tap on **[LCD test]**.
The entire display turns white.
- 2 ▪ Check the display for pixel errors and other irregularities.
 - Continue the test with [▶].
 Other colors and patterns will be displayed one after the other.
- 3 ▪ Repeat step 2 until the **System / Diagnosis** dialog is displayed again.
- 4 ▪ Contact your Metrohm representative in the event of faults or irregularities.

32.11.2 Formatting an external storage medium

System ► Diagnosis ► Format storage medium

With this function, you have the option of formatting an external storage medium directly on the 916 Ti-Touch (quick formatting).

The following file systems can be selected:

- **FAT**

- **FAT32**
- **ExFAT**



CAUTION

If you format the storage medium with the **ExFAT** file system, it is possible that it will no longer be recognized by a PC. Read the pertinent information from Microsoft Support at <http://support.microsoft.com>.

Do not use this file system unless you need to be able to save more than 999 files in a single group.

32.11.3 Removing an external storage medium

System ► Diagnosis ► Remove storage medium

As long as the stored data is not being accessed (reading/saving data), you can plug in and disconnect the storage medium without difficulty at any time. The **[Remove storage medium]** function offers additional protection. This ensures that the storage medium cannot be removed while any data is being transferred.

32.11.4 Adjusting the touch screen

System ► Diagnosis ► Touch adjustment

With time the touch-sensitive screen may not register touches at their precise position. You can readjust the screen in this dialog.

Proceed as follows:

1 In the main dialog, tap on **[System]**.

2 Tap on **[Diagnosis]**.

3 Tap on **[Touch adjustment]**.

A message in English and a cross hair will appear in the center of the screen.

4 Touch a stylus (a special pen-shaped instrument for operating devices with touch screens) to the center of the cross hair and maintain this position until the cross hair moves to a new position on the screen.

5 Repeat this process as long as the cross hair continues to be repositioned on the screen.

A message in English will appear.

6 Tap at any location on the screen to confirm the adjustment.

7 Press [] to finish the adjustment.

32.11.5 Testing the touch screen

System ► Diagnosis ► Touch screen test

You can use this test to check the function of the touch-sensitive monitor screen.

A calibration is not possible with this test. If the Touch screen does not work correctly anymore, it will need to be replaced. For that please contact your Metrohm representative.



NOTICE

You can use the [] fixed key to stop the test at any time.

Proceed as follows:


1 Tap on **[Touch screen test]**.


A red square is displayed at the upper left.

2 Tap on the square as close to the center as possible.

The actual detected point of contact will be displayed by a cross hair, and the deviation from the center will also be displayed on the lower edge of the image.

The next red square will also be displayed.

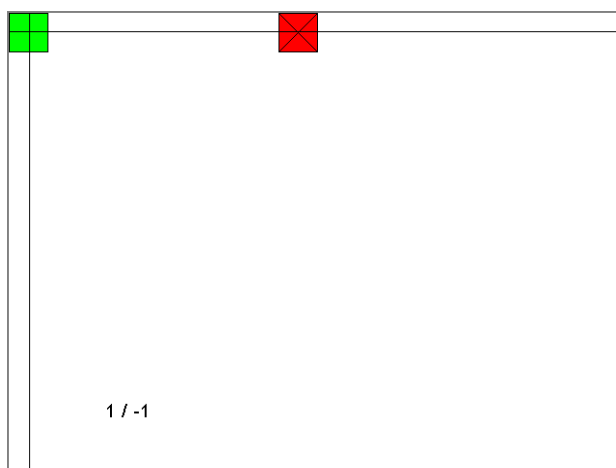

- 3 Repeat step 2 until all of the areas in the display have been tested.
A list with fixed keys will appear.
- 4 Tap on all of the fixed keys one after the other in any sequence.
The result is displayed in each case with a short information text, e.g. **[Print] key OK**.
- 5 Finish Touch screen test with [].

32.11.6 Software update (loading program versions and language files)

System ► Diagnosis ► Software update

New program versions or language files can be loaded from an external storage medium (e.g. USB flash drive). The corresponding files must be stored on the storage medium in the **916/SwUpdates** directory (see "Directory structure", page 141).



NOTICE

If you wish to update a connected Sample Processor together with the Ti-Touch, then do not fail to update the control software of the Ti-Touch first and then afterwards the device program (firmware) of the Sample Processor.



NOTICE

The memory will be reformatted when updating from software version 5.916.0043.

Before carrying out the software update, make a backup in order to save your data and the system settings (see chapter 12.3, page 142).

Use a separate USB flash drive for the software update. It must be removed from the instrument after the successful software update.

Program files

The files are instrument-specific. The file names are structured as follows:

- Control software for 916 Ti-Touch:
 - 5XXXyyyy.BIN**
 - 5XXXyyyy.MBIN**
 - XXX = instrument type (i.e. "916" for the 916 Ti-Touch)
 - yyyy = program version

- Firmware of the control instrument:
5XXXyyyy.BIN
 - XXX = instrument type (e.g. 814 for the 814 USB Sample Processor)
 - yyyy = program version

Language files

Language files may have the following content:

- one or more additional dialog languages
- the online help for one or more additional dialog languages
- supplements for existing dialog languages/online help documents

They can be recognized by means of the two-digit language code in the file name. The file name has the following structure:

- **5XXXyZZZML.BIN**
5XXXyZZZML.MBIN
 - XXX = instrument type (i.e. "916" for the 916 Ti-Touch)
 - y = version number of the language package
 - ZZZ = program version

32.11.6.1 Carrying out a software update



CAUTION

Make sure that the power supply is guaranteed during the entire update process. Otherwise there is the chance that the instrument will no longer be able to be switched on and that it will need to be sent in for repair.



NOTICE

Make sure that no USB/RS-232 adapter is connected!

Proceed as follows:


1 Copying files to an external storage medium



NOTICE

In the case of the program versions and language files for the Ti-Touch, it is imperative that both the BIN file and the associated MBIN file be copied.

- Copy the files to the "SwUpdates" directory.
 - Copy also the "MetrohmTools" folder with the "MUpdate-Tool.exe" file to the external storage medium.
- Make sure that the file is stored at the top level of the external storage medium. This also applies to the "916" folder.

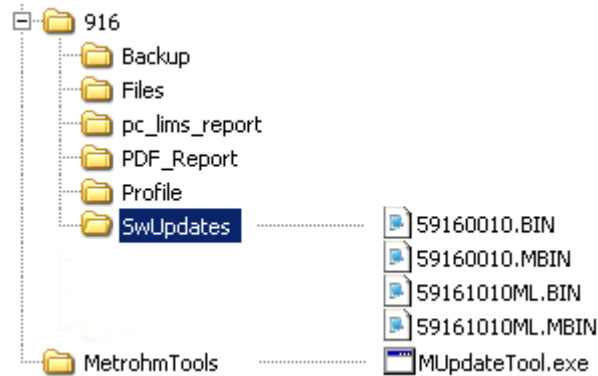

For the software update from version 5.916.0043, an additional folder that is on the same level as Metrohm-Tools with the name SRVPRF900 containing the file St900_SrPrf.cfg is required

If this structure is not maintained, then the files will not be found, because the directories are accessed directly during the update process.

- Plug in the external storage medium at the Ti-Touch.

2 Opening the update dialog

- Under **System ► Diagnosis**, tap on the button **Software update**.

3 Selecting the instrument

- Tap on the **Control device** list box and select the instrument to be updated.

4 Selecting the file

- Tap on the **Binary file** list box.
The selection list with the BIN files saved in the "SwUpdates" directory is opened.
- Select the required file.
- Tap on **[Select]**.
- Update of the Ti-Touch: Continue with Step 5.
Update of the Sample Processor: Continue with Step 6.

5 Starting the update (916 Ti-Touch)

- Tap on the **[Start]** button (NOT on the [▶] fixed key).
- The message **023-102 Program update** is displayed.
- Confirm the message with **[Yes]**.
The update process is started, it runs automatically. The device is automatically switched off and back on during this process, possibly several times. No user intervention is required.
- **After successfully updating the software, remove the external storage medium (USB flash drive) from the instrument.**

6 Starting the update (control instruments)

- Tap on the **[Start]** button (NOT on the [▶] fixed key).
If **Result = No errors** is displayed, then the update has been successful.
- **Remove the external storage medium (USB flash drive) from the instrument.**
- Switch the 916 Ti-Touch off and then back on again.

32.11.7 Service

System ► Diagnosis ► Service

Access to the service functions is password-protected and only accessible to Metrohm service technicians. We recommend that you have maintenance performed on the system regularly by the Metrohm Service Department. You can automatically monitor the time interval after which the next service visit is due (*see chapter 13.5.1, page 154*).

32.12 Arithmetic algorithms in the Ti-Touch

Some of the algorithms and numerical formats used in the device software are explained in this chapter.

Numerical format

All floating point calculations use the IEEE 754 (1985) numerical format, either in "single precision" (32 Bit) or "double precision" (64 Bit).

Rounding off process

Measured values and results are rounded to the defined number of decimal places (commercial rounding, in accordance with the US Pharmacopeia USP). If the digit at the first dropped decimal place is **1, 2, 3 or 4**, then it will be rounded off; if this digit is **5, 6, 7, 8 or 9**, then it will be rounded up. Negative digits will be rounded in accordance with their amount, i.e. away from zero.

Examples:

2.33 yields 2.3

2.35 yields 2.4

2.47 yields 2.5

-2.38 yields -2.4

-2.45 yields -2.5

Statistics

The arithmetic mean value and the absolute and relative standard deviations of results R, endpoints EP and variable C are calculated:

You can statistically evaluate a maximum of nine results ($1 \leq k \leq 9$) calculated in a determination. A statistical series can contain a maximum of 20 determinations ($1 \leq n \leq 20$).

The following convention applies to the subsequent formulas:

$1 \leq n \leq 20$ and $1 \leq k \leq 9$.

Mean value:

$$\bar{x}_k = \frac{1}{n} \cdot \sum_{i=1}^n R_{k,i}$$

Absolute standard deviation:

$$Sabs_k = +\sqrt{\frac{\sum_{i=1}^n (R_{k,i} - \bar{x}_k)^2}{n-1}}$$

Relative standard deviation (in %):

$$Srel_k = 100 \cdot \frac{Sabs_k}{\bar{x}_k}$$

Explanations

The individual values are incorporated in the statistics in the selected Precision (Round, Truncate or Full precision).

7 significant places are yielded when the 32 bit numerical format is applied for the floating point number in decimal presentation. For the 64 bit numerical format, there are 15 significant places.

The accuracy can be controlled by the selection of the prefix of the unit (milli, micro) and the number of decimal places.

Example (64-bit numerical format):

The result displayed, **1234.56789158763 mg/L**, has 15 significant places. It should be rounded off to three decimal places according to the above rounding-off process:

- **1234.568 mg/L.**

If the same result is expressed in "g/L" (**1.23456789158763 g/L**), and is also rounded off to three decimal place, this yields

- **1.235 g/L.**

I.e. you obtain the lowest losses in accuracy with rounding when you select the application and the numerical format in such a way that the numbers displayed have as many places before the decimal point as possible.

A complete recalculation of the statistics using a pocket calculator or PC calculation programs may exhibit deviations. This can be explained by the different binary numerical formats used by these computers.



NOTICE

The above losses of accuracy by rounding off in the range of significant places are only relevant theoretically. Most of the time they are lower by several orders of magnitude than – as an example – the uncertainties resulting from weighing out the sample.

33.2 Measuring inputs

The measuring cycle is 100 ms for all measuring modes.

33.2.1 Potentiometry

A high-ohm measuring input (**Ind.**) for pH, metal or ion-selective electrodes and a measuring input for separate reference electrodes (**Ref.**).

<i>Input resistance</i>	$> 1 \cdot 10^{12} \Omega$ (under reference conditions)
<i>Offset current</i>	$< 1 \cdot 10^{-12} \text{ A}$ (under reference conditions)
<i>Measuring mode</i>	
<i>pH</i>	
<i>Measuring range</i>	-13 to +20 pH
<i>Resolution</i>	0.001 pH
<i>Measuring accuracy</i>	$\pm 0.003 \text{ pH}$ (± 1 digit, without sensor error, under reference conditions)
<i>Measuring mode</i>	
<i>U</i>	
<i>Measuring range</i>	-1200 to +1200 mV
<i>Resolution</i>	0.1 mV
<i>Measuring accuracy</i>	$\pm 0.2 \text{ mV}$ (± 1 digit, without sensor error, under reference conditions)

33.2.2 Polarizer

One measuring input (**Pol.**) for polarizable electrodes.

<i>Measuring mode</i>	Determination with adjustable polarization current.
<i>I_{pol}</i>	
<i>Polarization current</i>	-120 to +120 μA (increment: 1 μA) -125 to -121 μA / +121 to +125 μA : non-guaranteed values, dependent on reference voltage +2.5 V
<i>Measuring range</i>	-1200 to +1200 mV
<i>Resolution</i>	0.1 mV
<i>Measuring accuracy</i>	$\pm 0.2 \text{ mV}$ (± 1 digit, without sensor error, under reference conditions)



<i>Measuring mode</i>	Determination with adjustable polarization voltage.
<i>Upol</i>	
<i>Polarization voltage</i>	–1200 to +1200 mV (increment: 10 mV) –1250 to –1210 mV / +1210 to +1250 mV: non-guaranteed values, dependent on reference voltage +2.5 V
<i>Measuring range</i>	–120 to +120 μ A
<i>Resolution</i>	0.01 μ A
<i>Measuring accuracy</i>	–

33.2.3 Temperature

A measuring input (**Temp.**) for temperature sensors of the Pt1000 or NTC type with automatic temperature compensation.

R (25 °C) and B value can be configured for NTC sensors.

<i>Measuring range</i>	
<i>Pt1000</i>	–150 to +250 °C
<i>NTC</i>	–5 to +250 °C (For an NTC sensor with R (25 °C) = 30'000 Ω and B (25/50) = 4'100 K)
<i>Resolution</i>	
<i>Pt1000</i>	0.1 °C
<i>NTC</i>	0.1 °C
<i>Measuring accuracy</i>	
<i>Pt1000</i>	\pm 0.2 °C (Applies for measuring range –20 to +150 °C)
<i>NTC</i>	\pm 0.6 °C (Applies for measuring range +10 to +40 °C)

33.3 Built-in magnetic stirrer (depending on the product version)

Rotational speed

Maximum rotational speed $\pm(1700\text{--}1900)$ rpm

Rotational speed settings ± 15 steps

Change in rotational speed per step $\pm(115\text{--}125)$ rpm

Power consumption 4 W

Fuse Electronic overload protection

33.4 Interfaces

USB connector Type A, for connecting USB devices.

MSB connector For connecting dosing devices, one additional stirrer or a Remote Box.

iConnect connector For connecting an 854 iConnect with iTrode.

Stirrer connector (depends on product version; only for instruments without built-in magnetic stirrer) For connecting an 802 Stirrer.

Ethernet connector For connection to a data network (LAN).



33.5 Power supply

Instrument

Input voltage 24 V DC

Max. power consumption 2.0 A

External power supply unit 6.2164.010, LPS conformity according to UL60950-1

Input current 1.5 A

Max. output current 2.7 A

33.6 Declarations of Conformity

Up-to-date information on the *EU conformity* and the *requirements for electromagnetic compatibility* for your product can be found on the Internet under <https://www.metrohm.com/en-us/products-overview/>.

Using the article number, you can download the *Declaration of conformity for ...* under the documents of the product.

33.7 Ambient temperature

Nominal function range +5 to +45 °C
at max. 80% relative humidity, non-condensing

Storage +5 to +45 °C

33.8 Reference conditions

Ambient temperature +25 °C (±3 °C)

Relative humidity ≤ 60%

Operating temperature status Instrument in operation at least 30 min

Validity of the data After adjustment

33.9 Dimensions

<i>Width</i>	193 mm
<i>Height</i>	
<i>without support rod</i>	135 mm
<i>with support rod</i>	430 mm
<i>Depth</i>	412 mm
<i>Weight (including power supply unit)</i>	4900 g
<i>Material</i>	
<i>Housing</i>	Poly(butylene terephthalate) (PBT)
<i>Base</i>	Steel, stainless, coated

33.10 Storage capacity

<i>Random access memory</i>	128 MB
<i>Internal memory</i>	1 GB of which 900 MB available.
<i>Required memory per method</i>	3–6 KB

34 Accessories

Up-to-date information on the scope of delivery and optional accessories for your product can be found on the Internet. You can download this information using the article number as follows:

Downloading the accessories list

- 1** Enter <https://www.metrohm.com/> into your Internet browser.
- 2** Enter the article number (e.g. **916**) into the search field.
The search result is displayed.
- 3** Click on the product.
Detailed information regarding the product is shown on various tabs.
- 4** On the **Included parts** tab, click on **Download the PDF**.
The PDF file with the accessories data is created.



NOTICE

Once you have received your new product, we recommend downloading the accessories list from the Internet, printing it out and keeping it together with the manual for reference purposes.

Index

800 Dosino	16
801 Stirrer	17
803 Ti Stand	17
804 Ti Stand	17
805 Dosimat	16

A

Acoustic signals	59
ADD	410
Audit Trail	57
Archive	472
Export	472
Filter	471
Open	469
AuditTrailViewer	
Installing	468
Automatic system test	145
Automation	423
Axial distance	110

B

Backup	142
Interval	154
Balance	19, 128, 445
Barcode reader	
Connect	21
Beaker radius	
Sample vessel	117
Special beaker	122
Beaker sensor	
Sample vessel	118
Special beaker	122
BEEP	434
Break point	
Evaluation	384
Buffer series	452

C

CAL	367
CALC	387
CALC LIVE	395
Calculation	387
Add	388
Delete	388
Edit	388
Load template	388
Note	392
Result options	393
Calibration	
History	90

Calibration buffer	452
Calibration buffers	175
Calibration data	88, 233
Limit values	92
Calibration interval	93
Calibration parameters	
CAL pH	367
Calibration position 1	118
Cell	
SET	333
Command	
ADD	410
BEEP	434
CAL	367
CALC	387
CALC LIVE	395
CONTROL RS	422
CTRL	420
DET	291
ELT	373
EMPTY	409
END	434
EVAL BREAK	384
EVAL FIX-EP	380
EVAL MIN/MAX	382
EVAL pK/HNP	381
EVAL RATE	386
Insert	186
LIFT	425
LQH	415
MAT	355
MEAS	362
MET	307
MOVE	423
PREP	408
PUMP	426
RACK	427
REPORT	403
REQUEST	433
SAMPLE	427
SCAN	418
SCAN RS	421
SET	324
SIGN	434
STAT	337
STIR	431
SUBSEQ	428
WAIT	432
Command list	185

Common variable	
Delete	156
Edit	156
Properties	158
Validity	159
Communication	418
Conditioning	
SET	331
Connect	
Balance	19
Barcode reader	21
Dosing device	16
MSB devices	15
PC keyboard	20
Power supply unit	11
Printer	18
Remote Box	17
Sample Processor	21
Stirrer or titration stand	17
USB devices	18
USB hub	18
Control	201
Control instrument	103
Parameters	
304, 320, 333, 350, 357, 364,	
371, 375, 411, 417	
Control parameters	
SET	325
STAT	338
Control Remote Box	97
CONTROL RS	422
CTRL	420
Curve options	243
Curves	242
Custom logo	177

D

Delete	
Signature	200
Statistics data	203
DET	1, 31, 291
Determination	
Calibration data	233
Cancel	222
Carry out	220
Data	229
Define filters	240
Delete signatures	200
Endpoint details	232

List of saved determinations	238
Load	237
Measuring point list	232
Message	233
Properties	234
Recalculate	244
Reevaluate	244
Save	242
Save automatically	197
Send PC/LIMS report	197
Sign	199
Titrant data	233
Determination run	220
Device	
Add	96
Balance	128
Delete	95
Edit	95
New	95
Printer	124
Sample Processor	107
Ti-Touch titrator	103
USB/RS-232 adapter	130
Device manager	95
Diagnosis	472
Dialog language	41
Dialog options	43
Dialog settings	
System-specific	43
Differential potentiometry	29
Directory structure	141
Dosing	269
Continuous	272
Fixed volume	273
Dosing curve	242
Dosing device	
Connect	16
Parameters	
306, 322, 335, 351, 358, 411,	
417	
Dosing key	442
Dosing parameters	
ADD	410
MAT	356
Dosing port	66
Dosing rate	
Maximum	443, 444
Dosing unit	
Data	65
GLP test	75
Preparing parameters	66
Replace reagent	277
Tubing parameters	68

E	
E-mail	98
Electrode	84
Parameters	
304, 320, 333, 350, 357, 365,	
371, 375	
Electrode slope	
Intervention limit	90
Warning limit	90
Electrode test parameters	374
Electrode type	
Custom	178
Electrostatic charge	7
ELT	373
EMPTY	409
Default parameters	443
Emptying	277
Default parameters	443
END	434
Endpoint details	232
ERC	304, 320
Ethernet connector	30
EVAL BREAK	384
EVAL FIX-EP	380
EVAL MIN/MAX	382
EVAL pK/HNP	381
EVAL RATE	386
Evaluation	
Break point	384
DET	303
Fixed endpoint	380
HNP	381
Maximum	382
MET	319
Minimum	382
pK value	381
Potentiometric	299, 315
Rate	386
Set windows	300, 316, 385
Evaluations	379
Exchange unit	
Data	71
GLP test	75
Preparing parameters	73
Tubing parameters	74
Expiry date	
Sensor	87
Titrant	64
External position	
Swing angle	114
F	
Favorite	205
Add	206

Delete	206
Edit	206
FDA Guidance 21 CFR Part 11	
.....	46, 57
File	
Copy	138
Properties	139
Rename	138
File manager	136
Filling	277
Filling rate	
Maximum	443, 444
Filter	
Determination	240
Result table	254
Fixed endpoint	
Evaluation	380
Fixed key	9
Format storage medium	473
Formula editor	397
G	
GLP manager	144
GLP test "Measurement"	146
GLP test "Titration"	146
System validation	150
Test tool	145
GLP test	
Dosing unit	75
Exchange unit	75
Sensor	90
H	
Half neutralization potential	
Evaluation	381
HNP	
Evaluation	381
Hysteresis	345
I	
ID profile	
Create	51
Identification profile	
Create	51
Initialization	435
Input line	
Template	170
Installation	
Set up instrument	11
Instrument	
Configure	96
Switch off	36
Switch on	35
Ti-Touch	97

- Intervention limit
 Electrode slope 90
 pH(0) 90
 Titer 79
- K**
- Key functions 132
 Keyboard
 Connect 20
- L**
- LCD test 473
 LIFT 425
 Lift rate 284
 Limit values
 Calibration data 92
 Sample size 194
 Liquid Handling parameters 415
 Live display 225
 Live modifications 223
 Live parameters 227
 Live result
 Note 396
 Load
 Determination 237
 Sample table 215
 Load result template 388
 Login
 Options 52
 Logout automatically 53
 LQH 415
- M**
- Magnetic stirrer
 Connect 17
 Manual control 265
 Continuous dosing 272
 Dosing 269
 Dosing fixed volume 273
 Emptying 277
 Filling 277
 Measuring 267
 Preparing 276
 Remote lines 280
 Sample Processor 281
 Stirring 278
 Troubleshooting 439
 Manual Dosing Controller 442
 MAT 2, 33, 355
 Max. swing angle 112
 Maximum evaluation 382
 Maximum stroke path 109
 MEAS 362
 Measured value
 Display resolution 59
- Measurement curve 242
 Measuring 267
 Measuring parameters
 MEAS 362
 Measuring point list 232
 Message
 Determination 233
 Send e-mail 98
 MET 2, 31, 307
 Method 182
 Command list 185
 Create 182
 Delete signatures 200
 Edit 185
 Insert command 186
 Load 184
 Note 196
 Options 187
 Properties 195
 Sample data 191
 Save 183
 Save as favorite 205
 Sign 199
 Start options 188
 Stop options 190
 Troubleshooting 437
 Metrohm Serial Bus MSB, see also
 "MSB" 15
 Min. beaker radius 109
 Minimum evaluation 382
 Miscellaneous commands 431
 Monitoring
 Calibration interval 93
 Limit values calibration data 92
 Limit values sample size 194
 STAT 344
 MOVE 423
 MSB
 Connect devices 15
- N**
- Network connection 30
 Note
 Method 196
 Numerical input 39
- O**
- Operation
 General 36
 Output line
 Template 172
- P**
- Paper size 128
- Parameters
 ADD 410
 BEEP 434
 CAL 367
 CALC 387
 CALC LIVE 395
 CONTROL RS 422
 CTRL 420
 DET 291
 ELT 373
 EMPTY 409
 END 434
 EVAL BREAK 384
 EVAL FIX-EP 380
 EVAL MIN/MAX 382
 EVAL pK/HNP 381
 EVAL RATE 386
 LIFT 425
 LQH 415
 MAT 355
 MEAS 362
 MET 307
 Modify live 227
 MOVE 423
 PREP 408
 PUMP 426
 RACK 427
 REPORT 403
 REQUEST 433
 SAMPLE 427
 SCAN 418
 SCAN RS 421
 SET 324
 SIGN 434
 STAT 337
 STIR 431
 SUBSEQ 428
 WAIT 432
- Password
 Forgotten 55
 Options 54
 PC keyboard
 Connect 20
 PC/LIMS report
 Save 99
 Send 197
 PDF settings 125
 Peripheral devices 95
 pH(0)
 Intervention limit 90
 Warning limit 90
 Pins 446
 pK value
 Evaluation 381

- Potentiometric evaluation
 DET 299
 MET 315
- Power supply unit
 Connect 11
- PREP 408
 Default parameters ... 443, 444
 Parameters for the dosing unit
 66
 Parameters for the exchange
 unit 73
- Preparing 276
 Default parameters ... 443, 444
 Parameters for the dosing unit
 66
 Parameters for the exchange
 unit 73
- Printer 18, 124
 PDF settings 125
- Printing 257
 Troubleshooting 438
- Program version
 Ti-Touch 97
- Properties
 Determination 234
 File 139
 Method 195
 Swing Head 110
 Tower 108
- PUMP 426
- R**
- Rack
 Copy 115
 Delete 115
 Edit 115
 Load 115
- RACK 427
- Rack adjustment 123
- Rack name 108
- Rack offset 118
- Rate evaluation 386
- Reagent 80
 Add 80
 Delete 80
 Edit 80
 Replace 277
- Reagent monitoring 81
- Reagent replacement 82
- Reason 56
 Modification option 56
- Recalculation 244
- Reevaluation 244
- Remote
 Input 446
 Interface 446
 Output 447
- Remote Box
 Connect 17
 Pin assignment 446
- Remote lines
 Manual control 280
- Report 257
 Print 257
- REPORT 403
- Report header 177
- REQUEST 433
- Resolution
 Measured value display 59
- Result 229
 Calculation formula 391
 Limits 392
 Save as a common variable
 160
 Template 166
- Result table 251
 Define filters 254
 Delete automatically 255
 Save 256
- Result variable 392
 As parameter setting 449
- Robotic arm 112
 Max. swing angle 112
 Rotation offset 113
 Swing direction 113
 Swing increment 112
 Swing offset 112
 Swing radius 113
 Swing rate 286, 288, 425
- Rotation increment 110
- Rotation offset 113
- Routine dialog 45
 Standard configuration 45
- S**
- Safety instructions 6
- SAMPLE 427
- Sample assignment table 162
- Sample data 209
 Modify live 223
 Settings 191
- Sample identification
 Modify designation 191
 Modify the designation 219
- Sample identification list 162
- Sample Processor 281
 Connect 21
- Sample rack 115
- Sample series
 Perform 221
- Sample size
 Fixed 191
 Limit value 194
 Limits 191
 Send from a balance 445
- Sample table 212
 Save as favorite 205
 Troubleshooting 437
- Save
 Determination 242
 Result table 256
 Sample table 214
- SCAN 418
- SCAN RS 421
- Sensor 84
 Add 85
 Calibration data 88
 Calibration interval 93
 Delete 85
 Edit 85
 Expiry date 87
 GLP test 90
 Parameters
 304, 320, 333, 350, 357, 365,
 371, 375
 Working life 87
- Sensor list 84
- Serial number
 Dosing cylinder 66, 72
 Dosing unit 65
 Exchange unit 72
- Service 6, 479
- SET 2, 32, 324
- Set the date 42
- Set the time 42
- Set up instrument 11
- Sign
 Determination 199
 Method 199
- SIGN 434
- Signature
 Delete 200
- Software update 476
- Special beakers
 Edit 120
- Spooler 128
- Start conditions
 DET 291
 MET 307
 SET 324
 STAT 337

- Start options 188
- STAT 2, 32, 337
- Statistics 203, 246
- Delete 203
 - Troubleshooting 438
- Status LED 442
- STIR 431
- Stirrer
- Connect 17
 - Parameters
 - 306, 323, 336, 354, 361, 366,
 - 372, 376, 414
 - Stirring 278
 - Stirring rate 444
 - Stop 442
 - Stop conditions
 - DET 298
 - MET 314
 - SET 330
 - STAT 342
 - Stop criterion
 - SET 328
 - Stop options 190
 - SUBSEQ 428
 - Subsequence 428
 - Load 429
 - Note 431
 - Options 430
 - Save 429 - Supply voltage 6
 - Swing direction 113
 - Swing Head
 - Properties 110
 - Swing increment 112
 - Swing offset 112
 - Swing radius 113
 - Swing rate 286, 288, 425
 - Switch off 36
 - Switch on 35
 - System initialization 435
 - System settings 41
 - System test 145
- System validation 150
- T**
- Tandem dosing 352, 359, 412
- Template
- Calibration buffers 175
 - Electrode type 178
 - Input line 170
 - Output line 172
 - Report header 177
 - Result template 166
 - Sample assignment table .. 162
 - Sample identification list ... 162
- Templates 162
- Text input 39
- Ti-Touch 97
- Time zone 42
- Titer
- History 78
 - Intervention limit 79
 - Properties 77
 - Warning limit 79
- Titrant 60
- Add 61
 - Delete 61
 - Edit 61
 - Expiry date 64
 - Working life 64
- Titrant data 233
- Titrant list 60
- Titration curve 242
- Titration mode
- DET 1
 - MAT 2
 - MET 2
 - SET 2
 - STAT 2
- Titration parameters
- DET 294
 - MET 310
 - SET 329
 - STAT 341
- Titration stand 17
- Connect 17
- Touch screen 36
- Test 475
- Tower
- Properties 108
- Tubing parameters
- Dosing unit 68
 - Exchange unit 74
- U**
- Uninterrupted dosing
- 352, 359, 412
- Update 476
- USB
- Connect devices 18
- USB hub
- Connect 18
- USB/RS-232 adapter 130
- User profile
- Create 51
- V**
- Validation interval 152
- Validity
- Calibration 93
 - Common variable 159
 - Titer 77
- Valve disk
- Shift direction 71
- W**
- WAIT 432
- Warning limit
- Electrode slope 90
 - pH(0) 90
 - Titer 79
- Working life
- Sensor 87
 - Titrant 64
- Write protection
- External memory 140