Agilent Cary UV Workstation

Software Registration Instructions

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Introduction

SubscribeNet allows users to manage software entitlements, order media or download software. We recommend users renew subscriptions annually to have full access to the newest updates.

After installation of Cary UV Workstation, you will have 60 days to complete software registration as described in these instructions. To connect to an instrument after the 60-day period, you must register your software.

You will need the following:

- Authorization code(s) – found in the lavender envelope(s) shipped with the software.
- The host name and MAC address of the computer.
- An internet connection to register your software.
 - Is your computer not connected to the internet?
Perform the steps to register your software on another computer with internet access.
 - USB drive (customer-supplied)– copy the License file generated during the registration process onto the USB and then onto the computer where the Cary UV Workstation software is installed.
- SubscribeNet login credentials if your company has a pre-existing account.
 - If you aren't sure, first check with your IT department or email.
subscribenet_support@agilent.com

Registration

The main steps to register your software are:

- 1 Get the computer name and MAC address:
 - For Cary UV Workstation or Cary UV Workstation Plus, this is the computer name and MAC address.
 - For Cary UV Networked Workstation, this is the Shared Services server name and MAC address.
- 2 Create a SubscribeNet account or login with existing SubscribeNet username and password.
- 3 Enter any additional authorization codes, if required.
- 4 Create the License file.
- 5 Import the License file into OpenLab Control Panel.

Obtaining the computer name and MAC address


To find the computer name and MAC address:

- 1 Double-click the OpenLab Control Panel icon on the computer desktop or click the Windows **Start** button > **Agilent Technologies** > **Control Panel**.

NOTE

For Cary UV Networked Workstation installations, perform this step on the server.

- 2 Enter administrator credentials and then click **Log In**.
- 3 Click **Licenses** from the **Administration** tab.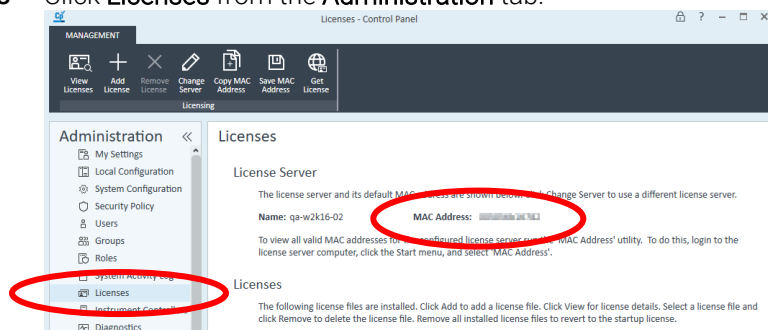

- 4 Take note of the **MAC Address**. You will need it later in this procedure.
- 5 Right-click the Microsoft Windows **Start** icon and then select **System**.
- 6 Take note of the 'Device name'. This is equivalent to the 'Host' name in the Control Panel application. You will need it later in this procedure.

NOTE

Do not use 'localhost' as the device name when creating the License.

Creating and logging into a SubscribeNet account

To create a SubscribeNet account or login with existing SubscribeNet credentials:

- 1 Go to the following website: <https://agilent.subscribe.net>.

NOTE

If your computer is not connected to the internet, print these instructions and take them, and the USB, to a computer that will be used to continue your software registration.

- If you already have a SubscribeNet login and password from previously registered software, enter in that information, click **Login** and then skip to [Register additional authorization codes](#) following the procedure on Page 4. Otherwise, go to Step 3.

NOTE

If you are generating additional Academic Licenses from an already registered Academic License authorization code go to [Generating the License File](#) on Page 4.

- Click '**CLICK HERE**' to create a Login ID for a new account.

Electronic Software and License Delivery

Please login. Your Login ID is your Email address.

Login ID

Password

☐ Remember my password until I logout

Login

If you have forgotten your login ID, password, or are not sure whether you have an account use our [Password Finder](#).

SubscribeNet new account registration.

Customers who have an authorization code from their Agilent product purchase may [CLICK HERE](#) to register and create a new SubscribeNet Account and Login ID.

- Enter in the required details. The Authorization Code(s) are printed on stickers in the lavender envelope(s) provided in the software box.

Electronic Software and License Delivery**Registration Information**

Registration information is used to create a new SubscribeNet Account and Login ID (email address) for the Account. An authorization code you received from your product purchase is required.

Important Email Notes

- Please ensure the email address is entered correctly or you will not receive a welcome email with your Login ID and Password.
- If the email entered has already been used for an account Login ID, you will not receive a welcome email. The new account will display as an additional selection after you log in to SubscribeNet.

* Indicates a required field

Authorization Code*

Email*

Re-enter Email to Verify*

Company*

Department*

First Name*

Last Name*

Job Title*

Phone*

Address 1*

Address 2*

Address 3*

City*

State/Province*

Postal Code*

Country*

Submit

- Press **Submit**. You will receive an email with your username and password. If you do not receive an email within 15 minutes first check your email Spam or Junk folder. Otherwise, contact subscribenet_support@agilent.com.
- Open the email and click the SubscribeNet link. Enter your Login ID (the email address you registered with) and password and log in.
- Follow the prompts to change the password.

Register additional authorization codes

Depending on which software package you have, you may have additional features with separate authorization codes that require registration. These include:

- Multizone add-on
- Universal connection License (for instrument connections when using OpenLab Server/OpenLab Basic Server/ECM XT systems)

To register additional authorization codes:

- 1 Click Register Authorization Code.

The screenshot shows the Agilent Technologies website interface. The header includes the Agilent logo and the tagline 'Our measure is your success.' Below the header, there's a navigation bar with 'Software' selected. Under 'Software', the 'Licenses' option is highlighted, and 'Register Authorization Code' is circled in red. The main content area is titled 'Electronic Software and License Delivery' and contains the 'Register Authorization Code' section. It explains that authorization codes represent license or subscription products and provides a text input field for the code, followed by a 'Submit' button.

- 2 Enter in the Authorization Code included in the lavender envelope provided in the Cary UV Workstation software box.
- 3 Press **Submit**.
- 4 Repeat until all Authorization codes are entered.

Generate the License file

NOTE

When registering an Academic License authorization code or additional Cary UV Workstation or WorkStation Plus or Multizone authorization codes, perform the following steps to generate the License file on *each* computer where the Cary UV Workstation software is installed.

For Cary UV Networked Workstation, when registering authorization codes, perform the following steps to generate the License file on the Shared Services server.

You can use the same SubscribeNet credentials for each registration. The MAC address found in Step 4 in the "Obtaining the computer name and MAC address" section, on Page 2, and the License file generated in this procedure must be specific to each computer where the Cary UV Workstation is installed or the Shared Services server.

To generate the License file:

- 1 Click **Generate** or **View Licenses**.

The screenshot shows the Agilent Technologies website interface. The left navigation bar has 'Licenses' selected, and 'Generate or View Licenses' is circled in red. The main content area is titled 'Electronic Software and License Delivery' and contains the 'Product Information' section. It provides instructions on how to maintain OpenLAB Licenses, including links to 'Subscriptions' and 'Licenses' in the License Pool, and a 'Generate or View Licenses' button.

- 2** Select the tick box of the License or Licenses to generate. Click **Generate**.

NOTE

You can select more than one active License to include in a generated License if all selections are all to be installed on the same MAC address.

- Software
- Product List
- Product Search
- Entitlements
- Licenses
- Register Authorization Code
- Generate or View Licenses
- View Licenses by Host
- Manage Licenses by Host
- Administration
- Account Members
- Change Password
- Email Preferences
- Product Preferences
- Your Profile
- Information
- Download Help
- FAQs
- User Manual

Electronic Software and License Delivery

Generate or View Licenses

For licenses in your License Pool, you may:

- Select one or more licenses to **Generate** by checking the **left** box or boxes.
- Select one or more generated licenses to **View** by clicking the 'View Details' link, and then checking the **right** box or boxes.

License Returns

Generated licenses may be returned to the License Pool by clicking the 'View Details' link, and then clicking the 'Return' button.

[More Help.](#)

Entitlement Number: License:

Entitlement Date: Sep 20, 2018

Generate

☐ Select All

Generate

View

☐ Select All

View

Agilent Cary 3500 Multizone Peltier UV-Vis

Total Number of Licenses: 1

<input type="checkbox"/>	Version	Description
<input type="checkbox"/>	1.0	Agilent Cary 3500 Multizone Peltier UV-Vis Software License
1 Available		

Generate

View

- On the SubscribeNet window, enter the MAC Address and computer name you noted in Step 4 in the "Obtaining the computer name and MAC address" section on Page 2. For Cary UV Networked Workstation, enter the MAC address and server name.

Enter the **MAC Address** into the **HostID** field and the **Name** into the **Hostname** field. Enter a Description of your computer or server.

- 4** Leave the 'Number of Licenses to Generate' box set as the default value of 1.

NOTE

When registering Academic License authorization codes, enter 1 in the 'Number of Licenses to Generate' box.

Software

Product List

Product Search

Entitlements

Licenses

Register Authorization Code

Generate or View Licenses

View Licenses by Host

Manage Licenses by Host

Administration

Account Members

Change Password

Email Preferences

Product Preferences

Your Profile

Information

Download Help

FAQs

User Manual

Support

Electronic Software and License Delivery

Generate Licenses

For your selected licenses, enter the Number of Licenses to Generate followed by the [License Host](#) information:

- Server Host ID**
This is the twelve-digit **MAC Address** of your computer. It may be entered as:
 - twelve digits
 - twelve digits separated by a colon (nn.nn.nn.nn.nn.nn)
 - twelve digits separated by a hyphen (nn-nn-nn-nn-nn-nn)
- Host Name**
This is your Computer Name without the Domain (For example: PC123).
- Host Description (Optional)**
User-friendly description when viewed within SubscribeNet.

[More Help](#)

Version	Description	Number of Licenses Available	Number of Licenses to Generate
1.0	Agilent Cary UV Workstation License	1	<input type="text" value="1"/>

Select an existing License Host from the list or enter a new one
[Existing hosts](#) | [Select an existing License Host](#)

Host ID *

MAC Address |

Hostname *

Description

Generate

- 5 Click **Generate** to generate the License file.
- 6 Click **Download License File**. A dialog box may appear. Follow the prompts to save the file or note where the file was automatically saved to, typically in your C:\\Users\\Username\\Downloads folder.

Agilent Cary UV Workstation Software Registration Instructions

Electronic Software and License Delivery

View Licenses by Host

Listed below are all licenses for the License Host(s) you selected. To save a license file for a Host, click on "Download License File".

[Back to Overview](#)

Host Name: DESKTOP-TTGSTG4
Host ID: 00155d016d08
Host Description:

[Download License File](#)

#Account Name: Agilent Technologies Pty Ltd,CSO
#Host: DESKTOP-TTGSTG4 (MAC-Address = 00155d016d08)
SERVER DESKTOP-TTGSTG4 00155d016d08
VENDOR AGTOL
USE_SERVER

NOTE

If you are performing this registration on a computer other than the one the Cary UV Workstation software was installed on, navigate to the saved License file and then save it to a USB drive.

Insert the USB drive into the computer that the Cary UV Workstation software is installed on before proceeding.

Import License files into Control Panel


To import the License file into OpenLab Control Panel:

- 1 If not already open, double-click the OpenLab **Control Panel** icon on the computer desktop or click the Windows **Start** button > **Agilent Technologies** > **Control Panel**.


NOTE

For Cary UV Networked Workstation, perform this step on the Shared Services server.

- 2 In the OpenLab Control Panel, click **Add License**.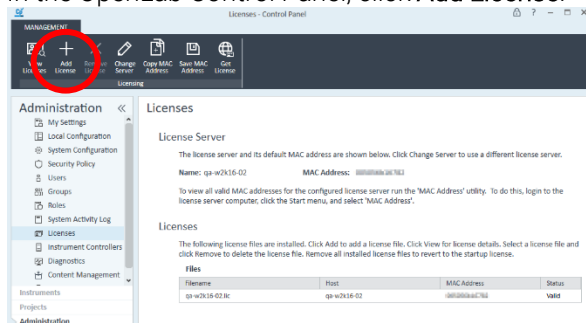

- 3 Click **Select a License file** and then click **Browse**.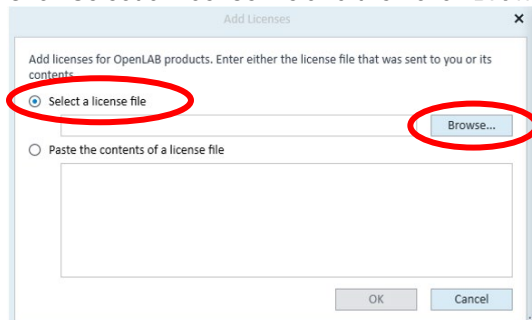

- 4 Navigate to and then select your previously saved License file on the computer or USB. Click **Open**.
- 5 Select **OK** to Add your License.
- 6 Your License will now appear in the list of Licenses in the OpenLab Control Panel and in SubscribeNet when you click 'Generate or View Licenses'.

- 7 Restart the computer to complete the process.

NOTE

When registering Academic License authorization codes, repeat the [Generate the License file](#) procedure on Page 4 and then [Import License files into Control Panel](#) on each computer the Cary UV Workstation software is installed on. You will need to log in to SubscribeNet using the same credentials as the first registration before continuing.

Return Cary UV Workstation and/or OpenLab Control Panel Licenses

NOTE

This procedure is only required if you are changing the computer the Cary UV Workstation and OpenLab Control panel software are currently installed on, or changing the server on which Cary UV Networked Workstation is installed.

- 1 Go to SubscribeNet: <http://agilent.subscribenet.com/>.
- 2 Enter the user name and password used when creating the original Licenses.
- 3 Click **Generate or view Licenses**.
- 4 Click **View Details**.
- 5 Click **Return** for the desired software.

NOTE

If “Return” is not available, then you have reached the maximum number of License returns. You only have two returns for each License product quantity in your License pool. For example, for a License product with a quantity of one, you will have two returns maximum.

A message appears indicating the License was successfully returned to the License pool.

NOTE

You must refresh your browser before attempting to generate a new License. Press F5 on your keyboard with the browser window open if you are using Microsoft Edge, Google Chrome, or Firefox.

- 6 Follow the previous registration instructions to generate and apply a new License
- 7 After completing the registration, restart your computer.

Upgrading License Files

Upgrade your Licenses in SubscribeNet before upgrading the Cary UV Workstation or Cary UV Networked Workstation software.

This involves the following steps:

- Obtain the MAC address of the affected computer or server in Control Panel
- Search for the host name in SubscribeNet
- Upgrade the License in SubscribeNet
- Enter the new License file on the affected computer or server in Control Panel


To upgrade and add the License file:

- 1 Open and login to Control Panel on the computer that Cary UV Workstation is installed on.

NOTE

For Cary UV Networked Workstation perform this step on the Shared Services server. If you are upgrading from Cary UV Networked Workstation version 1.3 applied to a computer, this License must be returned, regenerated with the server information, and then applied to the server.

- 2 Click the **Administration** tab.
- 3 Click **Licenses**.
- 4 Note down the host name. Either note down the MAC address or use the **Copy MAC Address** or **Save MAC Address** function to save it to the Clipboard.
- 5 Log into the Agilent Electronic Software and License Delivery (<https://agilent.subscribenet.com/>).
- 6 Navigate to **Manage Licenses by Host**.
- 7 In the Host ID field, enter or paste the previously noted MAC address, and then click **Search**.

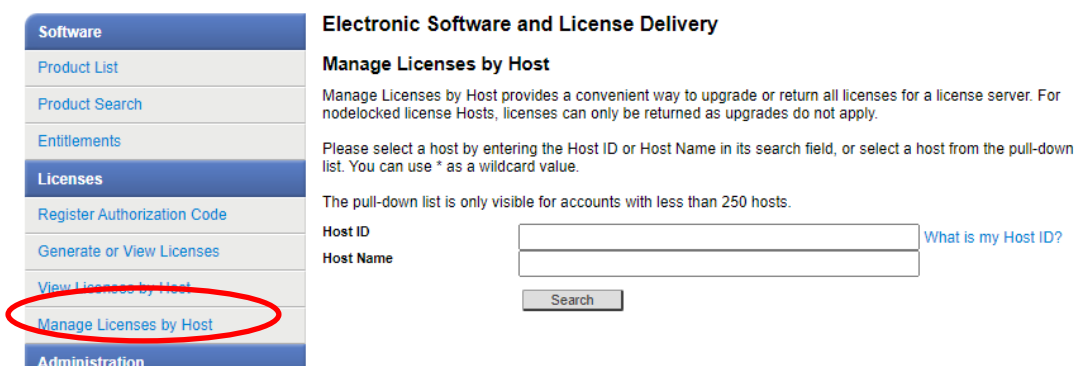The screenshot shows the 'Electronic Software and License Delivery' website. On the left is a navigation menu with the following items: Software, Product List, Product Search, Entitlements, Licenses (highlighted with a red circle), Register Authorization Code, Generate or View Licenses, View Licenses by Host, Manage Licenses by Host (also highlighted with a red circle), and Administration. The main content area is titled 'Manage Licenses by Host' and contains instructions on how to use the search functionality. It includes input fields for 'Host ID' and 'Host Name', a 'Search' button, and a link that says 'What is my Host ID?'. A note states: 'The pull-down list is only visible for accounts with less than 250 hosts.'

NOTE

If the relevant host name does not appear, you may be managing your Licenses in multiple SubscribeNet accounts. You will need to log into those accounts to upgrade those Licenses.

NOTE

If your License(s) are eligible for an upgrade, you will see the **Upgrade All** button. Otherwise you will need to contact your Agilent Sales Representative to renew your Software Maintenance Agreement.

- 8 Click **Upgrade All** to proceed with generating your upgrade License.
- 9 On the **Upgrade All Licenses for License Host** page, review the data, and confirm by clicking **Upgrade All**.
This upgrades the License file to the most current version. SubscribeNet will send you an email with a new License file.
- 10 Add the new License file on your system. See the next section for instructions.
If you have multiple standalone Workstations, repeat this step for each individual workstation.
Note that each workstation's MAC address is the file name. This helps identify the correct License file to import into the workstation's Control Panel.

Add an Upgraded License File to the System

If you have purchased new options or have upgraded a License, and regenerated your license in SubscribeNet, the upgraded License file must be re-applied to the system.

NOTE

For Cary UV Networked Workstation apply the updated License on the Shared Services server. If you are upgrading from Cary UV Networked Workstation version 1.3, this License must first be returned on the computer, and then apply the License generated in the previous section to the server.

-
- 1 Open and login to Control Panel on the computer that Cary UV Workstation is installed on.
 - 2 Navigate to **Administration > Licenses**.
 - 3 Select a License, then click **Remove License**.
 - 4 In the ribbon, click **Add License**.
 - 5 Browse to and open the License file saved from the License generation process in SubscribeNet.

NOTE

A full restart is required in order for any License to have an immediate effect.

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Generate Licenses

For your selected licenses, enter the Number of Licenses to Generate followed by the License Host information:

- Server Host ID**
This is the twelve-digit MAC Address of your computer. It may be entered as:
 - twelve digits
 - twelve digits separated by a colon (nn.nn.nn.nn.nn.nn)
 - twelve digits separated by a hyphen (nn-nn-nn-nn-nn-nn)
- Host Name**
This is your Computer Name without the Domain (For example: PC123).
- Host Description (Optional)**
User-friendly description when viewed within SubscribeNet.

[More Help...](#)

Version	Description	Number of Licenses Available	Number of Licenses to Generate
1.0	Agilent Cary 3500 Multizone Peltier UV-Vis Software License	1	1

Select an existing License Host from the list or enter a new one

Existing hosts: [Select an existing License Host...](#)

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View Licenses by Host

Listed below are all licenses generated for the License Host you selected. To save a license file, click on "Download License File".

[Back to Overview](#)

Host Name: localhost
Host ID:
Host Description: Instrument PC-1

Download License File

#Account Name: Agilent Technologies Pty Ltd,CSO
#Host: localhost (MAC-Address =)
SERVER localhost
VENDOR AGTOL
USE_SERVER

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Manage Licenses by Host

Manage Licenses by Host provides a convenient way to upgrade or return all licenses for a license server. For nodelocked license Hosts, licenses can only be returned as upgrades do not apply.

Please select a host by entering the Host ID or Host Name in its search field, or select a host from the pull-down list. You can use * as a wildcard value.

The pull-down list is only visible for accounts with less than 250 hosts.

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© Agilent Technologies, Inc. 2023
Agilent Technologies Australia [M] Pty Ltd