

Getting Started with Sample Scheduler

1 Select view

1 Start screen

2 Analyses

4 My Instruments

1 Start screen

1 My Instruments

4 Instruments (My Currently Used)

2 Analyses

3 Analyses (Pending)

3 Analyses (Unassigned)

2 Analyses toolbar

4 Analyses action toolbar

4 Create Analyses

3 View or Edit analysis parameters

3 Edit vial position

4 Process analyses, e.g., schedule, stop, delete, or restart an analysis

1 Click or tap elements on the **Start** Screen to view details about your laboratory's workload.

2 Manage your daily workload in the **Analyses** view, for example, create analyses, edit parameters in the **Details** pane, and process analyses or sequences.

3 Filter your daily workload by state. Click or tap on a segment of the half-circle radar or the corresponding category above to view analyses with a specific workload state. Click or tap on the **Unassigned** button to view all unassigned analyses.

4 Click or tap the **My Instruments** button to view your currently used instruments for analysis requests with the status **Scheduled** or **Pending**.