

# Development of an MRM based phospholipid profiling method in human plasma using an inert C18 column

Masaki Yamada, Naoko Nagano, Yohei Arai, Yutaka Umakoshi

Shimadzu Corporation, 1, Nishinokyo-Kuwabaracho Nakagyo-ku, Kyoto 604-8511, Japan

## 1. Overview

- The wider targeted monitoring method of human plasma phospholipids has been developed as a new tool for biomarker discovery.
- An LC/MS system combined with ultra fast triple quadrupole mass spectrometer LCMS-8060NX, Nexera™ UHPLC and a new inert C18 "Shim-pack Scepter™ Claris" column were used.
- Totally 662 MRM transitions were piled up on 20 minutes chromatographic condition. Over 170 phospholipids were detected in human plasma within CV of 50% without significant carryover.

## 2. Introduction

Plasma phospholipids are attractive targets to finding disease biomarker or pathophysiological research. Here we developed a phospholipid profiling method in human plasma based on multiple reaction monitoring, MRM, coupled with reversed phase chromatography using inert C18 column. Totally 170 phospholipids were detected including 130 diacyl-phospholipids, 28 lyso-phospholipids and 12 sphingomyelins in 20 minutes' analytical cycle. Fatty acid composition was determined for all the 130 diacyl-phospholipids by detecting corresponding fatty acid product ion in negative ion mode. Curved gradient was effective to differentiate fatty acid composition isomers such as phosphatidylcolins, PC 18:2\_20:4 and 18:1\_20:5, and 16:0\_22:6. Reproducibility was evaluated by 50 times repeated analyses of human plasma extract.

## 3. Methods and Materials

Pooled heparin human plasma was purchased from (Kojin-Bio Corp., Japan). An LC-MS system consisting of Nexera UHPLC and an LCMS-8060NX mass spectrometer (Shimadzu Corp.) was used. Twenty mM ammonium formate and acetonitrile/isopropanol (1/1, v/v) were used for mobile phase A and B, respectively. Shim-pack Scepter Claris C18, 2.1 x 100 mm, 1.9 μm, (Shimadzu Corp.) was used for chromatographic separation. Polarity switching time was 5 msec. Both dwell time and pause time were set at 1 msec for an MRM transition. One mL of methanol containing 0.1% formic acid was added to 20 μL of human plasma and vigorously mixed a few minutes. After centrifugation, 3 μL of supernatant was injected to LC-MS system.

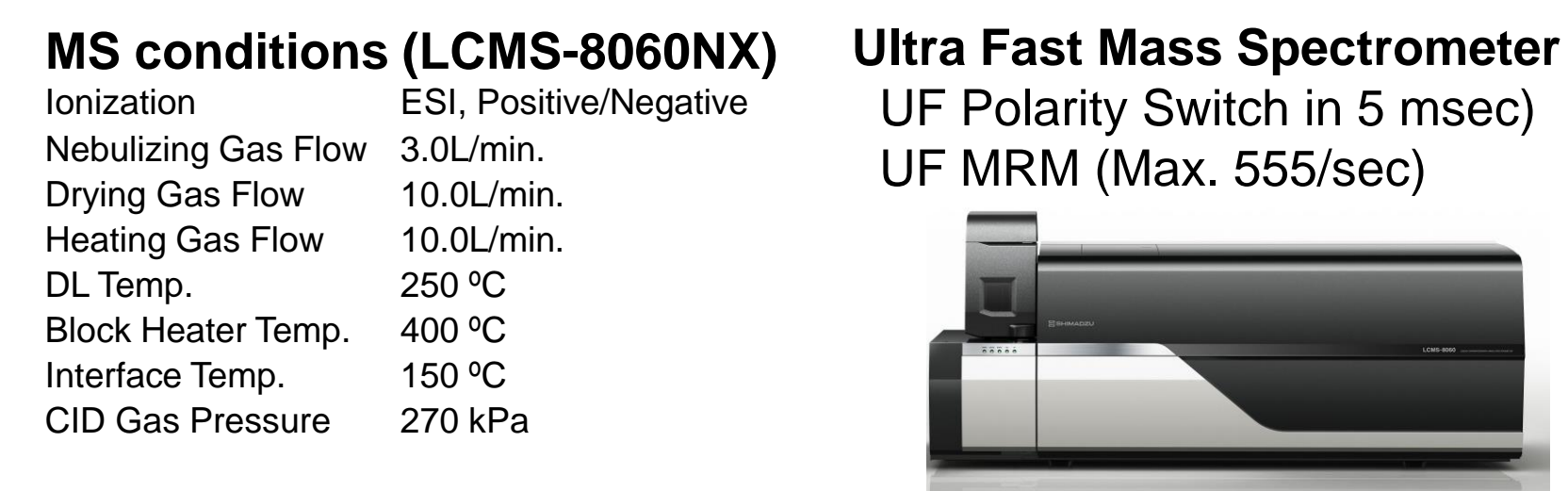


Figure 1 LCMS-8060NX triple quadrupole mass spectrometer

## 4. Results

### 4-1. Method development

For phospholipid profiling in human plasma, 662 MRM transitions were set on 20 minutes chromatographic condition using a new inert C18 column. An MRM to monitor polar head group such as colin derived *m/z* 184, and one or two MRM were set for a diacyl-phospholipid with the same fatty acids or different ones, respectively. As shown in Figure 2 for PC 38:6, 11 MRM were set including one for polar head group in positive ion mode and other 10 MRM for monitor fatty acids in negative ion mode. Three isomers PC 18:2\_20:4, 18:1\_20:5, and 16:0\_22:6 were assigned at 10.15 min, 10.25 min, and 10.45 min, respectively. The other isomers PC 16:1\_22:5 and 18:3\_20:3 were monitored but not detected.

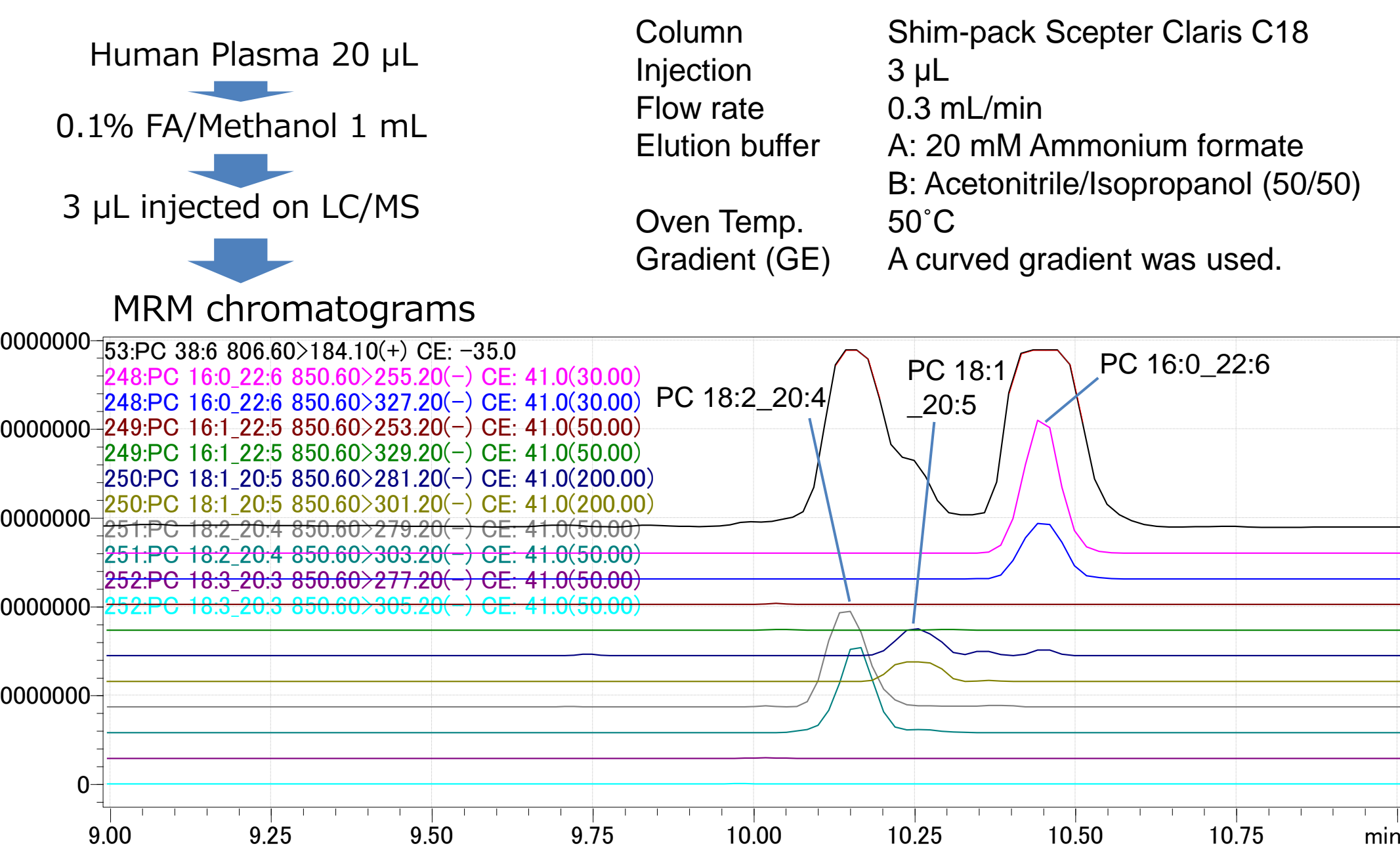


Figure 2 Pre-treatment workflow, chromatographic conditions, and MRM chromatograms for PC 38:6.

### 4-2. Phospholipids in human plasma

Table 1 170 phospholipids detected in human plasma. CV and mean area values for 50 repeated analyses. Each lipid class was listed in order of area.

Compound name	Mean area, N=50	CV%	Compound name	Mean area, N=50	CV%
1 LPC 16:0	12590891	5.5	69 PC 18:2_22:1	20967	40.1
2 LPC 18:0	10586633	6.5	70 PC 18:2_22:0	20693	28.3
3 LPC 18:2	5366848	6.8	71 PC 14:0_18:0	20314	38.9
4 LPC 18:1	4262403	7.3	72 PC 14:0_16:1	18351	42.5
5 LPC 20:4	599730	10.2	73 PC 14:0_14:0	16715	41.3
6 LPC 16:1	446275	12.4	74 PC 20:2_20:3	13224	48.8
7 LPC 17:0	371317	8.7	75 PC 20:1_20:3	12652	44.0
8 LPC 14:0	317125	8.2	76 PC 14:0_22:6	12605	46.1
9 LPC 20:3	314656	14.0	77 PC 18:3_20:4	11102	36.5
10 LPC 18:3	90463	11.9	78 PC 16:1_18:3	9065	41.0
11 LPC 20:2	89312	14.0	79 PC 20:0_20:3	8859	49.0
12 LPC 20:1	89018	13.2	80 PC 14:0_18:3	8356	42.2
13 LPC 15:0	86394	8.4			
14 LPC 17:1	55291	12.5	1 LPE 18:0	1535415	13.0
15 LPC 20:0	50618	15.1	2 LPE 18:2	1274312	12.6
16 LPC 22:5	22311	22.3	3 LPE 18:1	631482	8.5
17 LPC 20:5	19539	15.3	4 LPE 16:0	468721	8.7
18 LPC 22:6	15240	20.6	5 LPE 20:4	261270	12.1
19 LPC 22:4	12977	20.0	6 LPE 20:3	27132	18.7
			7 LPE 16:1	14670	18.9
			8 LPE 22:6	14210	19.9
1 PC 16:0_18:2	58965589	5.7	1 PE 18:0_20:4	1568647	8.7
2 PC 18:0_18:2	46587815	6.5	2 PE 18:0_18:2	1428584	10.5
3 PC 16:0_18:1	30741183	7.0	3 PE 18:1_18:1	568689	17.5
4 PC 16:0_20:4	12579677	6.2	4 PE 18:0_22:6	432873	12.3
5 PC 18:2_18:2	7641465	9.8	5 PE 16:0_18:2	408937	11.0
6 PC 18:0_20:4	6600442	8.0	6 PE 16:0_22:6	384514	17.6
7 PC 18:0_20:3	5237183	9.8	7 PE 18:1_18:2	374369	12.5
8 PC 16:0_20:3	4653924	8.3	8 PE 16:0_20:4	368134	11.0
9 PC 18:1_18:1	4268667	8.2	9 PE 18:0_18:1	339326	14.3
10 PC 16:0_18:1	4050703	9.2	10 PE 18:2_18:2	289988	15.7
11 PC 18:1_18:2	4005264	7.4	11 PE 18:1_20:4	159165	18.2
12 PC 16:0_16:0	2408516	9.7	12 PE 16:0_18:1	138010	22.2
13 PC 18:0_22:6	1989134	10.3	13 PE 18:0_20:3	121059	23.3
14 PC 16:0_22:6	1877485	9.0	14 PE 18:0_22:5	85505	25.9
15 PC 16:0_18:3	1568830	8.6	15 PE 18:2_20:4	68122	23.7
16 PC 16:1_18:2	1441655	10.9	16 PE 16:0_22:5	67460	16.3
17 PC 16:0_16:1	1378650	10.1	17 PE 16:0_20:3	66800	30.9
18 PC 16:0_20:2	1182801	10.1	18 PE 18:0_20:5	43813	31.2
19 PC 18:1_20:4	1140317	10.1	19 PE 16:0_22:4	32222	8.2
20 PC 18:1_20:3	1038014	11.9	20 PE 18:0_22:4	30734	44.5
21 PC 16:0_22:4	960176	11.9	21 PE 18:0_18:3	23283	44.4
22 PC 18:2_20:4	890457	10.6	22 PE 16:0_18:0	14284	20.0
23 PC 16:1_18:1	859248	12.4	23 PE 16:0_16:1	10187	39.8
24 PC 18:0_22:4	722544	12.1			
25 PC 16:0_20:5	698004	9.3	1 LPI 18:2	119082	11.2
26 PC 16:0_20:5	588343	12.5			
27 PC 16:0_18:0	559245	11.8	2 PI 18:0_20:4	8877499	9.3
28 PC 16:0_20:2	1182801	10.1	1 PI 18:0_18:2	4634368	10.9
29 PC 18:1_20:4	1140317	10.1	2 PI 18:0_18:1	113979	11.9
30 PC 18:1_20:3	1038014	11.9	3 PI 18:0_20:4	113159	11.1
31 PC 16:0_22:4	960176	11.9	4 PI 18:0_16:0	127246	13.5
32 PC 18:2_20:4	890457	10.6	5 PI 16:0_20:2	1132277	17.0
33 PC 16:1_18:1	859248	12.4	6 PI 16:0_20:4	850955	13.3
34 PC 18:0_22:4	722544	12.1	7 PI 18:1_18:2	681031	16.8
35 PC 16:0_22:5	698004	9.3	8 PI 18:1_18:1	657390	12.5
36 PC 16:0_20:5	588343	12.5	9 PI 16:0_18:1	558950	12.1
37 PC 16:0_18:0	559245	11.8	10 PI 18:0_22:6	350873	15.6
38 PC 18:0_18:0	523205	15.4	11 PI 18:1_20:4	289337	33.9
39 PC 18:0_20:5	487668	12.8	12 PI 18:0_22:5	176527	17.8
40 PC 18:2_20:2	387673	17.4	13 PI 16:0_20:3	150091	20.1
41 PC 17:0_18:2	362518	13.4	14 PI 16:0_16:1	128075	18.7
42 PC 18:0_22:5	362167	17.9	15 PI 16:0_20:3	127246	13.5
43 PC 14:0_16:1	348127	13.1	16 PI 16:0_16:0	121160	19.8
44 PC 18:0_20:2	315901	14.2	17 PI 16:0_20:4	100834	19.7
45 PC 18:0_18:3	314397	17.1	18 PI 18:0_22:4	92649	23.8
46 PC 18:2_20:1	274734	15.2	19 PI 16:0_22:5	87496	17.1
47 PC 18:2_20:3	262792	15.1	20 PI 18:0_20:2	85624	21.8
48 PC 16:0_20:1	251451	14.6	21 PI 18:2_18:2	60805	21.9
49 PC 18:2_20:0	249253	15.3	22 PI 18:0_20:5	43190	37.5
50 PC 14:0_18:2	246595	14.4	23 PI 16:0_18:3	43105	39.3
51 PC 18:1_18:3	206099	17.8	24 PI 16:0_22:4	35770	36.6
52 PC 16:1_18:0	188327	17.0	25 PI 18:0_20:3	31629	40.0
53 PC 14:0_16:0	166944	16.0	26 PI 16:1_18:1	28354	34.0
54 PC 18:1_22:6	153107	21.6	27 PI 16:0_20:2	25429	49.3
55 PC 18:2_18:3	147099	21.7			
56 PC 18:1_20:2	133449	25.1	1 SM 18:1;02/16:0	17142983	5.9
57 PC 17:0_20:4	132322	26.5	2 SM 18:1;02/22:0	6624463	5.8
58 PC 17:0_18:1	126394	18.2	3 SM 18:1;02/18:0	3366477	6.8
59 PC 18:1_20:1	106560	22.8	4 SM 18:1;02/22:1	3293764	7.5
60 PC 16:1_20:3	105760	28.1	5 SM 18:1;02/20:0	3213034	6.7
61 PC 16:1_20:4	103850	29.6	6 SM 18:1;02/16:1	3201541	7.5
62 PC 15:0_18:2	96415	22.3	7 SM 18:1;02/18:1	1921720	7.3
63 PC 20:4_20:4	81609	22.8	8 SM 18:1;02/14:0	141925	7.3
64 PC 18:2_22:6	66777	22.5	9 SM 18:1;02/20:1	899574	10.8
65 PC 14:0_20:4	66737	22.6	10 SM 18:1;02/18:2	119539	21.3
66 PC 18:1_20:5	64989	25.1	11 SM 18:1;02/20:2	30145	40.3
67 PC 20:0_20:4	60219	33.9	12 SM 18:1;02/18:3	10910	44.0
68 PC 18:1_22:4	59651	30.7			

- Phosphatidyl colins, LPC and PC were more than half in the 170 phospholipids as shown in Figure 3.
- Reproducibility was estimated by 50 times repeated analyses of human plasma. 107 of 170 phospholipids were within 20% CV (Figure 4).
- In the chromatograms for 22:5 including phospholipids, probable structural isomers were detected as arrowed peaks in Figure 5 and shown in red on Table 1.

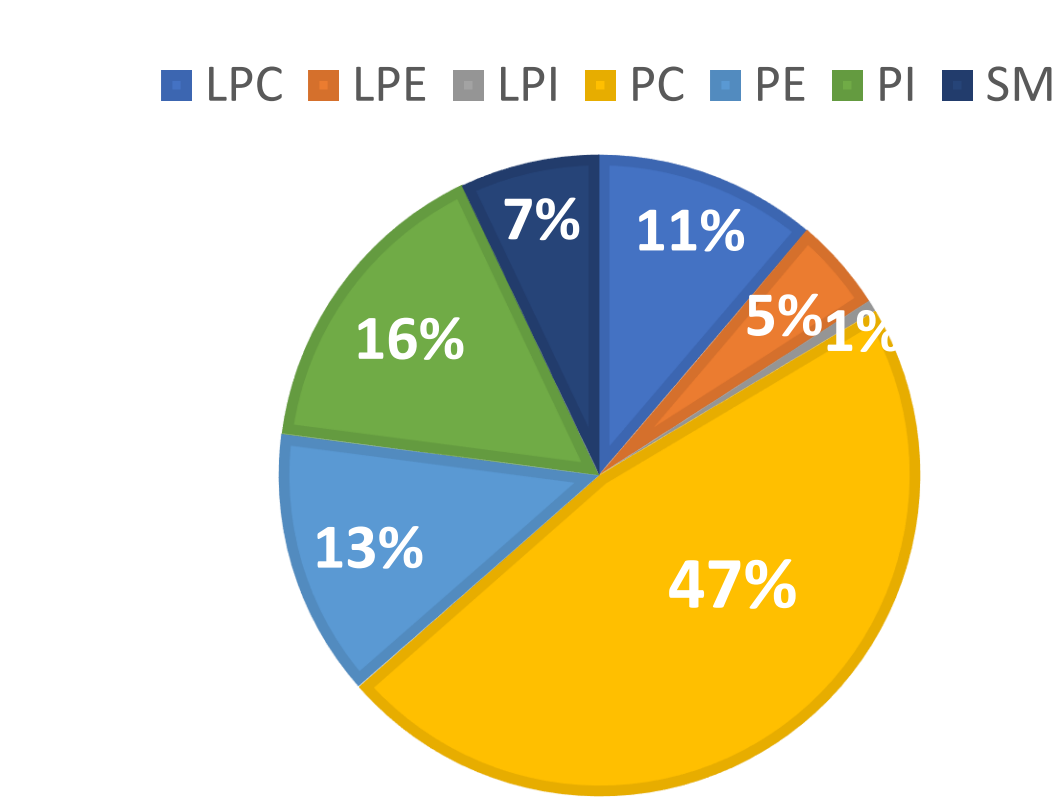


Figure 3 The percentage of 170 phospholipids were shown.

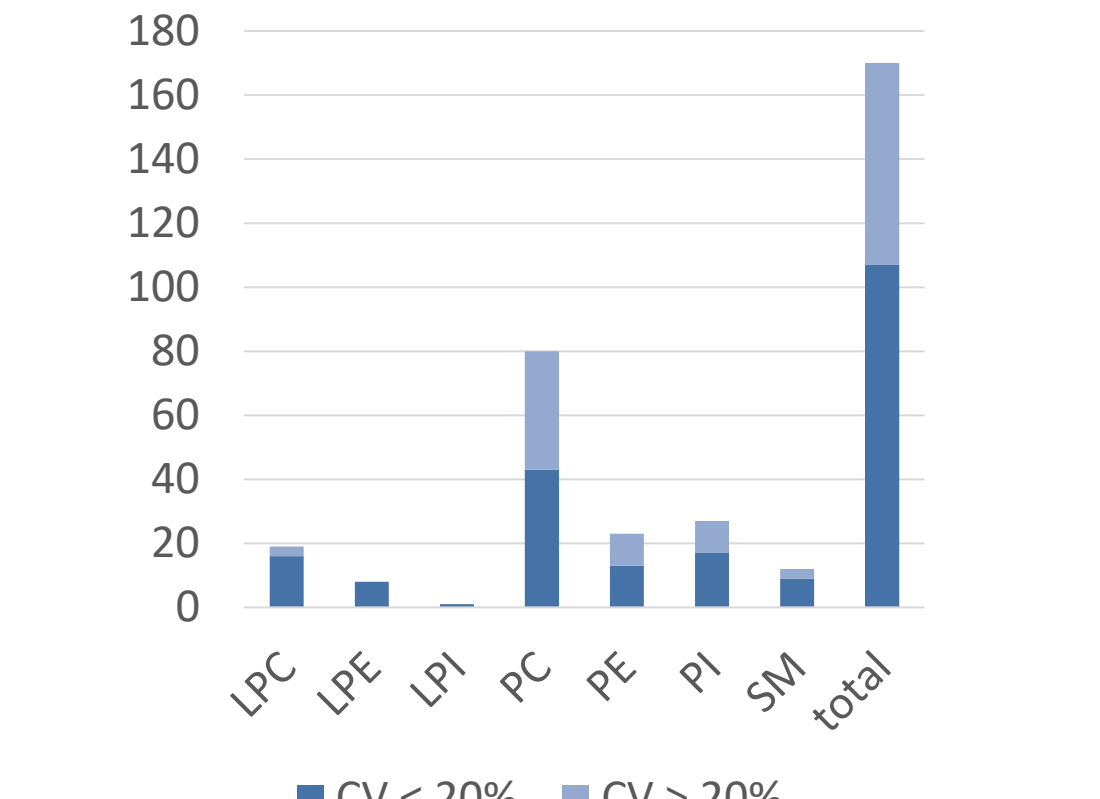


Figure 4 The number of phospholipids less or more than 20% of CV.

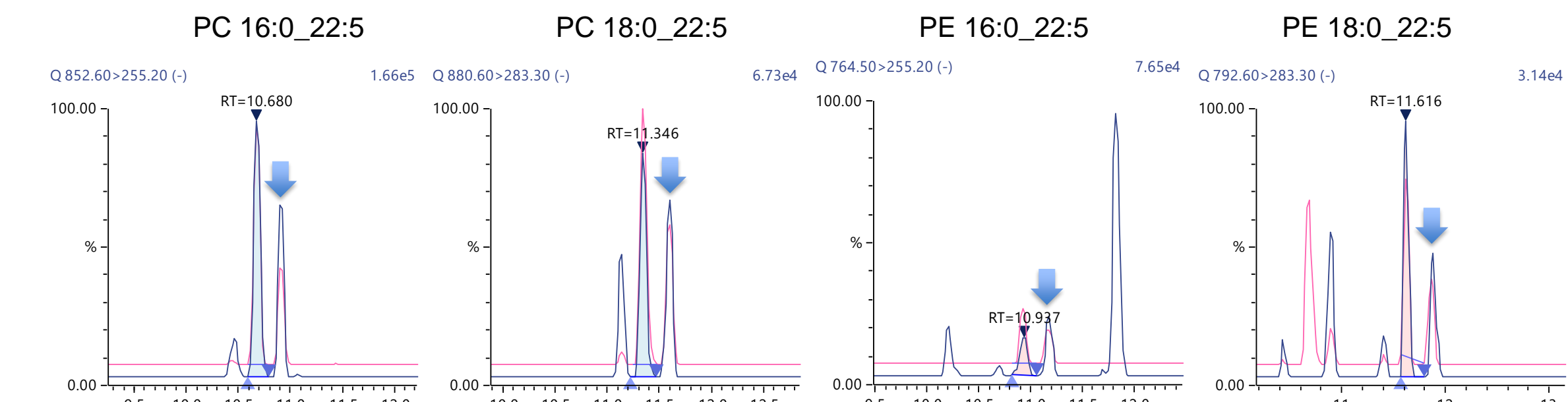


Figure 5 Probable structural isomers of fatty acid 22:5 including phospholipids were arrowed on MRM chromatograms for PC 16:0\_22:5, PC 18:0\_22:5, PE 16:0\_22:5, and PE 18:0\_22:5. The assigned peak was shown with retention time.

## 5. Conclusions

- We have developed a phospholipid profiling method with 20 minutes chromatographic condition. Totally 170 phospholipids were detected. Fatty acid combination was assigned for all 130 diacyl-phospholipids.
- Carryover could be estimated below 0.01% by most abundant peak area of PC 16:0\_18:2 in methanol blank analysis after plasma analysis (data not shown).
- Omega-3 fatty acids such as EPA 20:5 and DHA 22:6 bound to phospholipids were clearly detected as shown in blue on Table 1.
- We believe that this method is available for evaluation of "fatty acid balance" in human plasma sample.

## Acknowledgement

- We would like to thank to Prof. Yoshiya Oda, Dr. Yoshihiro Kita and Dr. Suzumi Tokuoaka at the University of Tokyo for their advice on the method development.

For Research Use Only. Not for use in diagnostic procedures. Nexera and Shim-pack Scepter are trademarks of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.