



# MALDI-nanochip based Screening of Exosomal Biomarkers: Application to Cancer Diagnostics

<u>Michael D. Nairn<sup>1</sup></u>; Michael Wuczkowski<sup>2</sup>; Jesús Jiménez<sup>3</sup>; Iris Prinz<sup>4</sup>; Marco Rissoglio<sup>5</sup>; Emanuele Barborini<sup>5,6</sup>; Gerald Stübiger<sup>2,7</sup> <sup>1</sup>Shimadzu, Manchester, United Kingdom; <sup>2</sup>Medical University of Vienna, Vienna, Austria; <sup>3</sup>Clover Bioanalytical Software, Granda, Spain; <sup>4</sup>Stratec Consumables, Salzburg, Austria; <sup>5</sup>Tethis, Milan, Italy; <sup>6</sup>Luxembourg Institute of Science and Technology; <sup>7</sup>Comprehensive Cancer Center, Vienna, Austria

### 1. Overview

We previously demonstrated we could rapidly distinguish fluorouraci resistant cancer sample groups based on protein profiling of extracellular vesicles using a linear benchtop MALDI TOF instrument [1]. The aim of this follow up work is to identify the proteins that are differentially expressed in the different sample groups in order to better understand the disease processes and to support the rapid screening approach developed previously

Here we present the results from this study using a high performance reflectron MS/MS MALDI-TOF platform (Fig. 1) for the comparative proteomic profiling of circulating extracellular vesicles (EV) extracted from plasma samples of patients with colorectal cancer, postoperative colorectal cancer patients, IBD patients and a healthy control group in view of liquid biopsy applications (as a potential application for liquid biopsy oncological diagnosis).

### 2. Introduction

Exosomes are small cell-derived vesicles (50-150 nm) which are increasingly recognised as a promising source of circulating biomarkers for non-invasive diagnostics from body fluids (liquid biopsy). MALDI-MS profiling of exosomal proteins was demonstrated as being capable to detect cancer-cell specific molecular signatures which can be used to differentiate between cancer types and stages as well as different grades of chemoresistance of cancer cells [1, 2]. This distinguishes MALDI-MS as promising tool for application in liquid biopsy based cancer diagnostics. However, the identification of the disease-related exosomal biomarkers represents a challenging task. Here we present a MALDI-nanochip platform in combination with bioinformatics data analysis for the

comparative profiling and detection of exosomal proteins as potential cancer biomarkers.

#### 3. Methods

Analysis (Clover Biosoft) and eMSTAT (Shimadzu) software.

### 4. Results

40 plasma samples from patients with colon cancer (CRC pre/post operative) inflammatory bowel disease (IBD) and healthy controls were used for evaluation. First, protein extracts were analysed by MALDI-MS in the range of m/z 2000-20000 which has previously been shown to contain most informative peaks of exosomes [2]. A comparison of the mass spectra showed distinct differences particularly in the range above m/z 8000 after exosome isolation. PLS-DA of the whole dataset recorded on the MALDI-nanochip showed a good clustering and separation of the samples belonging to the four study groups (Fig. 2). Next, tryptic digests of the individual samples were subjected to peptide mass fingerprinting (PMF) in order to identify the discriminatory peptide peaks between the study groups based on multivariate data analysis (**Fig. 3**) and MS/MS analysis



**Figure 1** – Overview of the sample preparation and analysis workflow.

Blood samples were prepared according to standard methods and exosomes were isolated using sequential (ultra)centrifugation. Proteins were solvent-extracted, dried under vacuum and stored at -80°C before analysis. Proteins were directly analysed and subsequently subjected to tryptic digestion after application to the MALDI-nanochips (Tethis) (Fig. 1). On-chip digests were dried, washed and covered with 0.5  $\mu$ L CHCA in ACN:2.5%TFA = 70:30 (v/v). Alternatively, samples were digested in-solution, desalted using C18-ZipTips and applied to FlexiMass-DS targets (Shimadzu). For protein profiling the AXIMA-Performance (Shimadzu) instrument was used. Protein identification was performed using the MALDI-7090 MALDI-TOF/TOF mass spectrometer (Shimadzu) with Mascot protein database search (Swiss/Uniprot). Statistical analysis was performed using Clover MS Data





(e.g. alpha-1-antitrypsin, immunoglobulin heavy constant alpha 1, haptoglobin, fibrinogen gamma chain, etc.) were identified in selected samples of the study groups. These proteins were also found in the MALDI-nanochip processed samples but they showed no differentiation between the study groups. In contrast, several

## **WP-031**

2. Stübiger G. et al. MALDI-MS Protein Profiling of Chemoresistance in Extracellular Vesicles of Cancer Cells. Anal Chem, 90, 13178-13182 (2018).