

Application Data Sheet

No. 1

GCMS

Gas Chromatograph Mass Spectrometer

Analysis of Amino Acids Contained in Alcohol

Amino acids contained in alcohol were treated with EZ:faast™ (Phenomenex, Inc.), which enables easy pretreatment, and then analyzed with a GC-MS system.

Experiment

Pretreatment

Three types of alcohol (Japanese *sake*, beer, and wine) were treated using EZ:faast. Norvaline was added as an internal standard.

Instrument

A GCMS-QP2010 Ultra (with high-power oven) was used for the measurements. The analysis conditions, shown in Table 1, were in conformity with the "Amino Acid Analysis Methods" in the "GC/MS Metabolic Components Database."

Table 1: Analysis Conditions (GC/MS Metabolic Components Database: Amino Acid Analysis Methods)

GC-MS	: GCMS-QP2010 Ultra (with high-power oven)	[MS]	
Column	: ZB-AAA (10 mL × 0.25 mm I.D.) (Phenomenex, Inc.)	Interface temperature	: 280°C
[GC]		Ion source temperature	: 200°C
Injection quantity	: 1 µL	Solvent elution time	: 0.4 min
Vaporization chamber temperature	: 280°C	Data sampling time	: 0.5 min to 7 min
Column oven temperature	: 110°C → (30 °C/min) → 320°C	Measurement mode	: Scan
Control mode	: Constant pressure (15 kPa)	Mass range	: m/z 45-450 (3,333u/sec)
Injection mode	: Split	Event time	: 0.15 sec
Split ratio	: 15		
Carrier gas	: Helium		

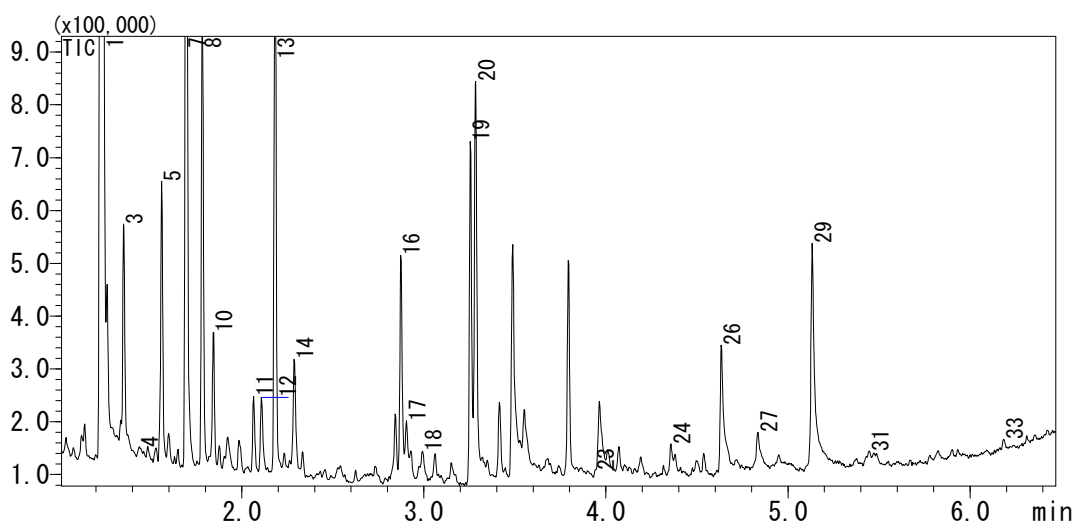


Fig. 1: Total Ion Current Chromatogram for Amino Acid Derivatives in Japanese Sake
The numbers for each component follow the serial numbers in the "GC/MS Metabolic Components Database."

1 Alanine	8 Leucine	14 Asparagine	20 Phenylalanine	29 Tyrosine
3 Glycine	10 Isoleucine	16 Aspartic acid	23 Glutamine	31 Tryptophan
4 alpha-aminobutyric acid	11 Threonine	17 Methionine	24 Ornithine	33 Cystine
5 Valine	12 Serine	18 4-Hydroxyproline	26 Lysine	
7 Norvaline(I.S.)	13 Proline	19 Glutamic acid	27 Histidine	

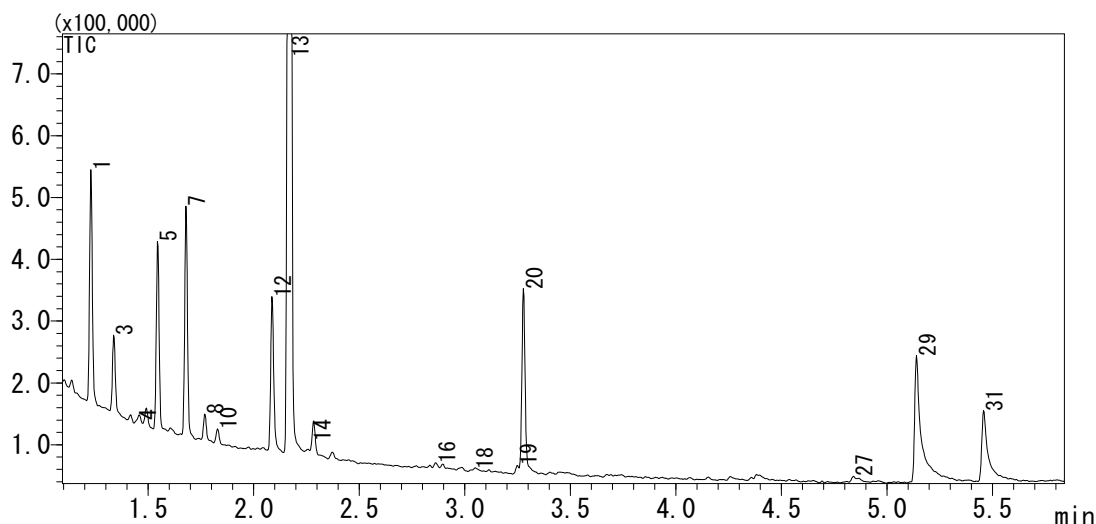


Fig. 2: Total Ion Current Chromatogram for Amino Acid Derivatives in Beer
The numbers for each component follow the serial numbers in the "GC/MS Metabolic Components Database."

1 Alanine	8 Leucine	16 Aspartic acid	29 Tyrosine
3 Glycine	10 Isoleucine	18 4-Hydroxyproline	31 Tryptophan
4 alpha-aminobutyric acid	12 Serine	19 Glutamic acid	
5 Valine	13 Proline	20 Phenylalanine	
7 Norvaline(I.S.)	14 Asparagine	27 Histidine	

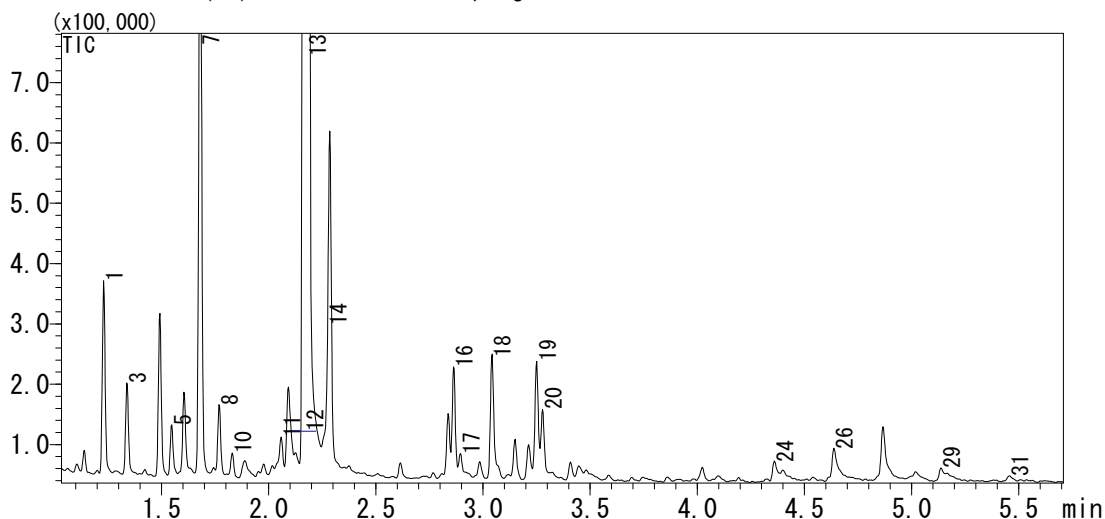


Fig. 3: Total Ion Current Chromatogram for Amino Acid Derivatives in Wine
The numbers for each component follow the serial numbers in the "GC/MS Metabolic Components Database."

1 Alanine	10 Isoleucine	16 Aspartic acid	24 Ornithine
3 Glycine	11 Threonine	17 Methionine	26 Lysine
5 Valine	12 Serine	18 4-Hydroxyproline	29 Tyrosine
7 Norvaline(I.S.)	13 Proline	19 Glutamic acid	31 Tryptophan
8 Leucine	14 Asparagine	20 Phenylalanine	

Summary

Pretreatment using the EZ:fast kit, following by analysis using the GCMS-QP2010 Ultra, which is equipped with a high-speed scanning function, enabled rapid analysis of amino acids. With this combination, it took only 15 minutes per sample from pretreatment to analysis.

(Reference: Shimadzu Application News No. M246 Analysis of Amino Acids Using Fast-GC/MS and Metabolite Database)

For Research Use Only. Not for use in diagnostic procedures.

Shimadzu Corporation ("Shimadzu") reserves all rights including copyright in this publication. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to, or arising out of the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

First Edition: September 2011



SHIMADZU CORPORATION. International Marketing Division

3, Kanda-Nishikicho 1-chome, Chiyoda-ku, Tokyo 101-8448, Japan

Phone: 81(3)3219-5641 Fax: 81(3)3219-5710

URL <http://www.shimadzu.com>