


GC CAPILLARY COLUMN APPLICATIONS: INDUSTRIAL SOLVENTS

Residual Solvents in Drugs USP Method 467

COLUMN: 007-1,
Dimethylpolysiloxane
30M. x 0.32mm I.D. x
10.0µm film
Cat. No.: 007-1-30W-10.0F
Temperature: 50° (4 min. hold)
(5°/min.) - 200°C
Injector: 220°C
Detector: 290°C, FID
Carrier Gas: 25 cm/sec., Helium

1. methanol
2. ethanol
3. acetonitrile
4. acetone
5. isopropanol
6. diethylether
7. methylene chloride
8. tert-butanol
9. propanol
10. methyl ethyl ketone
11. sec-butanol
12. ethyl acetate
13. chloroform & hexane
14. tetrahydrofuran
15. benzene
16. cyclohexane
17. 1,4-dioxane
18. trichloroethylene
19. heptane
20. pyridine
21. dimethylformamide
22. toluene

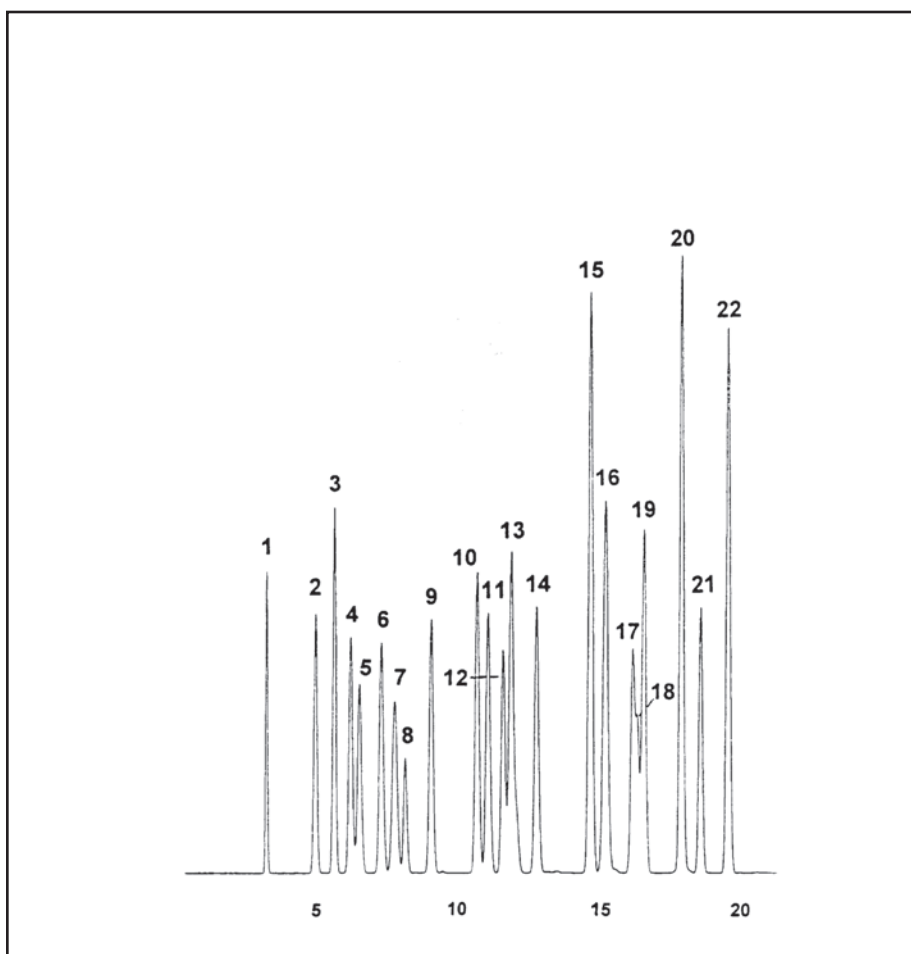

PO Box 3881, Woodbridge, CT 06525 • (203) 393-3112 • FAX (203) 393-0391
TOLLFREE: 800-275-7033 • E-MAIL: tech@quadrexcorp.com