



TwisTorr 305 Turbo Pumps Foreline Flange Installation Instructions


The TwisTorr 305 pumps (all models) are designed to be operated with KF16 and KF25 foreline flanges. While the KF16 is already installed on the pump, the KF25 can be installed following a few simple steps described below.

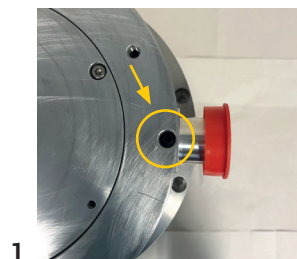


1 Turn the pump upside down and remove screw from its location

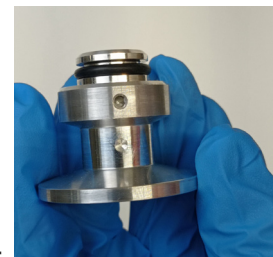


4 Make sure the o-ring is intact, and then install the provided foreline flange (PN X3513-68000) on the pump

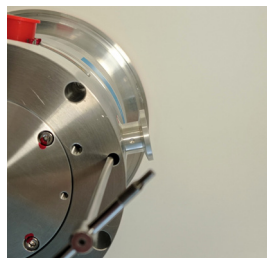


2 Remove foreline flange from the pump

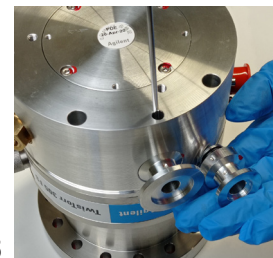

5 Place the screw back into place (use a 1.5 Nm torque)

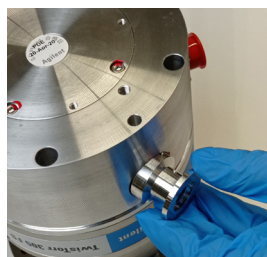

3

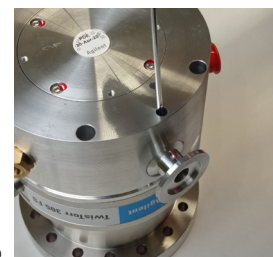

6

Visit the product web pages:



[TwisTorr 305 FS](#)



[TwisTorr 305-IC](#)



[TwisTorr 305 FSQ](#)



[TwisTorr 305-ICQ](#)



[TwisTorr 305 SF](#)



[TwisTorr 305-IC SF](#)

Please contact Agilent Customer Service for more information

[Click here](#)
to get your KF25 foreline flange
for TwisTorr 305 FS turbo pumps



Contact information

Americas

Agilent Technologies
121 Hartwell Avenue,
Lexington, MA 02421 USA
Toll free: +1 800 882 7426
vpl-customer@agilent.com

Europe, Middle East, Africa, India

Agilent Technologies Italia SpA
Via F.lli Varian 54,
10040 Leini (Torino), Italy
Tel: +39 011 9979 111
Toll free: 00 800 234 234 00
vpt-customer@agilent.com

China

Beijing Office Agilent Technologies (China) Co. Ltd.
No.3, Wang Jing Bei Lu,
Chao Yang District,
Beijing, 100102, China
Toll free: 800 820 6778
Contacts.vacuum@agilent.com

For more information, please contact
your Agilent representative or visit
www.agilent.com/chem/vacuum where
you can chat live with a vacuum expert.



DE18111894

This information is subject to change without notice.

© Agilent Technologies, Inc. 2023
Published in Italy, April 24th, 2023
5994-6003EN